

## 苏木素染色液

| 产品编号        | 产品名称   | 包装    |
|-------------|--------|-------|
| C0107-100ml | 苏木素染色液 | 100ml |
| C0107-500ml | 苏木素染色液 | 500ml |

### 产品简介:

- 碧云天生产的苏木素染色液(Hematoxylin Staining Solution)综合了多种经典的苏木素染色方法,例如Harris法、Mayer法等,简化了操作步骤,缩短了操作时间,并且染色液内不使用汞、甲醇等有毒试剂。可以用于组织切片或培养细胞的染色。
- 苏木素染色(hematoxylin staining or haematoxylin staining),也被称作苏木精染色,是最常用的组织和细胞的染色方法之一。无色的苏木素(hematoxylin)氧化后形成有醌环结构(quinoid ring)的氧化苏木素(hematein or haematein),从而可以和三价的铝离子、铁离子等形成有颜色的带正电荷的复合物(如hematein-Al<sup>3+</sup> complexes)。氧化苏木素(也称氧化苏木精)和铝离子等形成有色的复合物的过程也被称为Dye Lake Formation。细胞核内基因组DNA的双螺旋结构中,双链上的磷酸基团向外,带负电荷,呈酸性,很容易与带正电荷的氧化苏木素复合物结合,从而形成细胞核染色。苏木素染色液有多种不同的配制方法,不同的方法可以把细胞核染色成不同深浅的蓝色或蓝紫色[1-3]。
- 本苏木素染色液染色后细胞核呈蓝色,本产品用于石蜡切片染色的效果图参考图1。

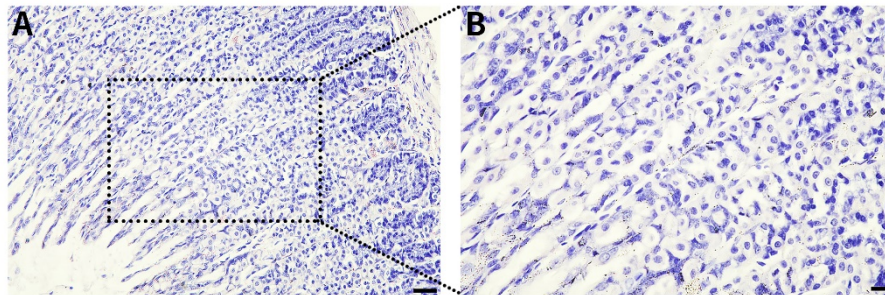


图1. 碧云天的苏木素染色液(C0107)用于大鼠胃部石蜡切片的染色效果图。B图为A图局部放大图片。图中可见非常清晰的细胞核着色。图A中的Scale bar = 100μm, 图B中的Scale bar = 10μm。本图仅作参考,不同的样品不同的检测条件,实际获得的结果可能有所差别。

- 本染色液可以和免疫荧光染色或免疫组化染色配合使用。一方面可以在本苏木素染色液染色后进行免疫荧光染色或其它染料的染色,另一方面也可以在免疫组化染色后再进行苏木素复染。
- 本染色液可以重复使用,直至认为效果不佳时再换用新的染色液。一个100ml包装的本染色液至少可以染色200个样品,一个500ml包装的本染色液至少可以染色1000个样品。

### 包装清单:

| 产品编号        | 产品名称   | 包装    |
|-------------|--------|-------|
| C0107-100ml | 苏木素染色液 | 100ml |
| C0107-500ml | 苏木素染色液 | 500ml |
| —           | 说明书    | 1份    |

### 保存条件:

室温避光保存,至少一年有效。

### 注意事项:

- 对于石蜡切片样品处理时间的长短视使用溶液的新鲜程度、温度及组织切片的大小而定,避免短时间脱蜡脱水不充分,影响后续染色。
- 染色后的分化为选做步骤,但分化后细胞核着色更清晰。
- 染色液可以重复使用多次,认为效果不佳时再更换新的染色液。
- 样品数量较多时,可以使用碧云天生产的染色架和染色缸,便于操作。
- 第一次使用本试剂盒时建议先取1-2个样品做预实验。
- 本产品仅限于专业人员的科学研究用,不得用于临床诊断或治疗,不得用于食品或药品,不得存放于普通住宅内。

➤ 为了您的安全和健康，请穿实验服并戴一次性手套操作。

## 使用说明：

### 1. 需要用户自己准备的试剂

- 固定液：碧云天的免疫染色固定液(P0098)，或4%多聚甲醛固定液(P0099)；
- 分化液(碧云天的盐酸乙醇分化液(C0161、C0163、C0165)或5%的乙酸溶液或0.5%的盐酸乙醇溶液)；
- 如果需要脱水、透明和封片处理，还需自备二甲苯，和封片剂(如碧云天的C0185 PVP封片液、P0126 抗荧光淬灭封片液或中性树胶等其它封片剂)，及80%乙醇、90%乙醇、无水乙醇；
- 70%乙醇。

### 2. 样品处理

#### a. 对于石蜡切片：

- 二甲苯中脱蜡5-10分钟。
- 换用新鲜的二甲苯，再脱蜡5-10分钟。
- 无水乙醇5分钟。
- 90%乙醇2分钟。
- 80%乙醇2分钟
- 70%乙醇2分钟。
- 蒸馏水2分钟。

#### b. 对于冰冻切片：

- 固定液固定10分钟以上。
- 蒸馏水2分钟。

#### c. 对于培养细胞：

- 固定液固定10分钟以上。
- 蒸馏水洗涤2分钟。
- 换用新鲜的蒸馏水，再洗涤2分钟。

### 3. 苏木素染色

对于上述处理好的样品：

- 苏木素染色液染色5-10分钟(可以根据染色结果和要求调整时间)。
- 浸蒸馏水或自来水中数秒，洗去多余的染色液。
- 选做：根据不同分化液的分化时间，分化约2-30秒。
- 自来水冲洗返蓝约10分钟。

此时，如果需要直接观察，可以用70%乙醇洗涤2次。如需脱水、透明后封片按后续步骤进行，70%乙醇洗涤后仍可按照后续步骤进行脱水、透明和封片处理。

注：如果用于免疫组化等染色后的复染，可以参考上述步骤在其它染色完成后直接进行苏木素染色。

### 4. 脱水、透明、封片或进行其它染色

#### a. 脱水、透明、封片：

- 70%乙醇10秒，80%乙醇10秒，90%乙醇10秒，无水乙醇10秒。二甲苯透明5分钟。
- 换用新鲜的二甲苯，再透明5分钟。
- 用中性树胶或其它封片剂封片。
- 显微镜下观察，细胞核呈蓝色。

#### b. 进行其它染色：

如果进行免疫荧光染色，或进行Hoechst等荧光染料的染色，在苏木素染色液染色后：

70%乙醇洗涤2次，每次2分钟。

PBS或生理盐水或TBS或TBST等用于免疫染色或荧光染料染色的溶液浸泡5分钟。

然后就可以进行免疫荧光染色或其它荧光染料的染色了。如果免疫荧光染色效果不佳，可能染料对抗体结合有影响，请单独染色。

## 参考文献：

- Smith C. MLO Med Lab Obs. 2006. 38(5):18, 20-2.
- Titford M. Biotech Histochem. 2005. 80(2):73-8.
- Fischer AH, Jacobson KA, Rose J, Zeller R. CSH Protoc. 2008. 2008:pdb.prot4986.

## 相关产品：

| 产品编号        | 产品名称           | 包装     |
|-------------|----------------|--------|
| C0105S      | 苏木素伊红(HE)染色试剂盒 | >200次  |
| C0105M      | 苏木素伊红(HE)染色试剂盒 | >1000次 |
| C0107-100ml | 苏木素染色液         | 100ml  |

|             |                 |       |
|-------------|-----------------|-------|
| C0107-500ml | 苏木素染色液          | 500ml |
| C0109       | 伊红染色液           | 100ml |
| C0115       | 甲基绿染色液          | 100ml |
| C0117       | 尼氏(Nissl)染色液    | 100ml |
| C0119       | 甲基绿-派洛宁染色液      | 100ml |
| C0121-100ml | 结晶紫染色液          | 100ml |
| C0121-500ml | 结晶紫染色液          | 500ml |
| C0123       | 中性红染色液          | 100ml |
| C0125       | 中性红染色液(活细胞染色用)  | 100ml |
| C0161S      | 盐酸乙醇慢速分化液       | 100ml |
| C0161M      | 盐酸乙醇慢速分化液       | 500ml |
| C0161L      | 盐酸乙醇慢速分化液(20X)  | 100ml |
| C0163S      | 盐酸乙醇快速分化液       | 100ml |
| C0163M      | 盐酸乙醇快速分化液       | 500ml |
| C0163L      | 盐酸乙醇快速分化液(20X)  | 100ml |
| C0165S      | 盐酸乙醇超快速分化液      | 100ml |
| C0165M      | 盐酸乙醇超快速分化液      | 500ml |
| C0165L      | 盐酸乙醇超快速分化液(20X) | 100ml |

### 使用本产品的文献：

- Jinliang Gao, Tao Luo, Jinke Wang. Gene interfered-ferroptosis therapy for cancers. *Nat Commun.* 2021 Sep 7;12(1):5311.
- Shasha Zhao, Yuelong Wang, Nian Yang, Min Mu, Zhiguo Wu, Hexian Li, Xin Tang, Kunhong Zhong, Zongliang Zhang, Cheng Huang, Ting Cao, Meijun Zheng, Guoqing Wang, Chunlai Nie, Hui Yang, Gang Guo, Liangxue Zhou, Xi Zheng, Aiping Tong. Genome-scale CRISPR-Cas9 screen reveals novel regulators of B7-H3 in tumor cells. *J Immunother Cancer.* 2022 Jun;10(6):e004875.
- Tengfei Huang, Tianqi Xu, Yangfan Wang, Yan Zhou, Dandan Yu, Zhiyuan Wang, Linfang He, Zhangpeng Chen, Yaliang Zhang, Don Davidson, Yuyuan Dai, Chunhua Hang, Xiangyu Liu, Chao Yan. Cannabidiol inhibits human glioma by induction of lethal mitophagy through activating TRPV4. *Autophagy.* 2021 Nov;17(11):3592-3606.
- Zhenzhen Chen, Chenxi Ouyang, Haizeng Zhang, Yuanrui Gu, Yue Deng, Congkuo Du, Changting Cui, Shuangyue Li, Wenjie Wang, Wei Kong, Jingzhou Chen, Jun Cai, Bin Geng. Vascular smooth muscle cell-derived hydrogen sulfide promotes atherosclerotic plaque stability via TFEB (transcription factor EB)-mediated autophagy. *Autophagy.* 2022 Oct;18(10):2270-2287.
- Xin Xu, Lifang Ma, Xiao Zhang, Susu Guo, Wanxin Guo, Yikun Wang, Shiyu Qiu, Xiaoting Tian, Yayou Miao, Yongchun Yu, Jiayi Wang. A positive feedback circuit between RN7SK snRNA and m6A readers is essential for tumorigenesis. *Mol Ther.* 2023 Jun 7;31(6):1615-1635.
- Liu H, Zhang C, Liang T, Song J, Hao J, Hou G. Inhibition of Pim2-prolonged skin allograft survival through the apoptosis regulation pathway. *Cell Mol Immunol.* 2012 Nov;9(6):503-10.
- Guangyu Ji, Wenjuan Zhou, Jingyi Du, Juan Zhou, Dong Wu, Man Zhao, Liping Yang, Aijun Hao. PCGF1 promotes epigenetic activation of stemness markers and colorectal cancer stem cell enrichment. *Cell Death Dis.* 2021 Jun 19;12(7):633.
- Yipu Fan, Yihan Zhang, Kunyu Shi, Shan Cheng, Duanqing Pei, Xiaodong Shu. Identification of a group of bisbenzylisoquinoline (BBIQ) compounds as ferroptosis inhibitors. *Cell Death Dis.* 2022 Nov 26;13(11):1000.
- Lingqiao Wang, Lixiong He, Hui Zeng, Wenjuan Fu, Jia Wang, Yao Tan, Chuanfen Zheng, Zhiqun Qiu, Jiaohua Luo, Chen Lv, Yujing Huang, Weiqun Shu. Low-dose microcystin-LR antagonizes aflatoxin B1 induced hepatocarcinogenesis through decreasing cytochrome P450 1A2 expression and aflatoxin B1-DNA adduct generation. *Chemosphere.* 2020 Jun;248:126036.
- Jingxin Yang, Yalin Zhang, Yiming Pan, Can Sun, Zuwang Liu, Ning Liu, Yu Fu, Xiaofeng Li, Ye Li, Juan Kong. The Protective Effect of 1,25(OH)2D3 on Myocardial Function is Mediated via Sirtuin 3-Regulated Fatty Acid Metabolism. *Front Cell Dev Biol.* 2021 Apr 26;9:627135.
- Jingjing Zhu, Shuangshuang Zheng, Hanbo Liu, Yiwei Wang, Zeren Jiao, Yi Nie, Hong Wang, Tianqing Liu, Kedong Song. Evaluation of anti-tumor effects of crocin on a novel 3D tissue-engineered tumor model based on sodium alginate/gelatin microbead. *Int J Biol Macromol.* 2021 Mar 31;174:339-351.
- Yiyun Deng, Zhe Zhang, Hui Yang, Jing Wang, Lijuan Feng, Yong Su, Dujuan Xu. Mycophenolic Acid Induces the Intestinal Epithelial Barrier Damage through Mitochondrial ROS. *Oxid Med Cell Longev.* 2022 Jul 5:2022:4195699.
- Luo J, Zhao X, Tan Z, Su Z, Meng F, Zhang M. Mesenchymal-like progenitors derived from human embryonic stem cells promote recovery from acute kidney injury via paracrine actions. *Cytotherapy.* 2013 Jun;15(6):649-62.
- Shi H, Tan B, Ji G, Lu L, Cao A, Shi S, Xie J. Zedoary oil (Ezhu You) inhibits proliferation of AGS cells. *Chinese Medical Journal.* 2013 Jun 28;8(1):13.
- Xi S, Yue L, Shi M, Peng Y, Xu Y, Wang X, Li Q, Kang Z, Li H, Wang Y. The Effects of Taoren-Honghua Herb Pair on Pathological Microvessel and Angiogenesis-Associated Signaling Pathway in Mice Model of CCl4-Induced Chronic Liver Disease. *EVID-BASED COMPL ALT.* 2016;2016:2974256.
- Zhou X, da Li W, Yuan BL, Niu LJ, Yang XY, Zhou ZB, Chen XH, Feng X. Lithium Treatment Prevents Apoptosis in Neonatal Rat Hippocampus Resulting from Sevoflurane Exposure. *Neurochem Res.* 2016 Aug;41(8):1993-2005.
- Wang B, Zhang Y, Cao W, Wei X, Chen J, Ying W. SIRT2 Plays Significant Roles in Lipopolysaccharides-Induced Neuroinflammation and Brain Injury in Mice. *Neurochem Res.* 2016 Sep;41(9):2490-500.

18. Zhang JH, Jiao LY, Li TJ, Zhu YY, Zhou JW, Tian J. GSK-3 $\beta$  suppresses HCC cell dissociation in vitro by upregulating epithelial junction proteins and inhibiting Wnt/ $\beta$ -catenin signaling pathway. *J Cancer*. 2017 Jun 3;8(9):1598-1608.
19. Chen YY, Sun ZW, Jiang JP, Kang XD, Wang LL, Shen YL, Xie XD, Zheng LR.  $\alpha$ -adrenoceptor-mediated enhanced inducibility of atrial fibrillation in a canine system inflammation model. *Mol Med Rep*. 2017 Jun;15(6):3767-3774.
20. He Z, Yu Y, Nong Y, Du L, Liu C, Cao Y, Bai L, Tang H. Hepatitis B virus X protein promotes hepatocellular carcinoma invasion and metastasis via upregulating thioredoxin interacting protein. *Oncol Lett*. 2017 Aug;14(2):1323-1332.
21. Wang X, Zhang W, Sun X, Lin Y, Chen W. Cancer-associated fibroblasts induce epithelial-mesenchymal transition through secreted cytokines in endometrial cancer cells. *Oncol Lett*. 2018 Apr;15(4):5694-5702.
22. Xia HB, Wang HJ, Fu LQ, Wang SB, Li L, Ru GQ, He XL, Tong XM, Mou XZ, Huang DS. Decreased CRHBP expression is predictive of poor prognosis in patients with hepatocellular carcinoma. *Oncol Lett*. 2018 Sep;16(3):3681-3689.
23. Han B, Feng D, Yu X, Liu Y, Yang M, Luo F, Zhou L, Liu F. MicroRNA-144 mediates chronic inflammation and tumorigenesis in colorectal cancer progression via regulating C-X-C motif chemokine ligand 11. *Exp Ther Med*. 2018 Sep;16(3):1935-1943.
24. Yin XF, Zhang Q, Chen ZY, Wang HF, Li X, Wang HX, Li HX, Kang CM, Chu S, Li KF, Li Y, Qiu YR. NLRP3 in human glioma is correlated with increased WHO grade, and regulates cellular proliferation, apoptosis and metastasis via epithelial-mesenchymal transition and the PTEN/AKT signaling pathway. *Int J Oncol*. 2018 Sep;53(3):973-986.
25. Peng H, Zhang K, Liu Z, Xu Q, You B, Li C, Cao J, Zhou H, Li X, Chen J, Cheng G, Shi R, Zhang G. VPO1 Modulates Vascular Smooth Muscle Cell Phenotypic Switch by Activating Extracellular Signal-regulated Kinase 1/2 (ERK 1/2) in Abdominal Aortic Aneurysms. *J Am Heart Assoc*. 2018 Sep 4;7(17):e010069. (IF 4.605)
26. Zhang C, Hao Y, Wu L, Dong X, Jiang N, Cong B, Liu J, Zhang W, Tang D, De Perrot M, Zhao X. Curcumin induces apoptosis and inhibits angiogenesis in murine malignant mesothelioma. *Int J Oncol*. 2018 Dec;53(6):2531-2541.
27. Xia YF, Chang J, Yang JF, Ouyang W, Pitt B, Billiar T, Zhang LM. Non-canonical Wnt signaling contributes to ventilator-induced lung injury through upregulation of WISP1 expression. *Int J Mol Med*. 2019 Mar 43(3):1217-1228.
28. Zhu Z, Yu J, Lin W, Tang H, Zhang W, Lu B. Molecular hydrogen accelerates the reversal of acute obstructive cholangitis-induced liver dysfunction by restoring gap and tight junctions. *Mol Med Rep*. 2019 Jun 19(6):5177-5184.
29. Chen L, Wu M, Zhang S, Tan W, Guan M, Feng L, Chen C, Tao J, Chen L, Qu L. Estrogen-related receptor  $\gamma$  regulates hepatic triglyceride metabolism through phospholipase A2 G12B. *FASEB J*. 2019 Jul 33(7):7942-7952.
30. Diao H, Cheng N, Zhao Y, Xu H, Dong H, Thamm DH, Zhang D, Lin D. Ivermectin inhibits canine mammary tumor growth by regulating cell cycle progression and WNT signaling. *BMC Vet Res*. 2019 Aug 2 15(1):276.
31. Wang Z, Zhang M, Shan R, Wang YJ, Chen J, Huang J, Sun LQ, Zhou WB. MTMR3 is upregulated in patients with breast cancer and regulates proliferation, cell cycle progression and autophagy in breast cancer cells. *Oncol Rep*. 2019 Nov 42(5):1915-1923.
32. Hu Z, Niu G, Ren J, Wang X, Chen L, Hong R, Ke C. TFAF5 promotes proliferation and migration in gastric cancer. *Mol Med Rep*. 2019 Nov 20(5):4477-4488.
33. Li J, Xiang X, Xu H, Shi Y. Cilostazol Promotes Angiogenesis and Increases Cell Proliferation After Myocardial Ischemia-Reperfusion Injury Through a cAMP-Dependent Mechanism. *Cardiovasc Eng Technol*. 2019 Dec 10(4):638-647.
34. Hai-Bing Xia, Hui-Ju Wang, Shu-Shu Song, Jun-Gang Zhang, Xiang-Lei He, Zhi-Ming Hu, Cheng-Wu Zhang, Dong-Sheng Huang, Xiao-Zhou Mou. Decreased DC-SIGNR expression in hepatocellular carcinoma predicts poor patient prognosis. *Oncol Lett*. 2020 Jan;19(1):69-76.
35. Yan Sun, Chuang Guo, Lin Yuan, Wen Li, Zhan-You Wang, Feng Yue, Jia-Yi Li. Cynomolgus Monkeys With Spontaneous Type-2-Diabetes-Mellitus-Like Pathology Develop Alpha-Synuclein Alterations Reminiscent of Prodromal Parkinson's Disease and Related Diseases. *FRONT NEUROSCI-SWITZ*. 2020 Feb 7;14:63.
36. Ning-Wei Zhu, Xiao-Lan Yin, Ren Lin, Xiao-Lan Fan, Shi-Jie Chen, Yuan-Ming Zhu, Xiao-Zhen Zhao. Possible mechanisms of lycopene amelioration of learning and memory impairment in rats with vascular dementia. *Neural Regen Res*. 2020 Feb;15(2):332-341.
37. Wang J, Zuo Y, Zhuang K, Luo K, Yan X, Li J, Zhang JH, Liu F. Recombinant Human Milk Fat Globule-Epidermal Growth Factor 8 Attenuates Microthrombosis after Subarachnoid Hemorrhage in Rats. *J STROKE CEREBROVASC*. 2020 Mar 29(3):104536.
38. Yuling Lian, Fangui Zhao, Wenjun Wang. Use of Bao Gui capsule in treatment of a polycystic ovary syndrome rat model. *Mol Med Rep*. 2020 Mar;21(3):1461-1470.
39. Z-Z Wang, S-C Chen, X-B Zou, L-L Tian, S-H Sui, N-Z Liu. Nesfatin-1 alleviates acute lung injury through reducing inflammation and oxidative stress via the regulation of HMGB1. *EUR REV MED PHARMACO*. 2020 May;24(9):5071-5081.
40. Xidan Zhu, Jia Feng, Wenguang Fu, Xiaojia Shu, Xue Wan, Jinbo Liu. Effects of cisplatin on the proliferation, invasion and apoptosis of breast cancer cells following  $\beta$ -catenin silencing. *Int J Mol Med*. 2020 Jun;45(6):1838-1850.
41. Xiuming Zhang, Zheng Kang, Xiaodong Xie, Wei Qiao, Lei Zhang, Zhen Gong, Yan Chen, Wenrong Shen. Silencing of HIF-1 $\alpha$  inhibited the expression of lncRNA NEAT1 to suppress development of hepatocellular carcinoma under hypoxia. *Am J Transl Res*. 2020 Jul 15;12(7):3871-3883.
42. Yiling Yang, Qianye Chen, Siru Zhou, Xinyi Gong, Hongyuan Xu, Yueyang Hong, Qinggang Dai, Lingyong Jiang. Skeletal Phenotype Analysis of a Conditional Stat3 Deletion Mouse Model. *JOVE-J VIS EXP*. 2020 Jul 3;(161).
43. Shubo Zhai, Baichao Sun, Yan Zhang, Lengyue Zhao, Li Zhang. IL-17 aggravates renal injury by promoting podocyte injury in children with primary nephrotic syndrome. *Exp Ther Med*. 2020 Jul;20(1):409-417.
44. Chengjie Ma, Peipei Peng, Yan Zhou, Tianya Liu, Lijuan Wang, Chen Lu. MicroRNA-93 promotes angiogenesis and attenuates remodeling via inactivation of the Hippo/Yap pathway by targeting Lats2 after myocardial infarction. *Mol Med Rep*. 2020 Jul;22(1):483-493.
45. Zhen Wang, Han Jiang, Lu-Yao Cai, Ning Ji, Xin Zeng, Yu Zhou, Ying-Qiang Shen, Qian-Ming Chen. Repurposing disulfiram to induce OSCC cell death by cristae dysfunction promoted autophagy. *Oral Dis*. 2020 Sep 28. doi: 10.1111/odi.13652.
46. Xiuhui Wu, Leng Liu, Hongxia Zhang. miR-802 inhibits the epithelial-mesenchymal transition, migration and invasion of cervical cancer by regulating BTF3. *Mol Med Rep*. 2020 Sep;22(3):1883-1891.
47. Bin Zhan, Linjin Huang, Yachun Chen, Wen Ye, Jingkun Li, Jianhui Chen, Sheng Yang, Wei Jiang. miR-196a-mediated downregulation of p27 kip1 protein promotes prostate cancer proliferation and relates to biochemical recurrence after radical prostatectomy. *Prostate*. 2020

- Sep;80(12):1024-1037.
48. Wenjing Wang, Wenxiao Ding, Hanpeng Huang, Yina Zhu, Ning Ding, Ganzhu Feng, Xilong Zhang. The role of mitophagy in the mechanism of genioglossal dysfunction caused by chronic intermittent hypoxia and the protective effect of adiponectin. *Sleep Breath*. 2020 Oct 8. doi: 10.1007/s11325-020-02211-0.
  49. Juan Tan, Ling Liu, Zhihua Zuo, Bin Song, Tingting Cai, Dafa Ding, Yibing Lu, Xiaolong Ye. Overexpression of novel long intergenic non-coding RNA LINC02454 is associated with a poor prognosis in papillary thyroid cancer. *Oncol Rep*. 2020 Oct;44(4):1489-1501.
  50. Yun Chen, Chaojin Hong, Xiaochen Chen, Zhiquan Qin. Demethoxycurcumin increases the sensitivity of cisplatin-resistant non-small lung cancer cells to cisplatin and induces apoptosis by activating the caspase signaling pathway. *Oncol Lett*. 2020 Nov;20(5):209.
  51. Hangshun Li, Haina Yang, Dandan Wang, Lixin Zhang, Tao Ma. Peroxiredoxin2 (Prdx2) Reduces Oxidative Stress and Apoptosis of Myocardial Cells Induced by Acute Myocardial Infarction by Inhibiting the TLR4/Nuclear Factor kappa B (NF- $\kappa$ B) Signaling Pathway. *MED SCI MONITOR*. 2020 Dec 3;26:e926281.
  52. Yingjie Lu, Dongliang Bian, Xuelin Zhang, Huibiao Zhang, Zhenghong Zhu. Inhibition of Bcl-2 and Bcl-xL overcomes the resistance to the third-generation EGFR tyrosine kinase inhibitor osimertinib in non-small cell lung cancer. *Mol Med Rep*. 2021 Jan;23(1):48.
  53. Zhi Yang, Weigang Wu, Pengcheng Ou, Minna Wu, Furong Zeng, Boping Zhou, Shipin Wu. MiR-122-5p knockdown protects against APAP-mediated liver injury through up-regulating NDRG3. *Mol Cell Biochem*. 2021 Feb;476(2):1257-1267.
  54. X-G Li, X Song, J-Y Wang, C-H Sun, Z-Q Li, L-L Meng, S-H Chi. Fibroblast growth factor 18 alleviates hyperoxia-induced lung injury in mice by adjusting oxidative stress and inflammation. *Eur Rev Med Pharmacol Sci*. 2021 Feb;25(3):1485-1494.
  55. Guangyu Ji, Wenjuan Zhou, Xian Li, Jingyi Du, Xinyue Li, Hongbo Hao. Melatonin inhibits proliferation and viability and promotes apoptosis in colorectal cancer cells via upregulation of the microRNA-34a/449a cluster. *Mol Med Rep*. 2021 Mar;23(3):187.
  56. Feifei Shao, Lingmin Zhou, Yu Zhang, Hongping Chen, Yu Zhang, Zhihui Guan. Gastrodin alleviates inflammatory injury of cardiomyocytes in septic shock mice via inhibiting NLRP3 expression. *In Vitro Cell Dev Biol Anim*. 2021 May;57(5):571-581.
  57. Xilin Zhang, Ying He, Yan Jiang, Ying Bao, Qiuqiang Chen, Dong Xie, Huanming Yu, Xiang Wang. TMM229A suppresses non-small cell lung cancer progression via inactivating the ERK pathway. *Oncol Rep*. 2021 Aug;46(2):176.
  58. Rui Wang, Jingjing Xu, Jinbo Wu, Shunheng Gao, Zhiping Wang. Angiotensin-converting enzyme 2 alleviates pulmonary artery hypertension through inhibition of focal adhesion kinase expression. *Exp Ther Med*. 2021 Oct;22(4):1165.
  59. Suna Zheng, Tianchi Zhuang, Yajun Tang, Ruihan Wu, Ting Xu, Tian Leng, Yao Wang, Zheng Lin, Minghui Ji. Leonurine protects against ulcerative colitis by alleviating inflammation and modulating intestinal microflora in mouse models. *Exp Ther Med*. 2021 Nov;22(5):1199.
  60. Bo He, Chao Wu, Weichao Sun, Yang Qiu, Jingyao Li, Zhihui Liu, Tao Jing, Haidong Wang, Yi Liao. miR-383 increases the cisplatin sensitivity of lung adenocarcinoma cells through inhibition of the RBM24-mediated NF- $\kappa$ B signaling pathway. *Int J Oncol*. 2021 Nov;59(5):87.
  61. Yu Tai, Yue Zhu, Dan Mei, Han Wang, Qianqian Yu, Chongyang Hong, Xiaoyu Cai, Li Xu, Jinru Ge, Faqin Liang, Chunru Jiang, Ziyang Xue, Ling Hu, Ruijin Liu, Tianjing Zhang, Pan Wang, Xianzheng Zhang, Feng Zhang, Wei Wei, Lingling Zhang. IgD promotes pannus formation by activating Wnt5A-Fzd5-CTHRC1-NF- $\kappa$ B signaling pathway in FLS of CIA rats and the regulation of IgD-Fc-Ig fusion protein. *Int Immunopharmacol*. 2021 Dec;101(Pt A):108261.
  62. Dehong Mao, Zhongmei He, Wei Xuan, Jiao Deng, Weichun Li, Xiaoying Fang, Linglong Li, Feng Zhang. Effect and mechanism of BDNF/TrkB signaling on vestibular compensation. *Bioengineered*. 2021 Dec;12(2):11823-11836.
  63. Huanhuan Zhu, Shaopeng Tong, Yan Cui, Xiaodong Wang, Mingyong Wang. Tanshinol alleviates ulcerative colitis by promoting the expression of VLDLR. *Drug Dev Res*. 2021 Dec;82(8):1258-1268.
  64. Jie Yang, Mengjia Sun, Ran Cheng, Hu Tan, Chuan Liu, Renzheng Chen, Jihang Zhang, Yuanqi Yang, Xubin Gao, Lan Huang. Pitavastatin activates mitophagy to protect EPC proliferation through a calcium-dependent CAMK1-PINK1 pathway in atherosclerotic mice. *Commun Biol*. 2022 Feb 10;5(1):124.
  65. Gang Yu, Wanjin Chen, Xianghua Li, Liang Yu, Yanyan Xu, Qiang Ruan, Yawei He, Yong Wang. TWIST1-EP300 Expedites Gastric Cancer Cell Resistance to Apatinib by Activating the Expression of COL1A2. *Anal Cell Pathol (Amst)*. 2022 Feb 22:2022:5374262.
  66. Yelin Song, Jie Yu, Lingling Li, Lei Wang, Liangle Dong, Guangmin Xi, Yun Jing Lu, Zuwei Li. Luteolin impacts deoxyribonucleic acid repair by modulating the mitogen-activated protein kinase pathway in colorectal cancer. *Bioengineered*. 2022 Apr;13(4):10998-11011.
  67. Zhen-Ye Zhang, Shi-Peng Dang, Shan-Shan Li, Ying Liu, Miao-Miao Qi, Ning Wang, Ling-Feng Miao, Ying Wu, Xiao-Yan Li, Chun-Xin Wang, Ling-Ling Qian, Ru-Xing Wang. Glucose Fluctuations Aggravate Myocardial Fibrosis via the Nuclear Factor- $\kappa$ B-Mediated Nucleotide-Binding Oligomerization Domain-Like Receptor Protein 3 Inflammasome Activation. *Front Cardiovasc Med*. 2022 May 3:9:748183.
  68. Li-Da Wu, Ying Liu, Feng Li, Jia-Yi Chen, Jie Zhang, Ling-Ling Qian, Ru-Xing Wang. Glucose fluctuation promotes cardiomyocyte apoptosis by triggering endoplasmic reticulum (ER) stress signaling pathway in vivo and in vitro. *Bioengineered*. 2022 May;13(5):13739-13751.
  69. Tao Luo, Yile Wang, Hailin Tang, Fei Zhou, Ying Chen, Bing Pei, Jinke Wang. An AAV-Based NF- $\kappa$ B-Targeting Gene Therapy (rAAV-DMP-miR533) to Inflammatory Diseases. *J Inflamm Res*. 2022 Jun 14:15:3447-3466.
  70. Jingxin Zhang, Hu Zhou, Pengfei Li, Huaxiang Shi, Xin Sui, Yongan Wang, Jingshan Shi, Liyun Wang. Hypothalamic response with PKA/CREB signaling is associated with direct cerebroventricular administration of bombesin-induced scratching. *Brain Res*. 2022 Aug 15:1789:147950.
  71. Xueli Tong, Chunqin Lao, Di Li, Junxi Du, Jingmian Chen, Weijie Xu, Lu Li, Huiling Ye, Xiaofeng Guo, Jiejing Li. An acetylated mannan isolated from Aloe vera induce colorectal cancer cells apoptosis via mitochondrial pathway. *Carbohydr Polym*. 2022 Sep 1:291:119464.
  72. Liping Qian, Lihua Mao, Weixing Mo, Rong Wang, Yunlong Zhang. Resveratrol Enhances the Radiosensitivity by Inducing DNA Damage and Antitumor Immunity in a Glioblastoma Rat Model under 3 T MRI Monitoring. *J Oncol*. 2022 Oct 13:2022:9672773.
  73. Zhang Junli, Wang Shuhan, Zhao Yajuan, Deng Xiaoling, Li Jiahuan, Xu Keshu. The Role and Mechanism of CREBH Regulating SIRT3 in Metabolic Associated Fatty Liver Disease. *Life Sci*. 2022 Oct 1:306:120838.
  74. Tao Luo, Yile Wang, Jinke Wang. Ferroptosis assassinate tumor. *J Nanobiotechnology*. 2022 Nov 3;20(1):467.

75. Lu Li, Weijie Xu, Yin Zhu Luo, Chunqin Lao, Xueli Tong, Junxi Du, Bihong Huang, Di Li, Jingmian Chen, Huiling Ye, Feng Cong, Xiaofeng Guo, Jiejing Li. Aloe polymeric acemannan inhibits the cytokine storm in mouse pneumonia models by modulating macrophage metabolism. *Carbohydr Polym.* 2022 Dec 1;297:120032.
76. Kunming Zhang, Xiwen Sun, Weikai Sun, Meng Wang, Fushi Han. Exosomal microRNA-506 inhibits biological activity of lung adenocarcinoma cells and increases sensitivity to cisplatin-based hyperthermia. *Cell Signal.* 2022 Dec;100:110469.
77. Qing Chen, Liping Shao, Yong Li, Mian Dai, He Liu, Nan Xiang, Hui Chen. Tanshinone IIA alleviates ovalbumin-induced allergic rhinitis symptoms by inhibiting Th2 cytokine production and mast cell histamine release in mice. *Pharm Biol.* 2022 Dec;60(1):326-333.
78. Qian Xie, Yi Wang, Guo-Liang Zou. Protective effects of lavender oil on sepsis-induced acute lung injury via regulation of the NF- $\kappa$ B pathway. *Pharm Biol.* 2022 Dec;60(1):968-978.
79. Huijun Cao, Zhiqiang Wang, Ying Wang, Lijuan Ye, Ruilei Li, Yuanbo Xue, Ke Li, Tiannan Di, Tao Li, Zonglin Fan, Yanyan Liu, Jiyin Guo, Hong Yao, Chunlei Ge. PPP1R14D promotes the proliferation, migration and invasion of lung adenocarcinoma via the PKC  $\alpha$  /BRAF/MEK/ERK signaling pathway. *Int J Oncol.* 2022 Dec;61(6):153.
80. Bin Liu, Jinchao Wang, Guangchuan Wang, Wanli Jiang, Zhen Li, Yongjun Shi, Junyong Zhang, Qingshan Pei, Guangjun Huang, Lifan Wang, Shengqiang Zhao, Lei Wu, Mingyan Zhang, Wenwen Wang, Xiao Li, Tong Mou, Chunqing Zhang, Qian Ding. Hepatocyte-derived exosomes deliver H2AFJ to hepatic stellate cells and promote liver fibrosis via the MAPK/STMN1 axis activation. *Int Immunopharmacol.* 2023 Feb;115:109605.
81. Jinfang Xu, Jianli Wang, Mingfei Zhao, Chenguang Li, Shen Hong, Jianmin Zhang. LncRNA LINC01018/miR-942-5p/KNG1 axis regulates the malignant development of glioma in vitro and in vivo. *CNS Neurosci Ther.* 2023 Feb;29(2):691-711.
82. Fanghua Chen, Yin Li, Ling Aye, Yingcheng Wu, Liangqing Dong, Zijian Yang, Qiang Gao, Shu Zhang. FUT8 is regulated by miR-122-5p and promotes malignancies in intrahepatic cholangiocarcinoma via PI3K/AKT signaling. *Cell Oncol (Dordr).* 2023 Feb;46(1):79-91.
83. Tao Li, Hong Ge, Qingyan Yang, Junmei Wang, Qian Yin, Hongbin Wang, Gaolei Hou. Oncogenic role of microRNA-19b-3p-mediated SOCS3 in glioma through activation of JAK-STAT pathway. *Metab Brain Dis.* 2023 Mar;38(3):945-960.
84. Kun Wu, Long Ma, Ting Xu, Jun Cao, Chengming Zhou, Xiangyou Yu, Yi Wang, Hao Wen. Transcription Factor YY1 Ameliorates Liver Ischemia-reperfusion Injury Through Modulating the miR-181a-5p/ESR1/ERBB2 Axis. *Transplantation.* 2023 Apr 1;107(4):878-889.
85. Shen Cheng, Chen Chen, Liling Wang. Gelsemine Exerts Neuroprotective Effects on Neonatal Mice with Hypoxic-Ischemic Brain Injury by Suppressing Inflammation and Oxidative Stress via Nrf2/HO-1 Pathway. *Neurochem Res.* 2023 May;48(5):1305-1319.

Version 2025.01.24