



碧云天生物技术/Beyotime Biotechnology
订货热线: 400-1683301或800-8283301
订货e-mail: order@beyotime.com
技术咨询: info@beyotime.com
网址: http://www.beyotime.com

Hoechst 33342

产品编号	产品名称	包装
C1022	Hoechst 33342	10mg

产品简介:

- Hoechst 33342, 也称bisBenzimide H 33342或HOE 33342。分子式为C₂₇H₂₈N₆O·3HCl·3H₂O, 分子量为615.99, CAS Number 23491-52-3。
- Hoechst 33342是一种可以穿透细胞膜的蓝色荧光染料, 对细胞的毒性较低。
- Hoechst 33342染色常用于细胞凋亡检测, 染色后用荧光显微镜观察或流式细胞仪检测。Hoechst 33342也常用于普通的细胞核染色, 或常规的DNA染色。
- Hoechst 33342的最大激发波长为346nm, 最大发射波长为460nm; Hoechst 33342和双链DNA结合后, 最大激发波长为350nm, 最大发射波长为461nm。
- Hoechst 33342溶于水, 溶解度可达20mg/ml。
- 用于细胞核染色时, 推荐的Hoechst 33342工作浓度为0.5-10μg/ml。

包装清单:

产品编号	产品名称	包装
C1022	Hoechst 33342	10mg
—	说明书	1份

保存条件:

-20°C避光保存, 一年有效。

注意事项:

- 荧光染料都存在淬灭的问题, 建议染色后尽量当天完成检测。
- 为减缓荧光淬灭可以使用抗荧光淬灭封片液。抗荧光淬灭封片液(P0126)可以向碧云天订购。
- Hoechst 33342对人体有害, 操作时请小心, 并注意有效防护以避免直接接触人体或吸入体内。
- 本产品仅限于专业人员的科学的研究用, 不得用于临床诊断或治疗, 不得用于食品或药品, 不得存放于普通住宅内。
- 为了您的安全和健康, 请穿实验服并戴一次性手套操作。

相关产品:

产品编号	产品名称	包装
C1002	DAPI	5mg/ml×0.2ml
C1005	DAPI染色液	10ml
C1006	DAPI染色液	50ml
C1011	Hoechst 33258	10mg
C1017	Hoechst 33258染色液	10ml
C1018	Hoechst 33258染色液	50ml
C1022	Hoechst 33342	10mg
C1025	Hoechst 33342染色液	10ml
C1026	Hoechst 33342染色液	50ml
C1027	Hoechst 33342活细胞染色液(100X)	0.1ml
C1028	Hoechst 33342活细胞染色液(100X)	0.5ml
C1029	Hoechst 33342活细胞染色液(100X)	3ml

使用本产品的文献:

1. Yao G, Ling L, Luan J, Ye D, Zhu P. Nonylphenol induces apoptosis of Jurkat cells by a caspase-8 dependent mechanism. Int Immunopharmacol. 2007 Apr;7(4):444-53.
2. Wang Q, Wang Y, Liang C, Song J, Chen X. Identification of a hydrophobic domain of HA2 essential to morphogenesis of Helicoverpa armigera nucleopolyhedrovirus. J Virol. 2008 Apr;82(8):4072-81.
3. Badraoui R, Blouin S, Moreau MF, Gallois Y, Rebai T, Sahnoun Z, Baslé M, Chappard D. Effect of alpha tocopherol acetate in Walker 256/B cells-induced

- oxidative damage in a rat model of breast cancer skeletal metastases. *CHEM-BIOL INTERACT.* 2009;182(2-3):98-105.
4. Han M, He CX, Fang QL, Yang XC, Diao YY, Xu DH, He QJ, Hu YZ, Liang WQ, Yang B, Gao JQ. A novel camptothecin derivative incorporated in nano-carrier induced distinguished improvement in solubility, stability and anti-tumor activity both in vitro and in vivo. *PHARM RES-DORDR.* 2009 Apr;26(4):926-35.
 5. Chen Z, Li Z, Peng G, Chen X, Yin W, Kotlikoff MI, Yuan ZQ, Ji G. Extracellular ATP-induced nuclear Ca²⁺ transient is mediated by inositol 1,4, 5-trisphosphate receptors in mouse pancreatic beta-cells. *BIOCHEM BIOPH RES CO.* 2009 May 1;382(2):381-4.
 6. Liu Y, Sheng Z, Liu H, Wen D, He Q, Wang S, Shao W, Jiang RJ, An S, Sun Y, Bendena WG, Wang J, Gilbert LI, Wilson TG, Song Q, Li S. Juvenile hormone counteracts the bHLH-PAS transcription factors MET and GCE to prevent caspase-dependent programmed cell death in *Drosophila*. *Development.* 2009 Jun;136(12):2015-25.
 7. Ma K, Hu M, Qi Y, Qiu L, Jin Y, Yu J, Li B. Structure-transfection activity relationships with glucocorticoid-polyethyl-enimine conjugate nuclear gene delivery systems. *Biomaterials.* 2009 Aug;30(22):3780-9.
 8. Zeng X, Nan F, Liang C, Song J, Wang Q, Vlak JM, Chen X. Functional analysis of the *Autographa californica* nucleopolyhedrovirus IAP1 and IAP2. *SCI CHINA LIFE SCI.* 2009 Aug;52(8):761-70.
 9. Ma K, Hu MX, Qi Y, Zou JH, Qiu LY, Jin Y, Ying XY, Sun HY. PAMAM-triamcinolone acetonide conjugate as a nucleus-targeting gene carrier for enhanced transfer activity. *Biomaterials.* 2009 Oct;30(30):6109-18.
 10. Gong Y, Wu J, Huang Y, Shen S, Han X. Nonylphenol induces apoptosis in rat testicular Sertoli cells via endoplasmic reticulum stress. *Toxicol Lett.* 2009;186(2):84-95.
 11. Yang M, Zhang M, Chen J, Mukherjee R, Zhang L, Lin S, Gu Y. Angiopoietin-1 inhibits mouse glomerular endothelial cell senescence via Tie2 receptor-modulated ERK1/2 signaling. *Am J Nephrol.* 2010;31(6):490-500.
 12. Chen YJ, Huang XB, Li ZX, Yin LL, Chen WQ, Li L. Tenuigenin protects cultured hippocampal neurons against methylglyoxal-induced neurotoxicity. *Eur J Pharmacol.* 2010;645(1-3):1-8.
 13. Zhang L, Li HZ, Gong X, Luo FL, Wang B, Hu N, Wang CD, Zhang Z, Wan JY. Protective effects of Asiaticoside on acute liver injury induced by lipopolysaccharide/D-galactosamine in mice. *Phytomedicine.* 2010;17(10):811-9.
 14. Jiang X, Zhang J, Xia M, Qiu W, Wang H, Zhao D, Wang Y. Role of activating transcription factor 3 (ATF3) in sublytic C5b-9-induced glomerular mesangial cell apoptosis. *Cell Mol Immunol.* 2010;7(2):143-51.
 15. Wu H, Lin Y, Shi Y, Qian W, Tian Z, Yu Y, Huo K. SIAH-1 interacts with mammalian polyhomeotic homologues HPH2 and affects its stability via the ubiquitin-proteasome pathway. *BIOCHEM BIOPH RES CO.* 2010;397(3):391-6.
 16. Guo XF, Zhu XF, Shang Y, Zhang SH, Zhen YS. A bispecific enediyne-energized fusion protein containing ligand-based and antibody-based oligopeptides against epidermal growth factor receptor and human epidermal growth factor receptor 2 shows potent antitumor activity. *Clin Cancer Res.* 2010 Apr 1;16(7):2085-94.
 17. Peng ZG, Fan B, Du NN, Wang YP, Gao LM, Li YH, Li YH, Liu F, You XF, Han YX, Zhao ZY, Cen S, Li JR, Song DQ, Jiang JD. Small molecular compounds that inhibit hepatitis C virus replication through destabilizing heatshock cognate 70 messenger RNA. *Hepatology.* 2010 Sep;52(3):845-53.
 18. Zhang YP, Wan P, Wang HQ, Zhao H, Xu YX, Yang R, Zhu CQ. Effect of neuronal excitotoxicity on Munc18-1 distribution in nuclei of rat hippocampal neuron and primary cultured neuron. *Neurosci Bull.* 2011 Jun;27(3):163-72.
 19. Peng ZG, Zhao ZY, Li YP, Wang YP, Hao LH, Fan B, Li YH, Wang YM, Shan YQ, Han YX, Zhu YP, Li JR, You XF, Li ZR, Jiang JD. Host apolipoprotein B messenger RNA-editing enzyme catalytic polypeptide-like 3G is an innate defensive factor and drug target against hepatitis C virus. *Hepatology.* 2011 Apr;53(4):1080-9.
 20. Hu XL, Cheng X, Cai L, Tan GH, Xu L, Feng XY, Lu TJ, Xiong H, Fei J, Xiong ZQ. Conditional deletion of NRSF in forebrain neurons accelerates epileptogenesis in the kindling model. *Cereb Cortex.* 2011 Sep;21(9):2158-65.
 21. He XN, Su F, Lou ZZ, Jia WZ, Song YL, Chang HY, Wu YH, Lan J, He XY, Zhang Y. Ipr1 gene mediates RAW 264.7 macrophage cell line resistance to *Mycobacterium bovis*. *Scand J Immunol.* 2011 Nov;74(5):438-44.
 22. Wu H, Shi Y, Lin Y, Qian W, Yu Y, Huo K. Eukaryotic translation elongation factor 1 delta inhibits the ubiquitin ligase activity of SIAH-1. *Mol Cell Biochem.* 2011 Nov;357(1-2):209-15.
 23. Yang H, Taylor HS, Lei C, Cheng C, Zhang W. Hormonal regulation of galectin 3 in trophoblasts and its effects on endometrium. *Reprod Sci.* 2011 Nov;18(11):1118-27.
 24. Wang H, Xie X, Zhang F, Zhou Q, Tao Q, Zou Y, Chen C, Zhou C, Yu S. Evaluation of cholesterol depletion as a marker of nephrotoxicity in vitro for novel β-cyclodextrin derivatives. *Food Chem Toxicol.* 2011 Jun;49(6):1387-93.
 25. Zhou J, Lin J, Zhou C, Deng X, Xia B. Cytotoxicity of red fluorescent protein DsRed is associated with the suppression of Bcl-xL translation. *FEBS Lett.* 2011 Mar 9;585(5):821-7.
 26. Liu Q, Kou JP, Yu BY. Ginsenoside Rg1 protects against hydrogen peroxide-induced cell death in PC12 cells via inhibiting NF-κB activation. *J NEUROCHEM.* 2011 Jan;58(1):119-25.
 27. Kong FL, Cheng W, Chen J, Liang Y. d-Ribose glycates β(2)-microglobulin to form aggregates with high cytotoxicity through a ROS-mediated pathway. *CHEM-BIOL INTERACT.* 2011 Oct 15;194(1):69-78.
 28. Ji SZ, Xiao SC, Luo PF, Huang GF, Wang GY, Zhu SH, Wu MJ, Xia ZF. An epidermal stem cells niche microenvironment created by engineered human amniotic membrane. *Biomaterials.* 2011 Nov;32(31):7801-11.
 29. Zhang X, Wang J, Zhou Q, Xu Y, Pu S, Wu J, Xue Y, Tian Y, Lu J, Jiang W, Du D. Brain-derived neurotrophic factor-activated astrocytes produce mechanical allodynia in neuropathic pain. *Neuroscience.* 2011 Dec 29;199:452-60.
 30. Li Z, Mi L, Wang PN, Chen JY. Study on the visible-light-induced photokilling effect of nitrogen-doped TiO₂ nanoparticles on cancer cells. *Nanoscale Res Lett.* 2011 Apr 21;6(1):356.
 31. Liu SL, Zhang ZL, Sun EZ, Peng J, Xie M, Tian ZQ, Lin Y, Pang DW. Visualizing the endocytic and exocytic processes of wheat germ agglutinin by quantum dot-based single-particle tracking. *Biomaterials.* 2011 Oct;32(30):7616-24.
 32. Bai D, Gao Q, Li C, Ge L, Gao Y, Wang H. A conserved TGFβ1/HuR feedback circuit regulates the fibrogenic response in fibroblasts. *Cell Signal.* 2012 Jul;24(7):1426-32.
 33. Li B, Xu H, Li Z, Yao M, Xie M, Shen H, Shen S, Wang X, Jin Y. Bypassing multidrug resistance in human breast cancer cells with lipid/polymer particle assemblies. *INT J NANOMED.* 2012;7:187-97.
 34. Wang S, Tan Y, Lei W, Wang Y, Zhou X, Jia X, Zhang K, Chu L, Liu X, Qian W. Complete eradication of xenograft hepatoma by oncolytic adenovirus ZD55 harboring TRAIL-IETD-Smac gene with broad antitumor effect. *Hum Gene Ther.* 2012 Sep;23(9):992-1002.
 35. Zhai W, He C, Wu L, Zhou Y, Chen H, Chang J, Zhang H. Degradation of hollow mesoporous silica nanoparticles in human umbilical vein endothelial cells. *J BIOMED MATER RES B.* 2012 Jul;100(5):1397-403.
 36. Zhang J, Zhou F, Wu X, Zhang X, Chen Y, Zha BS, Niu F, Lu M, Hao G, Sun Y, Sun J, Peng Y, Wang G. Cellular pharmacokinetic mechanisms of adriamycin resistance and its modulation by 20(S)-ginsenoside Rh2 in MCF-7/Adr cells. *BRIT J PHARMACOL.* 2012 Jan;165(1):120-34.
 37. Zhang KJ, Qian J, Wang SB, Yang Y. Targeting Gene-Viro-Therapy with AFP driving Apoptin gene shows potent antitumor effect in hepatocarcinoma. *J Biomed Sci.* 2012 Feb 9;19:20.
 38. He G, Lei W, Wang S, Xiao R, Guo K, Xia Y, Zhou X, Zhang K, Liu X, Wang Y. Overexpression of tumor suppressor TSLC1 by a survivin-regulated oncolytic adenovirus significantly inhibits hepatocellular carcinoma growth. *J CANCER RES CLIN.* 2012 Apr;138(4):657-70.
 39. Xiong XY, Guo L, Gong YC, Li ZL, Li YP, Liu ZY, Zhou M. In vitro & in vivo targeting behaviors of biotinylated Pluronic F127/poly(lactic acid) nanoparticlesthrough biotin-avidin interaction. *Eur J Pharm Sci.* 2012 Aug 15;46(5):537-44.
 40. Yin QQ, Dong CF, Dong SQ, Dong XL, Hong Y, Hou XY, Luo DZ, Pei JJ, Liu XP. AGEs Induce Cell Death via Oxidative and Endoplasmic Reticulum Stresses in

- Both Human SH-SY5Y Neuroblastoma Cells and Rat Cortical Neurons. *Cell Mol Neurobiol.* 2012 Nov;32(8):1299-309.
41. Li C, Nie H, Wang M, Su L, Li J, Yu B, Wei M, Ju J, Yu Y, Yan M, Gu Q, Zhu Z, Liu B. MicroRNA-409-3p regulates cell proliferation and apoptosis by targeting PHF10 in gastric cancer. *Cancer Lett.* 2012 Jul 28;320(2):189-97.
 42. Guo HF, Li ZS, Dong SW, Chen WJ, Deng L, Wang YF, Ying DJ. Piezoelectric PU/PVDF electrospun scaffolds for wound healing applications. *COLLOID SURFACE B.* 2012 Aug 1;96:29-36.
 43. Yin T, Wang P, Zheng R, Zheng B, Cheng D, Zhang X, Shuai X. Nanobubbles for enhanced ultrasound imaging of tumors. *INT J NANOMED.* 2012;7:895-904.
 44. Zheng J, Chen Z, Yin W, Miao L, Zhou Z, Ji G. Ryanodine receptors are involved in nuclear calcium oscillation in primary pancreatic β -cells. *BIOCHEM BIOPH RES CO.* 2012 Jun 29;423(2):207-11.
 45. Tu Z, Li H, Ma Y, Tang B, Tian J, Akers W, Achilefu S, Gu Y. The enhanced antiproliferative response to combined treatment of trichostatin A with raloxifene in MCF-7 breast cancer cells and its relevance to estrogen receptor β expression. *Mol Cell Biochem.* 2012 Jul;366(1-2):111-22.
 46. Zheng X, Tian J, Weng L, Wu L, Jin Q, Zhao J, Wang L. Cytotoxicity of cadmium-containing quantum dots based on a study using a microfluidic chip. *Nanotechnology.* 2012 Feb 10;23(5):055102.
 47. Tu Z, Ma Y, Tian J, Li H, Akers W, Achilefu S, Gu Y. Estrogen receptor β potentiates the antiproliferative effect of raloxifene and affects the cell migration and invasion in HCT-116 colon cancer cells. *J CANCER RES CLIN.* 2012 Jul;138(7):1091-103.
 48. Li S, Zhang Z, Xue J, Liu A, Zhang H. Cold-inducible RNA binding protein inhibits H2O2-induced apoptosis in rat cortical neurons. *Brain Res.* 2012 Mar 2;1441:47-52.
 49. Cheng J, Tang W, Su Z, Guo J, Tong L, Wei Q. Calcineurin subunit B promotes TNF-alpha-induced apoptosis by binding to mitochondria and causing mitochondrial Ca2+ overload. *Cancer Lett.* 2012 Aug 28;321(2):169-78.
 50. Chen LP, Li ZF, Ping M, Li R, Liu J, Xie XH, Song XJ, Guo L. Regulation of Olig2 during astroglial differentiation in the subventricular zone of a cuprizone-induced demyelination mouse model. *Neuroscience.* 2012 Sep 27;221:96-107.
 51. Yu J, Xie X, Zheng M, Yu L, Zhang L, Zhao J, Jiang D, Che X. Fabrication and characterization of nuclear localization signal-conjugated glycol chitosan micelles for improving the nuclear delivery of doxorubicin. *INT J NANOMED.* 2012;7:5079-90.
 52. Zhang KJ, Zhang J, Wu YM, Qian J, Liu XJ, Yan LC, Zhou XM, Xiao RJ, Wang YG, Cao X, Wei N, Liu XR, Tang B, Jiao XY, Chen K, Liu XY. Complete eradication of hepatomas using an oncolytic adenovirus containing AFP promoter controlling E1A and an E1B deletion to drive IL-24 expression. *Cancer Gene Ther.* 2012 Sep;19(9):619-29.
 53. Du J, Tang B, Wang J, Sui H, Jin X, Wang L, Wang Z. Antiproliferative effect of alpinetin in BxPC-3 pancreatic cancer cells. *Int J Mol Med.* 2012 Apr;29(4):607-12.
 54. Zhang B, Wang Y, Pang X. Enhanced radiosensitivity of EC109 cells by inhibition of HDAC1 expression. *Med Oncol.* 2012 Mar;29(1):340-8.
 55. Shen HJ, Shi H, Ma K, Xie M, Tang LL, Shen S, Li B, Wang XS, Jin Y. Polyelectrolyte capsules packaging BSA gels for pH-controlled drug loading and release and their antitumor activity. *Acta Biomater.* 2013 Apr;9(4):6123-33.
 56. Peng Y, Zhang J, Xu H, He J, Ying X, Wang Y. Neuroprotective effect of protease-activated receptor-2 in the hypoxia-induced apoptosis of rat RGC-5 cells. *J Mol Neurosci.* 2013 May;50(1):98-108.
 57. Li J, Cheng Y, Zhang X, Zheng L, Han Z, Li P, Xiao Y, Zhang Q, Wang F. The in vivo immunomodulatory and synergistic anti-tumor activity of thymosin α 1-thymopentin fusion peptide and its binding to TLR2. *Cancer Lett.* 2013 Sep 1;337(2):237-47.
 58. Shen H, Shi H, Xie M, Ma K, Li B, Shen S, Wang X, Jin Y. Biodegradable chitosan/alginate BSA-gel-capsules for pH-controlled loading and release of doxorubicin and treatment of pulmonary melanoma. *J MATER CHEM B.* 2013 May;32:3906-17.
 59. Lei W, Liu HB, Wang SB, Zhou XM, Zheng SD, Guo KN, Ma BY, Xia YL, Tan WS, Liu XY, Wang YG. Tumor suppressor in lung cancer-1 (TSLC1) mediated by dual-regulated oncolytic adenovirus exerts specific tumor actions in a mouse model. *Acta Pharmacol Sin.* 2013 Apr;34(4):531-40.
 60. Han F, Xu Y, Jiang D, Qin Y, Chen H. A BODIPY-derived fluorescent probe for cellular pH measurements. *Anal Biochem.* 2013 Apr 15;435(2):106-13.
 61. Liu Y, He G, Wang Y, Guan X, Pang X, Zhang B. MCM-2 is a therapeutic target of Trichostatin A in colon cancer cells. *Toxicol Lett.* 2013 Jul 31;221(1):23-30.
 62. Zhu K, Guo C, Xia Y, Lai H, Yang W, Wang Y, Song D, Wang C. Transplantation of novel vascular endothelial growth factor gene delivery system manipulated skeletal myoblasts promote myocardial repair. *Int J Cardiol.* 2013 Oct 3;168(3):2622-31.
 63. Wang J, Chen Y, Zhang W, Zheng G, Meng S, Che H, Ke T, Yang J, Chen J, Luo W. Akt Activation Protects Liver Cells from Apoptosis in Rats during Acute Cold Exposure. *Int J Biol Sci.* 2013 May 25;9(5):509-17.
 64. Chen Q, Sun J. Implication of Oxidative Stress in Dose-Dependent Toxicity of Silica Nanoparticles in Hepatocytes. *ADV MATER.* 2013;647:42-5.
 65. Xu C, Liu C, Huang W, Tu S, Wan F. Effect of Mst1 overexpression on the growth of human hepatocellular carcinoma HepG2 cells and the sensitivity to cisplatin in vitro. *ACTA BIOCH BIOPH SIN.* 2013 Apr;45(4):268-79.
 66. Yin T, Wang P, Li J, Zheng R, Zheng B, Cheng D, Li R, Lai J, Shuai X. Ultrasound-sensitive siRNA-loaded nanobubbles formed by hetero-assembly of polymeric micelles and liposomes and their therapeutic effect in gliomas. *Biomaterials.* 2013 Jun;34(18):4532-43.
 67. Wang B, Yu T, Dong X, Zhang Z, Song L, Xu Y, Zhang XH. Edwardsiella tarda invasion of fish cell lines and the activation of divergent cell death pathways. *Vet Microbiol.* 2013 May 3;163(3-4):282-9.
 68. Wang JJ, Liu YL, Sun YC, Ge W, Wang YY, Dyce PW, Hou R, Shen W. Basic Fibroblast Growth Factor Stimulates the Proliferation of Bone Marrow Mesenchymal Stem Cells in Giant Panda (*Ailuropoda melanoleuca*). *PLoS One.* 2015 Sep 16;10(9):e0137712.
 69. Hou L, Li Y, Song H, Zhang Z, Sun Y, Zhang X, Wu K. Protective Macroautophagy Is Involved in Vitamin E Succinate Effects on Human Gastric Carcinoma Cell Line SGC-7901 by Inhibiting mTOR Axis Phosphorylation. *PLoS One.* 2015 Jul 13;10(7):e0132829.
 70. Wu X, Zhao B, Cheng Y, Yang Y, Huang C, Meng X, Wu B, Zhang L, Lv X, Li J. Melittin induces PTCH1 expression by down-regulating MeCP2 in human hepatocellular carcinoma SMMC-7721 cells. *TOXICOL APPL PHARM.* 2015 Oct 1;288(1):74-83.
 71. Gao S, Yu Y, Ma ZY, Sun H, Zhang YL, Wang XT, Wang C, Fan WM, Zheng QY, Ma CL. NMDAR-Mediated Hippocampal Neuronal Death is Exacerbated by Activities of ASIC1a. *Neurotox Res.* 2015 Aug;28(2):122-37.
 72. Wang X, Ma Z, Fu Z, Gao S, Yang L, Jin Y, Sun H, Wang C, Fan W, Chen L, Zheng QY, Bi G, Ma CL. Hydroxyafflor Yellow A Protects Neurons From Excitotoxic Death through Inhibition of NMDARs. *ASN Neuro.* 2016 Apr 10;8(2). pii: 1759091416642345.
 73. Yang F, Yan G, Li Y, Han Z, Zhang L, Chen S, Feng C, Huang Q, Ding F, Yu Y, Bi C, Cai B, Yang L. Astragalus Polysaccharide Attenuated Iron Overload-Induced Dysfunction of Mesenchymal Stem Cells via Suppressing Mitochondrial ROS. *CELL PHYSIOL BIOCHEM.* 2016;39(4):1369-79.
 74. Li L, Liu JC, Lai FN, Liu HQ, Zhang XF, Dyce PW, Shen W, Chen H. Di (2-ethylhexyl) Phthalate Exposure Impairs Growth of Antral Follicle in Mice. *PLoS One.* 2016 Feb 4;11(2):e0148350.
 75. Wang Y, Qiu Y, Yin S, Zhang L, Shi K, Gao H, Zhang Z, He Q. A functional nanocarrier that copenetrates extracellular matrix and multiple layers of tumor cells for sequential and deep tumor autophagy inhibitor and chemotherapeutic delivery. *Autophagy.* 2016 Dec 2:1-12.
 76. Liu Y, Zhi D, Li M, Liu D, Wang X, Wu Z, Zhang Z, Fei D, Li Y, Zhu H, Xie Q, Yang H, Li H. Shengmai Formula suppressed over-activated Ras/MAPK pathway in *C. elegans* by opening mitochondrial permeability transition pore via regulating cyclophilin D. *SCI REP-UK.* 2016 Dec 16;6:38934.

77. You F, Li Q, Jin G, Zheng Y, Chen J, Yang H. Genistein protects against A β 25-35 induced apoptosis of PC12cells through JNK signaling and modulation of Bcl-2 family messengers. *BMC Neurosci.* 2017 Jan;12(1):12.
78. Yan X, Xun M, Dou X, Wu L, Han Y, Zheng J. Regulation of Na $^{+}$ -K $^{+}$ -ATPase effected high glucose-induced myocardial cell injury through c-Srcdependent NADPH oxidase/ROS pathway. *Exp Cell Res.* 2017 Aug 15;357(2):243-251.
79. Wang Z, Liao K, Zuo W, Liu X, Qiu Z, Gong Z, Liu C, Zeng Q, Qian Y, Jiang L, Bu Y, Hong S, Hu G. Depletion of NFBD1/MDC1 Induces Apoptosis in Nasopharyngeal Carcinoma Cells Through the p53-ROS-Mitochondrial Pathway. *Oncol Res.* 2017 Jan 2;25(1):123-136.
80. Hu Y, Xiong LL, Zhang P, Wang TH. Microarray expression profiles of genes in lung tissues of rats subjected to focal cerebral ischemia-induced lung injury following bone marrow-derived mesenchymal stem cell transplantation. *Int J Mol Med.* 2017 Jan;39(1):57-70.
81. Geng YD, Zhang C, Lei JL, Yu P, Xia YZ, Zhang H, Yang L, Kong LY. Walsuronoid B induces mitochondrial and lysosomal dysfunction leading to apoptotic rather than autophagic cell death via ROS/p53 signaling pathways in liver cancer. *Biochem Pharmacol.* 2017 Oct 15;142:71-86.
82. Fan Y, Liu J, Liu D, Zhou Z, Bao Y, Wang J, Zhao Q, Xu Y. NSCA-1-a novel N-substituted coumalamide derivative-increases Adriamycin sensitivity in HepG2/adriamycin cells through modulating Akt/GSK-3 β signaling and p53-dependant apoptotic pathway. *ENVIRON TOXICOL CHEM.* 2017 Jan;49:1-7.
83. Chen GH, Xu CS, Zhang J, Li Q, Cui HH, Li XD, Chang LP, Tang RJ, Xu JY, Tian XQ, Huang PS, Xu J, Jin C, Yang YJ. Inhibition of miR-128-3p by Tongxinluo Protects Human Cardiomyocytes from Ischemia/reperfusion Injury via Upregulation of p70s6k1/p-p70s6k1. *Front Pharmacol.* 2017 Oct 30;8:775.
84. Hao J, Yang Z, Wang L, Zhang Y, Shu Y, Jiang L, Hu Y, Lv W, Dong P, Liu Y. Downregulation of BRD4 inhibits gallbladder cancer proliferation and metastasis and induces apoptosis via PI3K/AKT pathway. *Int J Oncol.* 2017 Sep;51(3):823-831.
85. Liu W, Wang X, Liu Z, Wang Y, Yin B, Yu P, Duan X, Liao Z, Chen Y, Liu C, Li X, Dai Y, Tao Z. SGK1 inhibition induces autophagy-dependent apoptosis via the mTOR-Foxo3a pathway. *BRIT J CANCER.* 2017 Oct 10;117(8):1139-1153.
86. Liu X, Qiu Z, Wang Z, Zuo W, Gong Z, Liu C, Zeng Q, Qian Y, Jiang L, Li Y, Bu Y, Hu G. NFBD1/MDC1 participates inthe regulation of proliferation and apoptosis in human laryngealsquamous cell carcinoma. *Clin Transl Oncol.* 2017 Sep 18.
87. Liu JC, Lai FN, Li L, Sun XF, Cheng SF, Ge W, Wang YF, Li L, Zhang XF, De Felici M, Dyce PW, Shen W. Di (2-ethylhexyl) phthalate exposure impairs meiotic progression and DNA damage repair in fetalmouse oocytes in vitro. *Cell Death Dis.* 2017 Aug 3;8(8):e2966.
88. Ying C, Xiao BD, Qin Y, Wang BR, Liu XY, Wang RW, Fang L, Yan H, Zhou XM, Wang YG. GOLPH2-regulated oncolytic adenovirus, GD55, exerts strong killing effect on human prostatecancer stem-like cells in vitro and in vivo. *Acta Pharmacol Sin.* 2017 Sep 7.
89. Zhai Y1, Li Y2, Wang Y3, Cui J2, Feng K4, Kong X2, Chen L5. Psoralidin, a prenylated coumestan, as a novel anti-osteoporosis candidate to enhance boneformation of osteoblasts and decrease bone resorption of osteoclasts. *Eur J Pharmacol.* 2017 Apr 15;801:62-71.
90. Wang Y1, Qiu Y1, Yin S1, Zhang L1, Shi K1, Gao H1, Zhang Z1, He Q1. A functional nanocarrier that copenetrates extracellular matrix and multiple layers of tumor cells for sequential and deep tumor autophagy inhibitor and chemotherapeutic delivery. *Autophagy.* 2017 Feb;13(2):359-370.
91. Yang LL1, Zhao Y1, Luo SM1, Ma JY1, Ge ZJ1, Shen W1, Yin S2. Toxic Effects and Possible Mechanisms of HydrogenSulfide and/or Ammonia on Porcine OocyteMaturation in vitro. *Toxicol Lett.* 2017 Dec 29. pii: S0378-4274(17)31529-1.
92. Guo Y, Song Z, Zhou M, Yang Y, Zhao Y, Liu B, Zhang X. Infiltrating macrophages in diabetic nephropathy promote podocytes apoptosis via TNF- α -ROS-p38MAPK pathway. *ONCOTARGET.* 2017 Jun 7;8(32):53276-53287.
93. Sun L, Li Y, Jia X, Wang Q, Li Y, Hu M, Tian L, Yang J, Xing W, Zhang W, Wang J, Xu H, Wang L, Zhang D, Ren H. Neuroprotection byIFN- γ via astrocyte-secreted IL-6 in acute neuroinflammation. *ONCOTARGET.* 2017 Jun 20;8(25):40065-40078.
94. Zhao C, Yao X, Chen X, Wu W, Xi F, Yang G, Yu T. Knockdown of ubiquitin D inhibits adipogenesis duringthe differentiation of porcineintramuscular and subcutaneous preadipocytes. *CELL PROLIFERAT.* 2017 Nov 23.
95. Wu X, Gao H, Sun W, Yu J, Hu H, Xu Q, Chen X. Nepetoidin B, α Natural Product, Inhibits LPS-stimulated Nitric Oxide Production via Modulation of iNOS Mediated by NF- κ B/MKP-5 Pathways. *Phytother Res.* 2017 Jul;31(7):1072-1077.
96. Xiong LL, Liu F, Lu BT, Zhao WL, Dong XJ, Liu J, Zhang RP, Zhang P, Wang TH. Bone Marrow Mesenchymal Stem-Cell Transplantation Promotes Functional ImprovementAssociated with CNTF-STAT3 Activation after Hemi-Sectioned Spinal Cord Injury in Tree Shrews. *Front Cell Neurosci.* 2017 Jun 28;11:172.
97. Li Y, Sun Y, Cao T, Su Q, Li Z, Huang M, Ouyang R, Chang H, Zhang S, Miao Y. A cation-exchange controlled core-shell MnS@Bi2S3 theranostic platform for multimodal imaging guided radiation therapy with hyperthermia boost. *Nanoscale.* 2017 Oct 5;9(38):14364-14375.
98. Li X, Yan X, Wang Y, Wang J, Zhou F, Wang H, Xie W, Kong H. NLRP3 inflammasome inhibition attenuates silica-induced epithelial to mesenchymal transition (EMT) in human bronchial epithelial cells. *Exp Cell Res.* 2017 Dec 16. pii: S0014-4827(17)30668-7.
99. Chang S, Ruan WC, Xu YZ, Wang YJ, Pang J, Zhang LY, Liao H, Pang T. The natural product 4,10-aromadendranediol induces neuritogenesis in neuronal cells in vitro through activation of the ERK pathway. *Acta Pharmacol Sin.* 2017 Jan;38(1):29-40.
100. Sun Y, Wang C, Meng Q, Liu Z, Huo X, Sun P, Sun H, Ma X, Peng J, Liu K. Targeting P-glycoprotein and SORCIN: Dihydromyricetin strengthens anti-proliferative efficiency of adriamycin via MAPK/ERK and Ca $^{2+}$ -mediated apoptosis pathways in MCF-7/ADR and K562/ADR. *J Cell Physiol.* 2018 Apr;233(4):3066-3079.
101. Zhai QY, Ge W, Wang JJ, Sun XF, Ma JM, Liu JC, Zhao Y, Feng YZ, Dyce PW, De Felici M, Shen W. Exposure to Zinc oxide nanoparticles during pregnancy induces oocyte DNA damage and affects ovarian reserve of mouse offspring. *AGING-US.* 2018 Aug 28;10(8):2170-2189.
102. Zhou B, Zhang JY, Liu XS, Chen HZ, Ai YI, Cheng K, Sun RY, Zhou D, Han J, Wu Q. Tom20 senses iron-activated ROS signaling to promote melanoma cell pyroptosis. *Cell Res.* 2018 Dec;28(12):1171-1185.
103. Sun ZY, Zhang P, Wang JJ, Liu JC, Li L, Shen W, Zhai QY. Melatonin alleviates meiotic defects in fetal mouse oocytes induced by Di (2-ethylhexyl) phthalate in vitro. *AGING-US.* 2018 Dec 26;10(12):4175-4187.
104. Zhang TY, Wu RY, Zhao Y, Xu CS, Zhang WD, Ge W, Liu J, Sun ZY, Zou SH, Shen W. Ochratoxin A exposure decreased sperm motility via the AMPK and PTEN signaling pathways. *TOXICOL APPL PHARM.* 2018 Feb 1;340:49-57.
105. Tang X, Wu S, Shen L, Wei Y, Cao X, Wang Y, Long C, Zhou Y, Li D, Huang F, Liu B, Wei G. Di-(2-ethylhexyl) phthalate (DEHP)-induced testicular toxicity through Nrf2-mediated Notch1 signaling pathway in Sprague-Dawley rats. *Environ Toxicol.* 2018 Jul;33(7):720-728.
106. Zhang GL, Song JL, Zhou Y, Zhang RQ, Cheng SF, Sun XF, Qin GQ, Shen W, Li L. Differentiation of sow and mouse ovarian granulosa cells exposed to zearalenone in vitro using RNA-seq gene expression. *TOXICOL APPL PHARM.* 2018 Jul 1;350:78-90.
107. Zhang GL, Song JL, Ji CL, Feng YL, Yu J, Nyachoti CM, Yang GS. Zearalenone Exposure Enhanced the Expression of Tumorigenesis Genes in Donkey Granulosa Cells via the PTEN/PI3K/AKT Signaling Pathway. *Front Genet.* 2018 Jul 31;9:293.
108. Wang JJ, Yu XW, Wu RY, Sun XF, Cheng SF, Ge W, Liu JC, Li YP, Liu J, Zou SH, De Felici M, Shen W. Starvation during pregnancy impairs fetal oogenesis and folliculogenesis in offspring in the mouse. *Cell Death Dis.* 2018 May 1;9(5):452.
109. Zhao Y, Zhao R, Wu J, Wang Q, Pang K, Shi Q, Gao Q, Hu Y, Dong X, Zhang J, Sun J. Melatonin protects against A β -induced neurotoxicity in primary neurons via miR-132/PTEN/AKT/FOXO3a pathway. *Biofactors.* 2018 Nov;44(6):609-618.
110. Shen L, Tang X, Wei Y, Long C, Tan B, Wu S, Sun M, Zhou Y, Cao X, Wei G. Vitamin E and vitamin C attenuate Di-(2-ethylhexyl) phthalate-induced blood-testis barrier disruption by p38 MAPK in immature SD rats. *Reprod Toxicol.* 2018 Oct;81:17-27.

111. Zhang RQ, Sun XF, Wu RY, Cheng SF, Zhang GL, Zhai QY, Liu XL, Zhao Y, Shen W, Li L. Zearalenone exposure elevated the expression of tumorigenesis genes in mouse ovarian granulosa cells. *TOXICOL APPL PHARM*. 2018 Oct 1;356:191-203.
112. Zhang T, Zhou Y, Li L, Zhao Y, De Felici M, Reiter RJ, Shen W. Melatonin protects prepuberal testis from deleterious effects of bisphenol A or diethylhexyl phthalate by preserving H3K9 methylation. *J Pineal Res*. 2018 Sep;65(2):e12497.
113. Zhu P, Liu Z, Zhou J, Chen Y. Tanshinol inhibits the growth, migration and invasion of hepatocellular carcinoma cells via regulating the PI3K-AKT signaling pathway. *ONCOTARGETS THER*. 12:87-99.
114. Sun Y, Liu W, Wang C, Meng Q, Liu Z, Huo X, Yang X, Sun P, Sun H, Ma X, Peng J, Liu K. Combination of dihydromyricetin and ondansetron strengthens antiproliferative efficiency of adriamycin in K562/ADR through downregulation of SORCIN: A new strategy of inhibiting P-glycoprotein. *J Cell Physiol*. 234(4):3685-3696.
115. Huang Z, Li Q, Yuan Y, Zhang C, Wu L, Liu X, Cao W, Guo H, Duan S, Xu X, Zhang B, Xing C. Renalase attenuates mitochondrial fission in cisplatin-induced acute kidney injury via modulating sirtuin-3. *Life Sci*. 222:78-87.
116. Su W, Wang Y, Wang F, Sun S, Li M, Shen Y, Yang H. Hsa_circ_0005379 regulates malignant behavior of oral squamous cell carcinoma through the EGFR pathway. *BMC Cancer*. 19(1):400.
117. Zhu D, Huang R, Fu P, Chen L, Luo L, Chu P, He L, Li Y, Liao L, Zhu Z, Wang Y. Investigating the Role of BATF3 in Grass Carp (Ctenopharyngodon idella) Immune Modulation: A Fundamental Functional Analysis. *Int J Mol Sci*. 20(7). pii: E1687.
118. Wang JJ, Zhai QY, Zhang RQ, Ge W, Liu JC, Li L, Sun ZY, De Felici M, Shen W. Effects of activin A on the transcriptome of mouse oogenesis in vitro. *J Cell Physiol*. 234(8):14339-14350.
119. Yin W, Yang L, Kong D, Nie Y, Liang Y, Teng CB. Guanine-rich RNA binding protein GRSF1 inhibits myoblast differentiation through repressing mitochondrial ROS production. *Exp Cell Res*. 381(1):139-149.
120. Chen J, Chen X, Yao J, Li M, Yang X. The combination of Decitabine and EPZ-6438 effectively facilitate adipogenic differentiation of induced pluripotent stem cell-derived mesenchymal stem cells. *BIOCHEM BIOPH RES CO*. 516(1):307-312.
121. Jiang S, Lyu C, Zhao P, Li W, Kong W, Huang C, Genin GM, Du Y. Cryoprotectant enables structural control of porous scaffolds for exploration of cellular mechano-responsiveness in 3D. *Nat Commun*. 10(1):3491.
122. Duan H, Lei Z, Xu F, Pan T, Lu D, Ding P, Zhu C, Pan C, Zhang S. PARK2 Suppresses Proliferation and Tumorigenicity in Non-small Cell Lung Cancer. *Front Oncol*. 9:790.
123. Chu P, He L, Zhu D, Huang R, Liao L, Li Y, Zhu Z, Wang Y. Identification, expression and functional characterisation of CYP1A in grass carp (Ctenopharyngodon idella). *FISH SHELLFISH IMMUN*. 95:35-43.
124. Han X, Wang R, Xu J, Chen Q, Liang C, Chen J, Zhao J, Chu J, Fan Q, Archibong E, Jiang L, Wang C, Liu Z. In situ thermal ablation of tumors in combination with nano-adjuvant and immune checkpoint blockade to inhibit cancer metastasis and recurrence. *Biomaterials*. 224:119490.
125. Yu X, Hao M, Liu Y, Ma X, Lin W, Xu Q, Zhou H, Shao N, Kuang H. Liraglutide ameliorates non-alcoholic steatohepatitis by inhibiting NLRP3 inflammasome and pyroptosis activation via mitophagy. *Eur J Pharmacol*. 864:172715.
126. Guo Q, Li J, Zhong L, Cai H, Wang H, Yan H, Xia G, Kang Y, Wang C. Prostaglandin-E2 deficiency during late pregnancy and the associated increase in interleukin-1 β derived from periaortic lymph nodes lead to abortion. *Mol Hum Reprod*. 25(12):825-837.
127. Li T, Chen Q, Zheng Y, Zhang P, Chen X, Lu J, Lv Y, Sun S, Zeng W. PAMAM-cRGD mediating efficient siRNA delivery to spermatogonial stem cells. *Stem Cell Res Ther*. 10(1):399.
128. Zhang N, Zeng X, Sun C, Guo H, Wang T, Wei L, Zhang Y, Zhao J, Ma X. LncRNA LINC00963 Promotes Tumorigenesis and Radioresistance in Breast Cancer by Sponging miR-324-3p and Inducing ACK1 Expression. *MOL THER-NUCL ACIDS*. 18:871-881.
129. Jiang Y, Zhang T, Vallesi A, Yang X, Gao F. Time-course analysis of nuclear events during conjugation in the marine ciliate *Euplotes vannus* and comparison with other ciliates (Protozoa, Ciliophora). *Cell Cycle*. 18(3):288-298.
130. Cao P, Feng Y, Deng M, Li J, Cai H, Meng Q, Fang W, Li Y, Ke J, Long X. MiR-15b is a key regulator of proliferation and apoptosis of chondrocytes from patients with condylar hyperplasia by targeting IGF1, IGF1R and BCL2. *OSTEOARTHR CARTILAGE*. 27(2):336-346.
131. Yang LL, Cui YX, Ma JY, Ge ZJ, Shen W, Yin S. Tributyltin oxide exposure impairs mouse oocyte maturation and its possible mechanisms. *J Cell Biochem*. 120(1):715-726.
132. Chu P, He L, Zhu D, Chen L, Huang R, Liao L, Li Y, Zhu Z, Wang Y. Identification, characterisation and preliminary functional analysis of IRAK-M in grass carp (Ctenopharyngodon idella). *FISH SHELLFISH IMMUN*. 84:312-321.
133. Zhou R, Wang H, Yang Y, Zhang C, Dong X, Du J, Yan L, Zhang G, Gu Z, Zhao Y. Tumor microenvironment-manipulated radiocatalytic sensitizer based on bismuth heteropolytungstate for radiotherapy enhancement. *Biomaterials*. 189:11-22.
134. Han B, Xu K, Liu Z, Ge W, Shao S, Li P, Yan N, Li X, Zhang Z. Oral yeast-based DNA vaccine confers effective protection from *Aeromonas hydrophila* infection on *Carassius auratus*. *FISH SHELLFISH IMMUN*. 84:948-954.
135. Yang R, Xing L, Zheng X, Sun Y, Wang X, Chen J. The circRNA circAGFG1 acts as a sponge of miR-195-5p to promote triple-negative breast cancer progression through regulating CCNE1 expression. *Mol Cancer*. 18(1):4.
136. Liu S, Xing Y, Wang J, Pan R, Li G, Tang H, Chen G, Yan L, Guo L, Jiang M, Gong Z, Lin L, Dong J. The Dual Role of HIV-1 gp120 V3 Loop-Induced Autophagy in the Survival and Apoptosis of the Primary Rat Hippocampal Neurons. *Neurochem Res*. 44(7):1636-1652.
137. Zhu D, Li Y, Huang R, Luo L, Chen L, Fu P, He L, Li Y, Liao L, Zhu Z, Wang Y. Molecular characterization and functional activity of Prx1 in grass carp (Ctenopharyngodon idella). *FISH SHELLFISH IMMUN*. 90:395-403.
138. Zhu J, Ye Q, Xu S, Chang YX, Liu X, Ma Y, Zhu Y, Hua S. Shengmai injection alleviates H2O2-induced oxidative stress through activation of AKT and inhibition of ERK pathways in neonatal rat cardiomyocytes. *J Ethnopharmacol*. 239:111677.
139. Tan T, Hu H, Wang H, Li J, Wang Z, Wang J, Wang S, Zhang Z, Li Y. Bioinspired lipoproteins-mediated photothermia remodels tumor stroma to improve cancer cell accessibility of second nanoparticles. *Nat Commun*. 10(1):3322.
140. Zhang CY, Gao Y, Zhu RX, Qiao YN, Zhou JC, Zhang JZ, Li Y, Li SW, Fan SH, Lou HX. Prenylated Bibenzyls from the Chinese Liverwort *Radula constricta* and Their Mitochondria-Derived Parapoptotic Cytotoxic Activities. *J Nat Prod*. 82(7):1741-1751.
141. Deng JL, Zhang R, Zeng Y, Zhu YS, Wang G. Statins induce cell apoptosis through a modulation of AKT/FOXO1 pathway in prostate cancer cells. *Cancer Manag Res*. 11:7231-7242.
142. Shen Z, Chen J, Lu H. Osthole induced apoptosis in human normal liver cells by regulating cell proliferation and endoplasmic reticulum stress. *Environ Toxicol*. 34(6):768-776.
143. Zhang S, Jiang H, Huang S, Li P, Wang F. Curdlan sulfate/O-linked quaternized chitosan nanoparticles acting as potential adjuvants promote multiple arms of immune responses. *CARBOHYD POLYM*. 213:100-111.
144. Zheng Y, Tu C, Zhang J, Wang J. Inhibition of multiple myeloma-derived exosomes uptake suppresses the functional response in bone marrow stromal cell. *Int J Oncol*. 54(3):1061-1070.
145. Ruan W, Li J, Xu Y, Wang Y, Zhao F, Yang X, Jiang H, Zhang L, Saavedra JM, Shi L, Pang T. MALAT1 Up-Regulator Polydatin Protects Brain Microvascular Integrity and Ameliorates Stroke Through C/EBP β /MALAT1/CREB/PGC-1 α /PPAR γ Pathway. *Cell Mol Neurobiol*. 39(2):265-286.
146. Tian J, Rui K, Hong Y, Wang X, Xiao F, Lin X, Ma J, Guo H, Xu H, Ma K, Xu D, Liu D, Zhao Y, Lu L, Wang S. Increased GITRL Impairs the Function of Myeloid-Derived Suppressor Cells and Exacerbates Primary Sjögren Syndrome. *J Immunol*. 202(6):1693-1703.
147. Li WD, Yu S, Luo SM, Shen W, Yin S, Sun QY. Melatonin defends mouse oocyte quality from benzo[ghi]perylene-induced deterioration. *J Cell Physiol*.

234(5):6220-6229.

148. Zhou X, Li T, Chen Y, Zhang N, Wang P, Liang Y, Long M, Liu H, Mao J, Liu Q, Sun X, Chen H. Mesenchymal stem cell-derived extracellular vesicles promote the in vitro proliferation and migration of breast cancer cells through the activation of the ERK pathway. *Int J Oncol.* 54(5):1843-1852.
149. Chen C, Ai Q, Chu S, Zhang Z, Zhou X, Luo P, Liu Y, Chen N. IMM-H004 protects against oxygen-glucose deprivation/reperfusion injury to BV2 microglia partly by modulating CKLF1 involved in microglia polarization. *Int Immunopharmacol.* 70:69-79.
150. Li J, Xing F, Chen F, He L, So KF, Liu Y, Xiao J. Functional 3D Human Liver Bud Assembled from MSC-Derived Multiple Liver Cell Lineages. *Cell Transplant.* 28(5):510-521.
151. Chu P, He L, Yang C, Li Y, Huang R, Liao L, Li Y, Zhu Z, Wang Y. Characterisation and function of TRIM23 in grass carp (*Ctenopharyngodon idella*). *FISH SHELLFISH IMMUN.* 88:627-635.
152. Liu H, Zhao J, Fu R, Zhu C, Fan D. The ginsenoside Rk3 exerts anti-esophageal cancer activity in vitro and in vivo by mediating apoptosis and autophagy through regulation of the PI3K/Akt/mTOR pathway. *PLoS One.* 14(5):e0216759.
153. Ye M, Zhu X, Yan R, Hui J, Zhang J, Liu Y, Li X. Sodium demethylcantharidate induces apoptosis in hepatocellular carcinoma cells via ER stress. *Am J Transl Res.* 11(5):3150-3158. eCollection 2019.
154. Qiumge N, He Z, Qin J, Sun Y, Wang X, Yu T, Dong W, Yang G, Pang W. Overexpression of DNMT3A promotes proliferation and inhibits differentiation of porcine intramuscular preadipocytes by methylating p21 and PPAR γ promoters. *Gene.* 696:54-62.
155. Li N, Wang LJ, Xu WL, Liu S, Yu JY. MicroRNA-379-5p suppresses renal fibrosis by regulating the LIN28/let-7 axis in diabetic nephropathy. *Int J Mol Med.* 44(5):1619-1628.
156. Deng H, Zhang W, Li J, Li J, Hu L, Yan W, Liu S, He J, Weng S. A signal transducers and activators of transcription (STAT) gene from *Scylla paramamosain* is involved in resistance against mud crab reovirus. *FISH SHELLFISH IMMUN.* 94:580-591.
157. Zhu D, Fu P, Huang R, Xiong L, Wang Y, He L, Liao L, Li Y, Zhu Z, Wang Y. Molecular characterization, tissue distribution and functional analysis of galectin 1-like 2 in grass carp (*Ctenopharyngodon idella*). *FISH SHELLFISH IMMUN.* 94:455-463.
158. Ying J, Han Z, Pei S, Su L, Ruan D. Effects of Stromal Cell-Derived Factor-1 α Secreted in Degenerative Intervertebral Disc on Activation and Recruitment of Nucleus Pulposus-Derived Stem Cells. *Stem Cells Int.* 2019:9147835.
159. Jing ZF, Bi JB, Li Z, Liu X, Li J, Zhu Y, Zhang XT, Zhang Z, Li Z, Kong CZ. Inhibition of miR-34a-5p can rescue disruption of the p53-DAPK axis to suppress progression of clear cell renal cell carcinoma. *Mol Oncol.* 13(10):2079-2097.
160. Chu P, He L, Yang C, Zeng W, Huang R, Liao L, Li Y, Zhu Z, Wang Y. Grass carp ATG5 and ATG12 promote autophagy but down-regulate the transcriptional expression levels of IFN-I signaling pathway. *FISH SHELLFISH IMMUN.* 92:600-611.
161. Chen R, Hou R, Hong X, Yan S, Zha J. Organophosphate flame retardants (OPFRs) induce genotoxicity in vivo: A survey on apoptosis, DNA methylation, DNA oxidative damage, liver metabolites, and transcriptomics. *Environ Int.* 130:104914.
162. Hou DD, Zhang W, Gao YL, Sun YZ, Wang HX, Qi RQ, Chen HD, Gao XH. Anti-inflammatory effects of quercetin in a mouse model of MC903-induced atopic dermatitis. *Int Immunopharmacol.* 74:105676.
163. Zhu D, Huang R, Yang C, Fu P, Chen L, Jiang Y, He L, Li Y, Liao L, Zhu Z, Wang Y. Identification and molecular characterization of peroxiredoxin 2 in grass carp (*Ctenopharyngodon idella*). *FISH SHELLFISH IMMUN.* 92:570-582.
164. Pan JA, Tang Y, Yu JY, Zhang H, Zhang JF, Wang CQ, Gu J. miR-146a attenuates apoptosis and modulates autophagy by targeting TAF9b/P53 pathway in doxorubicin-induced cardiotoxicity. *Cell Death Dis.* 10(9):668.
165. Zhong B, Shi D, Wu F, Wang S, Hu H, Cheng C, Qing X, Huang X, Luo X, Zhang Z, Shao Z. Dynasore suppresses cell proliferation, migration, and invasion and enhances the antitumor capacity of cisplatin via STAT3 pathway in osteosarcoma. *Cell Death Dis.* 10(10):687.
166. Li Y, Ruan X, Xu X, Li C, Qiang T, Zhou H, Gao J, Wang X. Shengmai Injection Suppresses Angiotensin II-Induced Cardiomyocyte Hypertrophy and Apoptosis via Activation of the AMPK Signaling Pathway Through Energy-Dependent Mechanisms. *Front Pharmacol.* 10:1095.
167. Zhan W, Liao X, Li L, Chen Z, Tian T, Yu L, Chen Z. In vitro mitochondrial-targeted antioxidant peptide induces apoptosis in cancer cells. *ONCOTARGETS THER.* 12:7297-7306.
168. Yang L, Zhu Y, Kong D, Gong J, Yu W, Liang Y, Nie Y, Teng CB. EGF suppresses the expression of miR-124a in pancreatic β cell lines via ETS2 activation through the MEK and PI3K signaling pathways. *Int J Biol Sci.* 15(12):2561-2575.
169. Wang XL, Ma YX, Xu RJ, Ma JJ, Zhang HC, Qi SB, Xu JH, Qin XZ, Zhang HN, Liu CM, Chen JQ, Li B, Yang HL, Saijilafu. c-Myc controls the fate of neural progenitor cells during cerebral cortex development. *J Cell Physiol.* 235(4):4011-4021.
170. Zhang Y, Huang WJ, Yang QR, Zhang HD, Zhu XJ, Zeng M, Zhou X, Wang ZY, Li WJ, Jing HS, Zhang XB, Shi YP, Hu H, Yan HX, Li ZH, Zhai B. Cryopreserved biopsy tissues of rectal cancer liver metastasis for assessment of anticancer drug response in vitro and in vivo. *Oncol Rep.* 43(2):405-414.
171. Li J, Yang X, Liu F, Song Y, Liu Y. Activated PI3K/AKT Signaling Pathway Associates with Oxidative Stress and Impaired Developmental Potential of Vitrified-Thawed Oocytes. *Reprod Sci.* 27(1):404-410.
172. Zhang B, Zhu L, Dai Y, Li H, Huang K, Luo Y, Xu W. An in vitro attempt at precision toxicology reveals the involvement of DNA methylation alteration in ochratoxin A-induced G0/G1 phase arrest. *Epigenetics.* 15(1-2):199-214.
173. Zhu D, Huang R, Chu P, Chen L, Li Y, He L, Li Y, Liao L, Zhu Z, Wang Y. Characterization and expression of galectin-3 in grass carp (*Ctenopharyngodon idella*). *Dev Comp Immunol.* 104:103567.
174. Deng H, Hu L, Li J, Yan W, Song E, Kuang M, Liu S, He J, Weng S. The NF- κ B family member dorsal plays a role in immune response against Gram-positive bacterial infection in mud crab (*Scylla paramamosain*). *Dev Comp Immunol.* 106:103581.
175. Liang L, Hui K, Hu C, Wen Y, Yang S, Zhu P, Wang L, Xia Y, Qiao Y, Sun W, Fei J, Chen T, Zhao F, Yang B, Jiang X. Autophagy inhibition potentiates the anti-angiogenic property of multikinase inhibitor anlotinib through JAK2/STAT3/VEGFA signaling in non-small cell lung cancer cells. *J EXP CLIN CANC RES.* 38(1):71.
176. Tang Q, Chen Z, Zhao L, Xu H. Circular RNA hsa_circ_0000515 acts as a miR-326 sponge to promote cervical cancer progression through up-regulation of ELK1. *AGING-US.* 11(22):9982-9999.
177. Chen R, Huang H, Liu H, Xi J, Ning J, Zeng W, Shen C, Zhang T, Yu G, Xu Q, Chen X, Wang J, Lu F. Friend or Foe? Evidence Indicates Endogenous Exosomes Can Deliver Functional gRNA and Cas9 Protein. *Small.* 15(38):e1902686.
178. Siyu Li, Xiaomeng Ren, Yuandong Wang, Jiangning Hu, Haitao Wu, Shuang Song, Chunhong Yan. Fucoxanthin alleviates palmitate-induced inflammation in RAW 264.7 cells through improving lipid metabolism and attenuating mitochondrial dysfunction. *Food Funct.* doi: 10.1039/d0fo00442a.
179. Wen-Xiang Liu, Shao-Jing Tan, Yu-Feng Wang, Lan Li, Xiao-Feng Sun, Jing Liu, Francesca Gioia Klinger, Massimo De Felici, Wei Shen, Shun-Feng Cheng. Melatonin ameliorates murine fetal oocyte meiotic dysfunction in F1 and F2 offspring caused by nicotine exposure during pregnancy. *Environ Pollut.* doi: 10.1016/j.envpol.2020.114519.
180. Yi Li, Jing Lu, Xin Cao, Hongwei Zhao, Longfei Gao, Peng Xia, Gang Pei. A Newly Synthesized Rhamnoside Derivative Alleviates Alzheimer's Amyloid- β -Induced Oxidative Stress, Mitochondrial Dysfunction, and Cell Senescence through Upregulating SIRT3. *Oxid Med Cell Longev.* doi: 10.1155/2020/7698560.
181. He Ding, Yang Song, Xiaowan Huang, Liansheng Wang, Shanzi Luo, Hao Zhang, Hao Pan, Wenwei Jiang, Jing Qian, Guangyu Yao, Longping Wen, Yunjiao Zhang. mTORC1-dependent TFEB nucleus translocation and pro-survival autophagy induced by zeolitic imidazolate framework-8. *BIOMATER SCI-UK.* doi: 10.1039/d0bm00773k.

182. Yujing Zhang, Huanfeng Bian, Yu Ma, Yuanyuan Xiao, Fang Xiao. Cr(VI)-induced overactive mitophagy contributes to mitochondrial loss and cytotoxicity in L02 hepatocytes. *Biochem J.* doi: 10.1042/BCJ20200262.
183. Wei-Dong Li, Chuan-Jie Zang, Shen Yin, Wei Shen, Qing-Yuan Sun, Minghui Zhao. Metformin protects against mouse oocyte apoptosis defects induced by arecoline. *CELL PROLIFERAT.* doi: 10.1111/cpr.12809.
184. Xiaochen Qiu, Kaisheng Dong, Jingzhi Guan, JianMiao He. Hydrogen attenuates radiation-induced intestinal damage by reducing oxidative stress and inflammatory response. *Int Immunopharmacol.* doi: 10.1016/j.intimp.2020.106517.
185. Chunming Zhao, Linlin Li, Zhiwei Li, Jiawen Xu, Qing Yang, Peng Shi, Kaining Zhang, Rui Jiang. A Novel Circular RNA hsa_circRPPH1_015 Exerts an Oncogenic Role in Breast Cancer by Impairing miRNA-326-Mediated ELK1 Inhibition. *Front Oncol.* doi: 10.3389/fonc.2020.00906.
186. Tian-Yu Zhang, Li Kong, Jia-Xing Hao, Han Wang, Zi-Hui Yan, Xiao-Feng Sun, Wei Shen. Effects of Ochratoxin A exposure on DNA damage in porcine granulosa cells in vitro. *Toxicol Lett.* doi: 10.1016/j.toxlet.2020.05.011.
187. Chao Tian, Shuai Li, Lang He, Xiaobo Han, Feng Tang, Rongqi Huang, Zuoxian Lin, Sihao Deng, Junjie Xu, Hualin Huang, Huifang Zhao, Zhiyuan Li. Transient receptor potential ankyrin 1 contributes to the lysophosphatidylcholine-induced oxidative stress and cytotoxicity in OLN-93 oligodendrocyte. *CELL STRESS CHAPERON.* doi: 10.1007/s12192-020-01131-y.
188. Jing-Cai Liu, Zi-Hui Yan, Bo Li, Hong-Chen Yan, Massimo De Felici, Wei Shen. Di (2-ethylhexyl) phthalate impairs primordial follicle assembly by increasing PDE3A expression in oocytes. *Environ Pollut.* doi: 10.1016/j.envpol.2020.116088.
189. Yuhao Wu, Junke Wang, Tianxin Zhao, Yuexin Wei, Lindong Han, Lianju Shen, Chunlan Long, Shengde Wu, Guanghui Wei. LncRNAs activate longevity regulation pathway due to aging of Leydig cells caused by DEHP exposure: A transcriptome-based study. *ECOTOX ENVIRON SAFE.* doi: 10.1016/j.ecoenv.2020.111798.
190. Wenjun Yu, Busong Wang, Liang Zhou, Guoqiang Xu. Endoplasmic Reticulum Stress-Mediated p62 Downregulation Inhibits Apoptosis via c-Jun Upregulation. *BIOMOL THER.* doi: 10.4062/biomolther.2020.089.
191. Yanyi Li, Qifan Hu, Weishuo Li, Shijing Liu, Kaiming Li, Xiaoyu Li, Junjie Du, Zexuan Yu, Cong Wang, Can Zhang. Simultaneous blockage of contextual TGF- β by cyto-pharmaceuticals to suppress breast cancer metastasis. *J Control Release.* doi: 10.1016/j.jconrel.2021.06.012.
192. Jing-Cai Liu, Chun-Hua Xing, Yi Xu, Zhen-Nan Pan, Hao-Lin Zhang, Yu Zhang, Shao-Chen Sun. DEHP exposure to lactating mice affects ovarian hormone production and antral follicle development of offspring. *J Hazard Mater.* doi: 10.1016/j.jhazmat.2021.125862.
193. Ruowen Guo, Qia Xu, Liwei Liu, Hui Liu, Yun Liu, Wenmei Wei, Yide Qin. Bioactive Hexapeptide Reduced the Resistance of Ovarian Cancer Cells to DDP by Affecting HSF1/HSP70 Signaling Pathway. *J Cancer.* doi: 10.7150/jca.62285.
194. Di Chen, Qiaoxia Zheng, Long Sun, Mingming Ji, Yan Li, Hongyu Deng, Hong Zhang. ORF3a of SARS-CoV-2 promotes lysosomal exocytosis-mediated viral egress. *Dev Cell.* doi: 10.1016/j.devcel.2021.10.006.
195. Na Kong, Xiaying Chen, Jiao Feng, Ting Duan, Shuiping Liu, Xueni Sun, Peng Chen, Ting Pan, Lili Yan, Ting Jin, Yu Xiang, Quan Gao, Chengyong Wen, Weirui Ma, Wencheng Liu, Mingming Zhang, Zuyi Yang, Wengang Wang, Ruohan Zhang, Bi Chen, Tian Xie, Xinbing Sui, Wei Tao. Baicalin induces ferroptosis in bladder cancer cells by downregulating FTH1. *Acta Pharm Sin B.* doi: 10.1016/j.apsb.2021.03.036.
196. Yibin Cheng, Fengling Lai, Xin Wang, Dantong Shang, Juan Zou, Majing Luo, Xizhong Xia, Hanhua Cheng, Rongjia Zhou. Srag Regulates Autophagy via Integrating into a Preexisting Autophagy Pathway in Testis. *Mol Biol Evol.* doi: 10.1093/molbev/msaa195.
197. Jianbing Du, Zhuo Wan, Cong Wang, Fan Lu, Mengying Wei, Desheng Wang, Qiang Hao. Designer exosomes for targeted and efficient ferroptosis induction in cancer via chemo-photodynamic therapy. *Theranostics.* doi: 10.7150/thno.59121.
198. Jinhong Qi, Li Zhou, Dongqing Li, Jingyuan Yang, He Wang, Hufang Cao, Yunlan Huang, Zhiming Zhang, Linlin Chang, Chenhao Zhu, Junlong Zhan, Yong Yuan. Oncogenic role of ALX3 in cervical cancer cells through KDM2B-mediated histone demethylation of CDC25A. *BMC Cancer.* doi: 10.1186/s12885-021-08552-7.
199. Xiaoli Zhou, Ya-Na Zhou, Ashaq Ali, Cuiqin Liang, Zhiqin Ye, Xiaomin Chen, Qing Zhang, Lihua Deng, Xinyi Sun, Qian Zhang, Jihong Luo, Wei Li, Kun Zhou, Shanshan Cao, Xiaowei Zhang, Xiao-Dong Li, Xian-En Zhang, Zongqiang Cui, Dong Men. Case Report: A Re-Positive Case of SARS-CoV-2 Associated With Glaucoma. *Front Immunol.* doi: 10.3389/fimmu.2021.701295.
200. Dinglei Ma, Lihong Zhao, Li Zhang, Yali Li, Lan Zhang, Lin Li. Icarin Promotes Survival, Proliferation, and Differentiation of Neural Stem Cells In Vitro and in a Rat Model of Alzheimer's Disease. *Stem Cells Int.* doi: 10.1155/2021/9974625.
201. Jun-Jie Wang, Yu Tian, Ming-Hao Li, Yan-Qin Feng, Li Kong, Fa-Li Zhang, Wei Shen. Single-cell transcriptome dissection of the toxic impact of Di (2-ethylhexyl) phthalate on primordial follicle assembly. *Theranostics.* doi: 10.7150/thno.55006.
202. Yu Tian, Ming-Yu Zhang, Ai-Hong Zhao, Li Kong, Jun-Jie Wang, Wei Shen, Lan Li. Single-cell transcriptomic profiling provides insights into the toxic effects of Zearalenone exposure on primordial follicle assembly. *Theranostics.* doi: 10.7150/thno.58433.
203. Fengling Lai, Yibin Cheng, Juan Zou, Haoyu Wang, Wang Zhu, Xin Wang, Hanhua Cheng, Rongjia Zhou. Identification of Histone Modifications Reveals a Role of H2b Monoubiquitination in Transcriptional Regulation of dmrt1 in Monopterus albus. *Int J Biol Sci.* doi: 10.7150/ijbs.59347.
204. Yuexin Wei, Yu Zhou, Chunlan Long, Huan Wu, Yifan Hong, Yan Fu, Junke Wang, Yuhao Wu, Lianju Shen, Guanghui Wei. Polystyrene microplastics disrupt the blood-testis barrier integrity through ROS-Mediated imbalance of mTORC1 and mTORC2. *Environ Pollut.* doi: 10.1016/j.envpol.2021.117904.
205. Wen-Heng Ji, Dan-Dan Li, Dan-Ping Wei, Ai-Qin Gu, Ying Yang, Jing-Pian Peng. Cytochrome P450 26A1 Modulates the Polarization of Uterine Macrophages During the Peri-Implantation Period. *Front Immunol.* doi: 10.3389/fimmu.2021.763067.
206. Xiuxiu Liu, Wenjuan Pu, Lingjuan He, Yan Li, Huan Zhao, Yi Li, Kuo Liu, Xizhen Huang, Wendong Weng, Qing-Dong Wang, Linghong Shen, Tao Zhong, Kun Sun, Reza Ardehali, Ben He, Bin Zhou. Cell proliferation fate mapping reveals regional cardiomyocyte cell-cycle activity in subendocardial muscle of left ventricle. *Nat Commun.* doi: 10.1038/s41467-021-25933-5.
207. Bo Chen, Gengzhen Zhu, An Yan, Jing He, Yang Liu, Lin Li, Xuerui Yang, Chen Dong, Kehkooi Kee. IGSF11 is required for pericentric heterochromatin dissociation during meiotic diplotene. *PLoS Genet.* doi: 10.1371/journal.pgen.1009778.
208. Wen-Xiang Liu, Yan-Jie Zhang, Yu-Feng Wang, Francesca Gioia Klinger, Shao-Jing Tan, Donatella Farini, Massimo De Felici, Wei Shen, Shun-Feng Cheng. Protective Mechanism of Luteinizing Hormone and Follicle-Stimulating Hormone Against Nicotine-Induced Damage of Mouse Early Folliculogenesis. *Front Cell Dev Biol.* doi: 10.3389/fcell.2021.723388.
209. Yan Xu, Zhuowei Zhou, Xinmei Kang, Lijie Pan, Chang Liu, Xiaoqi Liang, Jiajie Chu, Shuai Dong, Yanli Li, Qiuli Liu, Yuetong Sun, Shanshan Yu, Qi Zhang. Mettl3-mediated mRNA m6A modification controls postnatal liver development by modulating the transcription factor Hnf4a. *Nat Commun.* doi: 10.1038/s41467-022-32169-4.
210. Xiaopeng Hao, Huixia Fan, Jian Yang, Jinfu Tang, Junhui Zhou, Yuyang Zhao, Luqi Huang, Yong Xia. Network Pharmacology Research and Dual-omic Analyses Reveal the Molecular Mechanism of Natural Product Nodosin Inhibiting Muscle-Invasive Bladder Cancer in Vitro and in Vivo. *J Nat Prod.* doi: 10.1021/acs.jnatprod.2c00400.
211. Jile Xie, Jiaying Li, Jinjin Ma, Meimei Li, Xingran Wang, Xinya Fu, Yanxia Ma, Huilin Yang, Bin Li, Saijilafu. Magnesium Oxide/Poly(l-lactide-co- ϵ -caprolactone) Scaffolds Loaded with Neural Morphogens Promote Spinal Cord Repair through Targeting the Calcium Influx and Neuronal Differentiation of Neural Stem Cells. *Adv Healthc Mater.* doi: 10.1002/adhm.202200386.
212. Chunlei Li, Sirui Zhang, Jianhua Zhu, Weijuan Huang, Yuanyuan Luo, Hui Shi,

- Dongbo Yu, Liguo Chen, Liyan Song, Rongmin Yu. A Novel Peptide Derived from Arca inflata Induces Apoptosis in Colorectal Cancer Cells through Mitochondria and the p38 MAPK Pathway. *Mar Drugs.* doi: 10.3390/md20020110.
213. Manting Xu, Robert Kopajtich, Matthias Elstner, Hua Li, Zhimei Liu, Junling Wang, Holger Prokisch, Fang Fang. Identification of a novel m.3955G > A variant in MT-ND1 associated with Leigh syndrome. *Mitochondrion.* doi: 10.1016/j.mito.2021.10.002.
214. Qihao Sun, Yiran Zhou, Minghao Xiong, Yuying Chen, Wen-Song Tan, Haibo Cai. MEK1 activation enhances the ex vivo proliferation of haematopoietic stem/progenitor cell. *Cell Biochem Funct.* doi: 10.1002/cbf.3677.
215. Yuhao Wu, Junke Wang, Tianxin Zhao, Jiadong Chen, Lian Kang, Yuexin Wei, Lindong Han, Lianju Shen, Chunlan Long, Shengde Wu, Guanghui Wei. Di-(2-ethylhexyl) phthalate exposure leads to ferroptosis via the HIF-1 α /HO-1 signaling pathway in mouse testes. *J Hazard Mater.* doi: 10.1016/j.jhazmat.2021.127807.
216. Qiyang Yan, Zhuocheng Fei, Mei Li, Jingwen Zhou, Guocheng Du, Xin Guan. Naringenin Promotes Myotube Formation and Maturation for Cultured Meat Production. *Foods.* doi: 10.3390/foods11233755.
217. Bochuan Hu, Xiaohui Xiao, Pin Chen, Jieying Qian, Guotao Yuan, Yayi Ye, Lingli Zeng, Suqin Zhong, Xiaoli Wang, Xiaohua Qin, Yuedong Yang, Yue Pan, Yunjiao Zhang. Enhancing anti-tumor effect of ultrasensitive bimetallic RuCu nanoparticles as radiosensitizers with dual enzyme-like activities. *Biomaterials.* doi: 10.1016/j.biomaterials.2022.121811.
218. Ting-Li Tang, Yan Yang, Lin Guo, Shuang Xia, Bikui Zhang, Miao Yan. Sunitinib induced hepatotoxicity in L02 cells via ROS-MAPKs signaling pathway. *Front Pharmacol.* doi: 10.3389/fphar.2022.1002142.
219. Shuya Gao, Qingchen Yang, Yue Peng, Weixian Kong, Zekun Liu, Zhe Li, Jiawen Chen, Mengmeng Bao, Xie Li, Yubin Zhang, Xiaohong Bian, Liang Jin, Hanwen Zhang, Yuexin Zhang, Daniel Sanchis, Fangrong Yan, Junmei Ye. SIRT6 regulates obesity-induced oxidative stress via ENDOG/SOD2 signaling in the heart. *Cell Biol Toxicol.* doi: 10.1007/s10565-022-09735-z.

注：更多使用本产品的文献请参考产品网页。

Version 2024.03.12