碧云天生物技术/Beyotime Biotechnology
 订货热线: 400-1683301或800-8283301
 订货e-mail: order@beyotime.com
 技术咨询: info@beyotime.com
 网址: http://www.beyotime.com

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1. Qian YF, Yao WB, Wang H, Gao XD. The Mechanism of Kaixin San Inhibiting Apoptosis in Hydrogen Peroxide2induced PC12 Cells. Chin J Nat Med. 2007 Sep;5(5):379-84.
2. Liu Z, Zhao Y, Li J, Xu S, Liu C, Zhu Y, Liang S. The venom of the spider Macrothele raveni induces apoptosis in the myelogenous leukemia K562 cell line. Leuk Res. 2012 Aug;36(8):1063-6.
3. Xu S, Zhang F, Wang H, Liu Y, Li D, Wu Z, Liu Z. The Venom of the Centipede Scolopendra subspinipes mutilans Inhibits the Growth of Myelogenous Leukemia Cell Lines. Letters in Drug Design & Discovery. 2013 Jun;10(5):390-5.
4. Fang Q, Jiang S, Li C. Evodiamine Selectively Inhibits Multiple Myeloma Cell Growth by Triggering Activation of Intrinsic Apoptosis Pathway. ONCOTARGETS THER. 2019 Dec 24;12:11383-11391.
5. Baigang Zhang, Chenghui Huang, Qikun Lu, Hairong Liang, Jinliang Li, Dongmei Xu. Involvement of caspase in patulin-induced hepatotoxicity in vitro and in vivo. Toxicol. 2022 Jan 30;206:64-73.
6. Mingzhu Chen, Yizi Zhu, Huajun Li, Yubo Zhang, Mei Han. A Quantitative Proteomic Approach Explores the Possible Mechanisms by Which the Small Molecule Stemazole Promotes the Survival of Human Neural Stem Cells. Brain Sci. 2022 May 25;12(6):690.
7. Baigang Zhang, Dongmei Xu, Lin Shao, Hairong Liang, Jinliang Li, Chenghui Huang. Toxicity mechanism of patulin on 293 T cells and correlation analysis of Caspase family. Toxicol Res (Camb). 2022 Aug 25;11(5):758-764.

Version 2024.03.12