



碧云天生物技术/Beyotime Biotechnology
订货热线: 400-1683301或800-8283301
订货e-mail: order@beyotime.com
技术咨询: info@beyotime.com
网址: http://www.beyotime.com

Caspase 3活性检测试剂盒

产品编号	产品名称	包装
C1115	Caspase 3活性检测试剂盒	20次

产品简介:

- Caspase 3活性检测试剂盒(Caspase 3 Activity Assay Kit)是采用分光光度法检测细胞或组织裂解液中caspase 3酶活性或纯化的caspase 3酶活性的试剂盒。
- Caspase (Cysteine-requiring Aspartate Protease)是一个在细胞凋亡过程中起重要作用的蛋白酶家族。Caspase 3也称CPP32、Yama或apopain, 有时被写作caspase-3或caspase 3, 属于caspase家族的CED-3亚家族(CED-3 subfamily), 是细胞凋亡过程中的一个关键酶。Caspase 3是哺乳动物细胞中研究最多的一个caspase。Caspase 3可以剪切procaspase 2、6、7和9, 并可以直接特异性剪切许多caspase底物, 包括PARP (poly(ADP-ribose) polymerase), ICAD (Inhibitor of caspase-activated deoxyribonuclease), gelsolin和fodrin等。这些由caspase 3介导的蛋白剪切是细胞凋亡分子机制的重要组成部分。另外, caspase 3在细胞核凋亡过程中也起到了关键作用, 包括染色质固缩(chromatin condensation), DNA片段化(DNA fragmentation)等。同时caspase 3对细胞起泡(cell blebbing)也起到关键作用。
- 本Caspase 3活性检测试剂盒是基于caspase 3可以催化底物Ac-DEVD-pNA (acetyl-Asp-Glu-Val-Asp *p*-nitroanilide)产生黄色的pNA (*p*-nitroaniline), 从而可以通过测定吸光度来检测caspase 3的活性。pNA在405nm附近有强吸收。
- 试剂盒中提供了caspase 3催化产生的黄色产物pNA, 可以作为定量caspase 3酶活性的标准品。
- 本试剂盒用酶标仪检测或容量不超过100μl的分光光度检测杯检测时, 除标准曲线外可以检测20个样品。

包装清单:

产品编号	产品名称	包装
C1115-1	裂解液	8ml
C1115-2	检测缓冲液	8ml
C1115-3	Ac-DEVD-pNA (2mM)	200μl
C1115-4	pNA (10mM)	200μl
—	说明书	1份

保存条件:

-20℃保存, Ac-DEVD-pNA和pNA需避光保存。

注意事项:

- 须自备可以测定A405或A400的酶标仪或容量不超过100μl的分光光度检测杯及相应分光光度计。优先考虑测定A405, 如有困难可以测定A400。
- Ac-DEVD-pNA需尽量避免反复冻融, 请注意适当分装。
- 测定蛋白浓度需Bradford蛋白浓度测定试剂盒(P0006), 可向碧云天订购。建议样品用水稀释1倍后再用Bradford法测定蛋白浓度, 以降低DTT对蛋白浓度测定的干扰。
- 有文献报道少数类型的细胞凋亡检测不到caspase 3的激活。
- pNA (中文名为4-硝基苯胺) 对人体有毒, 操作时请特别小心, 并注意有效防护以避免直接接触人体或吸入体内。pNA (10mM) 在4℃、冰浴等较低温度情况下会凝固而粘在离心管管底、管壁或管盖内, 可以20-25℃水浴温育片刻至全部融解后使用。
- 本试剂盒的裂解液可以和碧云天生产的其它caspase活性检测试剂盒的裂解液通用, 即本试剂盒裂解液制备的蛋白样品可以用于碧云天其它caspase活性检测试剂盒的检测。
- 本产品仅限于专业人员的科学研究用, 不得用于临床诊断或治疗, 不得用于食品或药品, 不得存放于普通住宅内。
- 为了您的安全和健康, 请穿实验服并戴一次性手套操作。

使用说明:

1. 准备工作:

- 裂解液溶解后混匀并置于冰浴上备用。
- 检测缓冲液溶解后混匀并置于冰浴上备用。

2. 测定pNA标准曲线:

- 标准品稀释液的配制: 按照每0.9ml检测缓冲液加入0.1ml裂解液的比例配制适量的标准品稀释液。

- b. 把试剂盒提供的pNA (10mM)用标准品稀释液稀释为0、10、20、50、100和200μM，作为标准品。
- c. 每个浓度取100μl用酶标仪进行检测，或取适当量用容量不超过100μl的分光光度检测杯进行检测，测定A405。
- d. 每一个标准品的A405减去不含pNA的空白对照的A405计算出实际的因pNA而导致的吸光度，并制作出pNA浓度相对于A405的标准曲线。pNA标准曲线可以参考图1，在0-200μM范围内存在良好的线性关系。

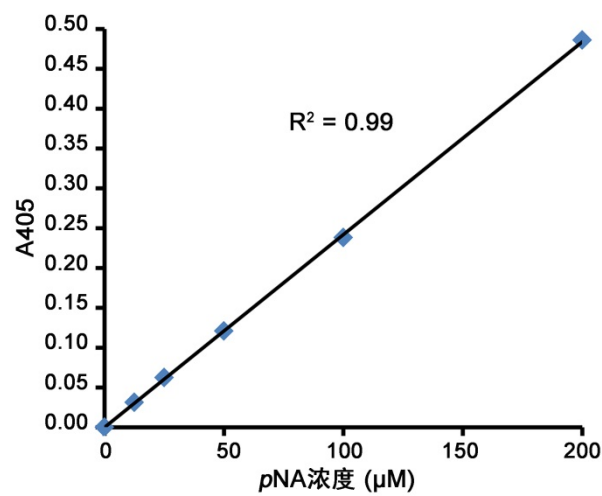


图1. pNA标准曲线。实测数据可能因实验条件、检测仪器等的不同而存在差异，图中数据仅供参考。

3. 样品的收集：

- a. **对于悬浮细胞：**把没有诱导凋亡的对照样品和诱导凋亡的样品，600g 4℃离心5分钟收集细胞，小心吸除上清，同时确保尽量没有细胞被吸除，PBS洗涤一次。同前吸尽上清后，按照每200万细胞加入100微升裂解液的比例加入裂解液（如果裂解不充分，可以把裂解液的用量提高至150或200微升），重悬沉淀，冰浴裂解15分钟。下转步骤3d。
- b. **对于贴壁细胞：**吸取细胞培养液，备用。用胰酶消化贴壁细胞，并收集至备用的细胞培养液中。600g 4℃离心5分钟收集细胞，小心吸除上清，同时确保尽量没有细胞被吸除，PBS洗涤一次。同前吸尽上清后，按照每200万细胞加入100微升裂解液的比例加入裂解液（如果裂解不充分，可以把裂解液的用量提高至150或200微升），重悬沉淀，冰浴裂解15分钟。下转步骤3d。
- c. **对于组织样品：**按照每3-10mg组织加入100微升裂解液的比例加入裂解液，在冰浴上用玻璃匀浆器匀浆。然后把匀浆液转移到1.5ml离心管中，冰浴再裂解5分钟。
- d. 4℃ 16,000-20,000g离心10-15分钟。
- e. 把上清转移到冰浴预冷的离心管中。
- f. 立即测定caspase 3的酶活性或-70℃保存样品。同时可以取少量样品用Bradford法测定蛋白浓度，尽量使蛋白浓度达到1-3mg/ml，相当于每10微升待测样品中至少含有10-30μg蛋白。如果细胞较小，可以适当增加细胞的用量。

4. Caspase 3酶活性的检测：

- a. 取出适量的Ac-DEVD-pNA (2mM)，置于冰浴上备用。
- b. 如下设置反应体系：

	空白对照	样品
检测缓冲液	40μl	40μl
待测样品	0μl	50μl
裂解液	50μl	0μl
Ac-DEVD-pNA (2mM)	10μl	10μl
总体积	100μl	100μl

注意：在设置反应体系时先加检测缓冲液，再加待测样品，适当混匀，注意避免在混匀时产生气泡。随后再加入10μl Ac-DEVD-pNA (2mM)。

- c. 加入Ac-DEVD-pNA (2mM)后混匀，注意避免在混匀时产生气泡。37℃孵育60-120分钟。发现颜色变化比较明显时即可测定A405。如果颜色变化不明显，可以适当延长孵育时间，甚至可以孵育过夜。
- d. 样品的A405扣除空白对照的A405，即为样品中caspase 3催化产生的pNA产生的吸光度。通过同步骤1中获得的标准曲线的对比就可以计算出样品中催化产生多少量的pNA。
- e. 参考Chemicon公司的caspase 3酶活力单位的定义：One unit is the amount of enzyme that will cleave 1.0nmol of the colorimetric substrate Ac-DEVD-pNA per hour at 37℃ under saturated substrate concentrations。即一个酶活力单位定义为当底物饱和时，在37℃一个小时内可以剪切1nmol Ac-DEVD-pNA产生1nmol pNA的caspase 3的酶量。这样就可以计算出样品中含有多少个酶活力单位的caspase 3。说明：在本试剂盒的检测体系中，底物的起始浓度为0.2mM，此时底物是饱和的，对于许多样品而言在37℃孵育2个小时以内底物都是饱和的；对于样品中caspase 3酶活力特别高的情况，须用裂解液适当稀释样品后再进行测定。
- f. 用Bradford法检测待测样品中的蛋白浓度(由于裂解液中含有较高浓度的DTT，不适合采用BCA法进行蛋白浓度测定)。这样就可以计算出一个样品单位重量蛋白中所含的caspase 3的酶活力单位。

常见问题：

1. 测定出的A405过低：

- a. 样品中蛋白含量太低，裂解样品时需设法使样品中的蛋白浓度至少达到1-3mg/ml。
- b. 样品中激活的caspase水平很低。首先确认凋亡现象是否明显，如果凋亡比较明显并且确认该caspase是可以被激活的，可以适当调节诱导细胞凋亡的时间，希望能找到一个caspase激活比较强的时间点，这样就可以检测出该caspase的激活。可以作一时间曲线，例如诱导凋亡0、2、4、8、16和24小时，或0、1、2、4、8和16小时，或0、1、2、4、6和8小时等。具体的诱导凋亡时间需根据具体情况而定。

2. 测定出的A405过高或者样品量不足：

测定出来的A405读数过高时，可以参考下表的反应体系适当减少样品的用量；样品量不足时也可以参考下表减少样品的用量。

	空白对照	样品
检测缓冲液	40μl	40μl
待测样品	0μl	xμl
裂解液	50μl	(50-x)μl
Ac-DEVD-pNA (2mM)	10μl	10μl
总体积	100μl	100μl

说明：其中x不超过50，其余检测方法同上面的使用说明所述。

相关产品：

产品编号	产品名称	包装
C1101	Caspase 1 活性检测试剂盒	20次
C1102	Caspase 1 活性检测试剂盒	100次
C1107	Caspase 2 活性检测试剂盒	20次
C1108	Caspase 2 活性检测试剂盒	100次
C1115	Caspase 3 活性检测试剂盒	20次
C1116	Caspase 3 活性检测试剂盒	100次
C1121	Caspase 4 活性检测试剂盒	20次
C1122	Caspase 4 活性检测试剂盒	100次
C1135	Caspase 6 活性检测试剂盒	20次
C1136	Caspase 6 活性检测试剂盒	100次
C1151	Caspase 8 活性检测试剂盒	20次
C1152	Caspase 8 活性检测试剂盒	100次
C1157	Caspase 9 活性检测试剂盒	20次
C1158	Caspase 9 活性检测试剂盒	100次
P0006	Bradford蛋白浓度测定试剂盒	1000次

使用本产品的文献：

1. Zhao CQ, Liu D, Li H, Jiang LS, Dai LY. Interleukin-1beta enhances the effect of serum deprivation on rat annular cell apoptosis. Apoptosis. 2007 Dec;12(12):2155-61.

2. Gan L, Zhu DX, Yang LP, Liu RS, Yan F, Zhang J. Involvement of transcription factor activator protein-2a in doxazosin-induced HeLa cell apoptosis1. Acta Pharmacol Sin. 2008 Apr; 29 (4): 465—472.

3. Yang Z, Mo X, Gong Q, Pan Q, Yang X, Cai W, Li C, Ma JX, He Y, Gao G. Critical effect of VEGF in the process of endothelial cell apoptosis induced by high glucose. Apoptosis. 2008 Nov;13(11):1331-43.

4. Li Y, Li Q, Wang Z, Liang D, Liang S, Tang X, Guo L, Zhang R, Zhu D. 15-HETE suppresses K(+) channel activity and inhibits apoptosis in pulmonary artery smooth muscle cells. Apoptosis. 2009 Jan;14(1):42-51.

5. Zhang SP, Zhou YJ, Liu Y, Cai YQ. Effect of liquiritigenin, a flavanone existed from Radix glycyrrhizae on pro-apoptotic in SMMC-7721 cells. Food Chem Toxicol. 2009 Apr;47(4):693-701.

6. Deng T, Zhang Y. 5-Aza-2'-deoxycytidine reactivates expression of RUNX3 by deletion of DNA methyltransferases leading to caspase independent apoptosis in colorectal cancer Lovo cells. Biomed Pharmacother. 2009 Aug; 63(7):492-500.

7. Song JQ, Teng X, Cai Y, Tang CS, Qi YF. Activation of Akt/GSK-3beta signaling pathway is involved in intermedin(1-53) protection. Apoptosis. 2009 Nov;14(11):1299-307.

8. Zhang L, Li HZ, Gong X, Luo FL, Wang B, Hu N, Wang CD, Zhang Z, Wan JY. Protective effects of Asiaticoside on acute liver injury induced by lipopolysaccharide/D-galactosamine in mice. Phytomedicine. 2010;17(10):811-9.

9. Zhu JR, Tao YF, Lou S, Wu ZM. Protective effects of ginsenoside Rb(3) on oxygen and glucose deprivation-induced ischemic injury in PC12 cells. Acta Pharmacol Sin. 2010;31(3):273-80.

10. Gong X, Luo FL, Zhang L, Li HZ, Wu MJ, Li XH, Wang B, Hu N, Wang CD, Yang JQ, Wan JY. Tetrandrine attenuates lipopolysaccharide-induced fulminant hepatic failure in D-galactosamine-sensitized mice. Int Immunopharmacol. 2010;10(3):357-63.

11. Xu AH, Hu ZM, Qu JB, Liu SM, Syed AK, Yuan HQ, Lou HX. Cyclic bisbibenzyls induce growth arrest and apoptosis of human prostate cancer PC3 cells. Acta Pharmacol Sin. 2010;31(5):609-15.

12. Wang Y, Gao J, Zhang D, Zhang J, Ma J, Jiang H. New insights into the antifibrotic effects of sorafenib on hepatic stellate cells and liver fibrosis. J Hepatol. 2010;53(1):132-44.

13. Shi X, Gu A, Ji G, Li Y, Di J, Jin J, Hu F, Long Y, Xia Y, Lu C, Song L, Wang S, Wang X. Developmental toxicity of cypermethrin in embryo-larval stages of zebrafish. Chemosphere. 2011 Oct;85(6):1010-6.

14. Huang B, Cao K, Li X, Guo S, Mao X, Wang Z, Zhuang J, Pan J, Mo C, Chen J, Qiu S. The expression and role of protein kinase C (PKC) epsilon in clear cell renal cellcarcinoma. J EXP CLIN CANC RES. 2011 Sep 28;30:88.
15. Feng Q, Cao HL, Xu W, Li XR, Ren YQ, Du LF. Apoptosis induced by genipin in human leukemia K562 cells: involvement of c-Jun N-terminal kinase in G2/M arrest. Acta Pharmacol Sin. 2011 Apr;32(4):519-27.
16. Zhang W, Liu J, Sun R, Zhao L, Du J, Ruan C, Dai K. Calpain activator dibucaine induces platelet apoptosis. Int J Mol Sci. 2011;12(4):2125-37.
17. Chen H, Mei L, Zhou L, Shen X, Guo C, Zheng Y, Zhu H, Zhu Y, Huang L. PTEN restoration and PIK3CB knockdown synergistically suppress glioblastoma growth invitroand in xenografts. NEURO-ONCOLOGY. 2011 Aug;104(1):155-67.
18. Tian LL, Yue W, Zhu F, Li S, Li W. Human mesenchymal stem cells play a dual role on tumor cell growth in vitro and in vivo. J Cell Physiol. 2011 Jul;226(7):1860-7.
19. Zhu S, Jin J, Wang Y, Ouyang Z, Xi C, Li J, Qiu Y, Wan J, Huang M, Huang Z. The endoplasmic reticulum stress response is involved in apoptosis induced by aloe-emodin inHK-2 cells. Food Chem Toxicol. 2012 Mar;50(3-4):1149-58.
20. Lv M, Zhang Y, Liang L, Wei M, Hu W, Li X, Huang Q. Effect of graphene oxide on undifferentiated and retinoic acid-differentiated SH-SY5Y cells line. Nanoscale. 2012 Jul 7;4(13):3861-6.
21. Ma J, Zhang L, Han W, Shen T, Ma C, Liu Y, Nie X, Liu M, Ran Y, Zhu D. Activation of JNK/c-Jun is required for the proliferation, survival, and angiogenesisinduced by EETin pulmonary artery endothelial cells. J Lipid Res. 2012 Jun;53(6):1093-105.
22. Tang J, Guo YS, Zhang Y, Yu XL, Li L, Huang W, Li Y, Chen B, Jiang JL, Chen ZN. CD147 induces UPR to inhibit apoptosis and chemosensitivity by increasing thetranscription ofBip in hepatocellular carcinoma. Cell Death Differ. 2012 Nov;19(11):1779-90.
23. Zhang L, Ma J, Shen T, Wang S, Ma C, Liu Y, Ran Y, Wang L, Liu L, Zhu D. Platelet-derived growth factor (PDGF) induces pulmonary vascular remodeling through 15-LO/15-HETE pathway under hypoxic condition. Cell Signal. 2012 Oct;24(10):1931-9.
24. Xu HL, Yu XF, Qu SC, Qu XR, Jiang YF, Sui da Y. Juglone, from Juglans mandshruica Maxim, inhibits growth and induces apoptosis inhumanleukemia cell HL-60 through a reactive oxygen species-dependent mechanism. Food Chem Toxicol. 2012 Mar;50(3-4):590-6.
25. Zhu X, Li Y, Luo X, Fei J. Inhibition of small GTPase RalA regulates growth and arsenic-induced apoptosis inchronicmyeloid leukemia (CML) cells. Cell Signal. 2012 Jun;24(6):1134-40.
26. Li Y, Liu L, Niu Y, Feng J, Sun Y, Kong X, Chen Y, Chen X, Gan H, Cao S, Mei Q. Modified apple polysaccharide prevents against tumorigenesis in a mouse model of colitis-associated colon cancer: role of galectin-3 and apoptosis in cancer prevention. Eur J Nutr. 2012 Feb;51(1):107-17.
27. Gao W, Ji L, Li L, Cui G, Xu K, Li P, Tang B. Bifunctional combined Au-Fe(2)O(3) nanoparticles for induction of cancer cell-specificapoptosis and real-time imaging. Biomaterials. 2012 May;33(14):3710-8.
28. Cheng J, Tang W, Su Z, Guo J, Tong L, Wei Q. Calcineurin subunit B promotes TNF-alpha-induced apoptosis by binding to mitochondria and causing mitochondrial Ca2+ overload. Cancer Lett. 2012 Aug 28;321(2):169-78.
29. Li Y, Wang H, Yang B, Yang J, Ruan X, Yang Y, Wakeland EK, Li Q, Fang X. Influence of carbon monoxide on growth and apoptosis of human umbilical artery smoothmuscle cells and veinendothelial cells. Int J Biol Sci. 2012;8(10):1431-46.
30. Guo C, Zeng X, Song J, Zhang M, Wang H, Xu X, Du F, Chen B. A Soluble Receptor for Advanced Glycation End-Products Inhibits Hypoxia/Reoxygenation-Induced Apoptosis inRat Cardiomyocytes via theMitochondrial Pathway. Int J Mol Sci. 2012;13(9):11923-40.
31. Liu J, Lu Y, Liang J. A novel fluorescence derivatization method combined with HPLC for determining theactivities of endogenouscaspase. Analyst. 2012 Nov 7;137(21):5097-104.
32. Yuan XH, Lu CL, Yao N, An LS, Yang BQ, Zhang CL, Ma X. Arsenic induced progesterone production in a caspase-3-dependent manner and changed redox status inpreovulatory granulosa cells. J Cell Physiol. 2012 Jan;227(1):194-203.
33. Zha YY, Yang B, Tang ML, Guo QC, Chen JT, Wen LP, Wang M. Concentration-dependent effects of fullereneol on cultured hippocampal neuron viability. INT J NANOMED. 2012;7:3099-109.
34. Gong K, Chen C, Zhan Y, Chen Y, Huang Z, Li W. Autophagy-related gene 7 (ATG7) and reactive oxygen species/extracellular signal-regulated kinase regulate tetrandrine-induced autophagy in human hepatocellularcarcinoma. J Biol Chem. 2012 Oct 12;287(42):35576-88.
35. Liu J, Pang Y, Chen J, Huang P, Huang W, Zhu X, Yan D. Hyperbranched polydiselenide as a self assembling broad spectrum anticancer agent. Biomaterials. 2012 Nov;33(31):7765-74.
36. hao WH, Gou BD, Zhang TL, Wang K. Lanthanum chloride bidirectionally influences calcification in bovine vascular smoothmuscle cells. J Cell Biochem. 2012 May;113(5):1776-86.
37. Wang H, Zhou X, Zhang Y, Zhu H, Zhao L, Fan L, Wang Y, Gang Y, Wu K, Liu Z, Fan D. Growth arrest-specific gene 1 is downregulated and inhibits tumor growth in gastriccancer. FEBS J. 2012 Oct;279(19):3652-64.
38. Yang G, Zhou X, Wang J, Zhang W, Zheng H, Lu W, Yuan J. MEHP-induced oxidative DNA damage and apoptosis in HepG2 cells correlates with p53-mediated mitochondria-dependent signaling pathway. Food Chem Toxicol. 2012 Jul;50(7):2424-31.
39. Li XY, Jing CQ, Lei WL, Li J, Wang JJ. Apoptosis caused by imidazolium-based ionic liquids in PC12 cells. ECOTOX ENVIRON SAFE. 2012 Sep;83:102-7.
40. Guo LD, Chen XJ, Hu YH, Yu ZJ, Wang D, Liu JZ. Curcumin inhibits proliferation and induces apoptosis of human colorectal cancer cellsby activating themitochondria apoptotic pathway. Phytother Res. 2013 Mar;27(3):422-30.
41. Li J, Du L, Yang Y, Wang C, Liu H, Wang L, Zhang X, Li W, Zheng G, Dong Z. MiR-429 is an independent prognostic factor in colorectal cancer and exerts its anti-apoptotic function bytargeting SOX2. Cancer Lett. 2013 Feb 1;329(1):84-90.
42. Wang J, Jiang Z, Ji J, Wang X, Wang T, Zhang Y, Tai T, Chen M, Sun L, Li X, Zhang L. Gene expression profiling and pathway analysis of hepatotoxicity induced by triptolidein Wistar rats. Food Chem Toxicol. 2013 Aug; 58:495-505.
43. Wu WY, Wang WY, Ma YL, Yan H, Wang XB, Qin Y, Su M, Chen T, Wang YP. Akt mediates sodium tanshinone IIA silate inhibition of oxygen-glucosedeprivation/reperfusion-induced myocardial cell apoptosis via suppression of the NF-κB/TNF-α pathway. BRIT J PHARMACOL. 2013 Mar 20.
44. Kun D, Xiang-Lin C, Ming Z, Qi L. Chlamydia inhibit host cell apoptosis by inducing Bag-1 via the MAPK/ERK survivalpathway. Apoptosis. 2013 Sep;18(9):1083-92.
45. Peng W, Chen J, Qin Y, Yang Z, Zhu YY. Long Double-Stranded Multiplex siRNAs for Dual Genes Silencing. Nucleic Acid Ther. 2013 August; 23(4):281-288.
46. Li ZL, Hu J, Li YL, Xue F, Zhang L, Xie JQ, Liu ZH, Li H, Yi DH, Liu JC, Wang SW. The effect of hyperoside on the functional recovery of the ischemic/reperfused isolatedrat heart: potential involvement of the extracellular signal-regulated kinase 1/2signaling pathway. FREE RADICAL BIO MED. 2013 Apr; 57:132-40.
47. Dan K, Yuehua G, Yanmei Z, Aodi L, Nannan D, Ying P, Yuan Y. The Biological Activity of H.Pylori SlyD in Vitro. Helicobacter. 2013 October;18(5): 347-355.
48. Jin M, Zhang L, Yu H, Meng J, Sun Z, Lu R. Protective effect of whey protein hydrolysates on H2O2-induced PC12 cells oxidativestress via a mitochondria-mediated pathway. Food Chemistry. 2013 Nov;141(2):847–52.
49. Li X, Chen N, Su Y, He Y, Yin M, Wei M, Wang L, Huang W, Fan C, Huang Q. Autophagy-sensitized cytotoxicity of quantum dots in PC12 cells. Adv Healthc Mater. 2014 Mar;3(3):354-9.
50. Sun C, Xu S, Guo J, Xu K, Wang C. The inhibitory and apoptotic effects of docetaxel-loaded mesoporous magnetic colloidalnanocrystal clusters on bladder cancer T24 cells in vitro. J BIOMED NANOTECHNOL. 2014 Mar;10(3):455-62.
51. Xu C, Hu Y, Hou L, Ju J, Li X, Du N, Guan X, Liu Z, Zhang T, Qin W, Shen N, Bilal MU, Lu Y, Zhang Y, Shan H. β-Blocker carvedilol protects cardiomyocytes against oxidative stress-induced apoptosisby up-regulating miR-133 expression. J Mol Cell Cardiol. 2014 Oct;75:111-21.
52. Qiu Q, Xiong W, Yang C, Dai X, Dan X, Yang Z, Jiao Y, Xiang Y, Liu G, Hardy P.

- Lymphocyte-derived microparticles induce apoptosis of airway epithelial cells through activation of p38 MAPK and production of arachidonic acid. *Apoptosis*. 2014 Jul;19(7):1113-27.
53. Zuo H, Lin T, Wang D, Peng R, Wang S, Gao Y, Xu X, Zhao L, Wang S, Su Z. RKIP Regulates Neural Cell Apoptosis Induced by Exposure to Microwave Radiation Partly Through the MEK/ERK/CREB Pathway. *Mol Neurobiol*. 2015;51(3):1520-9.
 54. C. Yu, Z. Zhou, J. Wang, J. Sun, W. Liu, Y. Sun, B. Kong, H. Yang and S. Yang. In depth analysis of apoptosis induced by silica coated manganese oxide nanoparticles in vitro. *Journal of Hazardous Materials*. 2015;283:519-28.
 55. D Zhu, H Wang, J Zhang, X Zhang, C Xin. Irisin improves endothelial function in type 2 diabetes through reducing oxidative/nitritative stresses. *Journal of Molecular and Cellular Cardiology*. 2015 Oct;87:138-47.
 56. J. Song, J. Li, F. Hou, X. Wang and B. Liu. Mangiferin inhibits endoplasmic reticulum stress-associated thioredoxin-interacting protein/NLRP3 inflammasome activation with regulation of AMPK in endothelial cells. *Metabolism*. 2015 Mar;64(3):428-37.
 57. J. Luo, N. Wu, B. Jiang, L. Wang, S. Wang, X. Li, B. Wang, C. Wang and D. Shi. Marine Bromophenol Derivative 3,4-Dibromo-5-(2-bromo-3,4-dihydroxy-6-isopropoxymethylbenzyl)benzene-1,2-diol Protects Hepatocytes from Lipid-Induced Cell Damage and Insulin Resistance via PTP1B Inhibition. *Mar Drugs*. 2015;13(7): 4452-69.
 58. Y. Yang, H. Meng, Q. Peng, X. Yang, R. Gan, L. Zhao, Z. Chen, J. Lu and Q. H. Meng. Downregulation of microRNA-21 expression restrains non-small cell lung cancer cell proliferation and migration through upregulation of programmed cell death 4. *Cancer Gene Ther*. 2015 Jan;22(1):23-9.
 59. J. Li, X. Li, W. Xu, S. Wang, Z. Hu, Q. Zhang, X. Deng, J. Wang, J. Zhang and C. Guo. Antifibrotic effects of luteolin on hepatic stellate cells and liver fibrosis by targeting AKT/mTOR/p70S6K and TGFβ/Smad signalling pathways. *Liver Int*. 2015 Apr;35(4): 1222-33.
 60. S. Yao, H. Tian, C. Miao, D. W. Zhang, L. Zhao, Y. Li, N. Yang, P. Jiao, H. Sang, S. Guo, Y. Wang and S. Qin. D4F alleviates macrophage-derived foam cell apoptosis by inhibiting CD36 expression and ER stress-CHOP pathway. *J Lipid Res*. 2015; 56(4): 836-47.
 61. Song YF, Liu DZ, Cheng Y, Liu M, Ye WL, Zhang BL, Liu XY, Zhou SY. Dual subcellular compartment delivery of doxorubicin to overcome drug resistant and enhance antitumor activity. *SCI REP-UK*. 2015 Nov 4;5:16125.
 62. Huang G, Mao J, Ji Z, Ailati A. Stachyose-induced apoptosis of Caco-2 cells via the caspase-dependent mitochondrial pathway. *Food Funct*. 2015 Mar;6(3):765-71.
 63. Xu Y, Zhu W, Wang Z, Yuan W, Sun Y, Liu H, Du Z. Combinatorial microRNAs suppress hypoxia-induced cardiomyocytes apoptosis. *CELL PHYSIOL BIOCHEM*. 2015;37(3):921-32.
 64. Song J, Li J, Hou F, Wang X, Liu B. Mangiferin inhibits endoplasmic reticulum stress-associated thioredoxin-interacting protein/NLRP3 inflammasome activation with regulation of AMPK in endothelial cells. *Metabolism*. 2015 Mar;64(3):428-37.
 65. Yu C, Zhou Z, Wang J, Sun J, Liu W, Sun Y, Kong B, Yang H, Yang S. In depth analysis of apoptosis induced by silica coated manganese oxide nanoparticles in vitro. *J Hazard Mater*. 2015;283:519-28.
 66. Sha WG, Shen L, Zhou L, Xu DY, Lu GY. Down-regulation of miR-186 contributes to podocytes apoptosis in membranous nephropathy. *Biomed Pharmacother*. 2015 Oct;75:179-84.
 67. Yang Y, Meng H, Peng Q, Yang X, Gan R, Zhao L, Chen Z, Lu J, Meng QH. Downregulation of microRNA-21 expression restrains non-small cell lung cancer cell proliferation and migration through upregulation of programmed cell death 4. *Cancer Gene Ther*. 2015 Jan;22(1):23-9.
 68. Li Y, Chen M, Xu Y, Yu X, Xiong T, Du M, Sun J, Liu L, Tang Y, Yao P. Iron-Mediated Lysosomal Membrane Permeabilization in Ethanol-Induced Hepatic Oxidative Damage and Apoptosis: Protective Effects of Quercetin. *Oxid Med Cell Longev*. 2016;2016:4147610.
 69. Wu H, Lin J, Liu P, Huang Z, Zhao P, Jin H, Ma J, Wen L, Gu N. Reactive oxygen species acts as executor in radiation enhancement and autophagy inducing by AgNPs. *Biomaterials*. 2016 Sep;101:1-9.
 70. Jiao S, Zhu H, He P, Teng J. Betulinic acid protects against cerebral ischemia/reperfusion injury by activating the PI3K/Akt signaling pathway. *Biomed Pharmacother*. 2016 Dec;84:1533-1537.
 71. Liu Q, Tao B, Liu G, Chen G, Zhu Q, Yu Y, Yu Y, Xiong H. Thromboxane A2 Receptor Inhibition Suppresses Multiple Myeloma Cell Proliferation by Inducing p38/c-Jun N-terminal Kinase (JNK) Mitogen-activated Protein Kinase (MAPK)-mediated G2/M Progression Delay and Cell Apoptosis. *J Biol Chem*. 2016 Feb 26;291(9):4779-92.
 72. Wang L, Shen M, Guo X, Wang B, Xia Y, Wang N, Zhang Q, Jia L, Wang X. Volume-sensitive outwardly rectifying chloride channel blockers protect against high glucose-induced apoptosis of cardiomyocytes via autophagy activation. *SCI REP-UK*. 2017 Mar 16;7:44265.
 73. Xu W, Chang J, Du X, Hou J. Long non-coding RNA PCAT-1 contributes to tumorigenesis by regulating FSCN1 via miR-145-5p in prostate cancer. *Biomed Pharmacother*. 2017 Nov;95:1112-1118.
 74. Liu T, Yu T, Hu H, He K. Knockdown of the long non-coding RNA HOTTIP inhibits colorectal cancer cell proliferation and migration and induces apoptosis by targeting SGK1. *Biomed Pharmacother*. 2017 Dec 20;98:286-296.
 75. Jiao C, Gao F, Ou L, Yu J, Li M, Wei P, Miao F. Tetrahydroxy stilbene glycoside (TSG) antagonizes Aβ-induced hippocampal neuron injury by suppressing mitochondrial dysfunction via Nrf2-dependent HO-1 pathway. *Biomed Pharmacother*. 2017 Dec;96:222-228.
 76. Han D, Li X, Li S, Su T, Fan L, Fan WS, Qiao HY, Chen JW, Fan MM, Li XJ, Wang YB, Ma S, Qiu Y, Tian ZH, Cao F. Reduced silent information regulator 1 signaling exacerbates sepsis-induced myocardial injury and mitigates the protective effect of a liver X receptor agonist. *FREE RADICAL BIO MED*. 2017 Dec;113:291-303.
 77. Liu Y, Yang Y, Wang T, Wang L, Wang X, Li T, Shi Y, Wang Y. Long non-coding RNA CCAL promotes hepatocellular carcinoma progression by regulating AP-2α and Wnt/β-catenin pathway. *Int J Biol Macromol*. 2017 Dec 22;109:424-434.
 78. Wang N, Zhang L, Lu Y, Zhang M, Zhang Z, Wang K, Lv J. Down-regulation of microRNA-142-5p attenuates oxygen-glucose deprivation and reoxygenation-induced neuron injury through up-regulating Nrf2/ARE signaling pathway. *Biomed Pharmacother*. 2017 May;89:1187-1195.
 79. Wu N, Hui H, Cui L, Yang F. GRIM-19 represses the proliferation and invasion of cutaneous squamous cell carcinoma cells associated with downregulation of STAT3 signaling. *Biomed Pharmacother*. 2017 Nov;95:1169-1176.
 80. Shen K, Feng X, Pan H, Zhang F, Xie H, Zheng S. Baicalin Ameliorates Experimental Liver Cholestasis in Mice by Modulation of Oxidative Stress, Inflammation, and NRF2 Transcription Factor. *Oxid Med Cell Longev*. 2017;2017:6169128.
 81. Du X, Chen L, Huang D, Peng Z, Zhao C, Zhang Y, Zhu Y, Wang Z, Li X, Liu G. Elevated Apoptosis in the Liver of Dairy Cows with Ketosis. *CELL PHYSIOL BIOCHEM*. 2017;43(2):568-578.
 82. Li R, Zhang Z, Wang J, Huang Y, Sun W, Xie R, Hu F, Lei T. Triptolide suppresses growth and hormone secretion in murine pituitary corticotroph tumor cells via NF-κB signaling pathway. *Biomed Pharmacother*. 2017 Nov;95:771-779.
 83. Song C, Zhao J, Fu B, Li D, Mao T, Peng W, Wu H, Zhang Y. Melatonin-mediated upregulation of Sirt3 attenuates sodium fluoride-induced hepatotoxicity by activating the MT1-PI3K/AKT-PGC-1α signaling pathway. *FREE RADICAL BIO MED*. 2017 Nov;112:616-630.
 84. Yu Y, Liu Q, Guo S, Zhang Q, Tang J, Liu G, Kong D, Li J, Yan S, Wang R, Wang P, Su X, Yu Y. 2, 3, 7, 8-Tetrachlorodibenzo-p-dioxin promotes endothelial cell apoptosis through activation of EP3/p38MAPK/Bcl-2 pathway. *J Cell Mol Med*. 2017 Dec;21(12):3540-3551.
 85. Wang Y, Xu Y, Liu Q, Zhang Y, Gao Z, Yin M, Jiang N, Cao G, Yu B, Cao Z, Kou J. Myosin IIA-related Actomyosin Contractility Mediates Oxidative Stress-induced Neuronal Apoptosis. *Front Mol Neurosci*. 2017 Mar 14;10:75.
 86. Yu J, Zhang W, Qian H, Tang H, Lin W, Lu B. SOCS1 regulates hepatic regenerative response and provides prognostic makers for acute obstructive cholangitis. *SCI REP-UK*. 2017 Aug 25;7(1):9482.

87. Sun L, Zhao M, Zhang J, Liu A, Ji W, Li Y, Yang X, Wu Z. MiR-144 promotes β -amyloid accumulation-induced cognitive impairments by targeting ADAM10 following traumatic brain injury. *ONCOTARGET*. 2017 Jul 22;8(35):59181-59203.
88. Yang Y, Tian Y, Hu S, Bi S, Li S, Hu Y, Kou J, Qi J, Yu B. Extract of Sheng-Mai-San Ameliorates Myocardial Ischemia-Induced Heart Failure by Modulating Ca²⁺-Calcineurin-Mediated Drp1 Signaling Pathways. *Int J Mol Sci*. 2017 Aug 25;18(9).
89. Zhang XT, Sun XQ, Wu C, Chen JL, Yuan JJ, Pang QF, Wang ZP. Heme oxygenase-1 induction by methylene blue protects RAW264.7 cells from hydrogen peroxide-induced injury. *Biochem Pharmacol*. 2018 Jan 5. pii: S0006-2952(18)30015-7.
90. Sun J, Shi X, Li S, Piao F. 2,5-hexanedione induces bone marrow mesenchymal stem cell apoptosis via inhibition of Akt/Bad signal pathway. *J Cell Biochem*. 2018 Apr;119(4):3732-3743.
91. Liu T, Yu T, Hu H, He K. Knockdown of the long non-coding RNA HOTTIP inhibits colorectal cancer cell proliferation and migration and induces apoptosis by targeting SGK1. *Biomed Pharmacother*. 2018 Feb;98:286-296.
92. Yin J, Ni B, Liao WG, Gao YQ. Hypoxia-induced apoptosis of mouse spermatocytes is mediated by HIF-1 α through a death receptor pathway and a mitochondrial pathway. *J Cell Physiol*. 2018 Feb;233(2):1146-1155.
93. Xiao G, Zhou J, Lu X, Huang R, Zhang H. Excessive UDPG resulting from the mutation of UAP1 causes programmed cell death by triggering reactive oxygen species accumulation and caspase-like activity in rice. *New Phytol*. 2018 Jan;217(1):332-343.
94. Dong RF, Zhang B, Tai LW, Liu HM, Shi FK, Liu NN. The Neuroprotective Role of MiR-124-3p in a 6-Hydroxydopamine-Induced Cell Model of Parkinson's Disease via the Regulation of ANAX5. *J Cell Biochem*. 2018 Jan;119(1):269-277.
95. Yang J, Pi C, Wang G. Inhibition of PI3K/Akt/mTOR pathway by apigenin induces apoptosis and autophagy in hepatocellular carcinoma cells. *Biomed Pharmacother*. 2018 Jul;103:699-707.
96. Xue R, Meng Q, Lu D, Liu X, Wang Y, Hao J. Mitofusin2 Induces Cell Autophagy of Pancreatic Cancer through Inhibiting the PI3K/Akt/mTOR Signaling Pathway. *Oxid Med Cell Longev*. 2018 Jun 26;2018:2798070.
97. Zhang J, Zhang Z, Shu B, Cui G, Zhong G. Cytotoxic and Apoptotic Activity of the Novel Harmine Derivative ZC-14 in Sf9 Cells. *Int J Mol Sci*. 2018 Mar 11;19(3). pii: E811.
98. Ding X, Sun W, Chen J. IL-2 augments the sorafenib-induced apoptosis in liver cancer by promoting mitochondrial fission and activating the JNK/TAZ pathway. *Cancer Cell Int*. 2018 Nov 9;18:176.
99. Li X, Ning L, Zhao X, Wan S. MicroRNA-543 promotes ovariectomy-induced osteoporosis through inhibition of AKT/p38 MAPK signaling pathway by targeting YAF2. *J Cell Biochem*. 2018 Dec 2.
100. Chen L, Li Q, Wang H, Chen Q, Wu Y, Shang Y. Paeoniflorin attenuated bupivacaine-induced neurotoxicity in SH-SY5Y cells via suppression of the p38 MAPK pathway. *J Cell Biochem*. 2018 Oct 28.
101. Cai W, Wu B, Li Z, He P, Wang B, Cai A, Zhang X. LncRNA NBR2 inhibits epithelial-mesenchymal transition by regulating Notch1 signaling in osteosarcoma cells. *J Cell Biochem*. 2018 Sep 6.
102. Liang B, Su J. Inducible Nitric Oxide Synthase (iNOS) Mediates Vascular Endothelial Cell Apoptosis in Grass Carp Reovirus (GCRV)-Induced Hemorrhage. *Int J Mol Sci*. 2019 Dec 16;20(24). pii: E6335.
103. Zhang L, Li S, Wang R, Chen C, Ma W, Cai H. Cytokine augments the sorafenib-induced apoptosis in Huh7 liver cancer cell by inducing mitochondrial fragmentation and activating MAPK-JNK signalling pathway. *Biomed Pharmacother*. 2019 Feb;110:213-223.
104. Wang J, Ullah SH, Li M, Zhang M, Zhang F, Zheng J, Yan X. DR region specific antibody ameliorated but ouabain worsened renal injury in nephrectomized rats through regulating Na,K-ATPase mediated signaling pathways. *AGING-US*. 2019 Feb 26;11(4):1151-1162.
105. Tong Z, Tang Y, Jiang B, Wu Y, Liu Y, Li Y, Xiao X. Phosphorylation of nucleolin is indispensable to upregulate miR-21 and inhibit apoptosis in cardiomyocytes. *J Cell Physiol*. 2019 Apr;234(4):4044-4053.
106. Li S, Chen Y, Zhang Y, Jiang X, Jiang Y, Qin X, Yang H, Wu C, Liu Y. Shear stress promotes anoikis resistance of cancer cells via caveolin-1-dependent extrinsic and intrinsic apoptotic pathways. *J Cell Physiol*. 2019 Apr;234(4):3730-3743.
107. Wu Y, Hu Y, Yuan Y, Luo Y, Lai D, Zhou H, Tong Z, Liu D. Gymnemic acid I triggers mechanistic target of rapamycin-mediated β cells cytoprotection through the promotion of autophagy under high glucose stress. *J Cell Physiol*. 2019 Jun;234(6):9370-9377.
108. Yang S, Zheng Y, Hou X. Lipoxin A4 restores oxidative stress-induced vascular endothelial cell injury and thrombosis-related factor expression by its receptor-mediated activation of Nrf2-HO-1 axis. *Cell Signal*. 2019 Aug;60:146-153.
109. Xue R, Yang J, Jia L, Zhu X, Wu J, Zhu Y, Meng Q. Mitofusin2, as a Protective Target in the Liver, Controls the Balance of Apoptosis and Autophagy in Acute-on-Chronic Liver Failure. *Front Pharmacol*. 2019 May 31;10:601.
110. Gao X, Wang Y, Li Y, Wang Y, Yan M, Sun H, Chen S, Pan X. Hupanpian, a traditional chinese medicine, inhibits liver cancer growth in vitro and in vivo by inducing autophagy and cell cycle arrest. *Biomed Pharmacother*. 2019 Dec;120:109469.
111. Ye Y, Wang W, Zhang W, Peng Y, Liu Y, Yu S, Chen Q, Geng L, Zhou L, Xie H, Lai M, Yu J, Zheng S. Galectin-1 attenuates hepatic ischemia reperfusion injury in mice. *Int Immunopharmacol*. 2019 Dec;77:105997.
112. Yu HF, Duan CC, Yang ZQ, Wang YS, Yue ZP, Guo B. HB-EGF Ameliorates Oxidative Stress-Mediated Uterine Decidualization Damage. *Oxid Med Cell Longev*. 2019 Dec 2;2019:6170936.
113. Zhang P, Ma J, Zhang Q, Jian S, Sun X, Liu B, Nie L, Liu M, Liang S, Zeng Y, Liu Z. Monosaccharide Analogues of Anticancer Peptide R-Lycosin-I: Role of Monosaccharide Conjugation in Complexation and the Potential of Lung Cancer Targeting and Therapy. *J Med Chem*. 2019 Sep 12;62(17):7857-7873.
114. Jia J, Nie L, Liu Y. Butyrate alleviates inflammatory response and NF- κ B activation in human degenerated intervertebral disc tissues. *Int Immunopharmacol*. 2020 Jan;78:106004.
115. Li R, Shi Y, Zhao S, Shi T, Zhang G. NF- κ B signaling and integrin- β 1 inhibition attenuates osteosarcoma metastasis via increased cell apoptosis. *Int J Biol Macromol*. 2019 Feb 15;123:1035-1043.
116. Wang K, Hu S, Wang B, Wang J, Wang X, Xu C. Genistein protects intervertebral discs from degeneration via Nrf2-mediated antioxidant defense system: An in vitro and in vivo study. *J Cell Physiol*. 2019 Feb 18.
117. Wang P, Wang H, Huang Q, Peng C, Yao L, Chen H, Qiu Z, Wu Y, Wang L, Chen W. Exosomes from M1-Polarized Macrophages Enhance Paclitaxel Antitumor Activity by Activating Macrophages-Mediated Inflammation. *Theranostics*. 2019 Feb 28;9(6):1714-1727.
118. Liu H, Huang H, Li R, Bi W, Feng L, E L, Hu M, Wen W. Mitophagy protects SH-SY5Y neuroblastoma cells against the TNF α -induced inflammatory injury: Involvement of microRNA-145 and Bnip3. *Biomed Pharmacother*. 2019 Jan;109:957-968.
119. Wang X, Zhang Y, Yang Y, Zhang W, Luo L, Han F, Guan H, Tao K, Hu D. Curcumin pretreatment protects against hypoxia/reoxygenation injury via improvement of mitochondrial function, destabilization of HIF-1 α and activation of Epac1-Akt pathway in rat bone marrow mesenchym. *Biomed Pharmacother*. 2019 Jan;109:1268-1275.
120. Xiang Y, Huang Y, Sun H, Pan Y, Wu M, Zhang J. Dereglulation of miR-520d-3p promotes hepatocellular carcinoma development via lncRNA MIAT regulation and EPHA2 signaling activation. *Biomed Pharmacother*. 2019 Jan;109:1630-1639.
121. Luo XY, Meng XJ, Cao DC, Wang W, Zhou K, Li L, Guo M, Wang P. Transplantation of bone marrow mesenchymal stromal cells attenuates liver fibrosis in mice by regulating macrophage subtypes. *Stem Cell Res Ther*. 2019 Jan 11;10(1):16.
122. Chen K, Cai H, Zhang H, Zhu H, Gu Z, Gong Q, Luo K. Stimuli-responsive polymer-doxorubicin conjugate: Antitumor mechanism and potential as

- p nano-prodrug.
- Acta Biomater.*
- 2019 Jan 15;84:339-355.
123. Yao Y, Zhou L, Liao W, Chen H, Du Z, Shao C, Wang P, Ding K. HH1-1, a novel Galectin-3 inhibitor, exerts anti-pancreatic cancer activity by blocking Galectin-3/EGFR/AKT/FOXO3 signaling pathway. *CARBOHYD POLYM.* 2019 Jan 15;204:111-123.
 124. Zhong W, Yang J, Li M, Li L, Li A. Long noncoding RNA NEAT1 promotes the growth of human retinoblastoma cells via regulation of miR-204/CXCR4 axis. *J Cell Physiol.* 2019 Jul;234(7):11567-11576.
 125. Zhong Y, Yu C, Qin W. LncRNA SNHG14 promotes inflammatory response induced by cerebral ischemia/reperfusion injury through regulating miR-136-5p /ROCK1. *Cancer Gene Ther.* 2019 Jul;26(7-8):234-247.
 126. Liang R, Cheng S, Dong Y, Ju H. Intracellular antioxidant activity and apoptosis inhibition capacity of PEF-treated KDHCH in HepG2 cells. *Food Res Int.* 2019 Jul;121:336-347.
 127. Zhang Q, Liu RX, Chan KW, Hu J, Zhang J, Wei L, Tan H, Yang X, Liu H. Exosomal transfer of p-STAT3 promotes acquired 5-FU resistance in colorectal cancer cells. *J EXP CLIN CANC RES.* 2019 Jul 19;38(1):320.
 128. Zheng WF, Zhang SY, Ma HF, Chang XW, Wang H. C1qTNF-related protein-6 protects against doxorubicin-induced cardiac injury. *J Cell Biochem.* 2019 Jun;120(6):10748-10755.
 129. Chang G, Chen Y, Zhang H, Zhou W. Trans sodium crocetinate alleviates ischemia/reperfusion-induced myocardial oxidative stress and apoptosis via the SIRT3/FOXO3a/SOD2 signaling pathway. *Int Immunopharmacol.* 2019 Jun;71:361-371.
 130. Ma D, Zhao Y, She J, Zhu Y, Zhao Y, Liu L, Zhang Y. NLRX1 alleviates lipopolysaccharide-induced apoptosis and inflammation in chondrocytes by suppressing the activation of NF- κ B signaling. *Int Immunopharmacol.* 2019 Jun;71:7-13.
 131. Huang H, Lai S, Luo Y, Wan Q, Wu Q, Wan L, Qi W, Liu J. Nutritional Preconditioning of Apigenin Alleviates Myocardial Ischemia/Reperfusion Injury via the Mitochondrial Pathway Mediated by Notch1/Hes1. *Oxid Med Cell Longev.* 2019 Mar 20;2019:7973098.
 132. Liu Z, Li T, Zhu F, Deng S, Li X, He Y. Regulatory roles of miR-22/Redd1-mediated mitochondrial ROS and cellular autophagy in ionizing radiation-induced BMSC injury. *Cell Death Dis.* 2019 Mar 7;10(3):227.
 133. Yu H, Zhang J, Ji Q, Yu K, Wang P, Song M, Cao Z, Zhang X, Li Y. Melatonin alleviates aluminium chloride-induced immunotoxicity by inhibiting oxidative stress and apoptosis associated with the activation of Nrf2 signaling pathway. *ECOTOX ENVIRON SAFE.* 2019 May 30;173:131-141.
 134. Yue L, Ailin W, Jinwei Z, Leng L, Jianan W, Li L, Haiming C, Ling H, Chuanjian L. PSORI-CM02 ameliorates psoriasis in vivo and in vitro by inducing autophagy via inhibition of the PI3K/Akt/mTOR pathway. *Phytomedicine.* 2019 Nov;64:153054.
 135. Deng PC, Chen WB, Cai HH, An Y, Wu XQ, Chen XM, Sun DL, Yang Y, Shi LQ, Yang Y. LncRNA SNHG14 potentiates pancreatic cancer progression via modulation of annexin A2 expression by acting as a competing endogenous RNA for miR-613. *J Cell Mol Med.* 2019 Nov;23(11):7222-7232.
 136. Du Y, Qian B, Gao L, Tan P, Chen H, Wang A, Zheng T, Pu S, Xia X, Fu W. Aloin Preconditioning Attenuates Hepatic Ischemia/Reperfusion Injury via Inhibiting TLR4/MyD88/NF- κ B Signal Pathway In Vivo and In Vitro. *Oxid Med Cell Longev.* 2019 Nov 20;2019:3765898.
 137. Li S, Zhao X, Lazarovici P, Zheng W. Artemether Activation of AMPK/GSK3 β (ser9)/Nrf2 Signaling Confers Neuroprotection towards β -Amyloid-Induced Neurotoxicity in 3xTg Alzheimer's Mouse Model. *Oxid Med Cell Longev.* 2019 Nov 21;2019:1862437.
 138. Ren H, Wang D, Zhang L, Kang X, Li Y, Zhou X, Yuan G. Catalpol induces autophagy and attenuates liver steatosis in ob/ob and high-fat diet-induced obese mice. *AGING-US.* 2019 Nov 7;11(21):9461-9477.
 139. Huang N, Guo W, Ren K, Li W, Jiang Y, Sun J, Dai W, Zhao W. LncRNA AFAP1-AS1 Suppresses miR-139-5p and Promotes Cell Proliferation and Chemotherapy Resistance of Non-small Cell Lung Cancer by Competitively Upregulating RRM2. *Front Oncol.* 2019 Oct 22;9:1103.
 140. Sun J, Yue F. Suppression of REDD1 attenuates oxygen glucose deprivation/reoxygenation-evoked ischemic injury in neuron by suppressing mTOR-mediated excessive autophagy. *J Cell Biochem.* 2019 Sep;120(9):14771-14779.
 141. Li P, Fu D, Sheng Q, Yu S, Bao X, Lv Z. TUDCA attenuates intestinal injury and inhibits endoplasmic reticulum stress-mediated intestinal cell apoptosis in necrotizing enterocolitis. *Int Immunopharmacol.* 2019 Sep;74:105665.
 142. Hao Y, Chen H, Xu W, Gao J, Yang Y, Zhang Y, Tao L. Roundup® confers cytotoxicity through DNA damage and Mitochondria-Associated apoptosis induction. *Environ Pollut.* 2019 Sep;252(Pt A):917-923.
 143. Yang Q, Qu J, Jin C, Feng Y, Xie S, Zhu J, Liu G, Xie H, Qiu H, Qi Y, Mu J, Huang J. Schistosoma japonicum Infection Promotes the Response of Tfh Cells Through Down-Regulation of Caspase-3-Mediating Apoptosis. *Front Immunol.* 2019 Sep 13;10:2154.
 144. Li X, Luo S, Zhang J, Yuan Y, Jiang W, Zhu H, Ding X, Zhan L, Wu H, Xie Y, Song R, Pan Z, Lu Y. lncRNA H19 Alleviated Myocardial I/RI via Suppressing miR-877-3p/Bcl-2-Mediated Mitochondrial Apoptosis. *MOL THER-NUCL ACIDS.* 2019 Sep 6;17:297-309.
 145. Lu QB, Du Q, Wang HP, Tang ZH, Wang YB, Sun HJ. Salusin- β mediates tubular cell apoptosis in acute kidney injury: Involvement of the PKC/ROS signaling pathway. *Redox Biol.* 2020 Feb;30:101411.
 146. Guo Y, Chen Y, Han P, Liu Y, Li W, Zhu F, Fu K, Chu M. Biocompatible chitosan-carbon nanocage hybrids for sustained drug release and highly efficient laser and microwave co-irradiation induced cancer therapy. *Acta Biomater.* 2020 Feb;103:237-246.
 147. Tang YX, Fu WW, Xi ZC, Yang JL, Zheng CW, Lu Y, Shen ZW, Xu HX. Xanthone derivatives from the leaves of Garcinia oligantha. *Eur J Med Chem.* 2019 Nov 1;181:111536.
 148. Hongchao Yao, Like Yang, Linli Tian, Yan Guo, Yushan Li. LncRNA MSC-AS1 aggravates nasopharyngeal carcinoma progression by targeting miR-524-5p/nuclear receptor subfamily 4 group A member 2 (NR4A2). *Cancer Cell Int.* 2020 Apr 28;20:138.
 149. Xue Bai, Tian-Yang Tan, Yun-Xin Li, Yue Li, Ya-Fei Chen, Ru Ma, Shu-Yan Wang, Qiang Li, Zhen-Quan Liu. The protective effect of cordyceps sinensis extract on cerebral ischemic injury via modulating the mitochondrial respiratory chain and inhibiting the mitochondrial apoptotic pathway. *Biomed Pharmacother.* 2020 Apr;124:109834.
 150. Fan Yuning, Chen Liang, Wang Tenghuan, Nan Zhenhua, Gong Shengkai. Knockdown of lincRNA PADNA promotes bupivacaine-induced neurotoxicity by miR-194/FBXW7 axis. *Mol Med.* 2020 Aug 13;26(1):79.
 151. Pingping Xia, Fan Zhang, Yajing Yuan, Cheng Chen, Yan Huang, Longyan Li, E Wang, Qulian Guo, Zhi Ye. ALDH 2 conferred neuroprotection on cerebral ischemic injury by alleviating mitochondria-related apoptosis through JNK/caspase-3 signaling pathway. *Int J Biol Sci.* 2020 Feb 17;16(8):1303-1323.
 152. Gaozhi Chen, Yuyan Bao, Qiaoyou Weng, Yingxin Zhao, Xiaoyao Lu, Lili Fu, Lingfeng Chen, Zhiguo Liu, Xiaomin Zhang, Guang Liang. Compound 15c, a Novel Dual Inhibitor of EGFR L858R/T790M and FGFR1, Efficiently Overcomes Epidermal Growth Factor Receptor-Tyrosine Kinase Inhibitor Resistance of Non-Small-Cell Lung Cancers. *Front Pharmacol.* 2020 Jan 10;10:1533.
 153. Dongdong Zhang, Yujiao Liu, Ziyi Luo, Yanling Chen, Anjie Xu, Yuxing Liang, Balu Wu, Xiqin Tong, Xiaoyan Liu, Hui Shen, Li Liu, Yongchang Wei, Haibing Zhou, Yi Liu, Fuling Zhou. The novel thioredoxin reductase inhibitor A-Z2 triggers intrinsic apoptosis and shows efficacy in the treatment of acute myeloid leukemia. *FREE RADICAL BIO MED.* 2020 Jan;146:275-286.
 154. Yujing Zhang, Huanfeng Bian, Yu Ma, Yuanyuan Xiao, Fang Xiao. Cr(VI)-induced overactive mitophagy contributes to mitochondrial loss and cytotoxicity in L02 hepatocytes. *Biochem J.* 2020 Jul 31;477(14):2607-2619.
 155. Mingyan Yao, Jing Zhang, Zhihong Li, Shuqin Guo, Xue Zhou, Wenjing Zhang. Marein protects human nucleus pulposus cells against high glucose-induced injury and extracellular matrix degradation at least partly by inhibition of ROS/NF- κ B

- pathway. *Int Immunopharmacol.* 2020 Mar;80:106126.
156. Hao Zhang, Yanan Chen, Yueping Chen, Peilu Jia, Shuli Ji, Jianxiong Xu, Yue Li, Tian Wang. Comparison of the effects of resveratrol and its derivative pterostilbene on hepatic oxidative stress and mitochondrial dysfunction in piglets challenged with diquat. *Food Funct.* 2020 May 1;11(5):4202-4215.
157. Kewei Deng, Yanling Li, Mi Xiao, Fanghui Wang, Ping Zhou, Wei Zhang, Axel Heep, Xiaoquan Li. Lycium ruthenicum Murr polysaccharide protects cortical neurons against oxygen-glucose deprivation/reperfusion in neonatal hypoxic-ischemic encephalopathy. *Int J Biol Macromol.* 2020 May 4;158:562-568.
158. Hui Ren, Zhi Li, Zhengjun Tang, Jun Li, Xiaou Lang. Long noncoding MAGI2-AS3 promotes colorectal cancer progression through regulating miR-3163/TMEM106B axis. *J Cell Physiol.* 2020 May;235(5):4824-4833.
159. Wensi Fan, Ran Zhang, Dong Han, Zhenhua Jiang, Shuang Li, Jibin Zhang, Yanhua Li, Yabin Wang, Feng Cao. Reduced Sirtuin1 signalling exacerbates diabetic mice hindlimb ischaemia injury and inhibits the protective effect of a liver X receptor agonist. *J Cell Mol Med.* 2020 May;24(10):5476-5490.
160. Dejun Xu, Lingbin Liu, Yongju Zhao, Li Yang, Jianyong Cheng, Rongmao Hua, Zelin Zhang, Qingwang Li. Melatonin protects mouse testes from palmitic acid-induced lipotoxicity by attenuating oxidative stress and DNA damage in a SIRT1-dependent manner. *J Pineal Res.* 2020 Nov;69(4):e12690.
161. Chao Li, Sheng Zhang, Liming Li, Qing Hu, Shen Ji. Ursodeoxycholic Acid Protects Against Arsenic Induced Hepatotoxicity by the Nrf2 Signaling Pathway. *Front Pharmacol.* 2020 Oct 16;11:594496.
162. Yuan Zhang, Xiaoyan Wang, Zhiling Ma, Bing Bai, Jie Liu, Lei Yang, Gaowu Qin, Erlin Zhang. A potential strategy for in-stent restenosis: Inhibition of migration and proliferation of vascular smooth muscle cells by Cu ion. *MAT SCI ENG C-MATER.* 2020 Oct;115:111090.
163. Meijuan Liu, Pengwei Wang, Xiaomei Sui, Fang Ding, Li Liu, Zengyan Gao, Zhaozhong Cheng. Circular RNA circABCC4 regulates lung adenocarcinoma progression via miR-3186-3p/TNRC6B axis. *J Cell Biochem.* 2020 Oct;121(10):4226-4238.
164. Huipeng Li, Huayu Zhang, Guohua Wang, Zhinian Chen, Youmin Pan. LncRNA LEX2-AS1 facilitates abdominal aortic aneurysm through miR-4685-5p/LBX2 feedback loop. *Biomed Pharmacother.* 2020 Sep;129:109904.
165. Ji Yu, Anneng Yang, Naixiang Wang, Haifeng Ling, Jiajun Song, Xi Chen, Yadong Lian, Zhishan Zhang, Feng Yan, Min Gu. Highly sensitive detection of caspase-3 activity based on peptide-modified organic electrochemical transistor biosensors. *Nanoscale.* 2021 Feb 7;13(5):2868-2874.
166. Wen-Bin Zhang, Xin Lai, Xu-Feng Guo. Activation of Nrf2 by miR-152 Inhibits Doxorubicin-Induced Cardiotoxicity via Attenuation of Oxidative Stress, Inflammation, and Apoptosis. *Oxid Med Cell Longev.* 2021 Jan 26;2021:8860883.
167. Xuliang Zhang, Yucong Wang, Xu Yang, Menglin Liu, Wanyue Huang, Jian Zhang, Miao Song, Bing Shao, Yanfei Li. The nephrotoxicity of T-2 toxin in mice caused by oxidative stress-mediated apoptosis is related to Nrf2 pathway. *Food Chem Toxicol.* 2021 Mar;149:112027.
168. Xiujing Dou, Di Yan, Siqi Liu, Lujia Gao, Anshan Shan. Thymol Alleviates LPS-Induced Liver Inflammation and Apoptosis by Inhibiting NLRP3 Inflammasome Activation and the AMPK-mTOR-Autophagy Pathway. *Nutrients.* 2022 Jul 8;14(14):2809.
169. Ying Su, Shan Lu, Chenjian Hou, Kehan Ren, Meili Wang, Xiaoli Liu, Shanyu Zhao, Xiuping Liu. Mitigation of liver fibrosis via hepatic stellate cells mitochondrial apoptosis induced by metformin. *Int Immunopharmacol.* 2022 Jul;108:108683.
170. Du Guo, Fangting Bai, Xiangjun Zhan, Wenting Zhang, Tong Jin, Yutang Wang, Xiaodong Xia, Chao Shi. Citral mitigates inflammation of Caco-2 cells induced by *Cronobacter sakazakii*. *Food Funct.* 2022 Mar 21;13(6):3540-3550.
171. Guangyue Li, Zixiang Jian, He Wang, Ling Xu, Tingwei Zhang, Jinlin Song. Irisin Promotes Osteogenesis by Modulating Oxidative Stress and Mitophagy through SIRT3 Signaling under Diabetic Conditions. *Oxid Med Cell Longev.* 2022 Oct 10:2022:3319056.

注：更多使用本产品的文献请参考产品网页。

Version 2024.03.12