碧云天生物技术/Beyotime Biotechnology  
订货热线: 400-1683301或800-8283301  
订货e-mail: order@beyotime.com  
技术咨询: info@beyotime.com  
网址: http://www.beyotime.com

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1. Liu C, Yu K, Shi X, Wang J, Lam P, Wu R, Zhou, B. Induction of oxidative stress and apoptosis by PFOS and PFOA in primary cultured hepatocytes of freshwater tilapia (Oreochromis niloticus). Aquatic Toxicology. 2007; 82(2): 135-43.

2. Qian YF, Wang H, Yao WB, Gao XD. Aqueous extract of the Chinese medicine, Danggui-Shaoyao-San, inhibits apoptosis in hydrogen peroxide-induced PC12 cells by preventing cytochrome c release and inactivating of caspase cascade. Cell Biol Int. 2008 Feb; 32(2):304-11.

3. Cai L, Wang H, Li Q, Qian Y, Yao W. Salidroside inhibits H2O2-induced apoptosis in PC12 cells by preventing cytochrome c release and inactivating of caspase cascade. Acta Biochim Biophys Sin (Shanghai). 2008 Sep; 40(9):796-802.

4. Shi DH, Wu JH, Ge HM, Tan RX. Protective effect of hopeahainol A, a novel acetylcholinesterase inhibitor, on hydrogen peroxide-induced injury in PC12 cells. Environmental Toxicology and Pharmacology. 2009; Jul. 28(1) 30–36.

5. Liu H, Dong A, Gao C, Tan C, Xie Z, Zu X, Qu L, Jiang Y. New synthetic flavone derivatives induce apoptosis of hepatocarcinoma cells. Bioorg Med Chem. 2010; 18(17):6322-8.

6. Wang GF, Guo YW, Feng B, Li L, Huang CG, Jiao BH. Tanghinigenin from seeds of Cerbera manghas L. induces apoptosis in human promyelocytic leukemia HL-60 cells. Environmental Toxicology and Pharmacology. 2010 Jul;30(1):31-6.

7. Zhu JR, Tao YF, Lou S, Wu ZM. Protective effects of ginsenoside

Rb(3) on oxygen and glucose deprivation-induced ischemic injury in PC12 cells. Acta Pharmacol Sin. 2010; 31(3):273-80.

8. Shi YQ, Wang YP, Song Y, Li HW, Liu CJ, Wu ZG, Yang KD. p,p'-DDE induces testicular apoptosis in prepubertal rats via the Fas/FasL pathway. Toxicol Lett. 2010; 193(1):79-85.

9. Li W, Xie L, Chen Z, Zhu Y, Sun Y, Miao Y, Xu Z, Han X. Cantharidin, a potent and selective PP2A inhibitor, induces an oxidative stress-independent growth inhibition of pancreatic cancer cells through G2/M cell-cycle arrest and apoptosis. Cancer Sci. 2010 May; 101(5):1226-33.

10. Zhao Q, Guo YW, Feng B, Li L, Huang CG, Jiao BH. Neriifolin from seeds of Cerbera manghas L. induces cell cycle arrest and apoptosis in human hepatocellular carcinoma HepG2 cells. Fitoterapia 2011 Jul;82(5):735–41.

11. Zhang Y, Li RJ, Ying X, Tian W, Yao HJ, Men Y, Yu Y, Zhang L, Ju RJ, Wang XX, Zhou J, Chen JX, Li N, Lu WL. Targeting therapy with mitosomal daunorubicin plus amlodipine has the potential to circumvent intrinsic resistant breast cancer. Mol Pharm. 2011 Feb 7; 8(1):162-75.

12. Zhang YH, Zhao CQ, Jiang LS, Dai LY. Cyclic stretch-induced apoptosis in rat annulus fibrosus cells is mediated in part by endoplasmic reticulum stress through nitric oxide production. Eur Spine J. 2011 Aug; 20(8):1233-43.

13. Jiang CP, Ding H, Shi DH, Wang YR, Li EG, Wu JH. Pro-apoptotic effects of tectorigenin on human hepatocellular carcinoma HepG2

- cells. *World J Gastroenterol*. 2012 Apr 21;18(15):1753-64.
14. Wang J, Wang Q, Li J, Shen Q, Wang F, Wang L. Cadmium induces hydrogen peroxide production and initiates hydrogen peroxide-dependent apoptosis in the gill of freshwater crab, *Sinopotamon henanense*. *Comp Biochem Physiol C Toxicol Pharmacol*. 2012 Nov;156(3-4):195-201.
15. Chen T, Zhang L, Yue JQ, Lv ZQ, Xia W, Wan YJ, Li YY, Xu SQ. Prenatal PFOS exposure induces oxidative stress and apoptosis in the lung of rat offspring. *Reprod Toxicol*. 2012 Jul; 33(4):538-45.
16. Ma EL, Zhao DM, Li YC, Cao H, Zhao QY, Li JC, Sun LX. Activation of ATM-Chk2 by 16-dehydropregnenolone induces G1 phase arrest and apoptosis in HeLa cells. *J Asian Nat Prod Res*. 2012; 14(9):817-25.
17. Zhou Y, Tao J, Yu H, Ni J, Zeng L, Teng Q, Kim KS, Zhao GP, Guo X, Yao Y. Hcp family proteins secreted via the type VI secretion system coordinately regulate *Escherichia coli* K1 interaction with human brain microvascular endothelial cells. *Infect Immun*. 2012 Mar; 80(3):1243-51.
18. Zhao D, Lin F, Wu X, Zhao Q, Zhao B, Lin P, Zhang Y, Yu X. Pseudolaric acid B induces apoptosis via proteasome-mediated Bcl-2 degradation in hormone-refractory prostate cancer DU145 cells. *Toxicol In Vitro*. 2012 Jun; 26(4):595-602.
19. Zhang L, Ji Q, Ni ZH, Sun J. Prohibitin induces apoptosis in BGC823 gastric cancer cells through the mitochondrial pathway. *Asian Pac J Cancer Prev*. 2012; 13(8):3803-7.
20. Zhang H, Shao D, Wu Y, Cai C, Hu C, Shou X, Dai B, Ye B, Wang M, Jia X. Apoptotic responses of *Carassius auratus* lymphocytes to nodularin exposure in vitro. *Fish Shellfish Immunol*. 2012 Dec; 33(6):1229-37.
21. Qin J, Shang L, Ping AS, Li J, Li XJ, Yu H, Magdalou J, Chen LB, Wang H. TNF/TNFR signal transduction pathway-mediated anti-apoptosis and anti-inflammatory effects of sodium ferulate on IL-1 $\beta$ -induced rat osteoarthritis chondrocytes in vitro. *Arthritis Res Ther*. 2012 Nov 7; 14(6):R242.
22. Liu J, Lu Y, Liang J. A novel fluorescence derivatization method combined with HPLC for determining the activities of endogenous caspase. *Analyst*. 2012 Nov 7;137(21):5097-104.
23. Chen XM, Liu J, Wang T, Shang J. Colchicine-induced apoptosis in human normal liver L-02 cells by mitochondrial mediated pathways. *Toxicol In Vitro*. 2012 Aug; 26(5):649-55.
24. Mao Z, Xia W, Wang J, Chen T, Zeng Q, Xu B, Li W, Chen X, Xu S. Perfluorooctane sulfonate induces apoptosis in lung cancer A549 cells through reactive oxygen species-mediated mitochondrion-dependent pathway. *J Appl Toxicol*. 2013 Nov;33(11):1268-76.
25. Hu M, Xu L, Yin L, Qi Y, Li H, Xu Y, Han X, Peng J, Wan X. Cytotoxicity of dioscin in human gastric carcinoma cells through death receptor and mitochondrial pathways. *J Appl Toxicol*. 2013 Aug;33(8):712-22.
26. Zhang JF, Zhang JG, Kuai XL, Zhang H, Jiang W, Ding WF, Li ZL, Zhu HJ, Mao ZB. Reactivation of the homeotic tumor suppressor gene CDX2 by 5-aza-2'-deoxycytidine-induced demethylation inhibits cell proliferation and induces caspase-independent apoptosis in gastric cancer cells. *Exp Ther Med*. 2013 Mar; 5(3):735-741.
27. Fang XY, Chen W, Fan JT, Song R, Wang L, Gu YH, Zeng GZ, Shen Y, Wu XF, Tan NH, Xu Q, Sun Y. Plant cyclopeptide RA-V kills human breast cancer cells by inducing mitochondria-mediated apoptosis through blocking PDK1-AKT interaction. *Toxicol Appl Pharmacol*. 2013 Feb 15; 267(1):95-103.
28. Shu G, Yang T, Wang C, Su H, Xiang M. Gastradin stimulates anticancer immune response and represses transplanted H22 hepatic ascitic tumor cell growth: Involvement of NF- $\kappa$ B signaling activation in CD4 $^{+}$  T cells. *Toxicol Appl Pharmacol*. 2013 Jun 15; 269(3):270-9.
29. Qiu L, Chen Y, Gao M, Zheng C, Zhao Q. Phagocytic uptake and ROS-mediated cytotoxicity in human hepatic cell line of amphiphilic polyphosphazene nanoparticles. *J Biomed Mater Res A*. 2013 Jan; 101(1):285-97.
30. Zhang J, Wu GQ, Zhang Y, Feng ZY, Zhu SM. Propofol induces apoptosis of hepatocellular carcinoma cells by upregulation of microRNA-199a expression. *Cell Biol Int*. 2013 Mar; 37(3):227-32.
31. Qiu Q, Xiong W, Yang C, Gagnon C, Hardy P. Lymphocyte-derived microparticles induce bronchial epithelial cells' pro-inflammatory cytokine production and apoptosis. *Mol Immunol*. 2013 Oct ; 55(3-4):220-30.
32. Wu WY, Wang WY, Ma YL, Yan H, Wang XB, Qin Y, Su M, Chen T, Wang YP. Sodium tanshinone IIA silicate inhibition of oxygen-glucose deprivation/reperfusion-induced myocardial cell apoptosis via suppression of the NF- $\kappa$ B/TNF- $\alpha$  pathway. *Br J Pharmacol*. 2013 Mar 20.
33. Zhang D, Wang HW, Yao CL. Molecular and acute temperature stress response characterizations of caspase-8 gene in two mussels, *Mytilus coruscus* and *Mytilus galloprovincialis*. *Comp Biochem Physiol B Biochem Mol Biol*. 2014 Nov-Dec; 177-178:10-20.
34. Liang Z, Guo YT, Yi YJ, Wang RC, Hu QL, Xiong XY. Ganoderma lucidum polysaccharides target a Fas/caspase dependent pathway to induce apoptosis in human colon cancer cells. *Asian Pac J Cancer Prev*. 2014; 15(9):3981-6.
35. Yang L, Liu X, Wu D, Zhang M, Ran G, Bi Y, Huang H. Growth inhibition and induction of apoptosis in SGC 7901 human gastric cancer cells by evodiamine. *Mol Med Rep*. 2014 Apr; 9(4):1147-52.
36. Yi S, Wen L, He J, Wang Y, Zhao F, Zhao J, Zhao Z, Cui G, Chen Y. Deguelin, a selective silencer of the NPM1 mutant, potentiates apoptosis and induces differentiation in AML cells carrying the NPM1 mutation. *Ann Hematol*. 2015 Feb; 94(2): 201-10.
37. Lu J, Li Y, Shen Z, Lu C, Lu L. TNF- $\alpha$  is involved in apoptosis triggered by grass carp reovirus infection in vitro. *Fish Shellfish Immunol*. 2016 Aug;55:559-67.
38. Tian F, Yao J, Yan M, Sun X, Wang W, Gao W, Tian Z, Guo S, Dong Z, Li B, Gao T, Shan P, Liu B, Wang H, Cheng J, Gao Q, Zhang Z, Cao W, Tian Y. 5-Aminolevulinic Acid-Mediated Sonodynamic Therapy Inhibits RIPK1/RIPK3-Dependent Necroptosis in THP-1-Derived Foam Cells. *Sci Rep*. 2016 Feb 25;6:21992.
39. Yang C, Wang H, Zhang B, Chen Y, Zhang Y, Sun X, Xiao G, Nan K, Ren H, Qin S. LCL161 increases paclitaxel-induced apoptosis by degrading cIAP1 and cIAP2 in NSCLC. *J Exp Clin Cancer Res*. 2016 Sep 30;35(1):158.
40. Wang H, Yu Y, Li J, Wu H, Sun J, Zhang Z, Geng L, Yu X, Liu Z. Cadmium stimulates mouse skin fibroblast apoptosis by affecting intracellular homeostasis. *Drug Chem Toxicol*. 2017 Jan;40(1):74-84.
41. Wang J, Li W, Huang X, Liu Y, Li Q, Zheng Z, Wang K. A polysaccharide from *Lentinus edodes* inhibits human colon cancer cell proliferation and suppresses tumor growth in athymic nude mice. *Oncotarget*. 2017 Jan 3;8(1):610-623.
42. Fang Y, Xu Z, Shi Y, Pei F, Yang W, Ma N, Kimatu BM, Liu K, Qiu W, Hu Q. Protection mechanism of Se-containing protein hydrolysates from Se-enriched rice on Pb $^{2+}$ -induced apoptosis in PC12 and RAW264.7 cells. *Food Chem*. 2017 Mar 15;219:391-398.
43. Tian Q, Wang Y, Zhang Q, Luo J, Jiang H, Zhang B, Mei M, Wu F, Wu Y, Peng J, Long T, Luo Y, Guo X. Phosphoprotein Gene Contributes to the Enhanced Apoptosis Induced by Wild-Type Rabies Virus GD-SH-01 In Vitro. *Front Microbiol*. 2017 Sep 5;8:1697.

44. Chen X,Hao K,Yu X,Huang A,Zhu B,Wang GX,Ling F. Magnolol protects Ctenopharyngodon idella kidney cells from apoptosis induced by grass carp reovirus. *Fish Shellfish Immunol.* 2018 Mar;74:426-435.
45. Yang S,Bao H,Wang H,Li Q. Anti-tumour Effect and Pharmacokinetics of an Active Ingredient Isolated from *Inonotus hispidus*. *Biol Pharm Bull.* 2019;42(1):10-17.
46. Batulu H,Du GJ,Li DZ,Sailike D,Fan YH,Geng D. Effect of poly-arginine R18 on neurocyte cell growth via autophagy in traumatic brain injury. *Exp Ther Med.* 2019 May;17(5):4109-4115.
47. Wang M,Wang Z,Zhu X,Guan S,Liu Z. NFKB1-miR-612-FAIM2 pathway regulates tumorigenesis in neurofibromatosis type 1. *IN VITRO CELL DEV-AN.* 2019 Aug;55(7):491-500.
48. Yang X,Zhang X,Zhang J,Ji Q,Huang W,Zhang X,Li Y. Spermatogenesis disorder caused by T-2 toxin is associated with germ cell apoptosis mediated by oxidative stress. *Environ Pollut.* 2019 Aug;251:372-379.
49. Wang P,Zhang SD,Jiao J,Wang WYu L,Zhao XL,Wang LT,Meng D,Fu YJ. ROS -mediated p53 activation by juglone enhances apoptosis and autophagy in vivo and in vitro. *TOXICOL APPL PHARM.* 2019 Sep 15;379:114647.
50. Yang Y,Shao R,Jiang R,Zhu M,Tang L,Li L,Zhang L.  $\beta$  -Hydroxybutyrate exacerbates lipopolysaccharide/ d-galactosamine-induced inflammatory response and hepatocyte apoptosis in mice. *J BIOCHEM MOL TOXIC.* 2019 Sep;33(9):e22372.
51. Wu S,Yi Z,Ling M,Liu S,Sun Z,Guo X. DR4-Associated Death Receptor Signal Promotes Cartilage Damage in Patients With Kashin-Beck Disease. *Cartilage.* 2019 Nov 24;19476035.
52. Fang Q,Jiang S,Li C. Evodiamine Selectively Inhibits Multiple Myeloma Cell Growth by Triggering Activation of Intrinsic Apoptosis Pathway. *ONCOTARGETS THER.* 2019 Dec 24;12:11383-11391.
53. Dongdong Zhang,Yujiao Liu,Ziyi Luo,Yanling Chen,Anjie Xu,Yuxing Liang,Balu Wu,Xiqin Tong,Xiaoyan Liu,Hui Shen,Li Liu,Yongchang Wei,Haibing Zhou,Yi Liu,Fuling Zhou. The novel thioredoxin reductase inhibitor A-Z2 triggers intrinsic apoptosis and shows efficacy in the treatment of acute myeloid leukemia. *FREE RADICAL BIO MED.* 2020 Jan;146:275-286.
54. Xue Bai,Tian-Yang Tan,Yun-Xin Li,Yue Li,Ya-Fei Chen,Ru Ma,Shu-Yan Wang,Qiang Li,Zhen-Quan Liu. The protective effect of cordyceps sinensis extract on cerebral ischemic injury via modulating the mitochondrial respiratory chain and inhibiting the mitochondrial apoptotic pathway. *Biomed Pharmacother.* 2020 Apr;124:109834.
55. Hao Zhang,Yanan Chen,Yueping Chen,Peilu Jia,Shuli Ji,Jianxiong Xu,Yue Li,Tian Wang. Comparison of the effects of resveratrol and its derivative pterostilbene on hepatic oxidative stress and mitochondrial dysfunction in piglets challenged with diquat. *Food Funct.* 2020 May 1;11(5):4202-4215.
56. Rongli Sun,Zhaodi Man,Jiahui Ji,Shuangbin Ji,Kai Xu,Yunqiu Pu,Linling Yu,Juan Zhang,Lihong Yin,Yuepu Pu. l-Carnitine protects against 1,4-benzoquinone-induced apoptosis and DNA damage by suppressing oxidative stress and promoting fatty acid oxidation in K562 cells. *Environ Toxicol.* 2020 Oct;35(10):1033-1042.
57. Ying Liu,jiawen Wang,Juhui Qiao,Shichao Liu,Siming Wang,Daqing Zhao,Xueyuan Bai,Meichen Liu. Ginsenoside Rh2 inhibits HeLa cell energy metabolism and induces apoptosis by upregulating voltage-dependent anion channel 1. *Int J Mol Med.* 2020 Nov;46(5):1695-1706.
58. Lu Liu,Li Zhang,Longjiang Li,Mengting Chen,Zhe Wang,Yi Shen,jiayi Huang,Ling Tang. Sleep deprivation aggravated lipopolysaccharide/D-galactosamine-induced acute liver injury by suppressing melatonin production. *Inflamm Res.* 2020 Nov;69(11):1133-1142.
59. Qing Lu,Weili Chen,Yajie Ji,Yu Liu,Xiaohong Xue. Ursolic Acid Enhances Cytotoxicity of Doxorubicin-Resistant Triple-Negative Breast Cancer Cells via ZEB1-AS1/ miR-186-5p/ ABCC1 Axis. *CANCER BIOTHER RADIO.* 2021 Jan 25;doi: 10.1089/cbr.2020.4147.
60. Wei Yang,Yunlei Li,Shuai Liu,Weimin Sun,Hualan Huang,Qiqing Zhang,Jie Yan. Inhibition of ULK1 promotes the death of leukemia cell in an autophagy irrelevant manner and exerts the antileukemia effect. *Clin Transl Med.* 2021 Jan;11(1):e282.
61. Zhiping Li,Hui Bi,Hongbo Jiang,Jingjing Song,Qingfan Meng,Yizhi Zhang,Xiaofang Fei. Neuroprotective effect of emodin against Alzheimer's disease via Nrf2 signaling in U251 cells and APP/PS1 mice. *Mol Med Rep.* 2021 Feb;23(2):108.
62. Yanting Zhu,Jintuo Zhou,Peiguang Niu,Huajiao Chen,Daohua Shi. Cardamonin inhibits cell proliferation by caspase-mediated cleavage of Raptor. *N-S ARCH PHARMACOL.* 2021 Apr;394(4):809-817.
63. Pei Zhang, Nan Wu, Zhi-Jun Song, Zheng-Fu Tai. Paeoniflorin Enhances the Sensitivity of ER-Positive Breast Cancer Cells to Tamoxifen through Promoting Sirtuin 4. *Evid Based Complement Alternat Med.* 2022 Jan 15:2022:6730559.
64. Kedi Dong, Yuan Zhu, Qifeng Deng, Lanqing Sun, Sidi Yang, Kai Huang, Yu Cao, Yuanyuan Li, Shuyan Wu, Rui Huang. Salmonella pSLT-encoded effector SpvB promotes RIPK3-dependent necroptosis in intestinal epithelial cells. *Cell Death Discov.* 2022 Feb 2;8(1):44.
65. Du Guo, Fangting Bai, Xiangjun Zhan, Wenting Zhang, Tong Jin, Yutang Wang, Xiaodong Xia, Chao Shi. Citral mitigates inflammation of Caco-2 cells induced by *Cronobacter sakazakii*. *Food Funct.* 2022 Mar 21;13(6):3540-3550.
66. Hongbin Liu, Qinyun Li, Xiaodan Zhang, Yun Shi, Jinyi Li. Effect of ginkgolide K on calcium channel activity in Alzheimer's disease. *Exp Ther Med.* 2022 Jun;23(6):426.
67. Ying Su, Shan Lu, Chenjian Hou, Kehan Ren, Meili Wang, Xiaoli Liu, Shanyu Zhao, Xiuping Liu. Mitigation of liver fibrosis via hepatic stellate cells mitochondrial apoptosis induced by metformin. *Int Immunopharmacol.* 2022 Jul;108:108683.
68. Baigang Zhang, Dongmei Xu, Lin Shao, Hairong Liang, Jinliang Li, Chenghui Huang. Toxicity mechanism of patulin on 293 T cells and correlation analysis of Caspase family. *Toxicol Res (Camb).* 2022 Aug 25;11(5):758-764.
69. Yang Li, Cheng Wen, Jialin Zhong, Junqi Ling, Qianzhou Jiang. *Enterococcus faecalis* OG1RF induces apoptosis in MG63 cells via caspase-3/-8/-9 without activation of caspase-1/GSDMD. *Oral Dis.* 2022 Oct;28(7):2026-2035.
70. Huazhong Xie, Pengfei Qiang, Yao Wang, Fan Xia, Peiqing Liu, Min Li. Discovery and mechanism studies of a novel ATG4B inhibitor Ebselen by drug repurposing and its anti-colorectal cancer effects in mice. *Cell Biosci.* 2022 Dec 21;12(1):206.
71. Xin Liang, Ruyi Qian, Yiqun Ou, Dan Wang, Xianyong Lin, Chengliang Sun. Lipid peroxide-derived short-chain aldehydes promote programmed cell death in wheat roots under aluminum stress. *J Hazard Mater.* 2023 Feb 5;443(Pt A):130142.