Calcein AM (钙黄绿素AM)

产品编号	产品名称	包装
C2012-0.1ml	Calcein AM (钙黄绿素AM)	2mM × 0.1ml
C2012-0.5ml	Calcein AM (钙黄绿素AM)	2mM × 0.5ml

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1. Jie Cheng,Tao Xu,Chuanhui Xun,Hailong Guo,Rui Cao,Shutao Gao,Weibin Sheng. Carnosic acid protects against ferroptosis in PC12 cells exposed to erastin through activation of Nrf2 pathway. Life Sci. 2021 Feb 1;266:118905.

2. Shan Lei, Jing Zhang, Nicholas Thomas Blum, Meng Li, Dong-Yang Zhang, Weimin Yin, Feng Zhao, Jing Lin, Peng Huang. In vivo three-dimensional multispectral photoacoustic imaging of dual enzyme-driven cyclic cascade reaction for tumor catalytic therapy. Nat Commun. 2022 Mar 11;13(1):1298.

3. Liwei Ying, Chengzhen Liang, Yujie Zhang, Jingkai Wang, Chenggui Wang, Kaishun Xia, Kesi Shi, Chao Yu, Biao Yang, Haibin Xu, Yuang Zhang, Jiawei Shu, Xianpeng Huang, Hongyuan Xing, Fangcai Li, Xiaopeng Zhou, Qixin Chen. Enhancement of nucleus pulposus repair by glycoengineered adipose-derived mesenchymal cells. Biomaterials. 2022 Apr;283:121463.

4. Huan Yao, Qian Xie, Qingman He, Lei Zeng, Jing Long, Yuanyuan Gong, Xueping Li, Xueping Li, Weiwei Liu, Zhiyi Xu, Huihui Wu, Chuan Zheng, Yongxiang Gao. Pretreatment with Panaxatriol Saponin Attenuates Mitochondrial Apoptosis and Oxidative Stress to Facilitate Treatment of Myocardial Ischemia-Reperfusion Injury via the Regulation of Keap1/Nrf2 Activit. Oxid Med Cell Longev. 2022 May 27;2022:9626703.

5. Ji-Cheng Huang, Cui-Cui Duan, Shan Jin, Chuan-Bo Sheng, Yu-Si Wang, Zhan-Peng Yue, Bin Guo. HB-EGF induces mitochondrial dysfunction via estrogen hypersecretion in granulosa cells dependent on cAMP-PKA-JNK/ERK-Ca2+-FOXO1 pathway. Int J Biol Sci. 2022 Feb 28;18(5):2047-2059.

6. Ruiyin Guo, Junfang Liu, Jinwei Chai, Yahua Gao, Mohamed A Abdel-Rahman, Xueqing Xu. Scorpion Peptide Smp24 Exhibits a Potent Antitumor Effect on Human Lung Cancer Cells by Damaging the Membrane and Cytoskeleton In Vivo and In Vitro. Toxins (Basel). 2022 Jun 28;14(7):438.

7. Hui Wang, Qi Wang, Jianyue Dong, Weiwei Jiang, Linghong Kong, Qiong Zhang, Hanping Liu. New perspective of ceria nanodots for precise tumor therapy via oxidative stress pathway. Heliyon. 2022 Aug 23;8(8):e10370.

8. Zhiying Miao, Zhiruo Miao, Xiaohua Teng, Shiwen Xu. Melatonin alleviates lead-induced fatty liver in the common carps (Cyprinus carpio) via gut-liver