碧云天生物技术/Beyotime Biotechnology
订货热线: 400-1683301或800-8283301
订货e-mail: order@beyotime.com
技术咨询: info@beyotime.com
网址: http://www.beyotime.com

Ad-mCherry-GFP-LC3B

产品编号	产品名称	包装
C3011-1ml	Ad-mCherry-GFP-LC3B	1ml
C3011-10ml	Ad-mCherry-GFP-LC3B	10ml

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Species	Cell line	Origin	Transduction efficiency
Human	293	Embryonic Kidney	95%
	MCF-10A	Mammary epithelium	51%
	MDA-MB-231	Malignant breast cancer cell	72%
	HepG2	Hepatocarcinoma	70%
	T24	Transitional cell carcinoma	37%
	TCCSUP	Transitional cell carcinoma	53%
	PC-3	Prostate cancer metastases site	67%
	DU145	Prostate cancer metastases site	86%
	LNCaP	Prostate cancer metastases site	69%
	Jurkat	T cell leukemia	2%
	HRT	Rectal tumor	5%
Murine	MT1A2	Mammary tumor transformed with SV40 T antigen	92%
	NIH 3T3	Mouse fibroblast	12%
	WEHI	Fibrosarcoma	10%
	AML12	Liver	28%
	MS-K68	Spleen	48%
	EL4	T cell	2%
Bovine	FBRT HE1	Fetal retina transformed with Had5 E1	90%
	MDBK	Kidney	70%
	FBK-34	Kidney transformed with BAd3 E1	25%
	BT-5705	Turbinate	28%
	EBL	Embryonic lung	8%
Porcine	FPRT HE1-5	Fetal retina transformed with HAd5 E1	91%
	PK-15	Kidney	76%
	PK-M12	Kidney cell transformed with SV40 T antigen	71%
	PT-K75	Turbinate	22%
	EPL	Embryonic lung	53%
	SBC-1765	Derived from buffy coat cells	73%

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BSL	Agents	Practices	Primary Barriers and Safety Equipment	Facilities (Secondary Barriers)
1	Not known to consistently cause diseases in healthy	Standard microbiological	■ No primary barriers required.	Laboratory bench and sink required

	adults	practices	■ PPE: laboratory coats and gloves; eye, face protection, as needed	
2	■ Agents associated with human disease ■ Routes of transmission include percutaneous injury, ingestion, mucous membrane exposure	BSL-1 practice plus: ■ Limited access ■ Biohazard warning signs ■ ‘Sharps’ precautions ■ Biosafety manual defining any needed waste decontamination or medical surveillance policies	Primary barriers: ■ BSCs or other physical containment devices used for all manipulations of agents that cause splashes or aerosols of infectious materials ■ PPE: Laboratory coats, gloves, face and eye protection, as needed	BSL-1 plus: ■ Autoclave available
3	Indigenous or exotic agents that may cause serious or potentially lethal disease through the inhalation route of exposure	BSL-2 practice plus: ■ Controlled access ■ Decontamination of all waste ■ Decontamination of laboratory clothing before laundering	Primary barriers: ■ BSCs or other physical containment devices used for all open manipulations of agents ■ PPE: Protective laboratory clothing, gloves, face, eye and respiratory protection, as needed	BSL-2 plus: ■ Physical separation from access corridors ■ Self-closing, double-door access ■ Exhausted air not recirculated ■ Negative airflow into laboratory ■ Entry through airlock or anteroom ■ Hand washing sink near laboratory exit
4	■ Dangerous/exotic agents which post high individual risk of aerosol-transmitted laboratory infections that are frequently fatal, for which there are no vaccines or treatments ■ Agents with a close or identical antigenic relationship to an agent requiring BSL-4 until data are available to redesignate the level ■ Related agents with unknown risk of transmission	BSL-3 practices plus: ■ Clothing change before entering ■ Shower on exit ■ All material decontaminated on exit from facility	Primary barriers: ■ All procedures conducted in Class III BSCs or Class I or II BSCs in combination with full-body, air-supplied, positive pressure suit	BSL-3 plus: ■ Separate building or isolated zone ■ Dedicated supply and exhaust, vacuum, and decontamination systems ■ Other requirements outlined in the text

BSL, biosafety level; PPE, personal protective equipment.

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产品编号	产品名称	包装
C3006-1ml	Ad-GFP-LC3B	1ml
C3006-10ml	Ad-GFP-LC3B	10ml
C3011-1ml	Ad-mCherry-GFP-LC3B	1ml
C3011-10ml	Ad-mCherry-GFP-LC3B	10ml
C3012-1ml	AdPlus-mCherry-GFP-LC3B	1ml
C3012-10ml	AdPlus-mCherry-GFP-LC3B	10ml

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1. Yu P, Zhang C, Gao CY, Ma T, Zhang H, Zhou MM, Yang YW, Yang L, Kong LY. Anti-proliferation of triple-negative breast cancer cells with physagulide P: ROS/JNK signaling pathway induces apoptosis and autophagic cell death. *ONCOTARGET*. 2017 Jul 17;8(38):64032-64049.
2. Peng Y, Qiu L, Xu D, Zhang L, Yu H, Ding Y, Deng L, Lin J. M4IDP, a zoledronic acid derivative, induces G1 arrest, apoptosis and autophagy in HCT116 coloncarcinoma cells via blocking PI3K/Akt/mTOR pathway. *Life Sci*. 2017 Sep 15;185:63-72.
3. Ma Z, Wang C, Liu C, Yan DY, Deng Y, Liu W, Yang TY, Xu ZF, Xu B. The role S-nitrosylation in manganese-induced autophagy dysregulation in SH-SY5Y cells. *Environ Toxicol*. 2017 Dec;32(12):2428-2439.
4. He X, Yuan W, Li Z, Hou Y, Liu F, Feng J. 6-Hydroxydopamine induces autophagic flux dysfunction by impairing transcription factor EB activation and lysosomal function in dopaminergic neurons and SH-SY5Y cells. *Toxicol Lett*. 2018 Feb;283:58-68.
5. He X, Yuan W, Li Z, Hou Y, Liu F, Feng J. 6-Hydroxydopamine induces autophagic flux dysfunction by impairing transcription factor EB activation and lysosomal function in dopaminergic neurons and SH-SY5Y cells. *Toxicol Lett*. 2018 Feb;283:58-68.
6. Xiao Y, Zhou Y, Lu Y, Zhou K, Cai W. PHB2 interacts with LC3 and SQSTM1 is required for bile acids-induced mitophagy in cholestatic liver. *Cell Death Dis*. 2018 Feb 7;9(2):160.
7. Wang J, Hussain T, Yue R, Liao Y, Li Q, Yao J, Song Y, Sun X, Wang N, Xu L, Sreevatsan S, Zhao D, Zhou X. MicroRNA-199a Inhibits Cellular Autophagy and Downregulates IFN- β Expression by Targeting TBK1 in Mycobacterium bovis Infected Cells. *FRONT CELL INFECT MI*. 2018 Jul 10;8:238.
8. Zhang Z, Wei Y, Liu B, Wu Y, Wang H, Xie X, Feng Z, Shao G, Xiong Q. Hsp90/Sec22b promotes unconventional secretion of mature-IL-1 β through an autophagosomal carrier in porcine alveolar macrophages during Mycoplasma hyopneumoniae infection. *Mol Immunol*. 2018 Sep;101:130-139.
9. Liu M, Zhao G, Zhang D, An W, Lai H, Li X, Cao S, Lin X. Active fraction of clove induces apoptosis via PI3K/Akt/mTOR-mediated autophagy in human colorectal cancer HCT-116 cells. *Int J Oncol*. 2018 Sep;53(3):1363-1373.
10. Yin J, Ni B, Yang YD, Tang ZW, Gao ZQ, Feng L, Liao WG, Gao YQ. Elevation of autophagy rescues spermatogenesis by inhibiting apoptosis of mouse spermatocytes. *Reproduction*. 2018 Sep 27. pii: REP-18-0243.
11. Yang Y, Gao J, Zhang Y, Xu W, Hao Y, Xu Z, Tao L. Natural pyrethrins induce autophagy of HepG2 cells through the activation of AMPK/mTOR pathway. *Environ Pollut*. 2018 Oct;241:1091-1097.
12. Wei R, Mao L, Xu P, Zheng X, Hackman RM, Mackenzie GG, Wang Y. Suppressing glucose metabolism with epigallocatechin-3-gallate (EGCG) reduces breast cancer cell growth in preclinical models. *Food Funct*. 2018 Nov 14;9(11):5682-5696.
13. Zhuo FF, Zhang C, Zhang H, Xia Y, Xue GM, Yang L, Kong LY. Chrysanthemulide A induces apoptosis through DR5 upregulation via JNK-mediated autophagosome accumulation in human osteosarcoma cells. *J Cell Physiol*. 2018 Dec 17.
14. Liu B, Ding L, Zhang L, Wang S, Wang Y, Wang B, Li L. Baicalein Induces Autophagy and Apoptosis through AMPK Pathway in Human Glioma Cells. *Am J Chin Med*. 2019;47(6):1405-1418.
15. Ma Z, Li F, Chen L, Gu T, Zhang Q, Qu Y, Xu M, Cai X, Lu L. Autophagy promotes hepatic differentiation of hepatic progenitor cells by regulating the Wnt/ β -catenin signaling pathway. *J Mol Histol*. 2019 Feb;50(1):75-90.
16. Yan D, Ma Z, Liu C, Wang C, Deng Y, Liu W, Xu B. Corynoxine B ameliorates HMGB1-dependent autophagy dysfunction during manganese exposure in SH-SY5Y human neuroblastoma cells. *Food Chem Toxicol*. 2019 Feb;124:336-348.
17. Yang X, Gao J, Guo J, Zhao Z, Zhang SL, He Y. Anti-lung cancer activity and inhibitory mechanisms of a novel Calothrixin A derivative. *Life Sci*. 2019 Feb 15;219:20-30.
18. Chu Q, Zhang S, Chen M, Han W, Jia R, Chen W, Zheng X. Cherry Anthocyanins Regulate NAFLD by Promoting Autophagy Pathway. *Oxid Med Cell Longev*. 2019 Feb 25;2019:4825949.
19. Luo Z, Xu X, Shao T, Zhang J, Xu W, Yao J, Xu J. ROS-induced autophagy regulates porcine trophectoderm cell apoptosis, proliferation, and differentiation. *AM J PHYSIOL-CELL PH*. 2019 Feb 1;316(2):C198-C209.
20. Ren Z, Xiao W, Zeng Y, Liu MH, Li GH, Tang ZH, Qu SL, Hao YM, Yuan HQ, Jiang ZS. Fibroblast growth factor-21 alleviates hypoxia/reoxygenation injury in H9c2 cardiomyocytes by promoting autophagic flux. *Int J Mol Med*. 2019 Mar;43(3):1321-1330.
21. Wang X, Xiao D, Ma C, Zhang L, Duan Q, Zheng X, Mao M, Zhu D, Li Q. The effect of honokiol on pulmonary artery endothelium cell autophagy mediated by cyclophilin A in hypoxic pulmonary arterial hypertension. *J Pharmacol Sci*. 2019 Mar;139(3):158-165.
22. Zeng H, Zhang W, Gong Y, Xie C. Radiotherapy activates autophagy to increase CD8 $^{+}$ T cell infiltration by modulating major histocompatibility complex class-I expression in non-small cell lung cancer. *J Int Med Res*. 2019 Aug;47(8):3818-3830.
23. Gao S, Li S, Li Q, Zhang F, Sun M, Wan Z, Wang S. Protective effects of salvianolic acid B against hydrogen peroxide-induced apoptosis of human umbilical vein endothelial cells and underlying mechanisms. *Int J Mol Med*. 2019 Aug;44(2):457-468.
24. Chu P, He L, Yang C, Zeng W, Huang R, Liao L, Li Y, Zhu Z, Wang Y. Grass carp ATG5 and ATG12 promote autophagy but down-regulate the transcriptional expression levels of IFN-I signaling pathway. *Fish Shellfish Immunol*. 2019 Sep;92:600-611.
25. Huang Y, Li Y, Qu Y, Zheng Y, Ouyang M, Zhang Y, Lai W, Xu Q. UVA-induced photoaging inhibits autophagic degradation by impairing lysosomal function in dermal fibroblasts. *Biochem Biophys Res Commun*. 2019 Oct 22;518(4):611-618.
26. Li R, Du J, Yao Y, Yao G, Wang X. Adiponectin inhibits high glucose-induced angiogenesis via inhibiting autophagy in RF/6A cells. *J Cell Physiol*. 2019 Nov;234(11):20566-20576.
27. Hong X, Min SN, Zhang YY, Lin YT, Wang F, Huang Y, Yu GY, Wu LL, Yang HY. TNF- α Suppresses Autophagic Flux in Acinar Cells in IgG4-Related Sialadenitis. *J Dent Res*. 2019 Nov;98(12):1386-1396.
28. He J, Shi XY, Li ZM, Pan XH, Li ZL, Chen Y, Yan SJ, Xiao L. Proton pump inhibitors can reverse the YAP mediated paclitaxel resistance in epithelial ovarian cancer. *BMC Mol Cell Biol*. 2019 Nov 12;20(1):49.
29. Zhang Y, Guo W, Chen H, Gao J, Xu Z, Tao L, Li Z, Xu W. Spinetoram confers its cytotoxic effects by inducing AMPK/mTOR-mediated autophagy and oxidative DNA damage. *Ecotoxicol Environ Saf*. 2019 Nov 15;183:109480.
30. Yan Z, Su G, Gao W, He J, Shen Y, Zeng Y, Liu X. Fluid shear stress induces cell migration and invasion via activating autophagy in HepG2 cells. *Cell Adh Migr*. 2019 Dec;13(1):152-163.
31. Zhuo Ma, Can Wang, Chang Liu, Dong-Ying Yan, Xuan Tan, Kuan Liu, Meng-Jiao Jing, Yu Deng, Wei Liu, Bin Xu. Manganese induces autophagy dysregulation: The role of S-nitrosylation in regulating autophagy related proteins in vivo and in vitro. *Sci Total Environ*. 2020 Jan 1;698:134294.
32. Chang Liu, Dong-Ying Yan, Can Wang, Zhuo Ma, Yu Deng, Wei Liu, Bin Xu. Manganese activates autophagy to alleviate endoplasmic reticulum

- stress-induced apoptosis via PERK pathway. *J Cell Mol Med.* 2020 Jan;24(1):328-341.
33. Su Li,Zhilei Bo,Ying Jiang,Xianmin Song,Chun Wang,Yin Tong. Homoharringtonine promotes BCR-ABL degradation through the p62-mediated autophagy pathway. *Oncol Rep.* 2020 Jan;43(1):113-120.
 34. Chang Liu,Dong-Ying Yan,Can Wang,Zhuo Ma,Yu Deng,Wei Liu,Bin Xu. IRE1 signaling pathway mediates protective autophagic response against manganese-induced neuronal apoptosis in vivo and in vitro. *Sci Total Environ.* 2020 Apr 10;712:136480.
 35. Na Geng,Xiaozhou Wang,Xiaohui Yu,Run Wang,Yiran Zhu,Meihua Zhang,Jianzhu Liu,Yongxia Liu. Staphylococcus aureus Avoids Autophagy Clearance of Bovine Mammary Epithelial Cells by Impairing Lysosomal Function. *Front Immunol.* 2020 May 5;11:746.
 36. Ying Wang,Xin Zhang,Wenjia Chen,Lei Gao, Jihe Li,Tao Song,Jinyu Chi,Xiaohui Zhang,Zhiyu Shi,Yanghong Dong,Xinhua Yin,Yue Liu. Cortistatin ameliorates Ang II-induced proliferation of vascular smooth muscle cells by inhibiting autophagy through SSTR3 and SSTR5. *Life Sci.* 2020 Jul 15;253:117726.
 37. Chunli Xu,Qingke Bai,Chen Wang,Qiuyu Meng,Yuming Gu,Qiwei Wang,Wenjia Xu,Ying Han,Yong Qin,Song Jia,Junfang Zhang,Jie Xu,Jiao Li,Miao Chen,Feng Wang. miR-433 Inhibits Neuronal Growth and Promotes Autophagy in Mouse Hippocampal HT-22 Cell Line. *Front Pharmacol.* 2020 Nov 23;11:536913.
 38. Yong Zhao,Minling Zhu,Wenjia Chen,Qing Chang,Li Shen,Runan Yan,Liuping Yang,Jiashu Li,Rui Xu,Yu Fu,Yue Liu. TFP1 α alleviated vascular endothelial cell injury by inhibiting autophagy and the class III PI3K/Beclin-1 pathway. *Thromb Res.* 2020 Nov;195:151-157.
 39. Lixin Zhang,Cui Ma,Xiaoying Wang,June Bai,Siyu He,Junting Zhang,Wei Xin,Yiying Li,Yuan Jiang,Jianshuang Li,Daling Zhu. MicroRNA-874-5p regulates autophagy and proliferation in pulmonary artery smooth muscle cells by targeting Sirtuin 3. *Eur J Pharmacol.* 2020 Dec 5;888:173485.
 40. Li Zhang,Shixia Zhou,Tiejun Zhou,Kaifeng Yuan,Xiaoming Li,Junling Tang. S100A8 promotes chemoresistance via augmenting autophagy in B-cell lymphoma cells. *Oncol Rep.* 2021 Jan;45(1):151-158.
 41. Li-Juan Gao,Ping Li,Tengyun Ma,Zhan-Qiong Zhong,Shi-Jun Xu. Ligustilide alleviates neurotoxicity in SH-SY5Y cells induced by A β 25-35 via regulating endoplasmic reticulum stress and autophagy. *Phytother Res.* 2021 Mar;35(3):1572-1584.
 42. Xiaomeng Guo, Qian Li, Junfeng Xiang, Meirong Liu, Aijiao Guan, Yalin Tang, Hongxia Sun. A hybrid aggregate FRET probe from the mixed assembly of cyanine dyes for highly specific monitoring of mitochondria autophagy. *Anal Chim Acta.* 2021 Jun 22;1165:338561.
 43. Heng Zhang, Junfeng Lu, Hao Liu, Lingling Guan, Shiqing Xu, Zai Wang, Yang Qiu, Honglin Liu, Liang Peng, Xiuli Men. Ajugol enhances TFEB-mediated lysosome biogenesis and lipophagy to alleviate non-alcoholic fatty liver disease. *Pharmacol Res.* 2021 Dec;174:105964.
 44. Qingqiong Luo, Xiaoyan Li, Guifang Gan, Meng Yang, Xu Chen, Fuxiang Chen. PPT1 Reduction Contributes to Erianin-Induced Growth Inhibition in Oral Squamous Carcinoma Cells. *Front Cell Dev Biol.* 2021 Dec 24;9:764263.
 45. Sunli Hu, Chenxi Zhang, Tianchen Qian, Yue Bai, Liang Chen, Jiaoxiang Chen, Chongan Huang, Chenglong Xie, Xiangyang Wang, Haiming Jin. Promoting Nrf2/Sirt3-Dependent Mitophagy Suppresses Apoptosis in Nucleus Pulposus Cells and Protects against Intervertebral Disc Degeneration. *Oxid Med Cell Longev.* 2021 Jun 9;2021:6694964.
 46. Xin He, Wei Yuan, Fei Liu, Juan Feng, Yanxia Guo. Acylated Ghrelin is Protective Against 6-OHDA-induced Neurotoxicity by Regulating Autophagic Flux. *Front Pharmacol.* 2021 Jan 27;11:586302.
 47. Ning-Fang Wang, Chun-Xue Bai. Bone marrow-derived mesenchymal stem cells modulate autophagy in RAW264.7 macrophages via the phosphoinositide 3-kinase/protein kinase B/heme oxygenase-1 signaling pathway under oxygen-glucose depriva. *Chin Med J (Engl).* 2021 Feb 17;134(6):699-707.
 48. Shuai Wang, Rong Guo, Yongxia Su, Chen Yang, Yazhou Guo, Chengjian Tan, Baoyu Zhao. Swainsonine promotes apoptosis by impairing lysosomal function and inhibiting autophagic degradation in rat primary renal tubular epithelial cells. *Chem Biol Interact.* 2021 Feb 25;336:109319.
 49. Qian Zhang, Hong-Song Li, Rong Li, Jun-Hui Du, Cong Jiao. Autophagy dysregulation mediates the damage of high glucose to retinal pigment epithelium cells. *Int J Ophthalmol.* 2021 Jun 18;14(6):805-811.
 50. Yan Huang, Chenxu Zhao, Yaoquan Liu, Yezi Kong, Panpan Tan, Siqi Liu, Fangyuan Zeng, Yang Yuan, Xinwei Li, Guowen Liu, Baoyu Zhao, Jianguo Wang. NEFA Promotes Autophagosome Formation through Modulating PERK Signaling Pathway in Bovine Hepatocytes. *Animals (Basel).* 2021 Nov 28;11(12):3400.
 51. Yuxiao Xing, Lumei Wang, Guanlong Xu, Shuhua Guo, Meihua Zhang, Guodong Cheng, Yongxia Liu, Jianzhu Liu. Platycodon grandiflorus polysaccharides inhibit Pseudorabies virus replication via downregulating virus-induced autophagy. *Res Vet Sci.* 2021 Nov;140:18-25.
 52. Rong Li, Hongsong Li, Qian Zhang. Procyanidin protects human retinal pigment epithelial cells from high glucose by inhibiting autophagy. *Environ Toxicol.* 2022 Feb;37(2):201-211.
 53. Lei Li, Ya-Juan Cui, Yu Liu, Hui-Xin Li, Yu-Dong Su, Sheng-Nan Li, Lan-Lan Wang, Yue-Wen Zhao, Shuang-Xi Wang, Feng Yan, Bo Dong. ATP6AP2 knockdown in cardiomyocyte deteriorates heart function via compromising autophagic flux and NLRP3 inflammasome activation. *Cell Death Discov.* 2022 Apr 4;8(1):161.

Version 2024.03.12