碧云天生物技术/Beyotime Biotechnology
订货热线: 400-1683301或800-8283301
订货e-mail: order@beyotime.com
技术咨询: info@beyotime.com
网址: http://www.beyotime.com

AdPlus-mCherry-GFP-LC3B

产品编号	产品名称	包装
C3012-1ml	AdPlus-mCherry-GFP-LC3B	1ml
C3012-10ml	AdPlus-mCherry-GFP-LC3B	10ml

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Table 1. Summary of Recommended Biosafety Levels for Infectious Agents

BSL	Agents	Practices	Primary Barriers and Safety Equipment	Facilities (Secondary Barriers)
1	Not known to consistently cause diseases in healthy adults	Standard microbiological practices	■ No primary barriers required. ■ PPE: laboratory coats and gloves; eye, face protection, as needed	Laboratory bench and sink required
2	■ Agents associated with human disease ■ Routes of transmission include percutaneous injury, ingestion, mucous membrane exposure	BSL-1 practice plus: ■ Limited access ■ Biohazard warning signs ■ ‘Sharps’ precautions ■ Biosafety manual defining any needed waste decontamination or medical surveillance policies	Primary barriers: ■ BSCs or other physical containment devices used for all manipulations of agents that cause splashes or aerosols of infectious materials ■ PPE: Laboratory coats, gloves, face and eye protection, as needed	BSL-1 plus: ■ Autoclave available
3	Indigenous or exotic agents that may cause serious or potentially lethal disease through the inhalation route of exposure	BSL-2 practice plus: ■ Controlled access ■ Decontamination of all waste ■ Decontamination of laboratory clothing before	Primary barriers: ■ BSCs or other physical containment devices used for all open manipulations of agents ■ PPE: Protective laboratory clothing, gloves, face, eye and respiratory protection, as needed	BSL-2 plus: ■ Physical separation from access corridors ■ Self-closing, double-door access ■ Exhausted air not recirculated ■ Negative airflow into laboratory ■ Entry through airlock or anteroom

		laundrying		■ Hand washing sink near laboratory exit
4	■ Dangerous/exotic agents which post high individual risk of aerosol-transmitted laboratory infections that are frequently fatal, for which there are no vaccines or treatments ■ Agents with a close or identical antigenic relationship to an agent requiring BSL-4 until data are available to redesignate the level ■ Related agents with unknown risk of transmission	BSL-3 practices plus: ■ Clothing change before entering ■ Shower on exit ■ All material decontaminated on exit from facility	Primary barriers: ■ All procedures conducted in Class III BSCs or Class I or II BSCs in combination with full-body, air-supplied, positive pressure suit	BSL-3 plus: ■ Separate building or isolated zone ■ Dedicated supply and exhaust, vacuum, and decontamination systems ■ Other requirements outlined in the text

BSL, biosafety level; PPE, personal protective equipment.

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产品编号	产品名称	包装
C3006-1ml	Ad-GFP-LC3B	1ml
C3006-10ml	Ad-GFP-LC3B	10ml
C3011-1ml	Ad-mCherry-GFP-LC3B	1ml
C3011-10ml	Ad-mCherry-GFP-LC3B	10ml
C3012-1ml	AdPlus-mCherry-GFP-LC3B	1ml
C3012-10ml	AdPlus-mCherry-GFP-LC3B	10ml

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1. Yiyu Cai,Xinzhu Li,Zihao Pan,Yue Zhu, Ji Tuo,Qinghai Meng,Guoliang Dai,Guangming Yang,Yang Pan. Anthocyanin ameliorates hypoxia and ischemia induced inflammation and apoptosis by increasing autophagic flux in SH-SY5Y cells. Eur J Pharmacol. 2020 Sep 15;883:173360.

2. Zhiping Yan, Danfeng Guo, Ruolin Tao, Xiao Yu, Jiacheng Zhang, Yuting He, Jiakai Zhang, Jie Li, Shuijun Zhang, Wenzhi Guo. Fluid shear stress induces cell migration via RhoA-YAP1-autophagy pathway in liver cancer stem cells. Cell Adh Migr. 2022 Dec;16(1):94-106.