碧云天生物技术/Beyotime Biotechnology
订货热线: 400-1683301或800-8283301
订货e-mail: order@beyotime.com
技术咨询: info@beyotime.com
网址: http://www.beyotime.com

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1. Jia Y, Liu X. Expression of Na⁺/K⁺-ATPase Was Affected by Salinity Change in Pacific abalone *Haliotis discus hannai*. *Front Physiol*. 2018 Sep 7;9:1244.

2. Yang L, Sun J, Liu Q, Zhu R, Yang Q, Hua J, Zheng L, Li K, Wang S, Li A. Synergetic Functional Nanocomposites Enhance Immunotherapy in Solid Tumors by Remodeling the Immunoenvironment. *Adv Sci (Weinh)*. 2019 Feb 16 6(8):1802012.

3. Ji X, Tang Z, Shuai W, Zhang Z, Li J, Chen L, Cao J, Yin W. Endogenous peptide LYENRL prevents the activation of hypertrophic scar-derived fibroblasts by inhibiting the TGF-β1/Smad pathway. *Life Sci*. 2019 Aug 15 231:116674.

4. Peng T, Deng X, Tian F, Li Z, Jiang P, Zhao X, Chen G, Chen Y, Zheng P, Li D, Wang S. The interaction of LOXL2 with GATA6 induces VEGFA expression and angiogenesis in cholangiocarcinoma. *Int J Oncol*. 2019 Sep 55(3):657-670.

5. Gu Z, Zhao L, Ge C, Liu S, Fang G, Chen SS, Yang Z, Zhou R. Facet-regulated adhesion of double-stranded DNA on palladium surfaces. *Nanoscale*. 2019 Jan 23 11(4):1827-1836.

6. Mengzhu Wang, Zhigang Wang, Bin Qiao, Jin Cao, Luya Quan, Yuanli Luo, Hanwen Qi, Xiaowen Zhong, Yubei He, Xianquan Zhang, Lan Hao. Inhibited Metastasis and Amplified Chemotherapeutic Effects by Epigene-Transfection Based on a Tumor-Targeting Nanoparticle. *INT J NANOMED*. 2020 Jun 23;15:4483-4500.

7. Hang Gao, Siyue Lou, Huanwu Hong, Qiufu Ge, Huajun Zhao. Autophagy Inhibition Potentiates the Anticancer Effects of a Bendamustine Derivative NL-101 in Acute T Lymphocytic Leukemia. *Biomed Res Int*. 2020 Feb 14;2020:1520651.

8. Liuting Mo, Danlian Liang, Wanqi He, Chan Yang, Weiying Lin. Ratiometric and amplified fluorescence nanosensor based on a DNA tetrahedron for miRNA imaging in living cells. *J Mater Chem B*. 2021 Oct 13;9(39):8341-8347.

9. Min Qing, Sheng Liang Chen, Jiao Zhou, Hong Qun Luo, Nian Bing Li. Spatially localized amplification reaction with accelerated target conversion for sensitive microRNA detection. *Talanta*. 2021 Sep 1;232:122422.

10. Yangji Xue, Qingming Kong, Haojie Ding, Chengzuo Xie, Bin Zheng, Xunhui Zhuo, Jianzu Ding, Qunbo Tong, Di Lou, Shaohong Lu, Hangjun Lv. A novel loop-mediated isothermal amplification-lateral-flow-dipstick (LAMP-LFD) device for rapid detection of *Toxoplasma gondii* in the blood of stray cats and dogs. *Parasite*. 2021;28:41.

11. Jin Liu, Hongzhou Gu. Biotechnological production of ssDNA with DNA-hydrolyzing deoxyribozymes. *STAR Protoc*. 2021 May 10;2(2):100531.

12. Haipeng Tan, Ya'nan Song, Jing Chen, Ning Zhang, Qiaozi Wang, Qiyu Li, Jinfeng Gao, Hongbo Yang, Zheng Dong, Xueyi Weng, Zhengmin Wang, Dili Sun, Wusiman Yakufu, Zhiqing Pang, Zheyong Huang, Junbo Ge. Platelet-Like Fusogenic Liposome-Mediated Targeting Delivery of miR-21 Improves Myocardial Remodeling by Reprogramming Macrophages Post Myocardial Ischemia-Reperfusion Injury. *Adv Sci (Weinh)*. 2021 Aug;8(15):e2100787.

13. Yu Guo, Huan Zhu, Xiangchen Li, Caiyun Ma, Yanan Li, Tingting Sun, Yuanyuan Wang, Chunjing Wang, Weijun Guan, Changqing Liu. RepSox effectively promotes the induced differentiation of sheep fibroblasts into adipocytes via the inhibition of the TGF-β1/Smad pathway. *Int J Mol Med*. 2021 Aug;48(2):148.

14. Yang Liu, Baozeng Sun, Jingyu Pan, Yuancai Feng, Wei Ye, Jiahao Xu, Mingfu Lan, Hao Sun, Xiyang Zhang, Yuanjie Sun, Shuya Yang, Jingqi Shi, Fanglin Zhang, Linfeng Cheng, Dongbo Jiang, Kun Yang. Construction and evaluation of DNA vaccine encoding Ebola virus glycoprotein fused with lysosome-associated membrane protein. *Antiviral Res*. 2021 Sep;193:105141.

15. Ya Fan, Jia Wang, Wen Jin, Yifei Sun, Yuemei Xu, Yipin Wang, Xiubin Liang, Dongming Su. CircNR3C2 promotes HRD1-mediated tumor-suppressive effect via sponging miR-513a-3p in triple-negative breast cancer. *Mol Cancer*. 2021 Feb 2;20(1):25.

16. Jin Qiu, Qianqian Fan, Sainan Xu, Dongmei Wang, Juntong Chen, Sainan Wang, Tianhui Hu, Xinran Ma, Yiyun Cheng, Lingyan Xu. A fluorinated peptide with high serum- and lipid-tolerance for the delivery of siRNA drugs to treat obesity and metabolic dysfunction. *Biomaterials*. 2022 Jun;285:121541.

17. Jiajia Xu, Yangyang Li, Huan Zhu, Wenyu Wu, Yumeng Liu, Yu Guo, Weijun Guan, Changqing Liu, Caiyun Ma. Therapeutic function of a novel rat induced pluripotent stem cell line in a 6-OHDA-induced rat model of Parkinson's disease. *Int J Mol Med*. 2022 Dec;50(6):140.