



碧云天生物技术/Beyotime Biotechnology
订货热线: 400-1683301或800-8283301
订货e-mail: order@beyotime.com
技术咨询: info@beyotime.com
网址: http://www.beyotime.com

SDS-PAGE凝胶配制试剂盒

产品编号	产品名称	包装
P0012A	SDS-PAGE 凝胶配制试剂盒	可制30-50块胶

产品简介:

- 碧云天生产的SDS-PAGE凝胶配制试剂盒提供了配制SDS-PAGE凝胶所需的各种试剂，用户只需自备制胶器具和蒸馏水，即可配制PAGE胶(即聚丙烯酰胺凝胶)。
- SDS-PAGE凝胶配制试剂盒不仅可用于配制SDS-PAGE凝胶，也可用于配制非变性(native)PAGE凝胶。
- 本试剂盒约可配制30-50块常规大小的PAGE胶。

包装清单:

产品编号	产品名称	包装
P0012A-1	30% Acr-Bis (29:1)	100ml
P0012A-2	1M Tris-HCl, pH8.8	100ml
P0012A-3	10% SDS	5ml
P0012A-4	凝胶聚合催化剂	0.5g
P0012A-5	TEMED Substitute	0.5ml
P0012A-6	1M Tris-HCl, pH6.8	15ml
—	说明书	1份

保存条件:

1M Tris-HCl, pH8.8、10% SDS、凝胶聚合催化剂和1M Tris-HCl, pH6.8室温保存。30% Acr-Bis (29:1)和TEMED Substitute 4°C避光保存。凝胶聚合催化剂用水配制成10%溶液后，分装成小管-20°C保存，通常半年内有效。

注意事项:

- 凝胶聚合催化剂用水配制成10%溶液后，应当-20°C保存。同时应尽量减少室温存放时间，以防失效。
- TEMED Substitute易挥发，使用后请盖紧瓶盖。另外凝胶凝聚的速度和温度及光照关系密切，可通过适当调节凝胶聚合催化剂和TEMED Substitute的用量，控制在不同的室内环境下凝胶凝聚的速度。
- TEMED Substitute易燃，有腐蚀性，操作时请小心，并注意有效防护以避免直接接触人体或腐蚀其他物品。
- 本产品仅限于专业人员的科学的研究用，不得用于临床诊断或治疗，不得用于食品或药品，不得存放于普通住宅内。
- 为了您的安全和健康，请穿实验服并戴一次性手套操作。

使用说明:

- 10%凝胶聚合催化剂的配制：例如称取0.1g凝胶聚合催化剂，用水溶解，并定容至1ml，即为10%凝胶聚合催化剂。
- 根据目的蛋白的分子量大小选择合适的凝胶浓度，再按照下面的表格配制SDS-PAGE的分离胶(即下层胶)：

不同浓度的SDS-PAGE分离胶的最佳分离范围：

SDS-PAGE分离胶浓度	最佳分离范围
6%胶	50-150kD
8%胶	30-90kD
10%胶	20-80kD
12%胶	12-60kD
15%胶	10-40kD

成分	配制不同体积SDS-PAGE分离胶所需各成分的体积(毫升)					
	5	10	15	20	30	50
6%胶	5	10	15	20	30	50
蒸馏水	2.0	4.0	6.0	8.0	12.0	20.0
30%Acr-Bis(29:1)	1.0	2.0	3.0	4.0	6.0	10.0
1M Tris, pH8.8	1.9	3.8	5.7	7.6	11.4	19.0
10%SDS	0.05	0.1	0.15	0.2	0.3	0.5

10%凝胶聚合催化剂	0.05	0.1	0.15	0.2	0.3	0.5
TEMED Substitute	0.004	0.008	0.012	0.016	0.024	0.04
成分	配制不同体积SDS-PAGE分离胶所需各成分的体积(毫升)					
8%胶	5	10	15	20	30	50
蒸馏水	1.7	3.3	5	6.7	10.0	16.7
30%Acr-Bis(29:1)	1.3	2.7	4.0	5.3	8.0	13.3
1M Tris, pH8.8	1.9	3.8	5.7	7.6	11.4	19.0
10%SDS	0.05	0.1	0.15	0.2	0.3	0.5
10%凝胶聚合催化剂	0.05	0.1	0.15	0.2	0.3	0.5
TEMED Substitute	0.003	0.006	0.009	0.012	0.018	0.03
成分	配制不同体积SDS-PAGE分离胶所需各成分的体积(毫升)					
10%胶	5	10	15	20	30	50
蒸馏水	1.3	2.7	4.0	5.3	8.0	13.3
30%Acr-Bis(29:1)	1.7	3.3	5.0	6.7	10.0	16.7
1M Tris, pH8.8	1.9	3.8	5.7	7.6	11.4	19.0
10%SDS	0.05	0.1	0.15	0.2	0.3	0.5
10%凝胶聚合催化剂	0.05	0.1	0.15	0.2	0.3	0.5
TEMED Substitute	0.002	0.004	0.006	0.008	0.012	0.02
成分	配制不同体积SDS-PAGE分离胶所需各成分的体积(毫升)					
12%胶	5	10	15	20	30	50
蒸馏水	1.0	2.0	3.0	4.0	6.0	10.0
30%Acr-Bis(29:1)	2.0	4.0	6.0	8.0	12.0	20.0
1M Tris, pH8.8	1.9	3.8	5.7	7.6	11.4	19.0
10%SDS	0.05	0.1	0.15	0.2	0.3	0.5
10%凝胶聚合催化剂	0.05	0.1	0.15	0.2	0.3	0.5
TEMED Substitute	0.002	0.004	0.006	0.008	0.012	0.02
成分	配制不同体积SDS-PAGE分离胶所需各成分的体积(毫升)					
15%胶	5	10	15	20	30	50
蒸馏水	0.5	1.0	1.5	2.0	3.0	5.0
30%Acr-Bis(29:1)	2.5	5.0	7.5	10.0	15.0	25.0
1M Tris, pH8.8	1.9	3.8	5.7	7.6	11.4	19.0
10%SDS	0.05	0.1	0.15	0.2	0.3	0.5
10%凝胶聚合催化剂	0.05	0.1	0.15	0.2	0.3	0.5
TEMED Substitute	0.002	0.004	0.006	0.008	0.012	0.02

注：如果配制非变性胶，参考上述配方，不加10%SDS即可配制成非变性PAGE胶。

3. 按照如下表格配制SDS-PAGE的浓缩胶(也称堆积胶、积层胶或上层胶)：

成分	配制不同体积SDS-PAGE浓缩胶所需各成分的体积(毫升)					
5%胶	2	3	4	6	8	10
蒸馏水	1.4	2.1	2.7	4.1	5.5	6.8
30%Acr-Bis(29:1)	0.33	0.5	0.67	1.0	1.3	1.7
1M Tris, pH6.8	0.25	0.38	0.5	0.75	1.0	1.25
10%SDS	0.02	0.03	0.04	0.06	0.08	0.1
10%凝胶聚合催化剂	0.02	0.03	0.04	0.06	0.08	0.1
TEMED Substitute	0.002	0.003	0.004	0.006	0.008	0.01

注：通常10-30分钟内胶会凝固。注：具体的凝固时间和温度及光照有关，说明书中10%凝胶聚合催化剂和TEMED Substitute的正常推荐用量是室温为25°C时的推荐用量。为达到与25°C时相近的凝固时间，当室温低于25°C时，可以适当同时加大10%凝胶聚合催化剂和TEMED Substitute的用量，例如20°C时建议用量是正常推荐用量的1.5倍，15°C时建议用量是正常推荐用量的2倍。

使用本产品的文献：

- Deng XQ, Chen LL, Li NX. The expression of SIRT1 in nonalcoholic fatty liver disease induced by high-fat diet in rats. Liver Int. 2007 Jun; 27(5):708-15.
- Wang PH, Gu ZH, Huang XD, Liu BD, Deng XX, Ai HS, Wang J, Yin ZX, Weng SP,

Yu XQ, He JG. An immune deficiency homolog from the white shrimp, Litopenaeus vannamei, activates antimicrobial peptide genes. Mol Immunol. 2009 May;46(8-9):1897-904.

3. Liang QL, Wang BR, Li GH. DcR3 and survivin are highly expressed in colorectal carcinoma and closely correlated to its clinicopathologic parameters. *J Zhejiang Univ Sci B*. 2009;10(9):675-82.
4. Deng XQ, Cheng JL, Zhang YP, Li NX, Chen LL. Effects of caloric restriction on SIRT1 expression and apoptosis of islet beta cells in type 2 diabetic rats. Springer Verlag. 2010 Dec;47 Suppl 1:177-85.
5. Cao X, Zhang Y, Zou L, Xiao H, Chu Y, Chu X. Persistent oxygen-glucose deprivation induces astrocytic death through two different pathways and calpain-mediated proteolysis of cytoskeletal proteins during astrocytic oncosis. *Neurosci Lett*. 2010;479(2):118-22.
6. Cao X, Xiao H, Zhang Y, Zou L, Chu Y, Chu X. 1, 5-Dicaffeoylquinic acid-mediated glutathione synthesis through activation of Nrf2 protects against OGD/reperfusion-induced oxidative stress in astrocytes. *Brain Res*. 2010;1347:142-8.
7. Huang L, Bi HC, Liu YH, Wang YT, Xue XP, Huang M. CAR-mediated up-regulation of CYP3A4 expression in LS174T cells by Chinese herbal compounds. *Drug Metab Pharmacokinet*. 2011;26(4):331-40.
8. Li Q, Lei RX, Zhou XD, Kolosov VP, Perelman JM. Regulation of PMA-induced MUC5AC expression by heparin in human bronchial epithelial cells. *Mol Cell Biochem*. 2012 Jan;360(1-2):383-91.
9. Luan HF, Zhao ZB, Zhao QH, Zhu P, Xiu MY. Hydrogen sulfide postconditioning protects isolated rat hearts against ischemia and reperfusion injury mediated by the JAK2/STAT3 survival pathway. *Braz J Med Biol Res*. 2012 Oct;45(10):898-905.
10. Chen Z, Qing J, Qin G, Hu L. Construction and characterization of bifunctional TIM-3-EGFP fusion proteins. *Protein Expr Purif*. 2012 Nov;86(1):1-6.
11. Huang L, Huang M, Li YH, Li RM, Zeng Y, Kuang SY, Zhang L, Wang YT, Bi HC. Up-regulation of CYP3A expression through pregnent X receptor by praeeruptorin D isolated from Peucedanumpraeeruptorum Dunn. *J Ethnopharmacol*. 2013 Jul 9;148(2):596-602.
12. Zhang H, Wang ZW, Wu HB, Li Z, Li LC, Hu XP, Ren ZL, Li BJ, Hu ZP. Transforming growth factor- β 1 induces matrix metalloproteinase-9 expression in rat vascular smooth muscle cells via ROS-dependent ERK-NF- κ B pathways. *Mol Cell Biochem*. 2013 Mar;375(1-2):11-21.
13. Liu R, Liu X, Zheng Y, Gu J, Xiong S, Jiang P, Jiang X, Huang E, Yang Y, Ge D, Chu Y. MicroRNA-7 sensitizes non-small cell lung cancer cells to paclitaxel. *Oncol Lett*. 2014 Nov;8(5):2193-2200.
14. Zhao W, Zhao J, Hou M, Wang Y, Zhang Y, Zhao X, Zhang C, Guo D, Hu R and TIA1/TIAL1 are involved in regulation of alternative splicing of SIRT1 pre-mRNA. *Int J Mol Sci*. 2014 Feb 20;15(2):2946-58.
15. Chen D, Tao X, Wang Y, Tian F, Wei Y, Chen G, Shen H, Wang Z, Yu Z, Li H, Chen G. Curcumin accelerates reendothelialization and ameliorates intimal hyperplasia in balloon-injured rat carotid artery via the upregulation of endothelial cell autophagy. *Int J Mol Med*. 2015 Dec;36(6):1563-71.
16. Liu X, Jiang X, Liu R, Wang L, Qian T, Zheng Y, Deng Y, Huang E, Xu F, Wang JY, Chu Y. B cells expressing CD11b effectively inhibit CD4+ T-cell responses and ameliorate experimental autoimmune hepatitis in mice. *Hepatology*. 2015 Nov;62(5):1563-75.
17. He B, Tao H, Liu S, Wei A. Protective effect of carboxymethylated chitosan on hydrogen peroxide-induced apoptosis in nucleus pulposus cells. *Mol Med Rep*. 2015 Mar;11(3):1629-38.
18. Gao CQ, Zhao YL, Li HC, Sui WG, Yan HC, Wang XQ. Heat stress inhibits proliferation, promotes growth, and induces apoptosis in cultured Lantang swine skeletal muscle satellite cells. *J Zhejiang Univ Sci B*. 2015 Jun;16(6):549-59.
19. Li F, Luo J, Wu Z, Xiao T, Zeng O, Li L, Li Y, Yang J. Hydrogen sulfide exhibits cardioprotective effects by decreasing endoplasmic reticulum stress in a diabetic cardiomyopathy rat model. *Mol Med Rep*. 2016 Jul;14(1):865-73.
20. Wang Y, Zhang X, Zhong M, Liu T, Zhang G, Liu S, Guo W, Wei M, He Q, Sun D, Hu S. Improvements of Glucose and Lipid Metabolism After Jejuno-ileal Circuit Procedure in a Non-obese Diabetic Rat Model. *Obes Surg*. 2016 Aug;26(8):1768-76.
21. Wang D, Pan J, Song G, Gao N, Zheng Y, Zhang Q, Li A. Abundance and Significance of Neuroligin-1 and Neurexin II in the Enteric Nervous System of Embryonic Rats. *Biomed Res Int*. 2017;2017:1209360.
22. Wu Y, Si F, Ji X, Jiang K, Song S, Yi Q. Cardiac Protection of Valsartan on Juvenile Rats with Heart Failure by Inhibiting Activity of CaMKII via Attenuating Phosphorylation. *Biomed Res Int*. 2017;2017:4150158.
23. Lin XL, Liu Y, Liu M, Hu H, Pan Y, Fan XJ, Hu XM, Zou WW. Inhibition of Hydrogen Peroxide-Induced Human Umbilical Vein Endothelial Cells Aging by Allicin Depends on Sirtuin1 Activation. *MED SCI MONITOR*. 2017 Jan 31;23:563-570.
24. Gu YH, Wang Y, Bai Y, Liu M, Wang HL. Endoplasmic reticulum stress and apoptosis via PERK-eIF2 α -CHOP signaling in the methamphetamine-induced chronic pulmonary injury. *ENVIRON TOXICOL CHEM*. 2017 Jan;49:194-201.
25. Sun MH, Yang M, Xie FY, Wang W, Zhang L, Shen W, Yin S, Ma JY. DNA Double-Strand Breaks Induce the Nuclear Actin Filaments Formation in Cumulus-Enclosed Oocytes but Not in Denuded Oocytes. *PLoS One*. 2017 Jan 18;12(1):e0170308.
26. Zhang M, Wang D, Zhu T, Yin R. RASSF4 Overexpression Inhibits the Proliferation, Invasion, EMT, and Wnt Signaling Pathway in Osteosarcoma Cells. *Oncol Res*. 2017 Jan 2;25(1):83-91.
27. Zhang HY, Liang F, Zhang JW, Wang F, Wang L, Kang XG. Effects of long noncoding RNA-ROR on tamoxifen resistance of breast cancer cells by regulating microRNA-205. *CANCER CHEMOTH PHARM*. 2017 Feb;79(2):327-337.
28. Li Y, Min D, Wang K, Yin S, Zheng H, Liu L. MicroRNA-152 inhibits cell proliferation, migration and invasion by directly targeting MAFF in nasopharyngeal carcinoma. *Mol Med Rep*. 2017 Feb;15(2):948-956.
29. Wang Q, Shen W, Tao GQ, Sun J, Shi LP. Study on the proliferation of human gastric cancer cell AGS by activation of EGFR in H2O2. *EUR REV MED PHARMACO*. 2017 Mar;21(5):1006-1012.
30. Li S, Zhang X, Zhou Z, Huang Z, Liu L, Huang Z. Downregulation of nucleophosmin expression inhibited proliferation and induced apoptosis in salivary gland adenoid cystic carcinoma. *J Oral Pathol Med*. 2017 Mar;46(3):175-181.
31. Zheng X, Wang C, Xing Y, Chen S, Meng T, You H, Ojima I, Dong Y. SB-T-121205, a next-generation taxane, enhances apoptosis and inhibits migration/invasion in MCF-7/PTX cells. *Int J Oncol*. 2017 Mar;50(3):893-902.
32. Yu L, Yang H, Xiang Y, Guo X, Liu Z, Guo R. Association between Cx37 rs1764390 polymorphism and susceptibility to sepsis in Chinese population. *Infect Genet Evol*. 2017 Mar;48:64-70.
33. Su WJ, Zhang Y, Chen Y, Gong H, Lian YJ, Peng W, Liu YZ, Wang YX, You ZL, Feng SJ, Zong Y, Lu GC, Jiang CL. NLRP3 gene knockout blocks NF- κ B and MAPK signaling pathway in CUMS-induced depression mouse model. *Behav Brain Res*. 2017 Mar 30;322(Pt A):1-8.
34. Dai F, Xuan Y, Jin JJ, Yu S, Long ZW, Cai H, Liu XW, Zhou Y, Wang YN, Chen Z, Huang H. CtBP2 overexpression promotes tumor cell proliferation and invasion in gastric cancer and is associated with poor prognosis. *ONCOTARGET*. 2017 Apr 25;8(17):28736-28749.
35. Lu C, Gao S, Xu G. Geraniin inhibits TNF- α -induced impairments of osteogenesis through NF- κ B and p38 MAPK signalling pathways in bone marrow stem cells. *Stroke Vasc Neurol*. 2017 Apr 21;2(2):47-52.
36. Wang F, Lv P, Gu Y, Li L, Ge X, Guo G. Galectin-1 knockdown improves drug sensitivity of breast cancer by reducing P-glycoprotein expression through inhibiting the Raf-1/AP-1 signaling pathway. *ONCOTARGET*. 2017 Apr 11;8(15):25097-25106.
37. Cui C, Geng L, Shi J, Zhu Y, Yang G, Wang Z, Wang J, Chen M. Structural and electrophysiological dysfunctions due to increased endoplasmic reticulum stress in a long-term pacing model using human induced pluripotent stem cell-derived ventricular cardiomyocytes. *Stem Cell Res Ther*. 2017 May 11;8(1):109.
38. Chen JR, Jia XH, Wang H, Yi YJ, Li YJ. With no interaction, knockdown of Apollon and MDR1 reverse the multidrug resistance of human chronic myelogenous leukemia K562/ADM cells. *Oncol Rep*. 2017 May;37(5):2735-2742.

39. Mu J, Zhu D, Shen Z, Ning S, Liu Y, Chen J, Li Y, Li Z. The repressive effect of miR-148a on Wnt/β-catenin signaling involved in Glabridin-induced anti-angiogenesis in human breast cancer cells. *BMC Cancer*. 2017 May;17(1):307.
40. Wang W, Zhang H, Tang M, Liu L, Zhou Z, Zhang S, Wang L. MicroRNA-592 targets IGF-1R to suppress cellular proliferation, migration and invasion in hepatocellular carcinoma. *Oncol Lett*. 2017 May;13(5):3522-3528.
41. Ding Y, Zhang C, Zhang J, Zhang N, Li T, Fang J, Zhang Y, Zuo F, Tao Z, Tang S, Zhu W, Chen H, Sun X. miR-145 inhibits proliferation and migration of breast cancer cells by directly or indirectly regulating TGF-β1 expression. *Int J Oncol*. 2017 May;50(5):1701-1710.
42. Liu C, Wang M, Sun W, Cai F, Geng S, Su X, Shi Y. PU.1 serves a critical role in the innate defense against Aspergillus fumigatus via dendritic cell-associated C-type lectin receptor-1 and toll-like receptors-2 and 4 in THP-1-derived macrophages. *Mol Med Rep*. 2017 Jun;15(6):4084-4092.
43. Wang N, Zhang Q, Ning B, Luo L, Fang Y. β-Asarone promotes Temozolomide's entry into glioma cells and decreases the expression of P-glycoprotein and MDR1. *Biomed Pharmacother*. 2017 Jun;90:368-374.
44. Lin XL, Hu HJ, Liu YB, Hu XM, Fan XJ, Zou WW, Pan YQ, Zhou WQ, Peng MW, Gu CH. Allicin induces the upregulation of ABCA1 expression via PPARγ/LXRA signaling in THP-1 macrophage-derived foam cells. *Int J Mol Med*. 2017 Jun;39(6):1452-1460.
45. Guo X, Xiang Y, Yang H, Yu L, Peng X, Guo R. Association of the LOX-1 rs1050283 Polymorphism with Risk for Atherosclerotic Cerebral Infarction and its Effect on sLOX-1 and LOX-1 Expression in a Chinese Population. *J Atheroscler Thromb*. 2017 Jun 1;24(6):572-582.
46. Yu Y, Lv F, Liang D, Yang Q, Zhang B, Lin H, Wang X, Qian G, Xu J, You W. HOTAIR may regulate proliferation, apoptosis, migration and invasion of MCF-7 cells through regulating the P53/Akt/JNK signaling pathway. *Biomed Pharmacother*. 2017 Jun;90:555-561.
47. Sun R, Zhu G, Wang J, Tong L, Zhai J. Indirect effects of X-irradiation on proliferation and osteogenic potential of bone marrow mesenchymal stem cells in a local irradiated rat model. *Mol Med Rep*. 2017 Jun;15(6):3706-3714.
48. Wang Q, Ding W, Ding Y, Ma J, Qian Z, Shao J, Li Y. Homoharringtonine suppresses imatinib resistance via the Bcl-6/p53 pathway in chronic myeloid leukemia cell lines. *ONCOTARGET*. 2017 Jun 6;8(23):37594-37604.
49. Du L, Chen X, Duan Z, Liu C, Zeng R, Chen Q, Li M. MiR-146a negatively regulates dectin-1-induced inflammatory responses. *ONCOTARGET*. 2017 Jun 6;8(23):37355-37366.
50. Liu Z, Zeng W, Wang S, Zhao X, Guo Y, Yu P, Yin X, Liu C, Huang T. A potential role for the Hippo pathway protein, YAP, in controlling proliferation, cell cycle progression, and autophagy in BCPAP and KI thyroid papillary carcinoma cells. *Am J Transl Res*. 2017 Jul 15;9(7):3212-3223.
51. Yang M, Bai L, Yu W, Sun X, Xu G, Guan R, Yang Y, Qiu M, Zhang Y, Tian J, Fang H. Expression of autophagy-associated proteins in papillary thyroid carcinoma. *Oncol Lett*. 2017 Jul;14(1):411-415.
52. Huang Q, Liu X, Wu Y, Liao Y, Huang Y, Wei X, Ma M. P38 MAPK pathway mediates cognitive damage in pentylenetetrazole-induced epilepsy via apoptosis cascade. *Epilepsy Res*. 2017 Jul;133:89-92.
53. Zou Q, Tang Q, Pan Y, Wang X, Dong X, Liang Z, Huang D. MicroRNA-22 inhibits cell growth and metastasis in breast cancer via targeting of SIRT1. *Exp Ther Med*. 2017 Aug;14(2):1009-1016.
54. Cheng P, Wang T, Li W, Muhammad I, Wang H, Sun X, Yang Y, Li J, Xiao T, Zhang X. Baicalin Alleviates Lipopolysaccharide-Induced Liver Inflammation in Chicken by Suppressing TLR4-Mediated NF-κB Pathway. *Front Pharmacol*. 2017 Aug 18;8:547.
55. Xu H, Yuan Y, Wu W, Zhou M, Jiang Q, Niu L, Ji J, Liu N, Zhang L, Wang X. Hypoxia stimulates invasion and migration of human cervical cancer cell lines HeLa/SiHa through the Rab11 trafficking of integrin avβ3/FAK/PI3K pathway-mediated Rac1 activation. *J BIOSCI BIOENG*. 2017 Sep;42(3):491-499.
56. Huang JL, Liu W, Tian LH, Chai TT, Liu Y, Zhang F, Fu HY, Zhou HR, Shen JZ. Upregulation of long non-coding RNA MALAT-1 confers poor prognosis and influences cell proliferation and apoptosis in acute monocytic leukemia. *Oncol Rep*. 2017 Sep;38(3):1353-1362.
57. Li H, Xiang Z, Liu Y, Xu B, Tang J. MicroRNA-133b Inhibits Proliferation, Cellular Migration, and Invasion via Targeting LASP1 in Hepatocarcinoma Cells. *Oncol Res*. 2017 Sep 21;25(8):1269-1282.
58. Yang M, Lv Y, Tian X, Lou J, An R, Zhang Q, Li M, Xu L, Dong Z. Neuroprotective Effect of β -Caryophyllene on Cerebral Ischemia-Reperfusion Injury via Regulation of Necroptotic Neuronal Death and Inflammation: In Vivo and in Vitro. *FRONT NEUROSCI-SWITZ*. 2017 Oct 26;11:583.
59. Huang L, Deng M, Zhang S, Lu S, Gui X, Fang Y. β-asarone and levodopa coadministration increases striatal levels of dopamine and levodopa and improves behavioral competence in Parkinson's rat by enhancing dopa decarboxylase activity. *Biomed Pharmacother*. 2017 Oct;94:666-678.
60. Deng L, Liu T, Zhang B, Wu H, Zhao J, Chen J. Forkhead box C1 is targeted by microRNA-133b and promotes cell proliferation and migration in osteosarcoma. *Exp Ther Med*. 2017 Oct;14(4):2823-2830.
61. Zhang J, Wen X, Liu N, Li YQ, Tang XR, Wang YQ, He QM, Yang XJ, Zhang PP, Ma J, Sun Y. Epigenetic mediated zinc finger protein 671 downregulation promotes cell proliferation and tumorigenicity in nasopharyngeal carcinoma by inhibiting cell cycle arrest. *J EXP CLIN CANC RES*. 2017 Oct 19;36(1):147.
62. Wu X, Zhou H, Yue B, Li M, Liu F, Qiu C, Chen B, Ma X. Upregulation of microRNA-25-3p inhibits proliferation, migration and invasion of osteosarcoma cells in vitro by directly targeting SOX4. *Mol Med Rep*. 2017 Oct;16(4):4293-4300.
63. Zhao Y, Yang L, Huang Z, Lin L, Zheng G. Synergistic effects of caffeine and catechins on lipid metabolism in chronically fed mice via the AMP-activated protein kinase signaling pathway. *Eur J Nutr*. 2017 Oct;56(7):2309-2318.
64. Zhao M, Tang S, Xin J, Liu D. Influence of reactive oxygen species on secretory component in the intestinal epithelium during hyperoxia. *Exp Ther Med*. 2017 Nov;14(5):4033-4040.
65. Fang D, He Y, Luan Z. Simvastatin augments activation of liver regeneration through attenuating transforming growth factor-β 1 induced-apoptosis in obstructive jaundice rats. *Exp Ther Med*. 2017 Nov;14(5):4839-4845.
66. Du WL, Fang Q, Chen Y, Teng JW, Xiao YS, Xie P, Jin B, Wang JQ. Effect of silencing the T-Box transcription factor TBX2 in prostate cancer PC3 and LNCaP cells. *Mol Med Rep*. 2017 Nov;16(5):6050-6058.
67. Tao B, Yi J, Huang C, Xu W, Qin C, Chen L, Chen J, Gao Y, Wang R. microRNA-96 regulates the proliferation of nucleus pulposus cells by targeting ARID2/AKT signaling. *Mol Med Rep*. 2017 Nov;16(5):7553-7560.
68. Wang TT, Yuan JH, Ma JZ, Yang WJ, Liu XN, Yin YP, Liu Y, Pan W, Sun SH. CTGF secreted by mesenchymal-like hepatocellular carcinoma cells plays a role in the polarization of macrophages in hepatocellular carcinoma progression. *Biomed Pharmacother*. 2017 Nov;95:111-119.
69. Huang B, Tian Y, Zhang W, Ma Y, Yuan Y, Liu C. Strontium doping promotes bioactivity of rhBMP-2 upon calcium phosphate cement via elevated recognition and expression of BMPR-IA. *COLLOID SURFACE B*. 2017 Nov 1;159:684-695.
70. He X, Li T, Kang N, Zeng H, Ren S, Zong D, Li J, Cai S, Chen P, Chen Y. The protective effect of PRMT6 overexpression on cigarette smoke extract-induced murine emphysema model. *INT J CHRONIC OBSTR*. 2017 Nov 3;12:3245-3254.
71. Liang B, Xiao T, Long J, Liu M, Li Z, Liu S, Yang J. Hydrogen sulfide alleviates myocardial fibrosis in mice with alcoholic cardiomyopathy by downregulating autophagy. *Int J Mol Med*. 2017 Dec;40(6):1781-1791.
72. Lin Y, Wang X, Yu Y, Liu W, Xie F, Ouyang X, Huang Q. Expression and prognostic significance of cyclin-dependent kinase inhibitor 1A in patients with resected gastric adenocarcinoma. *Oncol Lett*. 2017 Dec;14(6):7473-7482.
73. Zhang C, Yu P, Zhu L, Zhao Q, Lu X, Bo S. Blockade of α7 nicotinic acetylcholine receptors inhibit nicotine-induced tumor growth and vimentin expression in non-small cell lung cancer through MEK/ERK signaling way. *Oncol Rep*. 2017 Dec;38(6):3309-3318.
74. Chang M, Wu M, Li H. Curcumin combined with glycyrrhetic acid inhibits the

- development of hepatocellular carcinoma cells by down-regulating the PTEN/PI3K/AKT signalling pathway. *Am J Transl Res*. 2017 Dec 15;9(12):5567-5575.
75. Li E, Zhao Z, Ma B, Zhang J. Long noncoding RNA HOTAIR promotes the proliferation and metastasis of osteosarcoma cells through the AKT/mTOR signaling pathway. *Exp Ther Med*. 2017 Dec;14(6):5321-5328.
 76. Huang W, Cheng P, Yu K, Han Y, Song M, Li Y. Hyperforin attenuates aluminum-induced A β production and Tau phosphorylation via regulating Akt/GSK-3 β signaling pathway in PC12 cells. *Biomed Pharmacother*. 2017 Dec;96:1-6.
 77. Jiao S, Liu Y, Yao Y, Teng J. miR-124 promotes proliferation and differentiation of neuronal stem cells through inactivating Notch pathway. *Cell Biosci*. 2017 Dec 11;7:68.
 78. Wen L, Jiang X, Sun J, Li X, Li X, Tian L, Li Y, Bai W. Cyanidin-3-O-glucoside promotes the biosynthesis of progesterone through the protection of mitochondrial function in Pb-exposed rat Leydig cells. *Food Chem Toxicol*. 2018 Feb;112:427-434.
 79. Wang LM, Zhao N, Zhang J, Sun QF, Yang CZ, Yang PS. Tumor necrosis factor-alpha inhibits osteogenic differentiation of pre-osteoblasts by downregulation of EphB4 signaling via activated nuclear factor-kappaB signaling pathway. *J Periodontal Res*. 2018 Feb;53(1):66-72.
 80. Pu T, Liu Y, Xu R, Li F, Chen S, Sun K. Identification of ZFPM2 mutations in sporadic conotruncal heart defect patients. *Mol Genet Genomics*. 2018 Feb;293(1):217-223.
 81. Wang N, Zhang Q, Luo L, Ning B, Fang Y. β -asarone inhibited cell growth and promoted autophagy via P53/Bcl-2/Bcln-1 and P53/AMPK/mTOR pathways in Human Glioma U251 cells. *J Cell Physiol*. 2018 Mar;233(3):2434-2443.
 82. Ouyang G, Xiong L, Liu Z, Lam B, Bui B, Ma L, Chen X, Zhou P, Wang K, Zhang Z, Huang H, Miao X, Chen W, Wen Y. Inhibition of autophagy potentiates the apoptosis-inducing effects of photodynamic therapy on human colon cancer cells. *PHOTODIAGN PHOTODYN*. 2018 Mar;21:396-403.
 83. Xia PP, Li LJ, Qi RD, Shi JJ, Ju WZ, Chen ML. Electrical and histological remodeling of the pulmonary vein in 2K1C hypertensive rats: Indication of initiation and maintenance of atrial fibrillation. *Anatol J Cardiol*. 2018 Mar;19(3):169-175.
 84. Wang S, Zhao Z, Fan Y, Zhang M, Feng X, Lin J, Hu J, Cheng Z, Sun C, Liu T, Xiong Z, Yang Z, Wang H, Sun D. Mst1 inhibits Sirt3 expression and contributes to diabetic cardiomyopathy through inhibiting Parkin-dependent mitophagy. *BBA-MOL BASIS DIS*. 2018 Apr 17. pii: S0925-4439(18)30133-9.
 85. Cui S, Sun Y, Liu Y, Liu C, Wang J, Hao G, Sun Q. [Corrigendum] MicroRNA-137 has a suppressive role in liver cancer via targeting EZH2. *Mol Med Rep*. 2018 May;17(5):7460.
 86. Liang H, Sun H, Yang J, Yi C. miR-145-5p reduces proliferation and migration of hepatocellular carcinoma by targeting KLF5. *Mol Med Rep*. 2018 Jun;17(6):8332-8338.
 87. Yu G, Chen J, Li S, Pu P, Huang W, Zhao Y, Peng X, Wang R, Lei H. RING finger protein 10 prevents neointimal hyperplasia by promoting apoptosis in vitro and in vivo. *Life Sci*. 2018 Sep 1;208:325-332.
 88. Feng Y, Guo X, Huang X, Wu M, Li X, Wu S, Luo X. Metformin reverses stem cell-like HepG2 sphere formation and resistance to sorafenib by attenuating epithelial-mesenchymal transformation. *Mol Med Rep*. 2018 Oct;18(4):3866-3872.
 89. Wang N, Han Y, Luo L, Zhang Q, Ning B, Fang Y. β -asarone induces cell apoptosis, inhibits cell proliferation and decreases migration and invasion of glioma cells. *Biomed Pharmacother*. 2018 Oct;106:655-664.
 90. Qiu X, Guo Q, Liu X, Luo H, Fan D, Deng Y, Cui H, Lu C, Zhang G, He X, Lu A. Pien Tze Huang Alleviates Relapsing-Remitting Experimental Autoimmune Encephalomyelitis Mice by Regulating Th1 and Th17 Cells. *Front Pharmacol*. 2018 Oct 31;9:1237.
 91. Xiao LY, Wang XR, Yang JW, Ye Y, Zhu W, Cao Y, Ma SM, Liu CZ. Acupuncture Prevents the Impairment of Hippocampal LTP Through β 1-AR in Vascular Dementia Rats. *Mol Neurobiol*. 2018 Oct;55(10):7677-7690.
 92. Liu H, Li X, Yu WQ, Liu CX. Upregulated EFNB2 and EPHB4 promotes lung development in a nitrofen-induced congenital diaphragmatic hernia rat model. *Int J Mol Med*. 2018 Nov;42(5):2373-2382.
 93. Tian X, Wang F, Luo Y, Ma S, Zhang N, Sun Y, You C, Tang G, Li S, Gong Y, Xie C. Protective Role of Nuclear Factor-Erythroid 2-Related Factor 2 Against Radiation-Induced Lung Injury and Inflammation. *Front Oncol*. 2018 Nov 23;8:542.
 94. Li SS, Zhang MX, Wang Y, Wang W, Zhao CM, Sun XM, Dong GK, Li ZR, Yin WJ, Zhu B, Cai HX. Reduction of PGRN increased fibrosis during skin wound healing in mice. *Histo Histopathol*. 2018 Dec 18:18076.
 95. Gao Z, Zhao GS, Lv Y, Peng D, Tang X, Song H, Guo QN. Anoikis-resistant human osteosarcoma cells display significant angiogenesis by activating the Src kinase-mediated MAPK pathway. *Oncol Rep*. 2019 Jan;41(1):235-245.
 96. Zhou ZQ, Chen Y, Chai M, Tao R, Lei YH, Jia YQ, Shu J, Ren J, Li G, Wei WX, Han YD, Han Y. Adipose extracellular matrix promotes skin wound healing by inducing the differentiation of adipose-derived stem cells into fibroblasts. *Int J Mol Med*. 2019 Feb;43(2):890-900.
 97. Wu XB, Feng X, Chang QM, Zhang CW, Wang ZF, Liu J, Hu ZQ, Liu JZ, Wu WD, Zhang ZP, Liu XQ. Cross-talk among AFAP1-AS1, ACVR1 and microRNA-384 regulates the stemness of pancreatic cancer cells and tumorigenicity in nude mice. *J Exp Clin Cancer Res*. 2019 Feb 28;38(1):107.
 98. Huang X, Han X, Huang Z, Yu M, Zhang Y, Fan Y, Xu B, Zhou K, Song L, Wang X, Lu C, Xia Y. Maternal pentachlorophenol exposure induces developmental toxicity mediated by autophagy on pregnancy mice. *Ecotoxicol Environ Saf*. 2019 Mar;169:829-836.
 99. Wang C, Ren YL, Zhai J, Zhou XY, Wu J. Down-regulated LAMA4 inhibits oxidative stress-induced apoptosis of retinal ganglion cells through the MAPK signaling pathway in rats with glaucoma. *Cell Cycle*. 2019 May;18(9):932-948.
 100. Xing Y, Ren S, Ai L, Sun W, Zhao Z, Jiang F, Zhu Y, Piao D. ZNF692 promotes colon adenocarcinoma cell growth and metastasis by activating the PI3K/AKT pathway. *Int J Oncol*. 2019 May;54(5):1691-1703.
 101. Chen YZ, Sun DQ, Zheng Y, Zheng GK, Chen RQ, Lin M, Huang LF, Huang C, Song D, Wu BQ. WISP1 silencing confers protection against epithelial-mesenchymal transition of renal tubular epithelial cells in rats via inactivation of the wnt/ β -catenin signaling pathway in uremia. *J Cell Physiol*. 2019 Jun;234(6):9673-9686.
 102. Shen Y, Chen Q, Li L. Endostar regulates EMT, migration and invasion of lung cancer cells through the HGF-Met pathway. *Mol Cell Probes*. 2019 Jun;45:57-64.
 103. Zuo R, Wang Y, Li J, Wu J, Wang W, Li B, Sun C, Wang Z, Shi C, Zhou Y, Liu M, Zhang C. Rapamycin Induced Autophagy Inhibits Inflammation-Mediated Endplate Degeneration by Enhancing Nrf2/Keap1 Signaling of Cartilage Endplate Stem Cells. *Stem Cells*. 2019 Jun;37(6):828-840.
 104. Liu AH, Chu M, Wang YP. Up-Regulation of Trem2 Inhibits Hippocampal Neuronal Apoptosis and Alleviates Oxidative Stress in Epilepsy via the PI3K/Akt Pathway in Mice. *Neurosci Bull*. 2019 Jun;35(3):471-485.
 105. Liu B, Zhang Q, Wu X, Fu Y, Wang H, Guan Y, Yang S, Liu Y, Cao W, Wang J. Effect of Bletilla striata on the Prevention of Postoperative Peritoneal Adhesions in Abrasion-Induced Rat Model. *Evid Based Complement Alternat Med*. 2019 Jun 10;2019:9148754.
 106. Yang YH, Liu TH, Zhang LD, Chen ZY, Huang XS. Role of the PERK-eIF2 α -CHOP Signaling Pathway in the Effect of Needle Knife Therapy on Knee Joint Chondrocyte Apoptosis. *Evid Based Complement Alternat Med*. 2019 Jun 16;2019:7164916.
 107. Zhuang L, Qu H, Cong J, Dai H, Liu X. MiR-181c affects estrogen-dependent endometrial carcinoma cell growth by targeting PTEN. *Endocr J*. 2019 Jun 28;66(6):523-533.
 108. Wang D, Hu Y. Long Non-coding RNA PVT1 Competitively Binds MicroRNA-424-5p to Regulate CARM1 in Radiosensitivity of Non-Small-Cell Lung Cancer. *Mol Ther Nucleic Acids*. 2019 Jun 7;16:130-140.
 109. Ji Wei Li, Li Wei, Zhijun Han, Zhong Chen. Mesenchymal Stromal Cells-Derived Exosomes Alleviate Ischemia/Reperfusion Injury in Mouse Lung by Transporting Anti-Apoptotic miR-21-5p. *Eur J Pharmacol*. 2019 Jun 5;852:68-76; doi: 10.1016/j.ejphar.2019.01.022.

110. Yuan-Zhen Chen,Dan-Qin Sun,Yi Zheng,Guang-Kuai Zheng,Rong-Quan Chen,Mei Lin,Lian-Fang Huang,Cong Huang,Dan Song,Ben-Qing Wu. WISP1 Silencing Confers Protection Against Epithelial-Mesenchymal Transition of Renal Tubular Epithelial Cells in Rats via Inactivation of the Wnt/ β -Catenin Signaling Pathway in Uremia. *J Cell Physiol.* 2019 Jun;234(6):9673-9686.;doi: 10.1002/jcp.27654.
111. Tian X,An R,Luo Y,Li M,Xu L,Dong Z. Tamibarotene Improves Hippocampus Injury Induced by Focal Cerebral Ischemia-Reperfusion via Modulating PI3K/Akt Pathway in Rats. *J Stroke Cerebrovasc Dis.* 2019 Jul;28(7):1832-1840.
112. Chen JX,Xu X,Zhang S. Silence of long noncoding RNA NEAT1 exerts suppressive effects on immunity during sepsis by promoting microRNA-125-dependent MCEMP1 downregulation. *IUBMB Life.* 2019 Jul;71(7):956-968.
113. Yang QS,Li B,Xu G,Yang SQ,Wang PTang HH,Liu YY. Long noncoding RNA LINC00483/microRNA-144 regulates radiosensitivity and epithelial-mesenchymal transition in lung adenocarcinoma by interacting with HOXA10. *J Cell Physiol.* 2019 Jul;234(7):11805-11821.
114. Zhou GF,Jiang YH,Ma DF,Wang YC,Yang JL,Chen JY,Chi CY,Han XW,Li ZY,Li X. Xiao-Qing-Long Tang Prevents Cardiomyocyte Hypertrophy, Fibrosis, and the Development of Heart Failure with Preserved Ejection Fraction in Rats by Modulating the Composition of the Gut Microbiota. *Biomed Res Int.* 2019 Jul 18;2019:9637479.
115. Cao J,Zhang C,Jiang GQ,Jin SJ,Gao ZH,Wang Q,Yu DC,Ke AW,Fan YQ,Li DW,Wang AQ,Bai DS. Expression of GLS1 in intrahepatic cholangiocarcinoma and its clinical significance. *Mol Med Rep.* 2019 Aug;20(2):1915-1924.
116. Su J,Wang J,Luo J,Li H. Ultrasound-mediated destruction of vascular endothelial growth factor (VEGF) targeted and paclitaxel loaded microbubbles for inhibition of human breast cancer cell MCF-7 proliferation. *Mol Cell Probes.* 2019 Aug;46:101415.
117. Li S,Cao H,Shen D,Chen C,Xing S,Dou F,Jia Q. Efficacy of Shoushen granule on adenosine triphosphate binding cassette transporter A1, proprotein convertase subtilisin/kexin type 9 and toll-like receptor 4/nuclear factor kappa-B signaling pathway. *J Tradit Chin Med.* 2019 Aug;39(4):524-534.
118. Ni T,Gao F,Zhang J,Lin H,Luo H,Chi J,Guo H. Impaired autophagy mediates hyperhomocysteinemia-induced HA-VSMC phenotypic switching. *J Mol Histol.* 2019 Aug;50(4):305-314.
119. Sun Y,Lei B,Huang Q. SOX18 Affects Cell Viability, Migration, Invasiveness, and Apoptosis in Hepatocellular Carcinoma (HCC) Cells by Participating in Epithelial-to-Mesenchymal Transition (EMT) Progression and Adenosine Mo. *Med Sci Monit.* 2019 Aug 20;25:6244-6254.
120. Li Q,Liu B,Breitkopf-Heinlein K,Weng H,Jiang Q,Dong P,Dooley S,Xu K,Ding H. Adenovirus-mediated overexpression of bone morphogenetic protein-9 promotes methionine choline deficiency-induced non-alcoholic steatohepatitis in non-obese mice. *Mol Med Rep.* 2019 Sep;20(3):2743-2753.
121. Jia Q,Cao H,Shen D,Li S,Yan L,Chen C,Xing S,Dou F. Quercetin protects against atherosclerosis by regulating the expression of PCSK9, CD36, PPAR γ , LXRx and ABCA1. *Int J Mol Med.* 2019 Sep;44(3):893-902.
122. Ren L,Qi K,Zhang L,Bai Z,Ren C,Xu X,Zhang Z,Li X. Glutathione Might Attenuate Cadmium-Induced Liver Oxidative Stress and Hepatic Stellate Cell Activation. *Biol Trace Elem Res.* 2019 Oct;191(2):443-452.
123. Gao J,Yin X,Yu X,Dai C,Zhou F. Long noncoding LINC01551 promotes hepatocellular carcinoma cell proliferation, migration, and invasion by acting as a competing endogenous RNA of microRNA-122-5p to regulate ADAM10 expression. *J Cell Biochem.* 2019 Oct;120(10):16393-16407.
124. Cao H,Jia Q,Shen D,Yan L,Chen C,Xing S. Quercetin has a protective effect on atherosclerosis via enhancement of autophagy in ApoE-/- mice. *Exp Ther Med.* 2019 Oct;18(4):2451-2458.
125. Li YJ,Yang Z,Wang YY,Wang Y. Long noncoding RNA ZNF667-AS1 reduces tumor invasion and metastasis in cervical cancer by counteracting microRNA-93-3p-dependent PEG3 downregulation. *Mol Oncol.* 2019 Nov;13(11):2375-2392.
126. Yan R,Jiang Y,Lai B,Lin Y,Wen J. The positive feedback loop FOXO3/CASC11/miR-498 promotes the tumorigenesis of non-small cell lung cancer. *Biochem Biophys Res Commun.* 2019 Nov 12;519(3):518-524.
127. Huang D,Liu M,Jiang Y. Mitochondric acid-5 attenuates TNF- α -mediated neuronal inflammation via activating Parkin-related mitophagy and augmenting the AMPK-Sirt3 pathways. *J Cell Physiol.* 2019 Dec;234(12):22172-22182.
128. Hua X,Liu Z,Zhou M,Tian Y,Zhao PP,Pan WH,Li CX,Huang XX,Liao ZX,Xian Q,Chen B,Hu Y,Leng L,Fang XW,Yu LN. LSAMP-AS1 binds to microRNA-183-5p to suppress the progression of prostate cancer by up-regulating the tumor suppressor DCN. *EBioMedicine.* 2019 Dec;50:178-190.
129. Qin S,Wang H,Liu G,Mei H,Chen M. miR-21-5p ameliorates hyperoxic acute lung injury and decreases apoptosis of AEC II cells via PTEN/AKT signaling in rats. *Mol Med Rep.* 2019 Dec;20(6):4953-4962.
130. Jia Q,Cao H,Shen D,Yan L,Chen C,Xing S. Fisetin, via CKIP-1/REGy, limits oxidized LDL-induced lipid accumulation and senescence in RAW264.7 macrophage-derived foam cells. *Eur J Pharmacol.* 2019 Dec 15;865:172748.
131. Cao H,Jia Q,Yan L,Chen C,Xing S,Shen D. Quercetin Suppresses the Progression of Atherosclerosis by Regulating MST1-Mediated Autophagy in ox-LDL-Induced RAW264.7 Macrophage Foam Cells. *Int J Mol Sci.* 2019 Dec 3;20(23). pii: E6093.
132. Jie Luo,Huaiming Wang,Huanjie Chen,Guolian Gan,Yifeng Zheng. CLDN4 Silencing Promotes Proliferation and Reduces Chemotherapy Sensitivity of Gastric Cancer Cells Through Activation of the PI3K/Akt Signalling Pathway. *Exp Physiol.* 2019 Dec 19.;doi: 10.1113/EP088112.
133. Shan-Shan Li,Hui Cao,Ding-Zhu Shen,Chuan Chen,San-Li Xing,Fang-Fang Dou,Qing-Ling Jia. Effect of Quercetin on Atherosclerosis Based on Expressions of ABCA1, LXRx and PCSK9 in ApoE -/- Mice. *Chin J Integr Med.* 2020 Feb;26(2):114-121.;doi: 10.1007/s11655-019-2942-9.
134. Wang J,Zuo Y,Zhuang K,Luo K,Yan X,Li J,Zhang JH,Liu F. Recombinant Human Milk Fat Globule-Epidermal Growth Factor 8 Attenuates Microthrombosis after Subarachnoid Hemorrhage in Rats. *J Stroke Cerebrovasc Dis.* 2020 Mar;29(3):104536.
135. Wenfeng Xie, Lei Chen, Li Chen, Qiuye Kou. Silencing of long non-coding RNA MALAT1 suppresses inflammation in septic mice: role of microRNA-23a in the down-regulation of MCEMP1 expression. *Inflamm Res.* 2020 Feb;69(2):179-190.
136. Xicheng Wang, Xining Wang, Yunxiu Xu, Maolin Yan, Wenxin Li, Jie Chen, Tao Chen. Effect of nicastrin on hepatocellular carcinoma proliferation and apoptosis through PI3K/AKT signalling pathway modulation. *Cancer Cell Int.* 2020 Mar 24;20:91.
137. Yuanyuan Li, Jiapo Li, Yue Hou, Ling Huang, Yue Bian, Guiyu Song, Chong Qiao. Circadian clock gene Clock is involved in the pathogenesis of preeclampsia through hypoxia. *Life Sci.* 2020 Apr 15;247:117441.
138. Yanjun Li, Ruimin Guo, Mengying Zhang, Peng Chen, Jingxin Li, Yanni Sun. Protective effect of emodin on intestinal epithelial tight junction barrier integrity in rats with sepsis induced by cecal ligation and puncture. *Exp Ther Med.* 2020 Jun;19(6):3521-3530.
139. Xidan Zhu, Jia Feng, Wenguang Fu, Xiaojia Shu, Xue Wan, Jinbo Liu. Effects of cisplatin on the proliferation, invasion and apoptosis of breast cancer cells following β -catenin silencing. *Int J Mol Med.* 2020 Jun;45(6):1838-1850.
140. Ping Wang, Jiming Xu, Weijing You, Yongfeng Hou, Shuiliang Wang, Yujie Ma, Jianming Tan, Zengli Zhang, Wentao Hu, Bingyan Li. Knockdown of CYP2A1 Aggravates 1 α ,25(OH)2 D 3-Inhibited Migration and Invasion of Mouse Ovarian Epithelial Cells by Suppressing EMT. *Front Oncol.* 2020 Jul 29;10:1258.
141. Bing Liu, Lihong Yan, Xuefei Jiao, Xiaozhi Sun, Zonggang Zhao, Junwei Yan, Mingjin Guo, Yunjin Zang. Lycopene Alleviates Hepatic Hypoxia/Reoxygenation Injury Through Nrf2/HO-1 Pathway in AML12 Cell. *J Interferon Cytokine Res.* 2020 Aug;40(8):406-417.
142. Kun Ma, Chuan Zhang, Wuyin Li. TGF- β is associated with poor prognosis and promotes osteosarcoma progression via PI3K/Akt pathway activation. *Cell Cycle.* 2020 Sep;19(18):2327-2339.
143. Yuan Bian, Qiu Li, Qiaolian Li, Ruigen Pan. Silencing of BRF2 inhibits the growth and metastasis of lung cancer cells. *Mol Med Rep.* 2020 Sep;22(3):1767-1774.
144. Xiaofeng Yang, Yanli Song, Yuexi Sun, Mengmeng Wang, Yang Xiang. Down-

- regulation of miR-361-5p promotes the viability, migration and tube formation of endothelial progenitor cells via targeting FGF1. *Biosci Rep.* 2020 Oct 30;40(10):BSR20200557.
145. Yufeng Han, Nanbu Wang, Jian Kang, Yongqi Fang. β -Asarone improves learning and memory in A β 1-42-induced Alzheimer's disease rats by regulating PINK1-Parkin-mediated mitophagy. *Metab Brain Dis.* 2020 Oct;35(7):1109-1117.
146. Niannian Yang, Hao Cheng, Qiao Mo, Xiaobiao Zhou, Minqiang Xie. miR-155-5p downregulation inhibits epithelial-to-mesenchymal transition by targeting SIRT1 in human nasal epithelial cells. *Mol Med Rep.* 2020 Nov;22(5):3695-3704.
147. Guangyu Qian, Xiaoyan Jin, Louwei Zhang. LncRNA FENDRR Upregulation Promotes Hepatic Carcinoma Cells Apoptosis by Targeting miR-362-5p Via NPR3 and p38-MAPK Pathway. *Cancer Biother Radiopharm.* 2020 Nov;35(9):629-639.
148. Yong Yang, Yong Zhou, Yiyong Wang, Xianglong Wei, Lihao Wu, Tingzhong Wang, Aiqun Ma. Exendin-4 reverses high glucose-induced endothelial progenitor cell dysfunction via SDF-1 β /CXCR7-AMPK/p38-MAPK/IL-6 axis. *Acta Diabetol.* 2020 Nov;57(11):1315-1326.
149. Pengcheng Dou, Yu He, Bo Yu, Juan Duan. Downregulation of microRNA-29b by DNMT3B decelerates chondrocyte apoptosis and the progression of osteoarthritis via PTHLH/CDK4/RUNX2 axis. *Aging (Albany NY)*. 2020 Nov 7;13(5):7676-7690.
150. Xiaojie Li, Lihe Yuan, Jin Wang, Zun Zhang, Shaojing Fu, Shaobin Wang, Xinhui Li. MiR-1b up-regulation inhibits rat neuron proliferation and regeneration yet promotes apoptosis via targeting KLF7. *Folia Neuropathol.* 2021;59(1):67-80.
151. Jing Huo, Taibin Liu, Fenghua Li, Xinfeng Song, Xiuli Hou. MicroRNA-21-5p protects melanocytes via targeting STAT3 and modulating Treg/Teff balance to alleviate vitiligo. *Mol Med Rep.* 2021 Jan;23(1):51.
152. Cheng Long, Shiqiang Cen, Zhou Zhong, Chang Zhou, Gang Zhong. FOXO3 is targeted by miR-223-3p and promotes osteogenic differentiation of bone marrow mesenchymal stem cells by enhancing autophagy. *Hum Cell.* 2021 Jan;34(1):14-27.
153. Peng Xiao, Xu Zhu, Jinpeng Sun, Yuhang Zhang, Weijian Qiu, Jianqiang Li, Xuejian Wu. LncRNA NEAT1 regulates chondrocyte proliferation and apoptosis via targeting miR-543/PLA2G4A axis. *Hum Cell.* 2021 Jan;34(1):60-75.
154. Qing Zhang, Ting Chen, Yun Zhang, Lingxia Lyu, Bohuan Zhang, Chengchen Huang, Xuhao Zhou, Yutao Wu, Zhoubin Li. MiR-30c-5p regulates adventitial progenitor cells differentiation to vascular smooth muscle cells through targeting OPG. *Stem Cell Res Ther.* 2021 Jan 19;12(1):67.
155. Chengwei Chen, Yaozeng Xu. Long noncoding RNA LINC00671 exacerbates osteoarthritis by promoting ONECUT2-mediated Smurf2 expression and extracellular matrix degradation. *Int Immunopharmacol.* 2021 Jan;90:106846.
156. Chensheng Zhao, Fan Yang, Xiaona Wei, Jingwen Zhang. miR-139-5p upregulation alleviated spontaneous recurrent epileptiform discharge-induced oxidative stress and apoptosis in rat hippocampal neurons via regulating the Notch pathway. *Cell Biol Int.* 2021 Feb;45(2):463-476.
157. Hui Jiang, Kefeng Cao, Chang Fan, Xiaoya Cui, Yanzhen Ma, Jian Liu. Transcriptome-Wide High-Throughput m6A Sequencing of Differential m6A Methylation Patterns in the Human Rheumatoid Arthritis Fibroblast-Like Synoviocytes Cell Line MH7A. *J Inflamm Res.* 2021 Feb 25;14:575-586.
158. Dezhi Chen, Min Zhang. GASS5 regulates diabetic cardiomyopathy via miR-221-3p/p27 axis-associated autophagy. *Mol Med Rep.* 2021 Feb;23(2):135.
159. Yan Hao, Jie Miao, Wenjia Liu, Li Peng, Yue Chen, Qing Zhong. Formononetin protects against cisplatin-induced acute kidney injury through activation of the PPAR α /Nrf2/HO-1/NQO1 pathway. *Int J Mol Med.* 2021 Feb;47(2):511-522.
160. Liangping Zhao, Li Wang, Daimin Zhang, Yuqi Chen, Fulu Jin. Puerarin alleviates coronary heart disease via suppressing inflammation in a rat model. *Gene.* 2021 Mar 1;771:145354.
161. Linyan Wang, Juan Zhang. Long intergenic ncRNA 00473 improves the invasion of trophoblastic cells via miR-16-5p. *Pregnancy Hypertens.* 2021 Mar;23:174-184.
162. Yunlin Ge, Shuchen Ding, Jiangbiao Feng, Jiangdong Du, Zenghui Gu. Diosgenin inhibits Wnt/ β -catenin pathway to regulate the proliferation and differentiation of MG-63 cells. *Cytotechnology.* 2021 Apr;73(2):169-178.
163. Jiye Chen, Yongjian Zhang, Yongcheng Wang, Ping Jiang, Guofeng Zhou, Zhaoyu Li, Jinlong Yang, Xiao Li. Potential mechanisms of Guizhi decoction against hypertension based on network pharmacology and Dahl salt-sensitive rat model. *Chin Med.* 2021 Apr 27;16(1):34.
164. Juan Xiang, Ying Bian. PWAR6 interacts with miR-106a-5p to regulate the osteogenic differentiation of human periodontal ligament stem cells. *Mol Med Rep.* 2021 Apr;23(4):268.
165. Shiwen Lu, Lifei Yu, Hao Liu. Trimetazidine alleviates hypoxia/reoxygenation-induced apoptosis in neonatal mice cardiomyocytes via up-regulating HMGB1 expression to promote autophagy. *J Recept Signal Transduct Res.* 2021 Apr;41(2):170-179.
166. Hong Xue, Baosong Xie, Nengluan Xu, Hongru Li, Qianshun Chen, Weiping Xie, Hong Wang. Etanercept Protected Against Cigarette Smoke Extract-Induced Inflammation and Apoptosis of Human Pulmonary Artery Endothelial Cells via Regulating TNFR1. *Int J Chron Obstruct Pulmon Dis.* 2021 May 11;16:1329-1345.
167. Changzhao Gao, Fei Yin, Ran Li, Qing Ruan, Chunyang Meng, Kunchi Zhao, Qingsan Zhu. MicroRNA-145-Mediated KDM6A Downregulation Enhances Neural Repair after Spinal Cord Injury via the NOTCH2/Abcb1a Axis. *Oxid Med Cell Longev.* 2021 May 25;2021:2580619.
168. Dong Xie, Yun-Long Li, Gui-Fen Wang, Jian Jiang, Li-Rong Sun. Ultraviolet light-emitting diode irradiation induces reactive oxygen species production and mitochondrial membrane potential reduction in HL-60 cells. *J Int Med Res.* 2021 May;49(5):3000605211016623.
169. Wei Feng. Tectorigenin attenuates cognitive impairments in mice with chronic cerebral ischemia by inhibiting the TLR4/NF- κ B signaling pathway. *Biosci Biotechnol Biochem.* 2021 Jun 24;85(7):1665-1674.
170. Xiaohui Xu, Yong Wu, Ke Yi, Yan Hu, Weiqun Ding, Chungen Xing. IRF1 regulates the progression of colorectal cancer via interferon-induced proteins. *Int J Mol Med.* 2021 Jun;47(6):104.
171. Lijie Wang, Lu Wang, Qi Wang. Constitutive activation of the NEAT1/miR-22-3p/Ltb4r1 signaling pathway in mice with myocardial injury following acute myocardial infarction. *Aging (Albany NY)*. 2021 Jun 3;13(11):15307-15319.
172. Hong Chen, Yuan Zhao, Jiajia Zhang, Yan Xie, Mulan Jin. Promoting effects of MiR-135b on human multiple myeloma cells via regulation of the Wnt/ β -catenin/Versican signaling pathway. *Cytokine.* 2021 Jun;142:155495.
173. Jianzhang Zheng, Yunshuo Lin, Faqiang Tang, Huiling Guo, Laipeng Yan, Shiping Hu, Hong Wu. Promotive Role of CircATRNL1 on Chondrogenic Differentiation of BMSCs Mediated by miR-338-3p. *Arch Med Res.* 2021 Jul;52(5):514-522.
174. Xuan Ma, Yang Liu, Hao Tian, Bo Zhang, Meiling Wang, Xia Gao. LINC01272 Suppressed Cell Multiplication and Induced Apoptosis Via Regulating MiR-7-5p/CRLS1 Axis in Lung Cancer. *J Microbiol Biotechnol.* 2021 Jul 28;31(7):921-932.
175. Weiting Chen, Yun Guan, Fangfang Xu, Beizhan Jiang. 4-Methylumbelliflferone promotes the migration and odontogenetic differentiation of human dental pulp stem cells exposed to lipopolysaccharide in vitro. *Cell Biol Int.* 2021 Jul;45(7):1415-1422.
176. Bin Liu, Wei Gan, Zhang Jin, Meng Wang, Guopeng Cui, Hongyu Zhang, Huifu Wang. The Role of miR-34c-5p in Osteogenic Differentiation of Bone Marrow Mesenchymal Stem Cells. *Int J Stem Cells.* 2021 Aug 30;14(3):286-297.
177. Huiming Guo, Jianyuan Gao, Yuan Qian, Huawei Wang, Jiang Liu, Qingyan Peng, Yong Zhou, Kunhua Wang. miR-125b-5p inhibits cell proliferation by targeting ASCT2 and regulating the PI3K/AKT/mTOR pathway in an LPS-induced intestinal mucosa cell injury model. *Exp Ther Med.* 2021 Aug;22(2):838.
178. Xingli Liu, Wei Song, Xiangmin Zhang, Fangmin Long, Junkun Yin, Xiao He, Liang Lv. Downregulating LncRNA XIST attenuated contrast-induced nephropathy injury via regulating miR-133a-3p/NLRP3 axis. *J Thromb Thrombolysis.* 2021 Aug;52(2):440-453.
179. Wensi Wu, Xiaojun Zhang, Jiaxin Zhou, Hongmei Yang, Junjun Chen, Le Zhao, Junying Zhong, Wei-Jye Lin, Zhi Wang. Clemastine Ameliorates Perioperative Neurocognitive Disorder in Aged Mice Caused by Anesthesia and Surgery. *Front Pharmacol.* 2021 Aug 23;12:738590.
180. Kangjun Zhang, Taishi Fang, Yajie Shao, Yanhui Wu. TGF- β -MTA1-SMAD7-

- SMAD3-SOX4-EZH2 Signaling Axis Promotes Viability, Migration, Invasion and EMT of Hepatocellular Carcinoma Cells. *Cancer Manag Res.* 2021 Sep 10;13:7087-7099.
181. Lan Ma, Di Wu. MicroRNA-383-5p regulates osteogenic differentiation of human periodontal ligament stem cells by targeting histone deacetylase 9. *Arch Oral Biol.* 2021 Sep;129:105166.
182. Qian Xu, Yonghui Song, Lili Lu. Overexpression of let-7d explains down-regulated KDM3A and ENO2 in the pathogenesis of preeclampsia. *J Cell Mol Med.* 2021 Sep;25(17):8127-8139.
183. Zhan Wang, Rui Chen, Zheming Xu, Wei Ru, Hongjuan Tian, Fan Yang, Chang Tao. MiR-155-5p promotes renal interstitial fibrosis in obstructive nephropathy via inhibiting SIRT1 signaling pathway. *J Recept Signal Transduct Res.* 2021 Oct;41(5):466-475.
184. Di Zhao, Bo Yang, Chen Ye, Shaoyi Zhang, Xiaoqiong Lv, Qiyi Chen. Enteral nutrition ameliorates the symptoms of Crohn's disease in mice via activating special pro-resolving mediators through innate lymphoid cells. *Innate Immun.* 2021 Oct;27(7-8):533-542.
185. Zhihong Lu, Xiujuan Zhu, Yuhong Ye, Haidong Fu, Jianhua Mao. PP2A protects podocytes against Adriamycin-induced injury and epithelial-to-mesenchymal transition via suppressing JIP4/p38-MAPK pathway. *Cytotechnology.* 2021 Oct;73(5):697-713.
186. Chunju Zheng, Ting Lu, Zhimin Fan. miR-200b-3p alleviates TNF- α -induced apoptosis and inflammation of intestinal epithelial cells and ulcerative colitis progression in rats via negatively regulating KHDRBS1. *Cytotechnology.* 2021 Oct;73(5):727-743.
187. Luping Liu, Li Sun, Jing Zheng, Li Cui. Berberine modulates Keratin 17 to inhibit cervical cancer cell viability and metastasis. *J Recept Signal Transduct Res.* 2021 Dec;41(6):521-531.
188. Jiahuan Li, Xiaoling Deng, Shuhan Wang, Qianqian Jiang, Keshu Xu. Resolvin D1 attenuates CCl4 Induced Liver Fibrosis by Inhibiting Autophagy-Mediated HSC activation via AKT/mTOR Pathway. *Front Pharmacol.* 2021 Dec 20;12:792414.
189. Yong Lin, Yushan Chen, Rongqiang Shen, Dingzhu Chen, Yimin Lin. MicroRNA-148a-3p suppresses cell proliferation and migration of esophageal carcinoma by targeting CEP55. *Cell Mol Biol Lett.* 2021 Dec 24;26(1):54.
190. Na Li, Rui Hou, Caixia Liu, Tian Yang, Chong Qiao, Jun Wei. Integration of transcriptome and proteome profiles in placenta accreta reveals trophoblast overmigration as the underlying pathogenesis. *Clin Proteomics.* 2021 Dec 29;18(1):31.
191. Wangchao Dong, Shanshan Wang, Weiqing Qian, Suming Li, Peimin Wang. Cedrol alleviates the apoptosis and inflammatory response of IL-1 β -treated chondrocytes by promoting miR-542-5p expression. *In Vitro Cell Dev Biol Anim.* 2021 Dec;57(10):962-972.
192. Li Hua, Xiaohong Zhang. MALAT1 regulates osteogenic differentiation of human periodontal ligament stem cells through mediating miR-155-5p/ETS1 axis. *Tissue Cell.* 2021 Dec;73:101619.
193. Xu Li, Xin Zhang, Yajun Liu, Ruihan Pan, Xiaolong Liang, Lifang Huang, Chao Yang. Exosomes derived from mesenchymal stem cells ameliorate oxygen-glucose deprivation/reoxygenation-induced neuronal injury via transferring MicroRNA-194 and targeting Bach1. *Tissue Cell.* 2021 Dec;73:101651.
194. Mingfang Huang, Xuqiang Liao, Liang Li, Gao Li, Minbiao Chen. MiR-552-3p facilitated cell proliferation, migration and invasion by sponging Fibulin 5 in non-small cell lung cancer via activation of ERK/GSK3 β /catenin signaling pathway. *Tissue Cell.* 2021 Dec;73:101672.
195. Huanhuan Zhu, Shaopeng Tong, Yan Cui, Xiaodong Wang, Minying Wang. Tanshinol alleviates ulcerative colitis by promoting the expression of VLDLR. *Drug Dev Res.* 2021 Dec;82(8):1258-1268.
196. Qingyu Liang, Jianqi Wu, Xin Zhao, Shuai Shen, Chen Zhu, Tianqi Liu, Xiao Cui, Ling Chen, Chunmi Wei, Peng Cheng, Wen Cheng, Anhua Wu. Establishment of tumor inflammasome clusters with distinct immunogenomic landscape aids immunotherapy. *Theranostics.* 2021 Oct 17;11(20):9884-9903.
197. Huina Wang, Xiuli Yi, Sen Guo, Sijia Wang, Jinyuan Ma, Tao Zhao, Qiong Shi, Yangzi Tian, Hao Wang, Lintao Jia, Tianwen Gao, Chunying Li, Weinan Guo. The XBP1-MARCH5-MFN2 Axis Confers Endoplasmic Reticulum Stress Resistance by Coordinating Mitochondrial Fission and Mitophagy in Melanoma. *J Invest Dermatol.* 2021 Dec;141(12):2932-2943.e12.
198. Wenjin Zhang, Yaowei Liu, Zhibin Li, Shu Xu, Jie Zhang, Kasper Hettinga, Peng Zhou. Effects of microfiltration combined with ultrasonication on shelf life and bioactive protein of skim milk. *Ultrason Sonochem.* 2021 Sep;77:105668.
199. Na Fan, Kepan Chen, Rong Zhu, Zhongwei Zhang, Hai Huang, Shugang Qin, Qian Zheng, Zhongshan He, Xi He, Wen Xiao, Yupei Zhang, Yongjun Gu, Changchun Zhao, Yongmei Liu, Xin Jiang, Shuaicheng Li, Yuquan Wei, Xiangrong Song. Manganese-coordinated mRNA vaccines with enhanced mRNA expression and immunogenicity induce robust immune responses against SARS-CoV-2 variants. *Sci Adv.* 2022 Dec 23;8(51):eabq3500.
200. Mingyang Li, Xinyang Jin, Tao Liu, Feng Fan, Feng Gao, Shuang Chai, Lihua Yang. Nanoparticle elasticity affects systemic circulation lifetime by modulating adsorption of apolipoprotein A-I in corona formation. *Nat Commun.* 2022 Jul 16;13(1):4137.
201. Weili Wang, Huizhen Zheng, Jun Jiang, Zhi Li, Dongpeng Jiang, Xiangru Shi, Hui Wang, Jie Jiang, Qianqian Xie, Meng Gao, Jianhong Chu, Xiaoming Cai, Tian Xia, Ruibin Li. Engineering micro oxygen factories to slow tumour progression via hyperoxic microenvironments. *Nat Commun.* 2022 Aug 2;13(1):4495.
202. Huan Wang, Qi Cheng, Kaikai Chang, Lingjie Bao, Xiaofang Yi. Integrated Analysis of Ferroptosis-Related Biomarker Signatures to Improve the Diagnosis and Prognosis Prediction of Ovarian Cancer. *Front Cell Dev Biol.* 2022 Jan 5;9:807862.

注：更多使用本产品的文献请参考产品网页。

Version 2023.11.13