



碧云天生物技术/Beyotime Biotechnology
订货热线: 400-168-3301 或 800-8283301
订货 e-mail: order@beyotime.com
技术咨询: info@beyotime.com
网址: <http://www.beyotime.com>

Western二抗稀释液

产品编号	产品名称	包装
P0023D-100ml	Western二抗稀释液	100ml
P0023D-500ml	Western二抗稀释液	500ml

产品简介:

- 碧云天生产的Western二抗稀释液(Secondary Antibody Dilution Buffer)可以用于Western时二抗(secondary antibody)的稀释。经本Western二抗稀释液稀释的二抗可以在1-2周内重复使用3-5次。
- 本产品的有效成分为经过反复优化的适量BSA、适当的去垢剂和二抗稳定试剂等，一方面能有效减少二抗的非特异性结合，另一方面能有效提升稀释后二抗的稳定保存时间。
- 按照每个二抗稀释10毫升计算，一个100ml包装的Western二抗稀释液可以稀释10个或10次二抗，一个500ml包装的Western二抗稀释液可以稀释50个或50次二抗。

包装清单:

产品编号	产品名称	包装
P0023D-100ml	Western二抗稀释液	100ml
P0023D-500ml	Western二抗稀释液	500ml
—	说明书	1份

保存条件:

4°C保存，一年有效。长期不使用可以-20°C保存。

注意事项:

- 本Western二抗稀释液经过滤除菌处理，使用过程中应尽量避免细菌污染。一旦启用后建议在一个月内用完，如果发现有沉淀或细菌污染请弃用。
- 本产品仅限于专业人员的科学研究用，不得用于临床诊断或治疗，不得用于食品或药品，不得存放于普通住宅内。
- 为了您的安全和健康，请穿实验服并戴一次性手套操作。

使用说明:

- 参考所用的二抗的使用说明，以及一抗的质量和样品中目的蛋白的含量，按照适当比例例如1:1000、1:500等比例稀释二抗。二抗稀释后即可直接用于Western。一次Western结束后，可以回收稀释的二抗，4°C保存，以用于下次的Western，通常在1-2周内可以重复使用3-5次。
- 详细的Western操作可以参考我们的相关网页：<http://www.beyotime.com/support/western.htm>。

相关产品:

产品编号	产品名称	包装
P0023A-100ml	Western一抗稀释液	100ml
P0023A-500ml	Western一抗稀释液	500ml
P0023B-100ml	Western封闭液	100ml
P0023B-500ml	Western封闭液	500ml
P0023D-100ml	Western二抗稀释液	100ml
P0023D-500ml	Western二抗稀释液	500ml
P0220	Quickblock™封闭液(PBS)	100ml
P0222	Quickblock™封闭液(PBSTw)	100ml
P0226	Quickblock™封闭液(PBSTx)	100ml
P0228	Quickblock™封闭液(TBS)	100ml
P0231	Quickblock™封闭液(TBSTw)	100ml
P0233	Quickblock™封闭液(TBSTx)	100ml
P0235	Quickblock™封闭液(10X)	100ml
P0252-100ml	QuickBlock™ Western封闭液	100ml

P0252-500ml	QuickBlock™ Western封闭液	500ml
P0256-100ml	QuickBlock™ Western一抗稀释液	100ml
P0256-500ml	QuickBlock™ Western一抗稀释液	500ml
P0258-100ml	QuickBlock™ Western二抗稀释液	100ml
P0258-500ml	QuickBlock™ Western二抗稀释液	500ml

使用本产品的文献：

1. Zhang Z, Wang Q, Li P, Zhou Y, Li S, Yi W, Chen A, Kong P, Hu C. Overexpression of the Interferon regulatory factor 4-binding protein in human colorectal cancer and its clinical significance. *Cancer Epidemiol.* 2009 Aug;33(2):130-6.
2. Mao SS, Hua R, Zhao XP, Qin X, Sun ZQ, Zhang Y, Wu YQ, Jia MX, Cao JL, Zhang YM. Exogenous Administration of PACAP Alleviates Traumatic Brain Injury in Rats through a Mechanism Involving the TLR4/MyD88/NF-κB Pathway. *J Neurotrauma.* 2012 Jul 1; 29(10):1941-59.
3. Qin X, Sun ZQ, Dai XJ, Mao SS, Zhang JL, Jia MX, Zhang YM. Toll-like receptor 4 signaling is involved in PACAP-induced neuroprotection in BV2 microglial cells under OGD/reoxygenation. *Neurol Res.* 2012 May;34(4):379-89.
4. Jian CX, Yang MZ, Li P, Xiong J, Zhang ZJ, Li CJ, Chen A, Hu CM, Zhou JX, Li SH. Ectopically expressed IBP promotes cell proliferation in oral squamous cell carcinoma. *Cancer Invest.* 2012 Dec;30(10):748-56.
5. Mao SS, Hua R, Zhao XP, Qin X, Sun ZQ, Zhang Y, Wu YQ, Jia MX, Cao JL, Zhang YM. Exogenous administration of PACAP alleviates traumatic brain injury in rats through a mechanism involving the TLR4/MyD88/NF-κB pathway. *J Neurotrauma.* 2012 Jul 1; 29(10):1941-59.
6. Qin X, Sun ZQ, Zhang XW, Dai XJ, Mao SS, Zhang YM. TLR4 signaling is involved in the protective effect of propofol in BV2 microglia against OGD/reoxygenation. *J Physiol Biochem.* 2013 Dec; 69(4):707-18.
7. Ye HH, Hua R, Yu L, Wu KJ, Fei SJ, Qin X, Song Y, Cao JL, Zhang YM. Abnormal Expression of Toll-Like Receptor 4 Is Associated with Susceptibility to Ethanol-Induced GastricMucosal Injury in Mice. *Dig Dis Sci.* 2013 Oct;58(10):2826-39.
8. Ye HH, Wu KJ, Fei SJ, Zhang XW, Liu HX, Zhang JL, Zhang YM. Propofol participates in gastric mucosal protection through inhibiting the toll-like receptor-4/nuclear factor kappa-B signaling pathway. *Clin Res Hepatol Gastroenterol.* 2013 Feb;37(1):e3-15.
9. Yang Z, Liu J, Liu H, Qiu M, Liu Q, Zheng L, Pang M, Quan F, Zhang Y. Isolation and characterization of SSEA3(+) stem cells derived from goat skin fibroblasts. *Cell Reprogram.* 2013 Jun;15(3):195-205.
10. Zhang Y, Shi W, Tang S, Li J, Yin S, Gao X, Wang L, Zou L, Zhao J, Huang Y, Shan L, Gounni AS, Wu Y, Yuan F, Zhang J. The influence of cathelicidin LL37 in human anti-neutrophils cytoplasmic antibody (ANCA)-associated vasculitis. *Arthritis Res Ther.* 2013 Oct 24; 15(5):R161.
11. Wang ZQ, Wang J, Ling WH, Zhang XG, Shi Q. Effects of CD40 ligation combined with chemotherapy drugs on human breast cancer cell lines. *J Int Med Res.* 2013 Oct;41(5):1495-504.
12. Wang Z, Li Q, Zhang B, Lu Y, Yang Y, Ban D, Zhang H. Single nucleotide polymorphism scanning and expression of the FRZB gene in pig populations. *Gene.* 2014 Jun 15;543(2):198-203.
13. Ye M, Zheng J, Chen X, Chen X, Wu X, Lin X, Liu Y. Prednisone inhibits the focal adhesion kinase/receptor activator of NF-κB ligand/mitogen-activated protein kinase signaling pathway in rats with adriamycin-induced nephropathy. *Mol Med Rep.* 2015 Nov; 12(5):7471-8.
14. Dai XJ, Li N, Yu L, Chen ZY, Hua R, Qin X, Zhang YM. Activation of BV2 microglia by lipopolysaccharide triggers an inflammatory reaction in PC12 cell apoptosis through a toll-like receptor 4-dependent pathway. *Cell Stress Chaperones.* 2015 Mar;20(2):321-31.
15. Zhao WJ, Song Q, Zhang ZJ, Mao L, Zheng WJ, Hu X, Lian HZ. The Kinetic Response of the Proteome in A549 Cells Exposed to ZnSO4 Stress. *PLoS One.* 2015 Jul 21;10(7):e0133451.
16. Zhang G, Zhao BX, Hua R, Kang J, Shao BM, Carbonaro TM, Zhang YM. Hippocampal microglial activation and glucocorticoid receptor down-regulation precipitate visceral hypersensitivity induced by colorectal distension in rats. *Neuropharmacology.* 2015 Dec 2; 102:295-303.
17. Liang W, Zhuo Y, Xiong C, Zheng Y, Chai Y, Yuan R. Ultrasensitive Cytosensor Based on Self-Enhanced Electrochemiluminescent Ruthenium-Silica Composite Nanoparticles for Efficient Drug Screening with Cell Apoptosis Monitoring. *Anal Chem.* 2015 Dec 15; 87(24):12363-71.
18. Li PP, Lu K, Geng LY, Zhou XX, Li XY, Wang X. Bruton's tyrosine kinase inhibitor restrains Wnt signaling in chronic lymphocytic leukemia. *Mol Med Rep.* 2016 Jun;13(6):4934-8.
19. Chen Y, Huang W, Jiang W, Wu X, Ye B, Zhou X. HIV-1 Tat Regulates Occludin and Aβ Transfer Receptor Expression in Brain Endothelial Cells viaRho/ROCK Signaling Pathway. *Oxid Med Cell Longev.* 2016;2016:4196572.
20. Xie L, Xiang GH, Tang T, Tang Y, Zhao LY, Liu D, Zhang YR, Tang JT, Zhou S, Wu DH. Capsaicin and dihydrocapsaicin induce apoptosis in human glioma cells via ROS and Ca2+ mediated mitochondrial pathway. *Mol Med Rep.* 2016 Nov;14(5):4198-4208.
21. Zhang G, Yu L, Chen ZY, Zhu JS, Hua R, Qin X, Cao JL, Zhang YM. Activation of corticotropin-releasing factor neurons and microglia in paraventricular nucleus precipitates visceral hypersensitivity induced by colorectal distension in rats. *Brain Behav Immun.* 2016 Jul;55:93-104.
22. Zhang B, Shang P, Qiangba Y, Xu A, Wang Z, Zhang H. The association of NR1H3 gene with lipid deposition in the pig. *Lipids Health Dis.* 2016 May 26;15:99.
23. Zhang B, Qiangba Y, Shang P, Lu Y, Yang Y, Wang Z, Zhang H. Gene expression of vascular endothelial growth factor A and hypoxic adaptation in Tibetan pig. *J Anim Sci Biotechnol.* 2016 Apr 2;7:21.
24. Zhang G, Zhao BX, Hua R, Kang J, Shao BM, Carbonaro TM, Zhang YM. Hippocampal microglial activation and glucocorticoid receptor down-regulation precipitate visceral hypersensitivity induced by colorectal distension in rats. *Neuropharmacology.* 2016 Mar; 102:295-303.
25. Zhang F, Cao S, Zhu Z, Yang Y, Wen X, Chang YF, Huang X, Wu R, Wen Y, Yan Q, Huang Y, Ma X, Zhao Q. Immuno protective Efficacy of Six In vivo-Induced Antigens against *Actinobacillus pleuropneumoniae* as Potential Vaccine Candidates in Murine Model. *Front Microbiol.* 2016 Oct 21;7:1623.
26. Liu B, Li Y, Ji H, Lu H, Li H, Shi Y. Glutamine attenuates obstructive cholestasis in rats via farnesoid X receptor-mediated regulation of Bsep and Mrp2. *Can J Physiol Pharmacol.* 2017 Feb;95(2):215-223.
27. Ye W, Zhong Z, Zhu S, Zheng S, Xiao J, Song S, Yu H, Wu Q, Lin Z, Chen J. Advanced oxidation protein products induce chondrocyte death through a redox-dependent, poly (ADP-ribose) polymerase-1-mediated pathway. *Apoptosis.* 2017 Jan;22(1):86-97.

Version 2019.09.30