BeyoMag™ Anti-Flag Magnetic Beads (Anti-Flag磁珠)

产品编号	产品名称	包装
P2115-0.5ml	BeyoMag™ Anti-Flag Magnetic Beads (Anti-Flag磁珠)	0.5ml
P2115-2ml	BeyoMag™ Anti-Flag Magnetic Beads (Anti-Flag磁珠)	2ml

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Characteristics	Description
Product content	10mg/ml magnetic beads in specific protective buffer
Beads size	~200nm
Magnetization	Superparamagnetic
Coupled antibody	Anti-Flag mouse monoclonal antibody
Isotype	IgG2b
M.W. of antibody	Approximately 150kDa
Antibody concentration	≥0.6mg Flag antibody per ml beads
Binding capacity	≥ 0.6mg Flag - tagged fusion protein per ml beads
Specificity	Met-Flag-Protein, Flag-Protein, Protein-Flag
Elution method	Acid, peptide competitive or SDS - PAGE loading buffer elution. Note: If elute with SDS-PAGE loading buffer, the light (~25kDa) and heavy (~50kDa) chain of Flag antibody will be denatured and release from the beads.
Application	IP, Co-IP, Protein purification

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产品编号	产品名称	包装
P2115-0.5ml	BeyoMag™ Anti-Flag Magnetic Beads (Anti-Flag磁珠)	0.5ml
P2115-2ml	BeyoMag™ Anti-Flag Magnetic Beads (Anti-Flag磁珠)	2ml
—	说明书	1份

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Problem	Possible Causes	Solution
Very few or no Flag - tagged protein exists in the eluate.	Protein is not completely eluted.	Change elution methods.
	No target protein expressed.	Make sure the protein of interest contains the Flag-tag by Western blot or dot blot analyses.
	Very low protein expression level.	1. Use larger volume of cell lysate. 2. Optimize expression conditions to raise the protein expression level.
	Washes are too stringent.	Reduce the time and number of washes.
	Incubation times are inadequate.	Increase the incubation time.
	Interfering substance is present in sample.	Lysates containing high concentration of DTT, 2-mercaptoethanol, or other reducing agents may destroy antibody function, and must be avoided.
	Detection system is inadequate.	If Western blot detection is used: 1. Check primary and secondary antibodies using proper controls to confirm binding and reactivity. 2. Verify that the transfer was adequate by using prestained protein marker or staining the membrane with Ponceau S. 3. Use fresh detection substrate or try a different detection system.
Background is too high.	Proteins bind nonspecifically to the Anti-Flag monoclonal antibody, insufficient washing on magnetic beads, or the microcentrifuge tubes.	1. Pre-clear lysate with Mouse IgG Magnetic Beads (P2171) to remove nonspecific binding proteins. 2. After suspending beads for the final wash, transfer entire sample to a clean microcentrifuge tube before centrifugation.
	Washes are insufficient.	1. Increase the number of washes. 2. Prolong duration of the washes, incubating each wash for at least 15 minutes. 3. Increase the salt and/or detergent concentrations in the wash solutions. 4. Centrifuge at lower speed to avoid nonspecific trapping of denatured proteins.
Multiple protein bands found in the eluate.	The protein is not stable at room temperature.	Purify the target protein at lower temperature, such as 4 $^{\circ}$ C.
	Protein degradation due to proteases activity during purification process.	Add protease inhibitors to cell lysate.
	Non - specific binding.	1. Prepare cell lysate again. 2. Add additional wash steps.

Version 2024.01.10