RNase Inhibitor

产品编号	产品名称	包装
R0102-2kU	RNase Inhibitor	2000U
R0102-10kU	RNase Inhibitor	10000U
R0102-50kU	RNase Inhibitor	50000U

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1. Chao Yu, Yin-Li Zhang, Heng-Yu Fan. Selective Smad4 knockout in ovarian preovulatory follicles results in multiple defects in ovulation. *Mol Endocrinol*. 2013 Jun;27(6):966-78.
2. Wang K, Zhang K, Lv Z, Zhu X, Zhu L, Zhou F. Ultrasensitive detection of microRNA with isothermal amplification and a time-resolved fluorescence sensor. *Biosens Bioelectron*. 2014 Jul 15;57:91-5.
3. Wu Y, Dong X, Kadowaki T. Characterization of the Copy Number and Variants of Deformed Wing Virus (DWV) in the Pairs of Honey Bee Pupa and Infesting *Varroa destructor* or *Tropilaelaps mercedesae*. *Front Microbiol*. 2017 Aug 22;8:1558.
4. Wang X, Hu H, Liu H. RNA binding protein Lin28B confers gastric cancer cells stemness via directly binding to NRP-1. *Biomed Pharmacother*. 2018 Aug;104:383-389.
5. Wang H, Wang X, Li X, Wang Q, Qing S, Zhang Y, Gao MQ. A novel long non-coding RNA regulates the immune response in MAC-T cells and contributes to bovine mastitis. *FEBS J*. 2019 May;286(9):1780-1795.
6. Ya-Xin Tan, Ying Hong, Shui Jiang, Min-Nan Lu, Shan Li, Bo Chen, Li Zhang, Tao Hu, Rui Mao, Rong Mei, Yan-Bin Xiyang. MicroRNA-449a regulates the progression of brain aging by targeting SCN2B in SAMP8 mice. *Int J Mol Med*. 2020 Apr;45(4):1091-1102.
7. Song Xiang, Xiaolong Liang, Shuo Yin, Jian Liu, Zheng Xiang. N6-methyladenosine methyltransferase METTL3 promotes colorectal cancer cell proliferation through enhancing MYC expression. *Am J Transl Res*. 2020 May 15;12(5):1789-1806.
8. Yunfei Wu, Qiushi Liu, Benjamin Weiss, Martin Kaltenpoth, Tatsuhiko Kadowaki. Honey Bee Suppresses the Parasitic Mite *Vitellogenin* by Antimicrobial Peptide. *Front Microbiol*. 2020 May 25;11:1037.
9. Weijia Xie, Youhao Wang, Yao Zhang, Ying Xiang, Na Wu, Long Wu, Chengying Li, Tongjian Cai, Xiangyu Ma, Zubin Yu, Li Bai, Yafei Li. Single-nucleotide polymorphism rs4142441 and MYC co-modulated long non-coding RNA OSER1-AS1 suppresses non-small cell lung cancer by sequestering ELAVL1. *Cancer Sci*. 2020 Oct 28.;doi:10.1111/cas.14713.
10. Xiaolong Liu, Zhongyang Shen. LncRNA TMPO-AS1 Aggravates the Development of Hepatocellular Carcinoma via miR-429/GOT1 Axis. *Am J Med Sci*. 2020 Dec;360(6):711-720.
11. Jing Yan, Tao Fang, Ming Zhang, Quan Zhou. LINC00467 facilitates osteosarcoma progression by sponging miR-217 to regulate KPNA4 expression. *Int J Mol Med*. 2021 Mar;47(3):26.
12. Jingsheng Yuan, Tao Lv, Jian Yang, Zhenru Wu, Lvnan Yan, Jiayin Yang, Yujun Shi. HDLBP-stabilized lncFAL inhibits ferroptosis vulnerability by diminishing Trim69-dependent FSP1 degradation in hepatocellular carcinoma. *Redox Biol*. 2022 Dec;58:102546.
13. Shun Peng, Dongbei Guo, Yuan Guo, Heyu Zhao, Jun Mei, Yakun Han, Rui Guan, Tianhua Wang, Teng Song, Keke Sun, Yunhan Liu, Ting Mao, Huan Chang, Jingshi Xue, Yingfan Cai, Dong Chen, Shui Wang. CONSTITUTIVE EXPRESSION OF PATHOGENESIS-RELATED GENES 5 is an RNA-binding protein controlling plant immunity via an RNA processing complex. *Plant Cell*. 2022 Apr 26;34(5):1724-1744.
14. Ying Yang, Ziyu Wang, Juan Wang, Xiaoxia Dai, Changjun You. Next-Generation Sequencing-Based Analysis of the Effects of N1- and N6-Methyldeoxyadenosine Adducts on DNA Transcription. *Anal*

- Chem. 2022 Aug 16;94(32):11248-11254.
15. Shouyi Li, Xuetao Gong, Liwen Yin, Xiaolei Pan, Yongxin Jin, Fang Bai, Zhihui Cheng, Un-Hwan Ha, Weihui Wu. Acetylation of CspC Controls the Las Quorum-Sensing System through Translational Regulation of *rsaL* in *Pseudomonas aeruginosa*. *mBio*. 2022 Jun 28;13(3):e0054722.
 16. Zhaoyi Pan, Yujie Feng, Zhihui Wang, Zhengyang Lei, Qiuju Han, Jian Zhang. MERS-CoV nsp1 impairs the cellular metabolic processes by selectively downregulating mRNAs in a novel granules. *Virulence*. 2022 Dec;13(1):355-369.
 17. Yiting Wan, Cancan Gu, Jueying Kong, Jin Sui, Ling Zuo, Yanhua Song, Jing Chen. Long noncoding RNA ADAMTS9-AS1 represses ferroptosis of endometrial stromal cells by regulating the miR-6516-5p/GPX4 axis in endometriosis. *Sci Rep*. 2022 Feb 16;12(1):2618.
 18. Pengchao Feng, Zhijiao Xu, Jialin Chen, Meizhen Liu, Yu Zhao, Daqi Wang, Lei Han, Li Wang, Bo Wan, Xingshun Xu, Dali Li, Yilai Shu, Yimin Hua. Rescue of mis-splicing of a common SLC26A4 mutant associated with sensorineural hearing loss by antisense oligonucleotides. *Mol Ther Nucleic Acids*. 2022 Mar 19;28:280-292.
 19. Shuping Zhou, Shaoli Zhang, Hui Zhang, Junxia Ma, Huangzhen Dai, Lili Qu, Meixiang Zhou. Clinical Potential of lncRNA PPP1R26-AS1 in Breast Cancer and Its Contribution to Cancer Progression. *Mol Biotechnol*. 2022 Jun;64(6):660-669.
 20. Kahaerjiang Abuduwaili, Xiaodan Zhu, Yanli Shen, Suqiong Lu, Chunling Liu. *circ_0008797* attenuates non-small cell lung cancer proliferation, metastasis, and aerobic glycolysis by sponging miR-301a-3p/SOCS2. *Environ Toxicol*. 2022 Jul;37(7):1697-1710.
 21. Yanliang Wang, Yu Chen Feng, Yujin Gan, Liu Teng, Li Wang, Ting La, Peilin Wang, Yue Gu, Lei Yan, Na Li, Lina Zhang, Limeng Wang, Rick F Thorne, Xu Dong Zhang, Huixia Cao, Feng-Min Shao. lncRNA MILIP links YBX1 to translational activation of Snail and promotes metastasis in clear cell renal cell carcinoma. *J Exp Clin Cancer Res*. 2022 Aug 26;41(1):260.
 22. Xiaoyu Yue, Yinfei Zheng, Lianxin Li, Zhihao Yang, Zhigang Chen, Yu Wang, Zhiwei Wang, Deran Zhang, Erbao Bian, Bing Zhao. Integrative analysis of a novel 5 methylated snoRNA genes prognostic signature in patients with glioma. *Epigenomics*. 2022 Sep;14(18):1089-1104.
 23. Mi Zhang, Ling Wu, Xudong Wang, Jiang Chen. lncKRT16P6 promotes tongue squamous cell carcinoma progression by sponging miR-3180 and regulating GATAD2A expression. *Int J Oncol*. 2022 Sep;61(3):111.
 24. Mengyao Zhao, Yian Wang, Fenghua Tan, Lingyun Liu, Xiangchan Hou, Chunmei Fan, Le Tang, Yongzhen Mo, Yumin Wang, Qijia Yan, Zhaojian Gong, Zheng Li, Qianjin Liao, Can Guo, He Huang, Xi Zeng, Guiyuan Li, Zhaoyang Zeng, Wei Xiong, Fuyan Wang. Circular RNA circCCNB1 inhibits the migration and invasion of nasopharyngeal carcinoma through binding and stabilizing TJP1 mRNA. *Sci China Life Sci*. 2022 Nov;65(11):2233-2247.
 25. Yanzhi Lu, Min Long, Zhaowei Gao, Chong Liu, Ke Dong, Huizhong Zhang. Long non-coding RNA ENST00000469812 promotes Enterovirus type 71 replication via targeting the miR-4443/NUPR1 axis in rhabdomyosarcoma cells. *Arch Virol*. 2022 Dec;167(12):2601-2611.
 26. Qiao Sun, Tingting Zhao, Biao Li, Mengying Li, Ping Luo, Chen Zhang, Gang Chen, Zhengguo Cao, Yicun Li, Mingyuan Du, Hong He. FTO/RUNX2 signaling axis promotes cementoblast differentiation under normal and inflammatory condition. *Biochim Biophys Acta Mol Cell Res*. 2022 Dec;1869(12):119358.
 27. Yutian Zou, Shaoquan Zheng, Xinhua Xie, Feng Ye, Xiaoqian Hu, Zhi Tian, Shu-Mei Yan, Lu Yang, Yanan Kong, Yuhui Tang, Wenwen Tian, Jindong Xie, Xinpei Deng, Yan Zeng, Zhe-Sheng Chen, Hailin Tang, Xiaoming Xie. N6-methyladenosine regulated FGFR4 attenuates ferroptotic cell death in recalcitrant HER2-positive breast cancer. *Nat Commun*. 2022 May 13;13(1):2672.
 28. Ruirui Jia, Jiamei Lin, Jin You, Shi Li, Ge Shan, Chuan Huang. The DEAD-box helicase Hlc regulates basal transcription and chromatin opening of stress-responsive genes. *Nucleic Acids Res*. 2022 Sep 9;50(16):9175-9189.
 29. Zhenxing Song, Jiamei Lin, Rui Su, Yu Ji, Ruirui Jia, Shi Li, Ge Shan, Chuan Huang. eIF3j inhibits translation of a subset of circular RNAs in eukaryotic cells. *Nucleic Acids Res*. 2022 Nov 11;50(20):11529-11549.
 30. Chuan-Jian Shi, Min-Yi Lv, Li-Qiang Deng, Wei-Qiang Zeng, Wei-Ming Fu, Jin-Fang Zhang. Linc-ROR drive adriamycin resistance by targeting AP-2 α /Wnt/ β -catenin axis in hepatocellular carcinoma. *Cell Biol Toxicol*. 2023 Aug;39(4):1735-1752.
 31. Youjing Yang, Qianmin Li, Yi Ling, Linxin Leng, Yu Ma, Lian Xue, Guoyuan Lu, Yue Ding, Jianzhong Li, Shasha Tao. m6A eraser FTO modulates autophagy by targeting SQSTM1/P62 in the prevention of canagliflozin against renal fibrosis. *Front Immunol*. 2023 Jan 4;13:1094556.

Version 2024.10.30