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- Zhao H, Li J, Han B, Li X, Chen J. Improvement of oxidative stress tolerance in *Saccharomyces cerevisiae* through global transcription machinery engineering. *J IND MICROBIOL BIOT* 2014 May;41(5):869-78.
- Liu XW, Zi Y, Liu YE, Zhang YB, Xiang LB, Hou MX. Melatonin exerts protective effect on N2a cells under hypoxia conditions through Zip1/ERK pathway. *Neurosci Lett* 2015 May 19;595:74-80.
- Yao J, Zou Z, Wang X, Ji X, Yang J. Pinoresinol Diglucoside Alleviates oxLDL-Induced Dysfunction in Human Umbilical Vein Endothelial Cells. *EVID-BASED COMPL ALT* 2016;2016:3124519.
- Shang B, Shi H, Wang X, Guo X, Wang N, Wang Y, Dong L. Protective effect of melatonin on myenteric neuron damage in experimental colitis in rats. *FUND CLIN PHARMACOL* 2016 Apr;30(2):117-27.
- Kong F, Wang H, Guo J, Peng M, Ji H, Yang H, Liu B, Wang J, Zhang X, Li S. Hsp70 suppresses apoptosis of BRL cells by regulating the expression of Bcl-2, cytochrome C, and caspase 8/3. *IN VITRO CELL DEV-AN* 2016 May;52(5):568-75.
- Meng J, Lv Z, Qiao X, Li X, Li Y, Zhang Y, Chen C. The decay of Redox-stress Response Capacity is a substantive characteristic Redox Biol 2016 Dec 28;11:365-374.
- Zhu Z, Fan X, Pan Y, Lu Y, Zeng W. Trehalose improves rabbit sperm quality during cryopreservation. *Cryobiology* 2017 Apr;75:45-51.
- Meng J, Lv Z, Qiao X, Li X, Li Y, Zhang Y, Chen C. The decay of Redox-stress Response Capacity is a substantive characteristic of aging: Revising the redox theory of aging. *Redox Biol* 2017 Apr;11:365-374.
- He J, Jiang YF, Liang L, Wang DJ, Wei WX, Ji PP, Huang YC, Song H, Lu XL, Zhao YX. Targeting of AUF1 to vascular endothelial cells as a novel anti-aging therapy. *J Geriatr Cardiol* 2017 Aug;14(8):515-523.
- Li A, Li L, Song K, Wang W, Zhang G. Temperature, energy metabolism, and adaptive divergence in two oyster subspecies. *Ecol Evol* 2017 Jun 30;7(16):6151-6162.
- Chang W, Bai J, Tian S, Ma M, Li W, Yin Y, Deng R, Cui J, Li J, Wang G, Zhang P, Tao K. Autophagy protects gastric mucosal epithelial cells from ethanol-induced oxidative damage via mTOR signaling pathway. *EXP BIOL MED* 2017 May;242(10):1025-1033.
- Zeng X, Yang J, Hu O, Huang J, Ran L, Chen M, Zhang Y, Zhou X, Zhu J, Zhang Q, Yi L, Mi M. Dihydromyricetin Ameliorates Nonalcoholic Fatty Liver Disease by Improving Mitochondrial Respiratory Capacity and Redox Homeostasis Through Modulation of SIRT3 Signaling. *ANTIOXID REDOX SIGN* 2018 Feb 21.
- Yin X, Wang SL, Alolga RN, Mais E, Li P, Yang P, Komatsu S, Qi LW. Label-free proteomic analysis to characterize ginger from China and Ghana. *Food Chem* 2018 May 30;249:1-7.
- Ren M, Wang T, Huang L, Ye X, Han Z. Mesoporous Silica Nanoparticles Rescue H2O2-induced Inhibition of Cardiac Differentiation. *Cell Struct Funct* 2018;43(2):109-117.
- Lu QB, Wan MY, Wang PY, Zhang CX, Xu DY, Liao X, Sun HJ. Chicoric acid prevents PDGF-BB-induced VSMC dedifferentiation, proliferation and migration by suppressing ROS/NFκB/mTOR/P70S6K signaling cascade. *Redox Biol* 2018 Apr;14:656-668.
- Lei Q, Tan J, Yi S, Wu N, Wang Y, Wu H. Mitochondrial acid 5 activates the MAPK-ERK-yap signaling pathways to protect mouse microglial BV-2 cells against TNFα-induced apoptosis via increased Bnip3-related mitophagy. *Cell Mol Biol Lett* 2018 Apr 5;23:14.
- Wang XX, Zhang TY, Dao GH, Hu HY. Interaction between 1,2-benzisothiazol-3(2H)-one and microalgae: Growth inhibition and detoxification mechanism. *Aquat Toxicol* 2018 Dec;205:66-75.

18. Xu S, Zhao X, Liu F, Cao Y, Wang B, Wang X, Yin M, Wang Q, Feng X Crucial role of oxidative stress in bactericidal effect of parthenolide against *Xanthomonas oryzae* pv. *oryzae*. *Pest Manag Sci* 2018 Dec;74(12):2716-2723.
19. Zhang B, Zhang Z, Ji H, Shi H, Chen S, Yan D, Jiang X, Shi B Grape seed proanthocyanidin extract alleviates urethral dysfunction in diabetic rats through modulating the NO-cGMP pathway. *Exp Ther Med* 2018 Jan;15(1):1053-1061.
20. Chen Y, Wang B, Zhao H Thymoquinone reduces spinal cord injury by inhibiting inflammatory response, oxidative stress and apoptosis via PPAR- γ and PI3K/Akt pathways. *Exp Ther Med* 2018 Jun;15(6):4987-4994.
21. Wang J, Zhang Y Neuroprotective effect of berberine agonist against impairment of learning and memory skills in severe traumatic brain injury via Sirt1/p38 MAPK expression. *Mol Med Rep* 2018 May;17(5):6881-6886.
22. Liu S, Yang L, Mu S, Fu Q Epigallocatechin-3-Gallate Ameliorates Glucocorticoid-Induced Osteoporosis of Rats in Vivo and in Vitro. *Front Pharmacol* 2018 May 9;9:447.
23. Wang J, Nisar M, Huang C, Pan X, Lin D, Zheng G, Jin H, Chen D, Tian N, Huang Q, Duan Y, Yan Y, Wang K, Wu C, Hu J, Zhang X, Wang X Small molecule natural compound agonist of SIRT3 as a therapeutic target for the treatment of intervertebral disc degeneration. *Exp Mol Med* 2018 Nov 12;50(11):146.
24. Tian HP, Sun YH, He L, Yi YF, Gao X, Xu DL Single-Stranded DNA-Binding Protein 1 Abrogates Cardiac Fibroblast Proliferation and Collagen Expression Induced by Angiotensin II. *Int Heart J* 2018 Nov 28;59(6):1398-1408.
25. Wang XX, Zhang TY, Dao GH, Hu HY Tolerance and resistance characteristics of microalgae *Scenedesmus* sp. LX1 to methylisothiazolinone. *Environ Pollut* 2018 Oct;241:200-211.
26. Wang QL, Guo C, Qi J, Ma JH, Liu FY, Lin SQ, Zhang CY, Xie WD, Zhuang JJ, Li X Protective effects of 3 β -angeloyloxy-8 β , 10 β -dihydroxyeremophila-7(11)-en-12, 8 α -lactone on paraquat-induced oxidative injury in SH-SY5Y cells. *J Asian Nat Prod Res* 2019 Apr;21(4):364-376.
27. Quan XJ, Liang CL, Sun MZ, Zhang L, Li XL Overexpression of steroid receptor coactivators alleviates hyperglycemia-induced endothelial cell injury in rats through activating the PI3K/Akt pathway. *Acta Pharmacol Sin* 2019 May;40(5):648-657.
28. Xiao X, Cheng Y, Song D, Li X, Hu Y, Lu Z, Wang F, Wang Y Selenium-enriched *Bacillus paralicheniformis* SR14 attenuates H2O2-induced oxidative damage in porcine jejunum epithelial cells via the MAPK pathway. *APPL MICROBIOL BIOT* 103(15):6231-6243. 2019 Aug
29. Liu K, Jing MJ, Liu C, Yan DY, Ma Z, Wang C, Deng Y, Liu W, Xu B Effect of trehalose on manganese-induced mitochondrial dysfunction and neuronal cell damage in mice. *BASIC CLIN PHARMACOL* 125(6):536-547. 2019 Dec
30. Pan X, Song X, Wang C, Cheng T, Luan D, Xu K, Tang B H2Se Induces Reductive Stress in HepG2 Cells and Activates Cell Autophagy by Regulating the Redox of HMGB1 Protein under Hypoxia. *Theranostics* 9(6):1794-1808. 2019 Feb 28
31. Ding XM, Zhao LJ, Qiao HY, Wu SL, Wang XH Long non-coding RNA-p21 regulates MPP+ induced neuronal injury by targeting miR-625 and derepressing TRPM2 in SH-SY5Y cells. *CHEM-BIOL INTERACT* 307:73-81. 2019 Jul 1
32. Wu J, Pan W, Wang C, Dong H, Xing L, Hou J, Fang S, Li H, Yang F, Yu B H2S attenuates endoplasmic reticulum stress in hypoxia-induced pulmonary artery hypertension. *BIOSCIENCE REP* 39(7). pii: BSR20190304. 2019 Jul 8
33. Zhu Z, Li R, Lv Y, Zeng W Melatonin protects rabbit spermatozoa from cryo-damage via decreasing oxidative stress. *Cryobiology* 88:1-8. 2019 Jun
34. Zhu X, Wu S, Guo H Active Vitamin D and Vitamin D Receptor Help Prevent High Glucose Induced Oxidative Stress of Renal Tubular Cells via AKT/UCP2 Signaling Pathway. *Biomed Res Int* 2019:9013904. 2019 May 28
35. Chen B, Li Y, Liu Y, Xu Z circLRP6 regulates high glucose-induced proliferation, oxidative stress, ECM accumulation, and inflammation in mesangial cells. *J Cell Physiol* 234(11):21249-21259. 2019 Nov
36. Li Y, Zhou T, Su YF, Hu ZY, Wei JJ, Wang W, Liu CY, Zhao K, Zhang HP Prokineticin 2 overexpression induces spermatocyte apoptosis in varicocele in rats. *Asian J Androl* 2019 Nov 19
37. Xiong H, Wang C, Wang Z, Jiang Z, Zhou J, Yao J Intracellular cascade activated nanosystem for improving ER+ breast cancer therapy through attacking GSH-mediated metabolic vulnerability. *J Control Release* 309:145-157. 2019 Sep 10
38. Jie Zhang, Changling Ding, Shuping Zhang, Yangyang Xu Neuroprotective effects of astaxanthin against oxygen and glucose deprivation damage via the PI3K/Akt/GSK3 β /Nrf2 signalling pathway in vitro *J Cell Mol Med* 2020 Aug;24(16):8977-8985.
39. Hang Yu, Zhidian Wu, Xiaozhi Wang, Chang Gao, Run Liu, Fuxin Kang, Mingming Dai Protective effects of combined treatment with mild hypothermia and edaravone against cerebral ischemia/reperfusion injury via oxidative stress and Nrf2 pathway regulation *Int J Oncol* 2020 Aug;57(2):500-508.
40. Yangyang Chen, Jichun Tang, Yinghong Zhang, Juan Du, Yuanli Wang, Hui Yu, Yanling He Astaxanthin alleviates gestational diabetes mellitus in mice through suppression of oxidative stress *Naunyn Schmiedebergs Arch Pharmacol* 2020 Dec;393(12):2517-2527.
41. Xin Wang, Yanan Zhang, Shuai Han, Hongshen Chen, Chen Chen, Lingling Ji, Bihu Gao Overexpression of miR-30c-5p reduces cellular cytotoxicity and inhibits the formation of kidney stones through ATG5 *Int J Mol Med* 2020 Feb;45(2):375-384.
42. Hui Chen, Yao Wang, Fang-Zhou Jiao, Fan Yang, Xun Li, Lu-Wen Wang Sinomenine Attenuates Acetaminophen-Induced Acute Liver Injury by Decreasing Oxidative Stress and Inflammatory Response via Regulating TGF- β /Smad Pathway in vitro and in vivo *Drug Des Devel Ther* 2020 Jun 17;14:2393-2403.
43. Hao Zhang, Ming Gong, Xinle Luo Methoxytetrahydro-2H-pyran-2-yl)methyl benzoate inhibits spinal cord injury in the rat model via PPAR- γ /PI3K/p-Akt activation *Environ Toxicol* 2020 Jun;35(6):714-721.
44. Lufei Ouyang, Xiaoyu Mu, Junying Wang, Qifeng Li, Yalong Gao, Haile Liu, Si Sun, Qinjuan Ren, Ruijuan Yan, Jingya Wang, Qiang Liu, Yuanming Sun, Changlong Liu, Hua He, Wei Long, Xiao-Dong Zhang Carbon dot targeting to nitrogen signaling molecules for inhibiting neuronal death *J Mater Chem B* 2020 Mar 18;8(11):2321-2330.
45. Xiao-Xiong Wang, Tian-Yuan Zhang, Guo-Hua Dao, Zi-Bin Xu, Yin-Hu Wu, Hong-Ying Hu Assessment and mechanisms of microalgae growth inhibition by phosphonates: Effects of intrinsic toxicity and complexation *Water Res* 2020 Nov 1;186:116333.
46. Wei-Qi Li, Ting Qing, Cheng-Cheng Li, Feng Li, Fei Ge, Jun-Jie Fei, Willie J G M Peijnenburg Integration of subcellular partitioning and chemical forms to understand silver nanoparticles toxicity to lettuce (*Lactuca sativa* L.) under different exposure pathways *Chemosphere* 2020 Nov;258:127349.
47. Zi-Sheng Huang, Da-Qi Xie, Li-Jun Xu, Chang-Shun Huang, Min Zheng, Yi-Jun Chen, Yin Cao Tetramethylpyrazine Ameliorates

- Lipopolysaccharide-Induced Sepsis in Rats via Protecting Blood-Brain Barrier, Impairing Inflammation and Nitrous Oxide Systems *Front Pharmacol* 2020 Oct 2;11:562084.
48. Qiushi Xu, Yunhui Fan, Juan J Looor, Yusheng Liang, Xudong Sun, Hongdou Jia, Chenxu Zhao, Chuang Xu Adenosine 5'-monophosphate-activated protein kinase ameliorates bovine adipocyte oxidative stress by inducing antioxidant responses and autophagy *J Dairy Sci* 2021 Apr;104(4):4516-4528.
 49. Jianjun Wu, Fan Yang, Xin Zhang, Guanghua Chen, Jilong Zou, Li Yin, Dawei Yang Hydrogen sulfide inhibits endoplasmic reticulum stress through the GRP78/mTOR pathway in rat chondrocytes subjected to oxidative stress *Int J Mol Med* 2021 Apr;47(4):34.
 50. Li Ma, Chengxu Li, Shuai Lian, Bin Xu, Hongming Lv, Yanzhi Liu, Jingjing Lu, Hong Ji, Shize Li, Jingru Guo, Huanmin Yang Procyanidin B2 alleviates liver injury caused by cold stimulation through Sonic hedgehog signalling and autophagy *J Cell Mol Med* 2021 Aug;25(16):8015-8027.
 51. Xu Yang, Pengli Liu, Xuliang Zhang, Jian Zhang, Yilong Cui, Miao Song, Yanfei Li T-2 toxin causes dysfunction of Sertoli cells by inducing oxidative stress *Ecotoxicol Environ Saf* 2021 Dec 1;225:112702.
 52. Lei Yan, Lei Li, Jingan Lei Long noncoding RNA small nucleolar RNA host gene 12/microRNA-138-5p/nuclear factor I/B regulates neuronal apoptosis, inflammatory response, and oxidative stress in Parkinson's disease *Bioengineered* 2021 Dec;12(2):12867-12879.
 53. Yang Xie, Shujiang Zhang, Zhiyu Lv, Ting Long, Ying Luo, Zuoxiao Li SOX21-AS1 modulates neuronal injury of MMP+-treated SH-SY5Y cells via targeting miR-7-5p and inhibiting IRS2 *Neurosci Lett* 2021 Feb 16;746:135602.
 54. Zhenling Ma, Wenwen Zhang, Yaru Wu, Menghao Zhang, Lei Wang, Yihan Wang, Yi Wang, Wei Liu Cyclophilin A inhibits A549 cell oxidative stress and apoptosis by modulating the PI3K/Akt/mTOR signaling pathway *Biosci Rep* 2021 Jan 29;41(1):BSR20203219.
 55. Qian-Lan Xu, Jie Wu Effects of Txk-mediated activation of NF- κ B signaling pathway on neurological deficit and oxidative stress after ischemia-reperfusion in rats *Mol Med Rep* 2021 Jul;24(1):524.
 56. Xue-Jiao Sun, Wen-Qi Ma, Yi Zhu, Nai-Feng Liu POSTN promotes diabetic vascular calcification by interfering with autophagic flux *Cell Signal* 2021 Jul;83:109983.
 57. Qingyuan Jiang, Ming Lu, Jinyu Li, Zhongsheng Zhu Ginkgolide B Protects Cardiomyocytes from Angiotensin II-Induced Hypertrophy via Regulation of Autophagy through SIRT1-FoxO1 *Cardiovasc Ther* 2021 Jun 23;2021:5554569.
 58. Jing Wang, Yan Ding, Linwu Zhuang, Zhenzhong Wang, Wei Xiao, Jingbo Zhu Ginkgolide B-induced AMPK pathway activation protects astrocytes by regulating endoplasmic reticulum stress, oxidative stress and energy metabolism induced by A β 1-42 *Mol Med Rep* 2021 Jun;23(6):457.
 59. Xufeng Fu, Hang Han, Yuanyuan Li, Bo Xu, Wenjie Dai, Yaoxu Zhang, Feng Zhou, Huiming Ma, Xiuying Pei Di-(2-ethylhexyl) phthalate exposure induces female reproductive toxicity and alters the intestinal microbiota community structure and fecal metabolite profile in mice *Environ Toxicol* 2021 Jun;36(6):1226-1242.
 60. Jiao Guo, Haiyu Yang, Ya Liu, Wei Liu, Ruiying Zhao, He Li, Wei Long, Wenqing Xu, Meili Guo, Xiaodong Zhang Atomically precise silver clusterzymes protect mice from radiation damages *J Nanobiotechnology* 2021 Nov 19;19(1):377.
 61. Yuan Liu, Yanlun Qiu, Qi Yin, Xinglong Li, Qunhua Bai, Yingli Li, Hong Xiao iTRAQ-based quantitative proteomic reveals proteomic changes in *Serratia* sp. CM01 and mechanism of Cr(VI) resistance *Ecotoxicol Environ Saf* 2021 Nov 22;228:112899.
 62. Mei Li, Zhi Zheng Protective effect of parecoxib sodium against ischemia reperfusion-induced intestinal injury *Mol Med Rep* 2021 Nov;24(5):776.
 63. Yanhong Duo, Guanghong Luo, Zihuang Li, Zide Chen, Xianming Li, Zhenyou Jiang, Binlu Yu, Hao Huang, Zhengbo Sun, Xue-Feng Yu Photothermal and Enhanced Photocatalytic Therapies Conduce to Synergistic Anticancer Phototherapy with Biodegradable Titanium Diselenide Nanosheets *Small* 2021 Oct;17(40):e2103239.
 64. Yuqin Wu, Lin Fan, Yun Wang, Jing Ding, Rongfu Wang Isorhamnetin Alleviates High Glucose-Aggravated Inflammatory Response and Apoptosis in Oxygen-Glucose Deprivation and Reoxygenation-Induced HT22 Hippocampal Neurons Through Akt/SIRT1/Nrf2/HO-1 Signaling Inflammation 2021 Oct;44(5):1993-2005.
 65. Fang Wan, Gaolong Zhong, Shaofeng Wu, Xuanxuan Jiang, Jianzhao Liao, Xiaoyong Zhang, Hui Zhang, Khalid Mehmood, Zhaoxin Tang, Lianmei Hu Arsenic and antimony co-induced nephrotoxicity via autophagy and pyroptosis through ROS-mediated pathway in vivo and in vitro *Ecotoxicol Environ Saf* 2021 Sep 15;221:112442.
 66. Z Dong, L Bian, Y-L Wang, L-M Sun Gastrodin protects against high glucose-induced cardiomyocyte toxicity via GSK-3 β -mediated nuclear translocation of Nrf2 *Hum Exp Toxicol* 2021 Sep;40(9):1584-1597.
 67. Chen Zhang, Han Wang, Xinhui Yang, Zi Fu, Xiuru Ji, Yifan Shi, Jie Zhong, Weiguo Hu, Youqiong Ye, Zhengting Wang, Dalong Ni Oral zero-valent-molybdenum nanodots for inflammatory bowel disease therapy *Sci Adv* 2022 Sep 16;8(37):eabp9882.
 68. Ye Zhang, Yanze Li, Qiangmin Qiu, Zhiyuan Chen, Yang Du, Xiuheng Liu MITD1 Deficiency Suppresses Clear Cell Renal Cell Carcinoma Growth and Migration by Inducing Ferroptosis through the TAZ/SLC7A11 Pathway *Oxid Med Cell Longev* 2022 Aug 22:2022:7560569.
 69. Weilun Miao, Mengli Chen, Minglong Chen, Chang Cui, Yue Zhu, Xiping Luo, Bangwei Wu Nr2f2 Overexpression Aggravates Ferroptosis and Mitochondrial Dysfunction by Regulating the PGC-1 α Signaling in Diabetes-Induced Heart Failure Mice *Mediators Inflamm* 2022 Aug 30:2022:8373389.
 70. Ya Liu, He Li, Wei Liu, Jiao Guo, Haiyu Yang, Haikang Tang, Maoye Tian, Hongmei Nie, Xiaodong Zhang, Wei Long Design of Monovalent Cerium-Based Metal Organic Frameworks as Bioinspired Superoxide Dismutase Mimics for Ionizing Radiation Protection *ACS Appl Mater Interfaces* 2022 Dec 14;14(49):54587-54597.
 71. Jichen Zhang, Xuemei Zhao, Hongling Zhu, Jingnan Wang, Junhua Ma, Mingjun Gu Crocin protects the renal tubular epithelial cells against high glucose-induced injury and oxidative stress via regulation of the SIRT1/Nrf2 pathway *Iran J Basic Med Sci* 2022 Feb;25(2):193-197.
 72. Huan Wan, Yaping Wang, Qingyun Pan, Xia Chen, Sijun Chen, Xiaohui Li, Weiguo Yao Quercetin attenuates the proliferation, inflammation, and oxidative stress of high glucose-induced human mesangial cells by regulating the miR-485-5p/YAP1 pathway *Int J Immunopathol Pharmacol* 2022 Jan-Dec;36:20587384211066440.
 73. Hong-Jie Yang, Bin Kong, Wei Shuai, Jing-Jing Zhang, He Huang Shensong Yangxin attenuates metabolic syndrome-induced atrial fibrillation via inhibition of ferroportin-mediated intracellular iron overload *Phytomedicine* 2022 Jul;101:154086.
 74. Kaiming Yue, Xudong Pu, Juan J Looor, Qianming Jiang, Jihong Dong, Taiyu Shen, Guojin Li, Wenwen Gao, Lin Lei, Xiliang Du, Yuxiang Song, Guowen Liu, Xinwei Li Impaired autophagy aggravates oxidative stress in mammary gland of dairy cows with clinical ketosis *J Dairy Sci* 2022 Jul;105(7):6030-6040.

75. Dan Zhu, Tingting Fan, Yaohua Chen, Xingyue Huo, Yuping Li, Danyong Liu, Yin Cai, Chi Wai Cheung, Jing Tang, Jian Cui, Zhengyuan Xia CXCR4/CX43 Regulate Diabetic Neuropathic Pain via Intercellular Interactions between Activated Neurons and Dysfunctional Astrocytes during Late Phase of Diabetes in Rats and the Effects of Antioxidant Oxid Med Cell Longev 2022 Jun 28:2022:8547563.
76. Tonghui Li, Liuliu Shi, Wenqiang Liu, Xuhao Hu, Yuanjian Hui, Maojun Di, Shen Xue, Yan Zheng, Mengjuan Yao, Chen Li, Kun Meng Aloe-Emodin Induces Mitochondrial Dysfunction and Pyroptosis by Activation of the Caspase-9/3/Gasdermin E Axis in HeLa Cells Front Pharmacol 2022 May 17:13:854526.

Version2024.03.12