



碧云天网站



微信公众号


碧云天生物技术/Beyotime Biotechnology
订货热线: 400-1683301 或 800-8283301
订货 e-mail: order@beyotime.com
技术咨询: info@beyotime.com
网址: http://www.beyotime.com

Fluo-3 AM (钙离子荧光探针, 5mM)

产品编号	产品名称	包装
S1056	Fluo-3 AM (钙离子荧光探针, 5mM)	20微升

产品简介:

- Fluo-3 AM是最常用的检测细胞内钙离子浓度的荧光探针之一。分子式为 $C_{51}H_{50}Cl_2N_2O_{23}$ ，分子量为1129.85。
- Fluo-3和Fura-2相比，其优点是：一方面可以被氩离子激光(argon-ion laser)的488nm激发光激发，便于检测；另一方面，Fluo-3和钙离子结合后荧光变化更强，即对钙离子浓度变化的检测更加灵敏；同时，Fluo-3和钙离子的结合能力较弱，这样可以比Fura-2检测到细胞内更高浓度的钙离子水平；此外，对于细胞内的钙离子的即时变化反应得更加准确，减小了因为和钙离子解离速度慢而导致的荧光变化滞后。
- 本Fluo-3 AM (钙离子荧光探针) 是配制于无水DMSO (anhydrous DMSO)中的储存液，浓度为5mM。
- Fluo-3 AM是一种可以穿透细胞膜的荧光染料。Fluo-3 AM的荧光非常弱，其荧光不会随钙离子浓度升高而增强。Fluo-3 AM进入细胞后可以被细胞内的酯酶剪切形成Fluo-3，从而被滞留在细胞内。Fluo-3可以和钙离子结合，结合钙离子后可以产生较强的荧光，最大激发波长为506nm，最大发射波长为526nm。实际检测时推荐使用的激发波长为488nm左右，发射波长为525-530nm。Fluo-3的激发光谱和发射光谱参考下图。

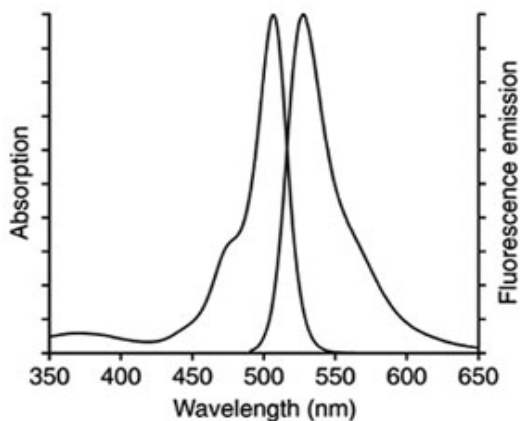

- 用于细胞内钙离子检测时，Fluo-3 AM的常用浓度为0.5-5 μ M。通常用含有0.5-5 μ M的Fluo-3 AM的适当溶液和细胞一起在20-37°C孵育15-60分钟进行荧光探针装载，随后适当洗涤，洗涤后可以适当再孵育20-30分钟以确保Fluo-3 AM在细胞内完全转变成Fluo-3。

包装清单:

产品编号	产品名称	包装
S1056	Fluo-3 AM (钙离子荧光探针, 5mM)	20微升
—	说明书	1份

保存条件:

-20°C避光保存，6个月有效。

注意事项:

- 本Fluo-3 AM在4°C、冰浴等较低温度情况下会凝固而粘在离心管管底、管壁或管盖内，可以20-25°C水浴温育片刻至全部融解后使用。
- 荧光染料均存在淬灭问题，请尽量注意避光，以减缓荧光淬灭。
- 本产品仅限于专业人员的科学研究用，不得用于临床诊断或治疗，不得用于食品或药品，不得存放于普通住宅内。
- 为了您的安全和健康，请穿实验服并戴一次性手套操作。

使用本产品的文献:

1. Kuang CY, Yu Y, Guo RW, Qian DH, Wang K, Den MY, Shi YK, Huang L. Silencing stromal interaction molecule 1 by RNA interference inhibits the proliferation and migration of endothelial progenitor cells. BIOCHEM BIOPH RES CO 2010;398(2):315-20.
2. Shi Y, Song M, Guo R, Wang H, Gao P, Shi W, Huang L. Knockdown of stromal interaction molecule 1 attenuates hepatocyte growth

- factor-induced endothelial progenitor cell proliferation. *EXP BIOL MED* 2010 Mar;235(3):317-25.
3. Jin H, Zhong X, Wang Z, Huang X, Ye H, Ma S, Chen Y, Cai J. Sonodynamic effects of hematoporphyrin monomethyl ether on CNE-2 cells detected by atomic force microscopy. *J Cell Biochem* 2011 Jan;112(1):169-78.
4. Li X, Ye X, Li X, Sun X, Liang Q, Tao L, Kang X, Chen J. Salidroside protects against MPP⁺-induced apoptosis in PC12 cells by inhibiting the NO pathway. *Brain Res* 2011 Mar 25;1382:9-18.
5. Wang M, Ruan Y, Chen Q, Li S, Wang Q, Cai J. Curcumin induced HepG2 cell apoptosis-associated mitochondrial membrane potential and intracellular free Ca²⁺ concentration. *Eur J Pharmacol* 2011 Jan 10;650(1):41-7.
6. Yang F, Zhang M, He W, Chen P, Cai X, Yang L, Gu N, Wu J. Controlled release of Fe₃O₄ nanoparticles in encapsulated microbubbles to tumor cells via sonoporation and associated cellular bioeffects. *Small* 2011 Apr 4;7(7):902-10.
7. Zhang J, Song W, Guo J, Zhang J, Sun Z, Ding F, Gao M. Toxic effect of different ZnO particles on mouse alveolar macrophages. *J Hazard Mater* 2012 Jun 15;219-220:148-55.
8. Li SP, Xie WL, Cai HH, Cai JY, Yang PH. Hydroxyl radical scavenging mechanism of human erythrocytes by quercetin-germanium (IV) complex. *Eur J Pharm Sci* 2012 Aug 30;47(1):28-34.
9. Jin H, Yang P, Cai J, Wang J, Liu M. Photothermal effects of folate-conjugated Au nanorods on HepG2 cells. *APPL MICROBIOL BIOT* 2012 Jun;94(5):1199-208.
10. Zhang HY, Wang WX, Yin H, Zhao XM, Du YG. Oligochitosan induces programmed cell death in tobacco suspension cells. *Carbohydrate Polymers* 2012 Feb;87(3):2270-8.
11. Li XY, Jing CQ, Zang XY, Yang S, Wang JJ. Toxic cytological alteration and mitochondrial dysfunction in PC12 cells induced by 1-octyl-3-methylimidazolium chloride. *Toxicol In Vitro* 2012 Oct;26(7):1087-92.
12. Che T, Sun H, Li J, Yu X, Zhu D, Xue B, Liu K, Zhang M, Kunze W, Liu C. Oxytocin hyperpolarizes cultured duodenum myenteric intrinsic primary afferent neurons by opening BK(Ca) channels through IP₃ pathway. *J Neurochem* 2012 May;121(4):516-25.
13. Zhang H, Shao D, Wu Y, Cai C, Hu C, Shou X, Dai B, Ye B, Wang M, Jia X. Apoptotic responses of *Carassius auratus* lymphocytes to nodularin exposure in vitro. *FISH SHELLFISH IMMUN* 2012 Dec;33(6):1229-37.
14. Lu N, Zhan M, Gao C, Wu G, Zhang H. I₄, a new synthetic sulfonylurea compound, inhibits the action of TXA₂ in vivo and in vitro on platelets and aorta vascular smooth muscle. *Thromb Res* 2012 Oct;130(4):e209-15.
15. Ouyang JS, Li YP, Li CY, Cai C, Chen CS, Chen SX, Chen YF, Yang L, Xie YP. Mitochondrial ROS-K⁺ channel signaling pathway regulated secretion of human pulmonary artery endothelial cells. *FREE RADICAL RES* 2012 Dec;46(12):1437-45.
16. Tan Q, Li Y, Wu J, Mei H, Zhao C, Zhang J. An optimized molecular inclusion complex of diferuloylmethane: enhanced physical properties and biological activity. *INT J NANOMED* 2012;5:5385-93.
17. Xu M, Lu N, Zhang H, Dai Q, Wei L, Li Z, You Q, Guo Q. Wogonin induced cytotoxicity in human hepatocellular carcinoma cells by activation of unfolded protein response and inactivation of AKT. *Hepatol Res* 2013 Aug;43(8):890-905.
18. Liang S, Su WW, Wang YG, Peng W, Nie YC, Li PB. Effect of quercetin 7-rhamnoside on glycochenodeoxycholic acid-induced L-02 human normal liver cell apoptosis. *Int J Mol Med* 2013 Aug;32(2):323-30.
19. Li Z, Pan X, Wang T, Wang PN, Chen JY, Mi L. Comparison of the killing effects between nitrogen-doped and pure TiO₂ on HeLa cells with visible light irradiation. *Nanoscale Res Lett* 2013 Feb 22;8(1):96.
20. Zhu K, Guo X, Guo X, Peng W, Zhou H. Protective effects of wheat germ protein isolate hydrolysates (WGPIH) against hydrogen peroxide-induced oxidative stress in PC12 cells. *Food Research International* 2013 Aug;53(1):297-303.
21. Chen G, Zhang P, Huang T, Yu W, Lin J, Li P, Chen K. Polysaccharides from *Rhizopus nigricans* mycelia induced apoptosis and G₂/M arrest in BGC-823 cells. *Carbohydrate Polymers* 2013 Sep 12;97(2):800-8.
22. Jiang L, Lan T, Chen Y, Sang J, Li Y, Wu M, Tao Y, Wang Y, Qian H, Gu L. PKG II inhibits EGF/EGFR-induced migration of gastric cancer cells. *PLoS One* 2013 Apr 16;8(4):e61674.
23. Yang D, Lü X, Hong Y, Xi T, Zhang D. The molecular mechanism of mediation of adsorbed serum proteins to endothelial cells adhesion and growth on biomaterials. *Biomaterials* 2013 Jul;34(23):5747-58.
24. Xin J, Zhan Y, Liu M, Hu H, Xia L, Nie Y, Wu K, Liang J, Tian J. ApoG₂ induces ER stress-dependent apoptosis in gastric cancer cells in vitro and its real-time evaluation by bioluminescence imaging in vivo. *Cancer Lett* 2013 Aug 19;336(2):260-9.
25. Chen Y, Chen JW, Zhai JH, Wang Y, Wang SL, Li X. Antitumor activity and toxicity relationship of annonaceous acetogenins. *Food Chem Toxicol* 2013 May 24;58C:394-400.
26. Pi J, Jin H, Liu R, Song B, Wu Q, Liu L, Jiang J, Yang F, Cai H, Cai J. Pathway of cytotoxicity induced by folic acid modified selenium nanoparticles in MCF-7 cells. *APPL MICROBIOL BIOT* 2013 Feb;97(3):1051-62.
27. Ma G, Zhou J, Tian C, Jiang D, Fang D, Chen H. Luminol electrochemiluminescence for the analysis of active cholesterol at the plasma membrane in single mammalian cells. *Anal Chem* 2013 Apr 16;85(8):3912-7.
28. Tan Q, Wu J, Li Y, Mei H, Zhao C, Zhang J. A supermolecular curcumin for enhanced antiproliferative and proapoptotic activities: molecular characteristics, computer modeling and in vivo pharmacokinetics. *Nanotechnology* 2013 Jan 25;24(3):035102.
29. Chen Z, Pan X, Georgakilas AG, Chen P, Hu H, Yang Y, Tian S, Xia L, Zhang J, Cai X, Ge J, Yu K, Zhuang J. Tetramethylpyrazine (TMP) protects cerebral neurocytes and inhibits glioma by down regulating chemokine receptor CXCR4 expression. *Cancer Lett* 2013 Aug 19;336(2):281-9.
30. Zhu XL, Qin XQ, Xiang Y, Tan YR, Qu XP, Liu HJ. Adipokine adiponectin is a potential protector to human bronchial epithelial cell for regulating proliferation, wound repair and apoptosis: comparison with leptin and resistin. *Peptides* 2013 Feb;40:34-41.
31. Tong J, Zhou XD, Perelman JM, Kolosov VP. Effect of epithelium ATP release on cyclic pressure-induced airway mucus secretion. *BIOSCIENCE REP* 2014 Feb 1;34(1). pii: e00088.
32. Xu SY, Hu YF, Li WP, Wu YM, Ji Z, Wang SN, Li K, Pan SY. Intermittent hypothermia is neuroprotective in an in vitro model of ischemic stroke. *Int J Biol Sci* 2014 Jul 29;10(8):873-81.
33. Guo Y, Wang W, Dong Y, Zhang Z, Zhou Y, Chen G. Carbon disulfide induces rat testicular injury via mitochondrial apoptotic pathway. *Chemosphere* 2014 Aug;108:367-75.
34. Tian J, Cheng J, Zhang J, Ye L, Zhang F, Dong Q, Wang H, Fu F. Protection of pyruvate against glutamate excitotoxicity is mediated by regulating DAPK1 protein complex. *PLoS One* 2014 Apr 22;9(4):e95777.
35. Li S, Zhang S, Jin X, Tan X, Lou J, Zhang X, Zhao Y. Singly protonated dehydronorcantharidin silver coordination polymer induces apoptosis of lung cancer cells via reactive oxygen species-mediated mitochondrial pathway. *Eur J Med Chem* 2014 Oct 30;86:1-11.
36. Wang Y, Yuan L, Yao C, Ding L, Li C, Fang J, Sui K, Liu Y, Wu M. A combined toxicity study of zinc oxide nanoparticles and vitamin C in food additives. *Nanoscale* 2014 Dec 21;6(24):15333-42.
37. Ding Y, Dai X, Zhang Z, Jiang Y, Ma X, Cai X, Li Y. Proanthocyanidins protect against early diabetic peripheral neuropathy by modulating endoplasmic reticulum stress. *J Nutr Biochem* 2014 Jul;25(7):765-72.
38. Duan Y, Wang Z, Zhang H, He Y, Fan R, Cheng Y, Sun G, Sun X. Extremely low frequency electromagnetic field exposure causes cognitive impairment associated with alteration of the glutamate level, MAPK pathway activation and decreased CREB phosphorylation in mice hippocampus: reversal by procyanidins extracted from the lotus seed pod. *Food Funct* 2014 Sep;5(9):2289-97.

39. Zhao Y, Zhou J, Xing D. Phytochrome B-mediated activation of lipoxygenase modulates an excess red light-induced defence response in Arabidopsis. *J Exp Bot* 2014 Sep;65(17):4907-18.
40. Yang Y, Li D, Katirai F, Zhang B, Xu Y, Xiong P, Gong F, Zheng F. Basophil activation through ASGM1 stimulation triggers PAF release and anaphylaxis-like shock in mice. *Eur J Immunol* 2014 Aug;44(8):2468-77.
41. Zhang S, Nie S, Huang D, Huang J, Feng Y, Xie M. Ganoderma atrum polysaccharide evokes antitumor activity via cAMP-PKA mediated apoptotic pathway and down-regulation of Ca(2+)/PKC signal pathway. *Food Chem Toxicol* 2014 Jun;68:239-46.
42. Li M, Wang N, Gong HQ, Li WZ, Liao XH, Yang XL, He HP, Cao DS, Zhang TC. Ca²⁺ signal-induced cardiomyocyte hypertrophy through activation of myocardin. *Gene* 2015 Feb 15;557(1):43-51.
43. Zeng Z, Leng T, Feng X, Sun H, Inoue K, Zhu L, Xiong ZG. Silencing TRPM7 in mouse cortical astrocytes impairs cell proliferation and migration via ERK and JNK signaling pathways. *PLoS One* 2015 Mar 23;10(3):e0119912.
44. Liu Z, Zhou Z, Wang L, Jiang S, Wang W, Zhang R, Song L. The immunomodulation mediated by a delta-opioid receptor for [Met(5)]-enkephalin in oyster *Crassostrea gigas*. *Dev Comp Immunol* 2015 Apr;49(2):217-24.
45. Hu XL, Guo LP, Song Q, Zhang Q, Chen Y, Wang J, Meng WH, Zhao QC. Kukoamine B, an amide alkaloid, protects against NMDA-induced neurotoxicity and potential mechanisms in vitro. *Neurochem Int* 2015 Aug;87:66-76.
46. Zhang J, Song J, Wu D, Wang J, Dong W. Hesperetin induces the apoptosis of hepatocellular carcinoma cells via mitochondrial pathway mediated by the increased intracellular reactive oxygen species, ATP and calcium. *Med Oncol* 2015 Apr;32(4):101.
47. Wang J, Ming H, Chen R, Ju JM, Peng WD, Zhang GX, Liu CF. CIH-induced neurocognitive impairments are associated with hippocampal Ca(2+) overload, apoptosis, and dephosphorylation of ERK1/2 and CREB that are mediated by overactivation of NMDARs. *Brain Res* 2015 Nov 2;1625:64-72.
48. Wang B, Li C, Huai R, Qu Z. Overexpression of ANO1/TMEM16A, an arterial Ca²⁺-activated Cl⁻ channel, contributes to spontaneous hypertension. *J Mol Cell Cardiol* 2015 May;82:22-32.
49. Hu XL, Gao LY, Niu YX, Tian X, Wang J, Meng WH, Zhang Q, Cui C, Han L, Zhao QC. Neuroprotection by Kukoamine A against oxidative stress may involve N-methyl-D-aspartate receptors. *BBA-BIOMEMBRANES* 2015 Feb;1850(2):287-98.
50. Li Z, Zhang Y, Peng D, Wang X, Peng Y, He X, Zhang X, Ma X, Huang L, Yan Y. Polyamine regulates tolerance to water stress in leaves of white clover associated with antioxidant defense and dehydrin genes via involvement in calcium messenger system and hydrogen peroxide signaling. *Front Physiol* 2015 Oct 12;6:280.
51. Liu Y, Wang XC, Hu D, Huang SR, Li QS, Li Z, Qu Y. Heat shock protein 70 protects PC12 cells against ischemia-hypoxia/reoxygenation by maintaining intracellular Ca(2+) homeostasis. *Neural Regen Res* 2016 Jul;11(7):1134-40.
52. Zhang J, Wang Y, Sun KM, Fang K, Tang X. A study of oxidative stress induced by two polybrominated diphenyl ethers in the rotifer *Brachionus plicatilis*. *Mar Pollut Bull* 2016 Dec 15;113(1-2):408-413.
53. Cai Q, Ma T, Li C, Tian Y, Li H. Catalpol Protects Pre-Myelinating Oligodendrocytes against Ischemia-induced Oxidative Injury through ERK1/2 Signaling Pathway. *Int J Biol Sci* 2016 Oct 26;12(12):1415-1426.
54. Chen C, Huang L, Zhang G, Li Y, Li L, Bai X, Liu W, Wang H, Li J. STK33 potentiates the malignancy of hypopharyngeal squamous carcinoma: Possible relation to calcium. *Cancer Biol Ther* 2016 Sep;17(9):976-84.
55. Zhou X, Wei Y, Qiu S, Xu Y, Zhang T, Zhang S. Propofol Decreases Endoplasmic Reticulum Stress-Mediated Apoptosis in Retinal Pigment Epithelial Cells. *PLoS One* 2016 Jun 16;11(6):e0157590.
56. Gong H, Liu L, Ni CX, Zhang Y, Su WJ, Lian YJ, Peng W, Zhang JP, Jiang CL. Dexamethasone rapidly inhibits glucose uptake via non-genomic mechanisms in contracting myotubes. *Arch Biochem Biophys* 2016 Aug 1;603:102-9.
57. Zhao F, Fu L, Yang W, Dong Y, Yang J, Sun S, Hou Y. Cardioprotective effects of baicalein on heart failure via modulation of Ca(2+) handling proteins in vivo and in vitro. *Life Sci* 2016 Jan 15;145:213-23.
58. Tao K, Wang B, Feng D, Zhang W, Lu F, Lai J, Huang L, Nie T, Yang Q. Salidroside Protects Against 6-Hydroxydopamine-Induced Cytotoxicity by Attenuating ER Stress. *Neurosci Bull* 2016 Feb;32(1):61-9.
59. Liu Z, Gan L, Wu T, Feng F, Luo D, Gu H, Liu S, Sun C. Adiponectin reduces ER stress-induced apoptosis through PPARα transcriptional regulation of ATF2 in mouse adipose. *Cell Death Dis* 2016 Nov 24;7(11):e2487.
60. Hu L, Sun J, Li H, Wang L, Wei Y, Wang Y, Zhu Y, Huo H, Tan Y. Differential mechanistic investigation of protective effects from imperatorin and sec-O-glucosylhamaudol against arsenic trioxide-induced cytotoxicity in vitro. *Toxicol In Vitro* 2016 Dec;37:97-105.
61. Jiang Q, Liu Y, Duan D, Gou M, Wang H, Wang J, Li Q, Xiao R. Anti-angiogenic activities of CRBGP from buccal glands of lampreys (*Lampetra japonica*). *Biochimie* 2016 Apr;123:7-19.
62. Yang L, Sui W, Li Y, Qi X, Wang Y, Zhou Q, Gao H. Substance P Inhibits Hyperosmotic Stress-Induced Apoptosis in Corneal Epithelial Cells through the Mechanism of Akt Activation and Reactive Oxygen Species Scavenging via the Neurokinin-1 Receptor. *PLoS One* 2016 Feb 22;11(2):e0149865.
63. Wu M, Wu D, Wang C, Guo Z, Li B, Zhuo Z. Hexabromocyclododecane exposure induces cardiac hypertrophy and arrhythmia by inhibiting miR-1 expression via up-regulation of the homeobox gene Nkx2.5. *J Hazard Mater* 2016 Jan 25;302:304-13.
64. Wang MD, Wu H, Huang S, Zhang HL, Qin CJ, Zhao LH, Fu GB, Zhou X, Wang XM, Tang L, Wen W, Yang W, Tang SH, Cao D, Guo LN, Zeng M, Wu MC, Yan HX, Wang HY. HBx regulates fatty acid oxidation to promote hepatocellular carcinoma survival during metabolic stress. *ONCOTARGET* 2016 Feb 9;7(6):6711-26.
65. Li M, He HP, Gong HQ, Zhang J, Ma WJ, Zhou H, Cao DS, Wang N, Zhang TC. NFATc4 and myocardin synergistically up-regulate the expression of LTCC α1C in ET-1-induced cardiomyocyte hypertrophy. *Life Sci* 2016 Jun 15;155:11-20.
66. Liu Z, Zhou Z, Wang L, Qiu L, Zhang H, Wang H, Song L. CgA1AR-1 acts as an alpha-1 adrenergic receptor in oyster *Crassostrea gigas* mediating both cellular and humoral immune response. *FISH SHELLFISH IMMUN* 2016 Nov;58:50-58.
67. Zhang C, Lu T, Wang GD, Ma C, Zhou YF. Costunolide, an active sesquiterpene lactone, induced apoptosis via ROS-mediated Biomed Pharmacother 2016 May;80:253-9.
68. Chen X, Sun FJ, Wei YJ, Wang LK, Zang ZL, Chen B, Li S, Liu SY, Yang H. Increased Expression of Transient Receptor Potential Vanilloid 4 in Cortical Lesions of Patients with Focal Cortical Dysplasia. *CNS Neurosci Ther* 2016 Apr;22(4):280-90.
69. Wang Y, Wang D, Zhang L, Ye F, Li M, Wen K. Role of JAK-STAT pathway in reducing cardiomyocytes hypoxia/reoxygenation injury induced by S1P postconditioning. *Eur J Pharmacol* 2016 Aug 5;784:129-36.
70. Chang R, Zhou R, Qi X, Wang J, Wu F, Yang W, Zhang W, Sun T, Li Y, Yu J. Protective effects of aloe on oxygen and glucose deprivation-induced injury in PC12 cells. *Brain Res Bull* 2016 Mar;121:75-83.
71. Lv S, Xu QY, Sun EC, Zhang JK, Wu DL. Dissection and integration of the autophagy signaling network initiated by bluetongue virus. *SCI REP-UK* 2016 Mar 15;6:23130.
72. Liu Z, Zhou Z, Wang L, Dong W, Qiu L, Song L. The cholinergic immune regulation mediated by a novel muscarinic acetylcholine receptor through TNF pathway in oyster *Crassostrea gigas*. *Dev Comp Immunol* 2016 Dec;65:139-48.
73. He C, Jiang S, Jin H, Chen S, Lin G, Yao H, Wang X, Mi P, Ji Z, Lin Y, Lin Z, Liu G. Mitochondrial electron transport chain identified as a novel molecular target of SPIO nanoparticles mediated cancer-specific cytotoxicity. *Biomaterials* 2016 Mar;83:102-14.
74. Meng X, Li X, Wang X, Gao J, Yang H, Chen H. Potential allergenicity response to structural modification of irradiated bovine α-lactalbumin. *Food Funct* 2016

- Jul 13;7(7):3102-10.
75. Liu M, Feng LX, Sun P, Liu W, Wu WY, Jiang BH, Yang M, Hu LH, Guo DA, Liu X. A Novel Bufalin Derivative Exhibited Stronger Apoptosis-Inducing Effect than Bufalin in A549 Lung Cancer Cells and Lower Acute Toxicity in Mice. *PLoS One* 2016 Jul 26;11(7):e0159789.
76. Liu Y, Zhi D, Li M, Liu D, Wang X, Wu Z, Zhang Z, Fei D, Li Y, Zhu H, Xie Q, Yang H, Li H. Shengmai Formula suppressed over-activated Ras/MAPK pathway in *C. elegans* by opening mitochondrial permeability transition pore via regulating cyclophilin D. *SCI REP-UK* 2016 Dec 16;6:38934.
77. Guo L, Lv G, Qiu L, Yang H, Zhang L, Yu H, Zou M, Lin J. Insights into anticancer activity and mechanism of action of a ruthenium(II) complex in human esophageal squamous carcinoma EC109 cells. *Eur J Pharmacol* 2016 Sep 5;786:60-71.
78. Cui X, Liang T, Liu C, Yuan Y, Qian J. Correlation of particle properties with cytotoxicity and cellular uptake of hydroxyapatite nanoparticles in human gastric cancer cells. *MAT SCI ENG C-MATER* 2016 Oct 1;67:453-60.
79. He B, Yuan L, Dai W, Gao W, Zhang H, Wang X, Fang W, Zhang Q. Dynamic bio-adhesion of polymer nanoparticles on MDCK epithelial cells and its impact on Nanoscale 2016 Mar 21;8(11):6129-45.
80. Zhao M, Li H, Liu X, Wei J, Ji J, Yang S, Hu Z, Wei S. Response of Human Osteoblast to n-HA/PEEK-Quantitative Proteomic Study of Bio-effects of Nano-Hydroxyapatite Composite. *SCI REP-UK* 2016 Mar 9;6:22832.
81. Sun T, Guo Z, Liu CJ, Li MR, Li TP, Wang X, Yuan DJ. Preservation of CGRP in myocardium attenuates development of cardiac dysfunction in diabetic rats. *Int J Cardiol* 2016 Oct 1;220:226-34.
82. Du C, Li Y, Guo Y, Han M, Zhang W, Qian H. The suppression of torulene and torularhodin treatment on the growth of PC-3 xenograft prostate tumors. *BIOCHEM BIOPH RES CO* 2016 Jan 22;469(4):1146-52.
83. Liu Q, Ji X, Ge Z, Diao H, Chang X, Wang L, Wu Q. Role of connexin 43 in cadmium-induced proliferation of human prostate epithelial cells. *J Appl Toxicol* 2017 Feb 8. doi: 10.1002/jat.3441.
84. Wang C, Qi S, Liu C, Yang A, Fu W, Quan C, Duan P, Yu T, Yang K. Mitochondrial Dysfunction and Ca²⁺ Overload in Injured Sertoli Cells Exposed to Bisphenol A. *Environ Toxicol* 2017 Mar;32(3):823-831.
85. Liu Z, Gu H, Gan L, Xu Y, Feng F, Saeed M, Sun C. Reducing Smad3/ATF4 was essential for Sirt1 inhibiting ER stress-induced apoptosis in mice brown adipose tissue. *ONCOTARGET* 2017 Feb 7;8(6):9267-9279.
86. Lin H, Ma X, Wang BC, Zhao L, Liu JX, Pu FF, Hu YQ, Hu HZ, Shao ZW. Edaravone ameliorates compression-induced damage in rat nucleus pulposus cells. *Life Sci* 2017 Nov 15;189:76-83.
87. Chen Y, Huang L, Zhang H, Diao X, Zhao S, Zhou W. Reduction in Autophagy by (-)-Epigallocatechin-3-Gallate (EGCG): a Potential Mechanism of Prevention of Mitochondrial Dysfunction After Subarachnoid Hemorrhage. *Mol Neurobiol* 2017 Jan;54(1):392-405.
88. Jianfei L, Min W, Chunlai M, Bicui C, Jiming Z, Bin W. The Ca²⁺/CaMKK2 axis mediates the telbivudine induced upregulation of creatine kinase: Implications for mechanism of antiviral nucleoside analogs' side effect. *Biochem Pharmacol* 2017 Dec 15;146:224-232.
89. Xu Z, Lu Y, Wang J, Ding X, Chen J, Miao C. The protective effect of propofol against TNF- α -induced apoptosis was mediated via inhibiting iNOS/NO production and maintaining intracellular Ca²⁺ homeostasis in mouse hippocampal HT22 cells. *Biomed Pharmacother* 2017 Jul;91:664-672.
90. Zhu G, Tao L, Wang R, Xue Y, Wang X, Yang S, Sun X, Gao G, Mao Z, Yang Q. Endoplasmic reticulum stress mediates distinct impacts of sevoflurane on different subfields of immature hippocampus. *J Neurochem* 2017 Jul;142(2):272-285.
91. Dai J, Wu S, Kong Y, Chi Z, Si L, Sheng X, Cui C, Fang J, Zhang J, Guo J. Nanosecond Pulsed Electric Fields Enhance the Anti-tumour Effects of the mTOR Inhibitor Everolimus against Melanoma. *SCI REP-UK* 2017 Jan 5;7:39597. doi: 10.1038/srep39597.
92. Wu XT, Sun LW, Yang X, Ding D, Han D, Fan YB. The potential role of spectrin network in the mechanotransduction of MLO-Y4 osteocytes. *SCI REP-UK* 2017 Jan 23;7:40940.
93. Zhang X, Xu L, Tian H, Wang C, Wang W, Ru S. Monocrotophos, an organophosphorus insecticide, disrupts the expression of HpNetrin and its receptor neogenin during early development in the sea urchin (*Hemicentrotus pulcherrimus*). *Neurotoxicology* 2017 Sep;62:130-137.
94. Wang X, Guo Q, Tao L, Zhao L, Chen Y, An T, Chen Z, Fu R. E platinum, a newly synthesized platinum compound, induces apoptosis through ROS-triggered ER stress in gastric carcinoma cells. *MOL CARCINOGEN* 2017 Jan;56(1):218-231.
95. Li N, Xu P, Jing WX, Hwang JS, Wang L. Toxic effects of Pb²⁺ entering sperm through Ca²⁺ channels in the freshwater crab *Sinopotamon henanense*. *Aquat Toxicol* 2017 Nov;192:24-29.
96. Qiu Y, Yu Q, Shi K, Zhang M, Zhou X, Yang Y, Liu Y, Tang J, Wang X, He Q. Cell-penetrating peptides induce apoptosis and necrosis through specific mechanism and cause impairment of Na⁺-K⁺-ATPase and mitochondria. *Amino Acids* 2017 Jan;49(1):75-88.
97. Li C, Li L, Lin B, Fang Y, Yang H, Liu H, Xi Z. Tris (1,3-dichloro-2-propyl) phosphate induces toxicity by stimulating CaMK2 in PC12 cells. *Environ Toxicol* 2017 Jun;32(6):1784-1791.
98. Hu Y, Zhang C, Li S, Jiao Y, Qi T, Wei G, Han G. Effects of Photodynamic Therapy Using Yellow LED-light with Concomitant Hypocrellin B on Apoptotic Signaling in Keloid Fibroblasts. *Int J Biol Sci* 2017 Feb 25;13(3):319-326.
99. Jiang Q, Li Q, Han J, Gou M, Zheng Y, Li B, Xiao R, Wang J. rLj-RGD3 induces apoptosis via the mitochondrial-dependent pathway and inhibits adhesion, migration and invasion of human HeyA8 cells via FAK pathway. *Int J Biol Macromol* 2017 Mar;96:652-668.
100. Liu S, Fei W, Shi Q, Li Q, Kuang Y, Wang C, He C, Hu X. CHAC2, downregulated in gastric and colorectal cancers, acted as a tumor suppressor inducing apoptosis and autophagy through unfolded protein response. *Cell Death Dis* 2017 Aug 24;8(8):e3009.
101. Huang YF, Zhu DJ, Chen XW, Chen QK, Luo ZT, Liu CC, Wang GX, Zhang WJ, Liao NZ. Curcumin enhances the effects of irinotecan on colorectal cancer cells through the generation of reactive oxygen species and activation of the endoplasmic reticulum stress pathway. *ONCOTARGET* 2017 Jun 20;8(25):40264-40275.
102. Li Y, Chen Y, Huang H, Shi M, Yang W, Kuang J, Yan J. Autophagy mediated by endoplasmic reticulum stress enhances the caffeine-induced apoptosis of hepatic stellate cells. *Int J Mol Med* 2017 Nov;40(5):1405-1414.
103. Wu DP, Lin TY, Bai LR, Huang JL, Zhou Y, Zhou N, Zhong SL, Gao S, Yin XX. Enhanced phototoxicity of photodynamic treatment by Cx26-composed GJIC via ROS, calcium- and lipid peroxide-mediated pathways. *J Biophotonics* 2017 Dec;10(12):1586-1596.
104. Zhao K, Li G, Yao Y, Zhou Y, Li Z, Guo Q, Lu N. Activation of phospholipase C- γ and translocation of phosphatidylinositol-3,4,5-trisphosphate 3-phosphatase contribute to GL-V9-induced apoptosis in human gastric cancer cells. *Exp Cell Res* 2017 Jul 1;356(1):8-19.
105. Chen LD, Liu ZH, Zhang LF, Yao JN, Wang CF. Sanggenon C induces apoptosis of colon cancer cells via inhibition of NO production, iNOS expression and ROS activation of the mitochondrial pathway. *Oncol Rep* 2017 Oct;38(4):2123-2131.
106. Qiu L, Ma Y, Luo Y, Cao Z, Lu H. Protective effects of isorhamnetin on N2a cell against endoplasmic reticulum stress-induced injury is mediated by PKC ϵ . *Biomed Pharmacother* 2017 Sep;93:830-836.
107. Lu Y, Gu Y, Ding X, Wang J, Chen J, Miao C. Intracellular Ca²⁺ homeostasis and JAK1/STAT3 pathway are involved in the protective effect of propofol on BV2 microglia against hypoxia-induced inflammation and apoptosis. *PLoS One* 2017 May 23;12(5):e0178098.
108. Wang P, Gao YM, Sun X, Guo N, Li J, Wang W, Yao LP, Fu YJ. Hepatoprotective

- effect of 2'-O-galloylhyperin against oxidative stress-induced liver damage through induction of Nrf2/ARE-mediated antioxidant pathway. *Food Chem Toxicol* 2017 Apr;102:129-142.
109. Ye W, Zhu S, Liao C, Xiao J, Wu Q, Lin Z, Chen J. Advanced oxidation protein products induce apoptosis of human chondrocyte through reactiveoxygen species-mediated mitochondrial dysfunction and endoplasmic reticulum stress pathways. *FUND CLIN PHARMACOL* 2017 Feb;31(1):64-74.
110. Wang XP, Yu X, Yan XJ, Lei F, Chai YS, Jiang JF, Yuan ZY, Xing DM, Du LJ. TRPM8 in the negative regulation of TNF α expression during cold stress. *SCI REP-UK* 2017 Mar 23;7:45155.
111. Xiang W, Liao W, Yi Z, He X, Ding Y. 25-Hydroxyvitamin D-1- α -hydroxylase in apolipoprotein E knockout mice: The role of protecting vascular smooth muscle cell from calcification. *Biomed Pharmacother* 2017 Apr;88:971-977.
112. Xu W, Guan R, Shi F, Du A, Hu S. Structural analysis and immunomodulatory effect of polysaccharide from *Atractylodismacrocephalae* Koidz. on bovine lymphocytes. *CARBOHYD POLYM* 2017 Oct 15;174:1213-1223.
113. Zhao P, Ma NT, Chang RY, Li YX, Hao YJ, Yang WL, Zheng J, Niu Y, Sun T, Yu JQ. Mechanism of *Lycium barbarum* polysaccharides on primary cultured rat hippocampal neurons. *Cell Tissue Res* 2017 Sep;369(3):455-465.
114. Yang K, Jiang X, Su Q, Wang J, Li C, Xia Y, Cheng S, Qin Q, Cao X, Chen C, Tu B. Disruption of glutamate neurotransmitter transmission is modulated by SNAP-25 in benzo[a]pyrene-induced neurotoxic effects. *Toxicology* 2017 Jun 1;384:11-22.
115. Xiang Y, Jing Z, Haixia W, Ruitao Y, Huaixiu W, Zenggen L, Lijuan M, Yiping W, Yanduo T. Antiproliferative Activity of Phenylpropanoids Isolated from *Lagotis brevifolia* Maxim. *Phytother Res* 2017 Oct;31(10):1509-1520.
116. Zhao T, Chang L, Zhang B, Lu M, Wang X, Orgah JO, Wang Y, Tian X, Yang J, Fan G, Zhang B, Zhu Y. Specific Combination of Salvianolic Acids As Core Active Ingredients of Danhong Injection for Treatment of Arterial Thrombosis and Its Derived Dry Gangrene. *Front Pharmacol* 2017 Jun 13;8:361.
117. Zhang Y, Wang L, Feng H, Hu G, Wang L, Liu J, Gao X, Shang J, Zhu T, Tang S, Jia G. Effects of 1,4-naphthoquinone aged carbon black particles on the cell membrane of human bronchial epithelium. *ENVIRON TOXICOL CHEM* 2017 Sep;36:21-27.
118. Sun N, Jin Z, Li D, Yin H, Lin S. An Exploration of the Calcium-Binding Mode of Egg White Peptide, Asp-His-Thr-Lys-Glu, and In Vitro Calcium Absorption Studies of Peptide-Calcium Complex. *J AGR FOOD CHEM* 2017 Nov 8;65(44):9782-9789.
119. Xiang QD, Yu Q, Wang H, Zhao MM, Liu SY, Nie SP, Xie MY. Immunomodulatory Activity of *Ganoderma atrum* Polysaccharide on Purified T Lymphocytes through Ca²⁺/CaN and Mitogen-Activated Protein Kinase Pathway Based on RNA Sequencing. *J AGR FOOD CHEM* 2017 Jul 5;65(26):5306-5315.
120. Liu L, Chang X, Zhang Y, Wu C, Li R, Tang L, Zhou Z. Fluorochloridone induces primary cultured Sertoli cells apoptosis: Involvement of ROS and intracellular calcium ions-mediated ERK1/2 activation. *Toxicol In Vitro* 2018 Mar;47:228-237.
121. Huang RZ, Huang XC, Zhang B, Jia HY, Liao ZX, Wang HS. 16-O-caffeoyl-16-hydroxylhexadecanoic acid, a medicinal plant-derived phenylpropanoid, induces apoptosis in human hepatocarcinoma cells through ROS-dependent endoplasmic reticulum stress. *Phytomedicine* 2018 Mar 1;41:33-44.
122. Zhao P, Chang RY, Liu N, Wang J, Zhou R, Qi X, Liu Y, Ma L, Niu Y, Sun T, Li YX, He YP, Yu JQ. Neuroprotective Effect of Oxysophocarpine by Modulation of MAPK Pathway in Rat Hippocampal Neurons Subject to Oxygen-Glucose Deprivation and Reperfusion. *Cell Mol Neurobiol* 2018 Mar;38(2):529-540.
123. Dong X, Gao X, Dai Y, Ran N, Yin H. Serum exosomes can restore cellular function in vitro and be used for diagnosis in dysferlinopathy. *Theranostics* 2018 Feb 2;8(5):1243-1255.
124. Maryam A, Mehmood T, Yan Q, Li Y, Khan M, Ma T. Proscillaridin A Promotes Oxidative Stress and ER Stress, Inhibits STAT3 Activation, and Induces Apoptosis in A549 Lung Adenocarcinoma Cells. *Oxid Med Cell Longev* 2018 Jan 11;2018:3853409.
125. Guo C, Ma R, Liu X, Chen T, Li Y, Yu Y, Duan J, Zhou X, Li Y, Sun Z. Silica nanoparticles promote oxLDL-induced macrophage lipid accumulation and apoptosis via endoplasmic reticulum stress signaling. *Sci Total Environ* 2018 Mar 10;631-632:570-579.
126. Ma K, Shi W, Xu M, Liu J, Zhang F. Genome-Wide Identification and Characterization of Long Non-Coding RNA in Wheat Roots in Response to Ca²⁺ Channel Blocker. *Front Plant Sci* 2018 Mar 6;9:244.
127. Yan D, Zhu D, Zhao X, Su J. SHP-2 restricts apoptosis induced by chemotherapeutic agents via Parkin-dependent autophagy in cervical cancer. *Cancer Cell Int* 2018 Jan 15;18:8.
128. Li Z, Tang Y, Zhu S, Li D, Han X, Gu G, Xing N, Ren J, Guo Z, Jiao W, Yan L, Xu Z, Zhang W. Ethanol extract of *Patrinia scabiosae* induces the death of human renal cell carcinoma 786-O cells via SIRT-1 and mTOR signaling-mediated metabolic disruptions. *Oncol Rep* 2018 Feb;39(2):764-772.
129. Shen XY, Luo T, Li S, Ting OY, He F, Xu J, Wang HQ. Quercetin inhibits okadaic acid-induced tau protein hyperphosphorylation through the Ca²⁺ calpain p25 CDK5 pathway in HT22 cells. *Int J Mol Med* 2018 Feb;41(2):1138-46.
130. Guo C, Wang J, Jing L, Ma R, Liu X, Gao L, Cao L, Duan J, Zhou X, Li Y, Sun Z. Mitochondrial dysfunction, perturbations of mitochondrial dynamics and biogenesis involved in endothelial injury induced by silica nanoparticles. *Environ Pollut* 2018 May;236:926-936.
131. Jia Y, Yang B, Dong W, Liu Z, Lv Z, Jia Z, Qiu L, Wang L, Song L. A serotonin receptor (Cg5-HTR-1) mediating immune response in oyster *Crassostrea gigas*. *Dev Comp Immunol* 2018 May;82:83-93.
132. Maryam A, Mehmood T, Yan Q, Li Y, Khan M, Ma T. Proscillaridin A Promotes Oxidative Stress and ER Stress, Inhibits STAT3 Activation, and Induces Apoptosis in A549 Lung Adenocarcinoma Cells. *Oxid Med Cell Longev* 2018 Jan 11;2018:3853409.
133. Liu Y, Wang H, Yang M, Liu N, Zhao Y, Qi X, Niu Y, Sun T, Li Y, Yu J. Cistanche deserticola polysaccharides protect PC12 cells against OGD/RP-induced injury. *Biomed Pharmacother* 2018 Mar;99:671-680.
134. Zhao P, Chang RY, Liu N, Wang J, Zhou R, Qi X, Liu Y, Ma L, Niu Y, Sun T, Li YX, He YP, Yu JQ. Neuroprotective Effect of Oxysophocarpine by Modulation of MAPK Pathway in Rat Hippocampal Neurons Subject to Oxygen-Glucose Deprivation and Reperfusion. *Cell Mol Neurobiol* 2018 Mar;38(2):529-540.
135. Zhang C, Xie L, Guan F, Cui Y. 3H-1,2-dithiole-3-thione protects PC12 cells against amyloid beta 1-42 (A β 1-42) induced apoptosis via activation of the ERK1/2 pathway. *Life Sci* 2018 Nov 15;213:74-81.
136. Li J, Ma L, Liao X, Liu D, Lu X, Chen S, Ye X, Ding T. Ultrasound-Induced *Escherichia coli* O157:H7 Cell Death Exhibits Physical Disruption and Biochemical Apoptosis. *Front Microbiol* 2018 Oct 16;9:2486.
137. Ma LQ, Yu Y, Chen H, Li M, Ihsan A, Tong HY, Huang XJ, Gao Y. Sweroside Alleviated Aconitine-Induced Cardiac Toxicity in H9c2 Cardiomyoblast Cell Line. *Front Pharmacol* 2018 Oct 25;9:1138.
138. Cheng X, Feng H, Wu H, Jin Z, Shen X, Kuang J, Huo Z, Chen X, Gao H, Ye F, Ji X, Jing X, Zhang Y, Zhang T, Qiu W, Zhao R. Targeting autophagy enhances apatinib-induced apoptosis via endoplasmic reticulum stress for human colorectal cancer. *Cancer Lett* 2018 Sep 1;431:105-114.
139. Chen X, Wu Q, Sun P, Zhao Y, Zhu M, Miao C. Propofol Disrupts Aerobic Glycolysis in Colorectal Cancer Cells via Inactivation of the NMDAR-CAMKII-ERK Pathway. *CELL PHYSIOL BIOCHEM* 46(2):492-504. 2018
140. Yang H, Li X, Liu Y, Li X, Li X, Wu M, Lv X, Chunhua C, Ding X, Zhang Y. Crocin Improves the Endothelial Function Regulated by Kca3.1 Through ERK and Akt Signaling Pathways. *CELL PHYSIOL BIOCHEM* 46(2):765-780. 2018
141. Du Y, Fu M, Wang YT, Dong Z. Neuroprotective Effects of Ginsenoside Rf on Amyloid- β -Induced Neurotoxicity in vitro and in vivo. *J Alzheimers Dis* 64(1):309-322. 2018
142. Zhu SY, Zhuang JS, Wu Q, Liu ZY, Liao CR, Luo SG, Chen JT, Zhong ZM. Advanced oxidation protein products induce pre-osteoblast apoptosis through a nicotinamide adenine dinucleotide phosphate oxidase-dependent,

mitogen-activated protein kinases-mediated intrinsic apopto Aging Cell 17(4):e12764. 2018 Aug

143. Guo C, Ma R, Liu X, Chen T, Li Y, Yu Y, Duan J, Zhou X, Li Y, Sun Z Silica nanoparticles promote oxLDL-induced macrophage lipid accumulation and apoptosis via endoplasmic reticulum stress signaling. Sci Total Environ 631-632:570-579. 2018 Aug 1

144. Guo X, Zheng M, Pan R, Zang B, Jin M Hydroxysafflor Yellow A Suppresses Platelet Activating Factor-Induced Activation of Human Small Airway Epithelial Cells. Front Pharmacol 9:859. 2018 Aug 3

145. Shen XY, Luo T, Li S, Ting OY, He F, Xu J, Wang HQ Quercetin inhibits okadaic acid-induced tau protein hyperphosphorylation through the Ca2+-calpain-p25-CDK5 pathway in HT22 cells. Int J Mol Med 41(2):1138-1146. 2018 Feb

146. Liao C, Xu D, Liu X, Fang Y, Yi J, Li X, Guo B Iridium (III) complex-loaded liposomes as a drug delivery system for lung cancer through mitochondrial dysfunction. INT J NANOMED 13:4417-4431. 2018 Jul 30

147. Jia Y, Miao Y, Yue M, Shu M, Wei Z, Dai Y Tetrandrine attenuates the bone erosion in collagen-induced arthritis rats by inhibiting osteoclastogenesis via spleen tyrosine kinase. FASEB J 32(6):3398-3410. 2018 Jun

148. Yuan XL, Zhao YP, Huang J, Liu JC, Mao WQ, Yin J, Peng BW, Liu WH, Han S, He XH A Kv1.3 channel-specific blocker alleviates neurological impairment through inhibiting T-cell activation in experimental autoimmune encephalomyelitis. CNS Neurosci Ther 24(10):967-977. 2018 Oct

149. Cui P, Lin S, Jin Z, Zhu B, Song L, Sun N In vitro digestion profile and calcium absorption studies of a sea cucumber ovum derived heptapeptide-calcium complex. Food Funct 9(9):4582-4592. 2018 Sep 19

150. Zhang Y, Feng T, Qu J, Sun N, Liu L Toxicity and haemolytic activity of a newly described dinoflagellate, *Heterocapsa bohainensis* to the rotifer *Brachionus plicatilis*. Harmful Algae 84:112-118. 2019 Apr

151. Zhong C, Mai Y, Gao H, Zhou W, Zhou D Mitochondrial targeting of TR3 is involved in TPA induced apoptosis in breast cancer cells. Gene 693:61-68. 2019 Apr 20

152. Dong G, Liang F, Sun B, Wang C, Liu Y, Guan X, Yang B, Xiu C, Yang N, Liu F, Lu T, Han W Presence and function of stress granules in atrial fibrillation. PLoS One 14(4):e0213769. 2019 Apr 3

153. La X, Zhang L, Yang Y, Li H, Song G, Li Z Tumor-secreted GRP78 facilitates the migration of macrophages into tumors by promoting cytoskeleton remodeling. Cell Signal 60:1-16. 2019 Aug

154. Zhang G, Wang Z, Chen W, Cao Y, Wu J, Qiang G, Ji A, Wu J, Jiang C Dual effects of gossypol on human hepatocellular carcinoma via endoplasmic reticulum stress and autophagy. INT J BIOCHEM CELL B 113:48-57. 2019 Aug

155. Fu S, Ding M, Liang Q, Yang Y, Chen M, Wei X, Wang A, Liao S, Ye J The key differentially expressed genes and proteins related to immune response in the spleen of pufferfish (*Takifugu obscurus*) infected by *Aeromonas hydrophila*. FISH SHELLFISH IMMUN 91:1-11. 2019 Aug

156. Yang Y, Cong H, Du N, Han X, Song L, Zhang W, Li C, Tien P Mitochondria Redistribution in Enterovirus A71 Infected Cells and Its Effect on Virus Replication. Virol Sin 34(4):397-411. 2019 Aug

157. Xiaoyu M, Xiuling D, Chunyu Z, Yi S, Jiangchao Q, Yuan Y, Changsheng L Polyglutamic acid-coordinated assembly of hydroxyapatite nanoparticles for synergistic tumor-specific therapy. Nanoscale 11(32):15312-15325. 2019 Aug 15

158. Li M, Xi P, Xu Y, Wang Z, Han X, Ren W, Phouthapane V, Miao J Taurine Attenuates Streptococcus uberis-Induced Bovine Mammary Epithelial Cells Inflammation via Phosphoinositides/Ca2+ Signaling. Front Immunol 10:1825. 2019 Aug 7

159. Feng Q, Liu C, Gao W, Geng XL, Dai N Salidroside-Mitigated Inflammatory Injury of Hepatocytes with Non-Alcoholic Fatty Liver Disease via Inhibition TRPM2 Ion Channel Activation. DIABET METAB SYND OB 12:2755-2763. 2019 Dec 27

160. Zhang X, Hong S, Qi S, Liu W, Zhang X, Shi Z, Chen W, Zhao M, Yin X NLRP3 Inflammasome Is Involved in Calcium-Sensing Receptor-Induced Aortic Remodeling in SHR. MEDIAT INFLAMM 2019:6847087. 2019 Feb 13

161. Cheng Q, Fang L, Feng D, Tang S, Yue S, Huang Y, Han J, Lan J, Liu W, Gao L, Luo Z Memantine ameliorates pulmonary inflammation in a mice model of COPD induced by cigarette smoke combined with LPS. Biomed Pharmacother 109:2005-2013. 2019 Jan

162. Hu Q, Niu Q, Song H, Wei S, Wang S, Yao L, Li YP Polysaccharides from *Portulaca oleracea* L. regulated insulin secretion in INS-1 cells through voltage-gated Na+ channel. Biomed Pharmacother 109:876-885. 2019 Jan

163. Tan Q, Zhang M, Geng L, Xia Z, Li C, Usman M, Du Y, Wei L, Bi H Hormesis of methylmercury-human serum albumin conjugate on N9 microglia via ERK/MAPKs and STAT3 signaling pathways. TOXICOL APPL PHARM 362:59-66. 2019 Jan 1

164. Wang C, Ming B, Wu X, Wu T, Cai S, Hu P, Tang J, Tan Z, Liu C, Zhong J, Zheng F, Dong L Sphingomyelin synthase 1 enhances BCR signaling to promote lupus-like autoimmune response. EBioMedicine 45:578-587. 2019 Jul

165. Bian X, Wu S, Yin X, Mu L, Yan F, Kong L, Guo Z, Wu L, Ye J Lyn is involved in host defense against *S. agalactiae* infection and BCR signaling in Nile tilapia (*Oreochromis niloticus*). Dev Comp Immunol 96:1-8. 2019 Jul

166. Zhao H, Chen G, Liang H Dual PI3K/mTOR Inhibitor, XL765, suppresses glioblastoma growth by inducing ER stress-dependent apoptosis. ONCOTARGETS THER 12:5415-5424. 2019 Jul 8

167. Wu DP, Bai LR, Lv YF, Zhou Y, Ding CH, Yang SM, Zhang F, Wang YY, Huang JL, Yin XX A Novel Role of Connexin 40-Formed Channels in the Enhanced Efficacy of Photodynamic Therapy. Front Oncol 9:595. 2019 Jul 9

168. Chen W, Ju XZ, Lu Y, Ding XW, Miao CH, Chen JW Propofol improved hypoxia-impaired integrity of blood-brain barrier via modulating the expression and phosphorylation of zonula occludens-1. CNS Neurosci Ther 25(6):704-713. 2019 Jun

169. Li S, Dai Z, Yang D, Li W, Dai H, Sun B, Liu X, Xie X, Xu R, Zhao X Targeting β 2 subunit of Na+/K+-ATPase induces glioblastoma cell apoptosis through elevation of intracellular Ca2. Am J Cancer Res 9(6):1293-1308. eCollection 2019. 2019 Jun 1

170. Li M, Fang XZ, Zheng YF, Xie YB, Ma XD, Liu XT, Xia Y, Shao DH Transient receptor potential vanilloid 4 is a critical mediator in LPS mediated inflammation by mediating calcineurin/NFATc3 signaling. BIOCHEM BIOPH RES CO 513(4):1005-1012.2019 Jun 11

171. Wang M, Wu Y, Yu Y, Fu Y, Yan H, Wang X, Li T, Peng W, Luo D Rutaecarpine prevented ox-LDL-induced VSMCs dysfunction through inhibiting overexpression of connexin 43. Eur J Pharmacol 853:84-92. 2019 Jun 15

172. Wang R, Wang R, Niu X, Cheng Y, Shang X, Li Y, Li S, Liu X, Shao J Role of astrocytes-derived d-serine in PFOS-induced neurotoxicity through NMDARs in the rat primary hippocampal neurons. Toxicology 422:14-24. 2019 Jun 15

173. Liu SM, Xiao ZF, Li X, Zhao YN, Wu XM, Han J, Chen B, Li JY, Fan CX, Xu B, Xue XY, Xue WW, Yang Y, Dai JW Vascular endothelial growth factor activates neural stem cells through epidermal growth factor receptor signal after spinal cord injury. CNS Neurosci Ther 25(3):375-385. 2019 Mar

174. Sun Z, Li Y, Zhou H, Cai M, Liu J, Gao S, Yang J, Tong L, Wang J, Zhou S, Hu Z, Wang Y, Wang K, Zhang L, Wang H, Zhang L, Shi F, Cao X, Zhang S, Ji Y, Zhao J Simulated microgravity reduces intracellular-free calcium concentration by inhibiting calcium channels in primary mouse osteoblasts. J Cell Biochem 120(3):4009-4020.2019 Mar

175. Wang Z, Wang L, Prabhakar N, Xing Y, Rosenholm JM, Zhang J, Cai K CaP coated mesoporous polydopamine nanoparticles with responsive membrane permeation ability for combined photothermal and siRNA therapy. Acta Biomater 86:416-428. 2019 Mar 1

176. Xu H, Cheng J, Wang X, Liu H, Wang S, Wu J, Xu B, Chen A, He F Resveratrol pretreatment alleviates myocardial ischemia/reperfusion injury by inhibiting STIM1-mediated intracellular calcium accumulation. J Physiol Biochem 75(4):607-618. 2019 Nov

177. Song J, Tang Z, Li H, Jiang H, Sun T, Lan R, Wang T, Wang S, Ye Z, Liu J Role of JAK2 in the Pathogenesis of Diabetic Erectile Dysfunction and an Intervention With Berberine. *J Sex Med* 16(11):1708-1720. 2019 Nov
178. Qi J, Wu Q, Zhu X, Zhang S, Chen X, Chen W, Sun Z, Zhu M, Miao C Propofol attenuates the adhesion of tumor and endothelial cells through inhibiting glycolysis in human umbilical vein endothelial cells. *ACTA BIOCH BIOPH SIN* 51(11):1114-1122.2019 Nov 18
179. Zhang L, Wu T, Olatunji OJ, Tang J, Wei Y, Ouyang Z N6-(2-hydroxyethyl)-adenosine from *Cordyceps cicadae* attenuates hydrogen peroxide induced oxidative toxicity in PC12 cells. *Metab Brain Dis* 34(5):1325-1334. 2019 Oct
180. Ding XW, Sun X, Shen XF, Lu Y, Wang JQ, Sun ZR, Miao CH, Chen JW Propofol attenuates TNF- α -induced MMP-9 expression in human cerebral microvascular endothelial cells by inhibiting Ca²⁺/CAMK II/ERK/NF- κ B signaling pathway. *Acta Pharmacol Sin* 40(10):1303-1313. 2019 Oct
181. Liu Y, Li J, Chen H, Cai Y, Sheng T, Wang P, Li Z, Yang F, Gu N Magnet-activatable nanoliposomes as intracellular bubble microreactors to enhance drug delivery efficacy and burst cancer cells. *Nanoscale* 11(40):18854-18865. 2019 Oct 28
182. Zhao C, He R, Shen M, Zhu F, Wang M, Liu Y, Chen H, Li X, Qin R PINK1/Parkin-Mediated Mitophagy Regulation by Reactive Oxygen Species Alleviates Rocaglamide A-Induced Apoptosis in Pancreatic Cancer Cells. *Front Pharmacol* 10:968. 2019 Sep 3
183. Deng Y, Wang J, Xie G, Zeng X, Li H Circ-HIPK3 Strengthens the Effects of Adrenaline in Heart Failure by MiR-17-3p - ADCY6 Axis. *Int J Biol Sci* 15(11):2484-2496.2019 Sep 7
184. Qian Yang, Xixi Cai, Muchen Huang, Lee Jia, Shaoyun Wang Immunomodulatory Effects of Pseudostellaria Heterophylla Peptide on Spleen Lymphocytes via a Ca²⁺/CaN/NFATc1/IFN- γ Pathway *Food Funct* 2019 Jun 19;10(6):3466-3476.
185. Yingshi Ji, Yang Hu, Jinghong Ren, Rajesh Khanna, Yuan Yao, Yang Chen, Qi Li, Li Sun CRMP2-derived Peptide ST2-104 (R9-CBD3) Protects SH-SY5Y Neuroblastoma Cells Against A β 25-35-induced Neurotoxicity by Inhibiting the pCRMP2/NMDAR2B Signaling Pathway *CHEM-BIOL INTERACT* 2019 May 25;305:28-39.
186. Zi-Qi Liu, Na Liu, Si-Si Huang, Miao-Miao Lin, Shu Qin, Jun-Chao Wu, Zhong-Qin Liang, Zheng-Hong Qin, Yan Wang NADPH protects against kainic acid-induced excitotoxicity via autophagy-lysosome pathway in rat striatum and primary cortical neurons *Toxicology* 2020 Apr 15;435:152408.
187. Jia-Peng Li, Wei Wei, Xiao-Xue Li, Ming Xu Regulation of NLRP3 inflammasome by CD38 through cADPR-mediated Ca²⁺ release in vascular smooth muscle cells in diabetic mice *Life Sci* 2020 Aug 15;255:117758.
188. Chunxia Chen, Wan Chen, Zhihuan Nong, Yichu Nie, Xiaoyu Chen, Xiaorong Pan, Ying Guo, Meicun Yao, Wenbin Deng Hyperbaric oxygen alleviated cognitive impairments in mice induced by repeated cerebral ischemia-reperfusion injury via inhibition of autophagy *Life Sci* 2020 Jan 15;241:117170.
189. Xiangnan Kong, Min Li, Kai Shao, Yinrong Yang, Qian Wang, Meijuan Cai Progesterone induces cell apoptosis via the CACNA2D3/Ca²⁺/p38 MAPK pathway in endometrial cancer *Oncol Rep* 2020 Jan;43(1):121-132.
190. Zi-Qiang Liu, Da-Yong Zhou, Yu-Xin Liu, Man-Man Yu, Bing Liu, Liang Song, Xiu-Ping Dong, Hang Qi, Fereidoon Shahidi Inhibitory effect of natural metal ion chelators on the autolysis of sea cucumber (*Stichopus japonicus*) and its mechanism *Food Res Int* 2020 Jul;133:109205.
191. Lin Che, Huan Yao, Chuan-Li Yang, Ni-Jun Guo, Jing Huang, Zi-Li Wu, Li-Yin Zhang, Yuan-Yuan Chen, Gang Liu, Zhong-Ning Lin, Yu-Chun Lin Cyclooxygenase-2 modulates ER-mitochondria crosstalk to mediate superparamagnetic iron oxide nanoparticles induced hepatotoxicity: an in vitro and in vivo study *Nanotoxicology* 2020 Mar;14(2):162-180.
192. Ke Wang, Yeliang Lou, Huang Xu, Xiaoming Zhong, Zhen Huang Harpagide from *Scrophularia* protects rat cortical neurons from oxygen-glucose deprivation and reoxygenation-induced injury by decreasing endoplasmic reticulum stress *J Ethnopharmacol* 2020 May 10;253:112614.
193. Fan Yang, Li Zou, Yong Wu, Zhihua Wu, Anshu Yang, Hongbing Chen, Xin Li Structure and allergenicity assessments of bovine β -lactoglobulin treated by sonication-assisted irradiation *J Dairy Sci* 2020 May;103(5):4109-4120.
194. Tao Ren, Yi-Jun Tang, Mei-Fang Wang, Han-Sheng Wang, Yan Liu, Xin Qian, Chan Chang, Ming-Wei Chen Triptolide induces apoptosis through the calcium/calmodulin-dependent protein kinase kinase β /AMP-activated protein kinase signaling pathway in non-small cell lung cancer cells *Oncol Rep* 2020 Nov;44(5):2288-2296.
195. Chunyuan Wang, Xianzong Meng, Meiling Meng, Min Shi, Wenping Sun, Xiaojing Li, Xiao Zhang, Ruihao Liu, Ying Fu, Laiyu Song Oxidative stress activates the TRPM2-Ca²⁺-NLRP3 axis to promote PM 2.5-induced lung injury of mice *Biomed Pharmacother* 2020 Oct;130:110481.
196. Jiaxin Li, Qian Zhou, Xin Zhou, Baodong Wei, Yingbo Zhao, Shujuan Ji Calcium Treatment Alleviates Pericarp Browning of 'Nanguo' Pears by Regulating the GABA Shunt After Cold Storage *Front Plant Sci* 2020 Sep 15;11:580986.
197. Li Luo, SiYi Liu, Dan Zhang, Fu Wei, NiNa Gu, Yan Zeng, XiaoYing Chen, Shan Xu, ShuKe Liu, Tao Xiang Chromogranin A (CGA)-derived polypeptide (CGA 47-66) inhibits TNF- α -induced vascular endothelial hyper-permeability through SOC-related Ca²⁺ signaling *Peptides* 2020 Sep;131:170297.
198. Jiheng Zhan, Xing Li, Dan Luo, Wanying Yan, Yonghui Hou, Yu Hou, Shudong Chen, Jiyao Luan, Qing Zhang, Dingkun Lin Polydatin Attenuates OGD/R-Induced Neuronal Injury and Spinal Cord Ischemia/Reperfusion Injury by Protecting Mitochondrial Function via Nrf2/ARE Signaling Pathway *Oxid Med Cell Longev* 2021 Apr 20;2021:6687212.
199. Bin Zeng, Lei Liu, Xiaoting Liao, Caixia Zhang Cardiomyocyte protective effects of thyroid hormone during hypoxia/reoxygenation injury through activating of IGF-1-mediated PI3K/Akt signalling *J Cell Mol Med* 2021 Apr;25(7):3205-3215.
200. Yuan Yao, Yingshi Ji, Jinghong Ren, Huanyu Liu, Rajesh Khanna, Li Sun Inhibition of autophagy by CRMP2-derived peptide ST2-104 (R9-CBD3) via a CaMKK β /AMPK/mTOR pathway contributes to ischemic postconditioning-induced neuroprotection against cerebral ischemia-reperfusion *Mol Brain* 2021 Aug 6;14(1):123.
201. Yujie Zhong, Chengni Jin, Jiahui Han, Jiachang Zhu, Qi Liu, Dianjun Sun, Xiaodong Xia, Yu Zhang, Xiaoli Peng Diosgenin Protects Against Kidney Injury and Mitochondrial Apoptosis Induced by 3-MCPD Through the Regulation of ER Stress, Ca²⁺ Homeostasis, and Bcl2 Expression *Mol Nutr Food Res* 2021 Aug;65(15):e2001202.
202. Jia Cheng, Jv Zhang, Jingyue Yang, Bing Yi, Gang Liu, Man Zhou, John P Kastelic, Bo Han, Jian Gao *Klebsiella pneumoniae* infection causes mitochondrial damage and dysfunction in bovine mammary epithelial cells *Vet Res* 2021 Feb 10;52(1):17.
203. Imran Rashid Rajput, Summra Yaqoob, Sun Yajing, Edmond Sanganyado, Liu Wenhua Polybrominated diphenyl ethers exert genotoxic effects in pantropic spotted dolphin fibroblast cell lines *Environ Pollut* 2021 Feb 15;271:116131.
204. Xiao-Yang Xie, Yu-Yun Li, Wen-Hui Ma, Ai-Fang Chen, Yu-Tong Sun, Ji Youn Lee, Aladdin Riad, Dao-Hua Xu, Robert H Mach, Yun-Sheng Huang Synthesis, binding, and functional properties of tetrahydroisoquinolino-2-alkyl phenones as selective σ 2R/TMEM97 ligands *Eur J Med Chem* 2021 Jan 1;209:112906.
205. Lin Che, Chuan-Li Yang, Yu Chen, Zi-Li Wu, Ze-Bang Du, Jia-Shen Wu, Cong-Ling Gan, Si-Ping Yan, Jing Huang, Ni-Jun Guo, Yu-Chun Lin, Zhong-Ning Lin Mitochondrial redox-driven mitofusin 2 S-glutathionylation promotes neuronal necroptosis via disrupting ER-mitochondria crosstalk in cadmium-induced neurotoxicity *Chemosphere* 2021 Jan;262:127878.
206. Jialin He, Jianyang Liu, Yan Huang, Xiangqi Tang, Han Xiao, Zuo Liu, Zheng Jiang, Liuwang Zeng, Zhiping Hu, Ming Lu OM-MSCs Alleviate the Golgi Apparatus Stress Response following Cerebral Ischemia/Reperfusion Injury via the PEDF-PI3K/Akt/mTOR Signaling Pathway *Oxid Med Cell Longev* 2021 Nov 13;2021:4805040.
207. Zhihao Liu, Xiyun Bian, Wenbo Gao, Jing Su, Chuanrui Ma, Xiaolin Xiao, Tian Yu,

- Han Zhang, Xiaozhi Liu, Guanwei Fan Rg3 promotes the SUMOylation of SERCA2a and corrects cardiac dysfunction in heart failure Pharmacol Res 2021 Oct;172:105843.
208. Yujie Zhong, Chengni Jin, Jiahui Han, Jiachang Zhu, Qi Liu, Dianjun Sun, Xiaodong Xia, Xiaoli Peng Inhibition of ER stress attenuates kidney injury and apoptosis induced by 3-MCPD via regulating mitochondrial fission/fusion and Ca2+ homeostasis Cell Biol Toxicol 2021 Oct;37(5):795-809.
209. Ruihong Ning, Yang Li, Zhou Du, Tianyu Li, Qinglin Sun, Lisen Lin, Qing Xu, Junchao Duan, Zhiwei Sun The mitochondria-targeted antioxidant MitoQ attenuated PM2.5-induced vascular fibrosis via regulating mitophagy Redox Biol 2021 Oct;46:102113.
210. Mingzhou Ye, Yi Zhao, Yuyuan Wang, Ruosen Xie, Yao Tong, John-Demian Sauer, Shaoqin Gong NAD(H)-loaded nanoparticles for efficient sepsis therapy via modulating immune and vascular homeostasis Nat Nanotechnol 2022 Aug;17(8):880-890.
211. Zequn Zhang, Chen Meng, Kun Hou, Zhigong Wang, Yan Huang, Xiaoying Lü The cytological and electrophysiological effects of silver nanoparticles on neuron-like PC12 cells PLoS One 2022 Dec 13;17(12):e0277942.
212. Jing Dai, Qi Li, Qinwei Zhou, Shen Zhang, Junru Chen, Yize Wang, Jing Guo, Yifan Gu, Fei Gong, Yueqiu Tan, Guangxiu Lu, Wei Zheng, Ge Lin IQCN disruption causes fertilization failure and male infertility due to manchette assembly defect EMBO Mol Med 2022 Dec 7;14(12):e16501.
213. Zhe Chen, Yucheng Jiao, Ying Zhang, Qingfeng Wang, Wenjian Wu, Jiancheng Zheng, Jitian Li G-Protein Coupled Receptor 35 Induces Intervertebral Disc Degeneration by Mediating the Influx of Calcium Ions and Upregulating Reactive Oxygen Species Oxid Med Cell Longev 2022 Jan 18;2022:5469220.
214. Qi Liu, Pengyue Du, Yue Zhu, Xintong Zhang, Jingzeng Cai, Ziwei Zhang Thioredoxin reductase 3 suppression promotes colitis and carcinogenesis via activating pyroptosis and necrosis Cell Mol Life Sci 2022 Jan 30;79(2):106.
215. Wenlong Zhang, Xin He, Haoli Yin, Wenmin Cao, Tingsheng Lin, Wei Chen, Wenli Diao, Meng Ding, Hao Hu, Wenjing Mo, Qing Zhang, Hongqian Guo Allosteric activation of the metabolic enzyme GPD1 inhibits bladder cancer growth via the lysoPC-PAFR-TRPV2 axis J Hematol Oncol 2022 Jul 14;15(1):93.
216. Jing Dai, Yongzhe Chen, Qi Li, Tianlei Zhang, Qinwei Zhou, Fei Gong, Guangxiu Lu, Wei Zheng, Ge Lin Pathogenic variant in ACTL7A causes severe teratozoospermia characterized by bubble-shaped acrosomes and male infertility Mol Hum Reprod 2022 Jul 29;28(8):gaac028.
217. Lishan Huang, Zhou Chen, Ruiyu Chen, Lu Lin, Lingjia Ren, Meilian Zhang, Libin Liu Increased fatty acid metabolism attenuates cardiac resistance to β -adrenoceptor activation via mitochondrial reactive oxygen species: A potential mechanism of hypoglycemia-induced myocardial injury in Redox Biol 2022 Jun;52:102320.
218. Lili Yang, Zhengjun Li, Yao Chen, Feiyan Chen, Haopeng Sun, Min Zhao, Yanqi Chen, Yali Wang, Wenlin Li, Li Zeng, Yaoyao Bian Elucidating the Novel Mechanism of Ligustrazine in Preventing Postoperative Peritoneal Adhesion Formation Oxid Med Cell Longev 2022 Mar 10;2022:9226022.
219. Peng Li, Xinhai Zhu, Guangchao Cao, Ruan Wu, Ke Li, Wenhui Yuan, Biyun Chen, Guodong Sun, Xichun Xia, Hua Zhang, Xiao Wang, Zhinan Yin, Ligong Lu, Yunfei Gao 1 α ,25(OH)2D3 reverses exhaustion and enhances antitumor immunity of human cytotoxic T cells J Immunother Cancer 2022 Mar;10(3):e003477.
220. Jinrong Bai, Na Xie, Ya Hou, Xiaorui Chen, Yao Hu, Yi Zhang, Xianli Meng, Xiaobo Wang, Ce Tang The enhanced mitochondrial dysfunction by cantleyoside confines inflammatory response and promotes apoptosis of human HFSL-RA cell line via AMPK/Sirt 1/NF- κ B pathway activation Biomed Pharmacother 2022 May;149:112847.

Version 2024.03.12